RANGER COUNT 3000

**2<sup>nd</sup> Scale Interface**  
Instruction Manual

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## INTRODUCTION

This Second Scale interface Kit is for use with the Ohaus product Ranger Count 3000. Through this kit you can connect additional platform with Ranger Count 3000.


## KIT CONTENTS

- 2<sup>nd</sup> Scale PC Board
- Pressure Button
- Screws (4)
- Instruction Manual


## INTERFACE INSTALLATION

Install the Interface module on the underside of the balance as shown.


1. Remove the red interface options cover. Break off the plastic blocker from the option cover.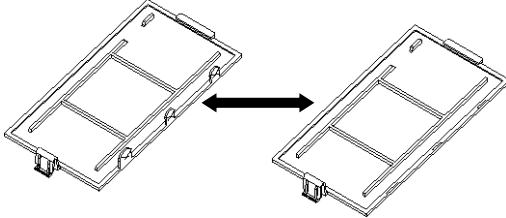

2. Install the Interface PCB. Push it through the pins and secure it with the 4 screws.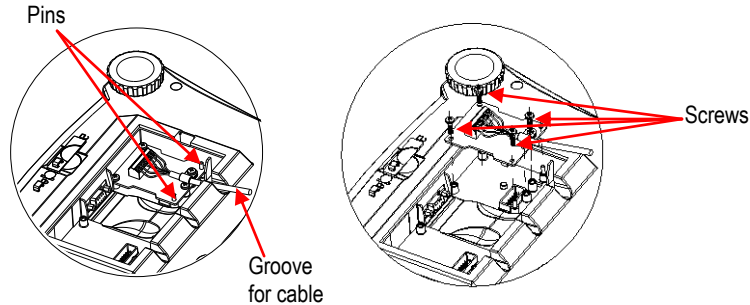


3. Connect a cable.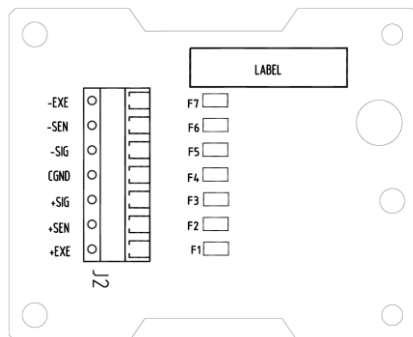

4. Place the cord in the groove and put the red interface option cover back on.

## SETUP

After the main platform is connected to the 2<sup>nd</sup> weighing platform, switch on the Dual Scale (D.Scale) menu item in the SETUP menu (see menu structure below).

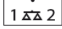Then set the CAP2 and Grad2 in the SETUP menu and do a calibration. Press and hold the  key to switch between the main platform and the secondary platform.

### CAL

Zero2  
Span1  
Span2  
Lin1  
Lin2  
GEO  
End

### SETUP

Reset  
**D.Scale**  
**Cap2**  
**Grad2**  
Pwr. Un  
Zero  
A.Opt  
A.Tare  
Bp.Opt  
Bp.Sig  
Bp.Key  
Accum  
End

**Note:** Make sure to perform a calibration of the scale before using the 2<sup>nd</sup> weighing platform. Make sure the weighing unit is in kg or lb when using high capacity platform. (If display shows error8.6, long press [**Units**] key to switch unit to kg or lb.)

## ACCESSORIES

For a complete listing of Ohaus printers and other accessories, contact Ohaus Corporation or visit [www.Ohaus.com](http://www.Ohaus.com)

## COMPLIANCE

This accessory has been tested and complies with the approvals listed in the applicable Instruction Manual.

### Disposal

In conformance with the European Directive 2002/96 EC on Waste Electrical and Electronic Equipment (WEEE) this device may not be disposed of in domestic waste. This also applies to countries outside the EU, per their specific requirements.

Please dispose of this product in accordance with local regulations at the collecting point specified for electrical and electronic equipment. If you have any questions, please contact the responsible authority or the distributor from which you purchased this device.

Should this device be passed on to other parties (for private or professional use), the content of this regulation must also be related.

For disposal instructions in Europe, refer to [www.ohaus.com/weee](http://www.ohaus.com/weee).

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P/N 30064481B © 2013 Ohaus Corporation, all rights reserved/ tous droits réservés/ todos los derechos reservados/ Alle Rechte vorbehalten/tutti i diritti riservati
OHAUS Corporation