

		<u>DNA Stain Type</u>	
		Ethidium Bromide	SmartGlow™ SYBR® Green
<b>Excitation Source</b>	SmartDoc blue illumination base	<b>Not recommended</b>	<b>E5001-ORANGE Orange filter</b>
	Blue light transilluminator	E5001-ORANGE Orange PMMA Filter	E5001-ORANGE Orange PMMA Filter
	UV transilluminator	E5001-590 590nm Filter	E5001-535 535nm Filter

**For optimum image quality:**

- Remove the phone's case
- Line up the camera lens with the photo filter

Place the photo filter into the recess on the top platform

Remove the extender for close ups or small gels



E5001-ORANGE Orange PMMA Filter (use with blue light)



E5001-535 535nm Filter (use with UV)



E5001-590 590nm Filter (use with UV)