



9935
LogWare II
Real-Time Data Acquisition Software
User's Guide

Fluke Corporation, Hart Scientific Division (Hart) warrants this product to be free from defects in material and workmanship under normal use and service for a period as stated in our current product catalog from the date of shipment. This warranty extends only to the original purchaser and shall not apply to any product which, in Hart's sole opinion, has been subject to misuse, alteration, abuse or abnormal conditions of operation or handling.

Software is warranted to operate in accordance with its programmed instructions on appropriate Hart products. It is not warranted to be error free.

Hart's obligation under this warranty is limited to repair or replacement of a product which is returned to Hart within the warranty period and is determined, upon examination by Hart, to be defective. If Hart determines that the defect or malfunction has been caused by misuse, alteration, abuse or abnormal conditions or operation or handling, Hart will repair the product and bill the purchaser for the reasonable cost of repair.

To exercise this warranty, the purchaser must forward the product after calling or writing Hart for authorization. Hart assumes NO risk for in-transit damage.

For service or assistance, please contact the manufacturer.

THE FOREGOING WARRANTY IS PURCHASER'S SOLE AND EXCLUSIVE REMEDY AND IS IN LIEU OF ALL OTHER WARRANTIES, EXPRESS OR IMPLIED, INCLUDING BUT NOT LIMITED TO ANY IMPLIED WARRANTY OR MERCHANTABILITY, OR FITNESS FOR ANY PARTICULAR PURPOSE OR USE. HART SHALL NOT BE LIABLE FOR ANY SPECIAL, INDIRECT, INCIDENTAL, OR CONSEQUENTIAL DAMAGES OR LOSS WHETHER IN CONTRACT, TORT, OR OTHERWISE.

Microsoft, MS-DOS, Windows, Window for Workgroups, and Windows 95, 98, ME, NT, 2000, XP, and VISTA are either registered trademarks or trademarks of Microsoft Corporation.

Fluke Corporation
Hart Scientific Division
799 E. Utah Valley Drive American Fork, UT 84003-9775 USA
Phone: +1.801.763.1600
Telefax: +1.801.763.1010
Email: support@hartscientific.com

www.hartscientific.com
Subject to change without notice. Copyright © 2008 Printed in USA

Table of Contents

1	Introduction	1
1.1	What is <i>LogWare</i>	1
1.2	License agreement	2
1.3	Features	2
1.4	Requirements	4
1.4.1	Communications requirements	4
1.4.2	Computer hardware requirements	4
1.4.3	Computer software requirements	5
1.4.4	Firmware requirements	5
1.5	Installation	6
1.6	Running <i>LogWare</i>	6
1.6.1	Getting started	6
1.7	Using <i>LogWare</i>	8
1.7.1	<i>LogWare</i> main display	9
1.8	<i>LogWare</i> options	11
2	Real-time logging	13
2.1	File menu	13
2.1.1	New	13
2.1.2	Open	13
2.1.3	Save as	14
2.2	Datalog menu	15
2.2.1	Setup	15
2.2.2	Start	15
2.2.3	Stop	15
2.2.4	Close	16
2.2.5	Mute all alarms	16
2.3	Log session settings dialog	16
2.3.1	Thermometer tab	19
2.3.2	Log Session tab	22
2.3.3	Channels tab	24
2.3.4	Delayed Start Date/Time dialog	26
2.3.5	Alarm Settings dialog	27
2.3.6	Statistics Settings dialog	30
2.4	Real-time log session	31
2.4.1	Channels grid	33
2.4.2	Statistics list	35
2.4.3	Graph	35
2.4.4	Current Alarm Settings dialog	37
2.4.5	Output file	37

2.4.6	Alarm log file	37
2.4.7	Trace color dialog	38
2.4.8	File Viewer window	39
2.4.9	Statistics dialog	40
2.5	Previous log session	40
2.5.1	Channels	42
2.5.2	Statistics list	43
2.5.3	Graph	44
2.5.4	Graph zooming	45
3	Downloading logged data	47
3.1	Download menu	47
3.1.1	Download logged data	47
3.1.2	Log status	47
3.2	Select Model dialog	48
3.3	Download options dialogs	50
3.3.1	1523 Handheld download options dialog	50
3.3.2	1524 Handheld download options dialog	51
3.3.3	1529 Chub-E4 download options dialog	53
3.3.4	1560 Black Stack download options dialog	55
3.3.5	1575A Super-Thermometer download options dialog	57
3.3.6	1590 Super-Thermometer II download options	58
3.4	Log status dialogs	60
3.4.1	1523 Handheld log status dialog	60
3.4.2	1524 Handheld log status dialog	61
3.4.3	1529 Chub-E4 log status dialog	62
3.4.4	1560 Black Stack log status dialog	62
4	Tools	65
4.1	Tools menu	65
4.1.1	Synchronize date/time	65
4.1.2	Logging labels	66
4.1.3	Thermometer settings	66
4.2	Select Model dialog	67
4.3	Logging labels dialogs	69
4.3.1	1523 Handheld logging label dialog	69
4.3.2	1524 Handheld logging labels dialog	70
4.3.3	1529 Chub-E4 logging labels	71
4.4	Thermometer settings dialogs	72
4.4.1	1523/1524 Handheld Settings	72
4.4.1.1	Sampling tab	74
4.4.1.2	Channels tab	75
4.4.1.3	Calibration tab	76
4.4.1.4	Misc tab	77
4.4.1.5	1523/1524 PRT/thermistor channel settings dialog	78
4.4.1.5.1	Probe tab	79
4.4.1.5.2	Misc tab	80
4.4.1.6	1523/1524 thermocouple channel settings dialog	81

	4.4.1.6.1	Probe tab	82
	4.4.1.6.2	Misc tab	84
4.4.2		1529 Chub-E4 settings dialog	84
4.4.2.1		Filtering tab	86
4.4.2.2		Sampling tab	87
4.4.2.3		Measurement tab	88
4.4.2.4		Channels Tab	88
4.4.2.5		Misc tab	89
4.4.2.6		1529 Chub-E4 PRT/thermistor channel settings dialog	90
	4.4.2.6.1	Probe tab	91
	4.4.2.6.2	Calibration tab	93
4.4.2.7		1529 Chub-E4 thermocouple channel settings	94
	4.4.2.7.1	Probe tab	95
	4.4.2.7.2	Calibration tab	96
4.4.3		1560 Black Stack settings dialog	96
4.4.3.1		Filtering tab	98
4.4.3.2		Sampling tab	99
4.4.3.3		Measurement tab	100
4.4.3.4		Modules tab	101
4.4.3.5		2560 SPRT Module Settings Dialog	101
	4.4.3.5.1	General tab	103
	4.4.3.5.2	Channels tab	103
	4.4.3.5.3	Calibration tab	104
4.4.3.6		2561 HPRT Module Settings dialog	105
	4.4.3.6.1	General tab	106
	4.4.3.6.2	Channels tab	106
	4.4.3.6.3	Calibration tab	107
4.4.3.7		2562 PRTS Module Settings dialog	108
	4.4.3.7.1	General tab	109
	4.4.3.7.2	Channels tab	110
	4.4.3.7.3	Calibration tab	110
4.4.3.8		2563 STHR Module Settings dialog	111
	4.4.3.8.1	General tab	113
	4.4.3.8.2	Channels tab	113
	4.4.3.8.3	Calibration tab	114
4.4.3.9		2564 THRS Module Settings dialog	115
	4.4.3.9.1	General tab	116
	4.4.3.9.2	Channels tab	117
	4.4.3.9.3	Calibration tab	117
4.4.3.10		2565 PTC Module Settings dialog	118
	4.4.3.10.1	Channels tab	120
	4.4.3.10.2	Calibration tab	121
4.4.3.11		2566 TCS Module Settings dialog	122
	4.4.3.11.1	Channels tab	123
	4.4.3.11.2	Calibration tab	124
		4.4.3.11.2.1 Inputs 1 to 6 tab	125
		4.4.3.11.2.2 Inputs 7 to 12 tab	126
4.4.3.12		2567 RTD Module Settings dialog	126
	4.4.3.12.1	General tab	128
	4.4.3.12.2	Channels tab	128
	4.4.3.12.3	Calibration tab	129
4.4.3.13		2568 RTDS Module Settings	130
	4.4.3.13.1	General tab	131
	4.4.3.13.2	Channels tab	132
	4.4.3.13.3	Calibration tab	132
4.4.3.14		1560 Black Stack PRT/Thermistor Channel Settings dialog	133
	4.4.3.14.1	Probe tab	134
4.4.3.15		1560 Black Stack Thermocouple Channel Settings dialog	135
	4.4.3.15.1	Probe tab	136
4.4.4		1575A Super-Thermometer Settings	137
4.4.4.1		Filtering tab	139
4.4.4.2		Sampling tab	140
4.4.4.3		Measurement tab	142
4.4.4.4		Channels tab	143
4.4.4.5		Calibration tab	145
4.4.4.6		1575A Super Thermometer/2575 Mighty Mux Channel Settings	146
	4.4.4.6.1	Probe tab	147
4.4.5		1590 Super-Thermometer II Settings dialog	148
4.4.5.1		Filtering tab	150
4.4.5.2		Sampling tab	151

4.4.5.3	Measurement tab	153
4.4.5.4	Channels tab	154
4.4.5.5	Calibration tab	156
4.4.5.6	1590 Super-Thermometer II/2590 'Mighty Mux II' Channel Settings dialog	157
4.4.5.6.1	Probe tab	158
4.4.6	1594A/1595A Hart Super-Thermometer Settings dialog	159
4.4.6.1	Filtering tab	161
4.4.6.2	Display tab	162
4.4.6.3	4.4.6.3 Timing tab	163
4.4.6.4	Scanning tab	164
4.4.6.5	Channel Setup tab	165
4.4.6.6	1594A/1595A Hart Super-Thermometer Probe Settings dialog	166
4.4.6.7	1594A/1595A Hart Super-Thermometer Channel Settings dialog	168
4.4.6.8	1594A/1595A Hart Super-Thermometer Rs Settings dialog	170

5 Menus 173

5.1	File menu	173
5.1.1	New	173
5.1.2	Open	173
5.1.3	Save as	174
5.1.4	Print	175
5.1.4.1	Readings	175
5.1.4.2	Graph	175
5.1.5	Exit	175
5.2	Datalog menu	175
5.2.1	Setup	175
5.2.2	Start	176
5.2.3	Stop	176
5.2.4	Close	176
5.2.5	Mute all alarms	176
5.3	Download menu	177
5.3.1	Download logged data	177
5.3.2	Log status	177
5.4	View menu	178
5.4.1	Toolbar	178
5.4.2	Status bar	178
5.4.3	Alarm settings	178
5.4.4	Show trace	179
5.4.5	Alarm log	179
5.4.6	Readings history	179
5.4.7	Open previously logged data file	179
5.4.8	MRU list	179
5.5	Graph menu	180
5.5.1	Graph settings dialog	180
5.5.1.1	Titles tab	180
5.5.1.2	Axis tab	183
5.5.1.3	Fonts tab	185
5.5.1.4	Trends tab	186
5.5.1.5	Background tab	189
5.5.1.6	Legend tab	190
5.5.2	Restore default settings	191
5.5.3	Trace color	191
5.5.4	Zoom in	191

5.5.5	Restore	191
5.5.6	Print graph	192
5.6	Tools menu	192
5.6.1	Synchronize date/time	192
5.6.2	Logging labels	192
5.6.3	Thermometer settings	193
5.6.4	Options	193
5.6.4.1	General tab	194
5.6.4.2	Folders tab	197
5.6.4.3	Tools tab	198
5.6.4.4	Admin. Passcode tab	199
5.7	Window menu	202
5.7.1	Cascade	202
5.7.2	Tile horizontal	202
5.7.3	Tile vertical	202
5.7.4	Arrange icons	202
5.7.5	Window list	202
5.8	Help menu	203
5.8.1	LogWare help	203
5.8.2	Technical support	203
5.8.3	About LogWare	203
6	How to	205
6.1	Setting up and starting a real-time log session	205
6.2	Starting a real-time log session unattended (delayed start)	205
6.3	Stopping a real-time log session automatically	206
6.4	Opening a configuration file automatically on start-up	207
6.5	Changing thermometer and probe settings	209
6.6	Importing logged data into a spreadsheet	210
6.7	Synchronizing the thermometer readout's clock to the computer	214
6.8	Protecting calibration parameters and probe conversion settings	215
6.9	Setting up an administrator passcode	216
6.10	Viewing remaining memory for logging data	218
6.11	Downloading logged data from the memory of a thermometer readout	218
6.12	Saving probe coefficients to a configuration file	219
6.13	Using an IR dongle with LogWare	220
6.14	Changing logging label names	221
6.15	Unlocking a demo version of LogWare	222
6.16	Setting up LogWare to read multiple scales	223
7	Files	227
7.1	Configuration files	227

7.1.1	Log session configuration file	227
7.1.2	Alarm configuration file	228
7.1.3	Conversion configuration files.	228
7.2	Data files	229
7.2.1	Output file	229
7.2.2	Download file	231
7.2.3	Alarm Log File	231
8	Download process	233
8.1	1523/1524 Handheld and 1529 Chub-E4 download process	233
8.2	1560 Black Stack download process.	234
8.3	1575A Super-Thermometer download process	235
8.4	1590 Super-Thermometer II download process	236
9	Conversion settings dialogs	239
9.1	ITS-90 Settings dialog	239
9.1.1	Low Range tab.	240
9.1.2	High Range tab	240
9.2	IPTS-68 settings dialog	242
9.3	Callendar-Van Dusen Settings dialog	243
9.4	PT100 Settings dialog	244
9.5	High-Temperature PRT Settings dialog	245
9.6	Polynomial Settings dialog.	246
9.7	W(T90) Settings dialog	247
9.8	Thermistor Settings dialog.	247
9.8.1	Thermistor T(R) Settings dialog.	248
9.8.2	Thermistor R(T) Settings dialog.	249
9.9	Thermocouple Polynomial Settings dialog	250
9.10	Thermocouple Table Settings dialog.	251
9.11	Thermocouple Type Settings dialog	252
9.12	Thermocouple Type Polynomial Settings dialog	254
10	Troubleshooting	257
10.1	Technical support	257
10.2	Troubleshooting common problems	259
10.3	Communication error	261

Figures and Tables

Figure 1	9935 LogWare II Main Display	9
Figure 2	File Toolbar	10
Figure 3	Datalog Toolbar	10
Figure 4	Graph Toolbar	10
Figure 5	Tools Toolbar	11
Figure 6	Help Toolbar	11
Figure 7	File Menu	13
Figure 8	Datalog Menu	15
Figure 9	Log Session Settings dialog	17
Figure 10	Log Session Settings - Thermometer tab.	19
Figure 11	Log Session Settings - Log Session tab	22
Figure 12	Log Session Settings - Channels Tab	24
Figure 13	Delayed Start Date/Time Dialog	26
Figure 14	Alarm Settings dialog	27
Figure 15	Statistics Settings dialog	30
Figure 16	Real-time log session window	31
Figure 17	Channels popup menu	34
Figure 18	Graph popup menu	36
Figure 19	Current Alarm Settings dialog	37
Figure 20	Trace color dialog	38
Figure 21	File Viewer window	39
Figure 22	Statistics dialog.	40
Figure 23	Previous Log Session window	41
Figure 24	Channels popup menu	43
Figure 25	Graph popup menu	44
Figure 26	Graph zooming	45
Figure 27	Download menu	47
Figure 28	Select Model dialog	48
Figure 29	1523 Download Options dialog	50
Figure 30	1524 Download Options dialog	52
Figure 31	1529 Chub-E4 Download Options dialog	54
Figure 32	1560 Black Stack Download Options dialog.	55
Figure 33	1575A ‘Super Thermometer’ Download Options dialog	57
Figure 34	1590 Super-Thermometer II Download Options dialog	59
Figure 35	1523 Log Status dialog.	60
Figure 36	1524 Log Status dialog.	61
Figure 37	1529 Chub-E4 Log Status dialog	62
Figure 38	1560 Black Stack Log Status dialog	63
Figure 39	Tools menu	65
Figure 40	Select Model dialog	67

Figure 41	1523 Logging Label dialog	69
Figure 42	1524 Logging Labels dialog	70
Figure 43	1529 Chub-E4 Logging Labels dialog	71
Figure 44	1523/1524 Settings dialog	73
Figure 45	1523/1524 Settings – Sampling tab	74
Figure 46	1523/1524 Settings – Channels tab	75
Figure 47	1523/1524 Settings – Calibration tab	76
Figure 48	1523/1524 Settings – Misc tab	77
Figure 49	1523/1524 PRT/Thermistor Channel Settings dialog	78
Figure 50	1523/1524 PRT/Thermistor Channel Settings – Probe tab.	79
Figure 51	1523/1524 PRT/Thermistor Channel Settings – Misc tab	80
Figure 52	1523/1524 Thermocouple Channel Settings dialog	81
Figure 53	1523/1524 Thermocouple Channel Settings – Probe tab	82
Figure 54	1523/1524 Thermocouple Channel Settings – Misc tab	84
Figure 55	1529 Chub-E4 Settings dialog	85
Figure 56	1529 Chub-E4 Settings - Filtering tab	86
Figure 57	1529 Chub-E4 Settings - Sampling tab	87
Figure 58	1529 Chub-E4 Settings - Measurement Tab	88
Figure 59	1529 Chub-E4 Settings - Misc tab	89
Figure 60	1529 Chub-E4 Settings - Channels tab.	89
Figure 61	1529 Chub-E4 PRT/thermistor channel settings dialog	90
Figure 62	1529 Chub-E4 PRT/thermistor channel settings - Probe tab.	91
Figure 63	1529 Chub-E4 PRT/thermistor channel settings - Calibration tab.	93
Figure 64	1529 Chub-E4 Thermocouple channel settings dialog.	94
Figure 65	1529 Chub-E4 thermocouple channel settings - Probe tab	95
Figure 66	1529 Chub-E4 thermocouple channel settings - Probe tab	96
Figure 67	1560 Black Stack Settings dialog	97
Figure 68	1560 Black Stack settings - Filtering tab.	98
Figure 69	1560 Black Stack Settings – Sampling tab.	99
Figure 70	1560 Black Stack settings - Measurement tab	100
Figure 71	1560 Black Stack settings - Modules tab	101
Figure 72	2560 SPRT Module Settings dialog	102
Figure 73	2560 SPRT Module Settings - General tab	103
Figure 74	2560 SPRT Module Settings - Channels tab	103
Figure 75	2560 SPRT Module Settings - Calibration tab	104
Figure 76	2561 HPRT Module Settings dialog	105
Figure 77	2561 HPRT Module Settings - General tab.	106
Figure 78	2561 HPRT Module Settings - Channels tab	106
Figure 79	2561 HPRT Module Settings - Calibration tab	107
Figure 80	2562 ‘PRTS’ Module Settings dialog.	108
Figure 81	2562 PRTS Module Settings - General tab	109
Figure 82	2562 ‘PRTS’ Module Settings - Channels tab	110
Figure 83	2562 ‘PRTS’ Module Settings - Calibration tab	110
Figure 84	2563 STHR Module Settings dialog	111
Figure 85	2563 STHR Module Settings - General tab.	113

Figure 86	2563 STHR Module Settings - Channels tab	113
Figure 87	2563 STHR Module Settings - Calibration tab	114
Figure 88	2564 THRS Module Settings dialog	115
Figure 89	2564 THRS Module Settings - General tab	116
Figure 90	2564 THRS Module Settings - Channels tab	117
Figure 91	2564 THRS Module Settings - Calibration tab	117
Figure 92	2565 PTC Module Settings dialog	119
Figure 93	2565 PTC Module Settings - Channels tab	120
Figure 94	2565 PTC Module Settings - Calibration tab	121
Figure 95	2566 TCS Module Settings dialog	122
Figure 96	2566 TCS Module Settings - Channels tab	123
Figure 97	2566 TCS Module Settings - Calibration tab	124
Figure 98	2566 TCS Module Settings - Inputs 1 to 6 tab	125
Figure 99	2566 TCS Module Settings - Inputs 7 to 12 tab	126
Figure 100	2567 RTD Module Settings dialog	127
Figure 101	2567 RTD Module Settings - General tab	128
Figure 102	2567 RTD Module Settings - Channels tab	128
Figure 103	2567 RTD Module Settings - Calibration tab	129
Figure 104	2568 RTDS Module Settings dialog	130
Figure 105	2568 RTDS Module Settings - General tab	131
Figure 106	2568 RTDS Module Settings - Channels tab	132
Figure 107	2568 RTDS Module Settings - Calibration tab	132
Figure 108	1560 Black Stack PRT/Thermistor Channel Settings dialog	133
Figure 109	1560 Black Stack PRT/Thermistor Channel Settings - Probe tab	134
Figure 110	1560 Black Stack Thermocouple Channel Settings dialog	135
Figure 111	1560 Black Stack PRT/Thermistor Channel Settings - Probe tab	136
Figure 112	1575A Super-Thermometer Settings dialog	137
Figure 113	1575A Super-Thermometer Settings - Filtering tab	139
Figure 114	1575A Super-Thermometer Settings - Sampling tab	140
Figure 115	1575A Super-Thermometer Settings - Measurement tab	142
Figure 116	1575A Super-Thermometer Settings - Channels tab	144
Figure 117	1575A Super-Thermometer Settings - Calibration tab	145
Figure 118	1575A Super Thermometer/2575 Mighty Mux Channel Settings dialog	146
Figure 119	1575A Super-Thermometer/2575 Mighty Mux Channel Settings - Probe tab	147
Figure 120	1590 Super-Thermometer II Settings dialog	149
Figure 121	1590 Super-Thermometer II Settings - Filtering tab	150
Figure 122	1590 Super-Thermometer II Settings - Sampling tab	151
Figure 123	1590 Super-Thermometer II Settings - Measurement tab	153
Figure 124	1590 Super-Thermometer II Settings - Channels tab	155
Figure 125	1590 Super-Thermometer II Settings - Calibration tab	156
Figure 126	1590 Super-Thermometer II/2590 'Mighty Mux II' Channel Settings dialog	157

Figure 127	1590 Super-Thermometer II/2590 ‘Mighty Mux II’ Channel Settings - Probe tab	158
Figure 128	1594A/1595A Hart Super-Thermometer settings	160
Figure 129	1594A/1595A Hart Super-Thermometer Filtering tab	161
Figure 130	1594A/1595A Hart Super-Thermometer Display tab	162
Figure 131	1594A/1595A Hart Super-Thermometer Timing tab	163
Figure 132	1594A/1595A Hart Super-Thermometer Scanning tab.	164
Figure 133	1594A/1595A Hart Super-Thermometer Channel setup tab	165
Figure 134	1594A/1595A Hart Super-Thermometer Probe settings dialog	166
Figure 135	1594A/1595A Hart Super-Thermometer Channel settings	168
Figure 136	1594A/1595A Hart Super-Thermometer Rs1 settings	170
Figure 137	File menu	173
Figure 138	Datalog menu	175
Figure 139	Download Menu	177
Figure 140	View Menu	178
Figure 141	Graph Menu.	180
Figure 142	Graph Settings - Titles tab	181
Figure 143	Graph Settings - Axis tab	183
Figure 144	Graph Settings - Fonts tab	185
Figure 145	Graph Settings - Trends tab	186
Figure 146	Graph Settings - Background tab	189
Figure 147	Graph Settings - Legend tab	190
Figure 148	Tools Menu	192
Figure 149	LogWare Options Dialog	193
Figure 150	LogWare Options - General tab.	194
Figure 151	LogWare Options - Folders tab	197
Figure 152	LogWare Options - Tools tab	198
Figure 153	LogWare Options - Admin. Passcode tab.	199
Figure 154	Enter Passcode dialog.	200
Figure 155	Change Passcode dialog	201
Figure 156	Window menu	202
Figure 157	Help menu	203
Figure 158	About LogWare dialog	203
Figure 159	Excel XP Text Import Wizard – Step 1 of 3	211
Figure 160	Excel XP data.	212
Figure 161	Excel XP Text Import Wizard – Step 2 of 3	212
Figure 162	Excel XP Chart Wizard – Step 1 of 4 – Chart Type	213
Figure 163	Excel XP graph	214
Figure 164	IR dongle	220
Figure 165	Unlock Code dialog.	223
Figure 166	Output file example.	230
Figure 167	Download file example	231
Figure 168	Alarm log file example	232
Figure 169	Downloaded Files dialog	234
Figure 170	ITS-90 Settings dialog	239

Figure 171 ITS-90 Settings – Low Range tab	240
Figure 172 ITS-90 Settings – High Range tab	241
Figure 173 IPTS-68 Settings dialog.	242
Figure 174 Callendar-Van Dusen Settings dialog.	243
Figure 175 PT100 Settings dialog	244
Figure 176 High-Temp PRT Settings dialog	245
Figure 177 Polynomial Settings dialog	246
Figure 178 W(T90) Settings dialog	247
Figure 179 Thermistor T(R) Settings dialog	248
Figure 180 Thermistor R(T) Settings dialog	249
Figure 181 Thermocouple Polynomial Settings dialog	250
Figure 182 Thermocouple Table Settings dialog	251
Figure 183 Thermocouple Polynomial Settings dialog	252
Figure 184 Thermocouple Type Polynomial Settings dialog	254
Table 1 Troubleshooting Common Problems	259

1 Introduction

1.1 What is LogWare

The 9935 LogWare II Software (LogWare) is a real-time data acquisition, storage and analysis program designed specifically for use with the following Hart Scientific's multi-channel digital thermometer readouts:

- 1523 and 1524 Handheld Thermometer readouts
- 1529, 1529-R, 1529-T Chub-E4 Thermometer readouts
- 1560 *Black Stack* Thermometer readout
 - ♦ 2560 SPRT Module
 - ♦ 2561 HTPRT Module
 - ♦ 2562 PRTS Module
 - ♦ 2563 STHR Module
 - ♦ 2564 THRS Module
 - ♦ 2565 PTC Module
 - ♦ 2566 TCS Module
 - ♦ 2567 RTD Module
 - ♦ 2568 RTDS Module
 - ♦ 3560 Communication Module
- 1575 Super Thermometer readout (**real-time logging only!**)
 - ♦ 2575 Mighty Mux Multiplexer
- 1575A Super Thermometer readout
 - ♦ 2575 Mighty Mux Multiplexer
- 1590 Super Thermometer II readout
 - ♦ Up to five 2590 Mighty Mux II Multiplexers
- 1594A/1595A Super-Thermometer Readout
 - ♦ Up to 2 Model 2590 Mighty Mux II Multiplexers

LogWare also provides utilities for:

- Downloading data that has been logged to the thermometer readout's logs or memory.
- Setting logging labels in the thermometer readout (for thermometer readouts that support this feature).
- Synchronizing thermometer readout's clock to the computer's clock.

- Viewing and editing many thermometer readout, module and probe settings including probe coefficients.

This manual is intended to document the use and features of LogWare. *To ensure correct results, verify that the thermometer readout is properly configured before collecting data. Refer to the individual thermometer readout's User's Guide for specifics pertaining to the instrument. Follow all warnings and cautions documented in the readout's User's Guide.*

1.2 License agreement

For information and details regarding the software End User License Agreement (EULA), refer to the online help file's License Agreement topic.

1.3 Features

LogWare provides an easy-to-use Windows® interface to acquire, store and analyze data from Hart Scientific's multi-channel digital thermometer readouts. LogWare also provides a method to download readings that are stored in the logs or memory of thermometer readouts. After being acquired or downloaded, the readings are written to a file that can be opened by LogWare, a spreadsheet or other analysis software.

LogWare's data logging features include:

- Quick and easy configuration
- Ability to save and open configuration files
- Selectable sample interval from 1 second to 60 hours (depending on thermometer readout)
- Selectable scan sequence interval from 1 minute to 60 hours
- Selectable delayed start date/time for off-hours logging
- Selectable end log options (to stop session after specified time or number of readings)
- Acquires readings to ASCII text files which can be opened, graphed and printed
- Graphs readings as they are being acquired
- Ability to acquire data in multiple scales simultaneously
- Maintains statistical data for all channels
- Selectable high and low alarm settings with user defined events
 - ♦ Can be configured to play .WAV file or launch an application (such as pager software) when an alarm trips
- User-customizable graph settings including titles, legend, grid lines and statistical and trend lines

LogWare's data analysis features include:

- Can open both real-time log files and downloaded files
- Spreadsheet-style grid to view all acquired readings
- Graph zooming capability
- User-customizable graph settings including titles, legend, grid lines and statistical and trend lines
- Printing graph and acquired readings
- Can have multiple windows open simultaneously
- Can superimpose up to 10 sets of data on the same graph

LogWare also provides utilities that allow access to the following digital thermometer readout features (as applicable):

- Downloading data stored in the thermometer readout's logs and/or memory
- Viewing log status and clearing log data
- Synchronizing thermometer readout's date/time with computer's clock
- Changing thermometer readout-specific parameters (filtering, sampling, measurement, calibration constants, etc.)
- Changing probe-specific parameters (serial numbers, conversion type and coefficients, calibration dates, cold junction settings, etc.)

New Features in Version 1.3

Version 1.3 has added support for Models 1594A and 1595A Super-Thermometer readouts. Features supported by these new thermometer readouts include:

- Real-Time logging
- Thermometer settings
- Synchronize Date/Time

Version 1.2

Version 1.2 has added support for Models 1523 and 1524 Handheld Thermometer readouts. Also, previously logged data files containing more than one data set (up to a maximum of 10 data sets) can now be opened.

Version 1.1

A list of features and enhancements in Version 1.1 follows

- **Simultaneous Logging of All Channels in Real Time** - A feature has been added to allow readings to be taken on all enabled channels of a real-time log session simultaneously (as fast as possible).

- **Download Logged Data from Serial Port on 3560 Communication Module** - A feature has been added to allow logged readings to be downloaded to the serial port on the Model 3560 Communication Module on a Model 1560 Black Stack thermometer readout.
- **Write All Real-Time Data to a Single File** - A new feature has been added to allow data from all channels of a real-time log session to be written to the same output file. In these output files, a new column has been added for the channel number.



Note: Although LogWare is still able to open these files, it does not separate the data for each channel. All readings are plotted on the graph as if all data was from a single channel.

- **Branding/Logo change** - The logos and graphics displayed in LogWare have been updated.

Also, various features have been added to the software to aid in troubleshooting issues related to real-time log sessions and downloading logged data.

1.4 Requirements

LogWare is designed to work with Hart Scientific's multi-channel digital thermometer readouts. Refer to Section 1.1, What is LogWare, for a complete list of supported thermometer readouts.

1.4.1 Communications requirements

The thermometer readout being used with LogWare must be connected to an RS-232 serial port (COM port) on a computer using a "null modem" cable. For more information on this type of cable, refer to the thermometer readout's User's Guide. The baud rate of the thermometer readout must be set to match the baud rate selected in LogWare. LogWare supports a wide choice of baud rate settings, depending on the thermometer readout being used.

Selected features of LogWare can also utilize an IR dongle with thermometer readouts that have an IR port for wireless communications.

1.4.2 Computer hardware requirements

The following minimum computer configuration is required to use LogWare:

- IBM Compatible 486 PC or better with 8MB RAM (Pentium® class processor with 16MB RAM or more recommended)
- VGA monitor or better
- CD-ROM drive for installation
- 1 - RS-232 serial (COM) port
- **Optional** - IR dongle

- Minimum of about 25MB of disk space for installation (additional space is required as data is acquired or downloaded)

1.4.3 Computer software requirements



Note: To use LogWare, Windows login accounts must be granted full access to the folder to which the software is installed and all subfolders. If the folder has read-only access, errors will occur when LogWare attempts to access its configuration files.

The LogWare Software requires one of the following operating systems:

- Windows® 95/98/ME
- Windows® NT with Service Pack 4 or later installed
- Windows® 2000
- Windows® XP
- Windows® Vista

1.4.4 Firmware requirements

The thermometer readouts supported by LogWare must have a certain firmware version installed to use the Thermometer Settings option in the Tools menu and the Download Logged Data option in the Download menu. Prior versions of firmware do not fully support these features. Real-time logging can be performed with any firmware version. The required firmware versions are indicated below:

- Model 1523 and 1524 – Requires firmware version 1.0 or later
- Model 1529, 1529-R and 1529-T Chub-E4 – Requires firmware version 1.35 or later
- Model 1560 Black Stack – Requires firmware version 1.31 or later
- Model 1575 Super-Thermometer – **NOT SUPPORTED BY TOOLS FEATURES!**
- Model 1575A Super-Thermometer – Requires firmware version 5.03n or 5.03o or later
- Model 1590 Super-Thermometer II – Requires firmware version 2.03n or 2.03o or later
- Model 1594A, 1595A – Requires firmware version 1.0 or later

1.5 Installation



Note: Before installing LogWare, be sure to read the License Agreement and Warranty.

A backup should always be made of your hard disk drive and registry before installing any software, and all running applications should be closed.



Note: To successfully install LogWare, you must be logged into Windows as an Administrator.

Installing LogWare

1. Insert the LogWare CD-ROM into your CD-ROM drive. The Setup program should run automatically.
2. Follow the on-screen instructions to install LogWare.
3. When prompted, enter the product serial number. The serial number is provided on a sticker typically located on the CD pouch on the inside back cover of the LogWare User's Guide.
4. After all files have been installed, a program group is created with icons for the software, the Help file, the README.TXT file, and an icon to uninstall the software. The installation process is complete.
5. The README.TXT file can be displayed by checking the appropriate check box before the setup program finishes. Read this file for important information that was not available when the User Guide was printed.

Uninstalling LogWare

To uninstall the software, use the Windows® Add/Remove Programs icon in the Control Panel. Follow the instructions given. The program files that were installed to your system are removed. Any files that were created or saved by users will not be removed from the system. These files, along with the folders that these files are in, must be manually deleted using Windows® Explorer to remove them from your system.

1.6 Running LogWare

The following sections describe how to get LogWare up and running quickly.

1.6.1 Getting started

The following information is intended to provide a brief overview on how to get LogWare up and running quickly.

To use this software:

1. Click the Start button, then locate and click the icon created by the setup process.
2. The software will start with the Log Session Settings dialog displayed. If the Log Session Settings dialog is not displayed when the software starts, select the Setup option in the Datalog menu to display the Log Session Settings dialog.



Note: In order to make the Log Session Settings dialog appear automatically every time LogWare starts, check the “Show Log Session Settings dialog on start-up” check box on the General tab of the Options dialog. The Options dialog can be displayed by selecting the Options option in the Tools menu. Please refer to Section 5.6.4, Options, for more details.

3. To start a log session, fill in the information on the Log Session Settings dialog. The following information must be entered before a log session can start:

Thermometer:

- Select the model of the thermometer readout to use, enter the serial number and select the session temperature scale from the available choices. The reading resolution may also be selected.
- Select the COM port that the thermometer readout is connected to and choose the baud rate to use.
- Power up the thermometer readout and make sure it is connected to the selected COM port and set to the selected baud rate. Click the Test Communication button to make sure LogWare can communicate with the thermometer readout.

Log session:

- Select the interval between readings.
- **Optional** - Enable and select the interval between scan sequences.
- To set a delayed start date/time, click the Delayed Start button and choose the appropriate date and time for the log session to start. Otherwise, the log session will begin when the Start button is clicked.
- Choose the appropriate end log session option and set the other end log settings as required.

Channels:

- Query the channels from the thermometer readout by clicking the Query Channels button.

- Enable or disable each channel as needed by selecting each channel in the Channels list and checking or unchecking the Acquire data from this channel box.
- **Optional** - To customize the statistical data for each channel, click the Statistics Settings button.
- **Optional** - To configure alarms for each channel, click the Alarm Settings button.
- Enter the channel name, output filename, and select the trace color for each channel.

These settings can be saved to a configuration file by clicking the Save As button. To open the current configuration file every time the software runs, click the Set as Default button.

After entering this information, click the Start button to start the real-time log session.

1.7 Using LogWare

The following sections describe some of the main features of LogWare.

1.7.1 LogWare main display

The main display consists of the menus, toolbar, a workspace area and a status bar.

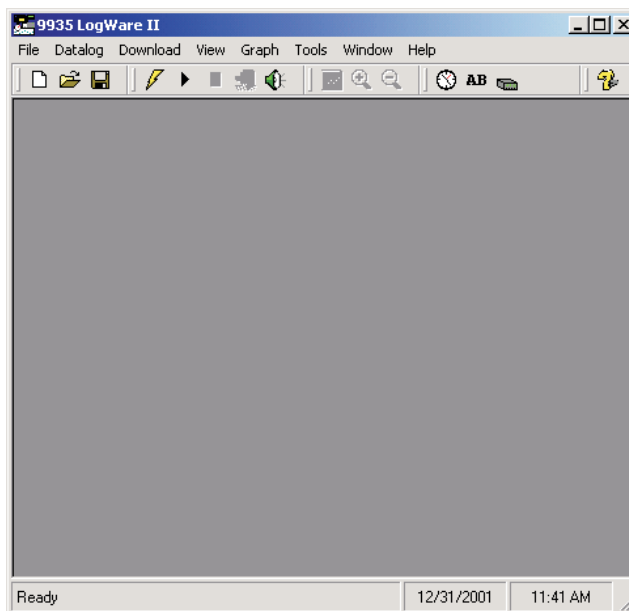


Figure 1 9935 LogWare II Main Display

The menus allow access to the various features of the software. A general description of the options in each menu is given below. These menus are described in more detail in Section 5, Menus:

- **File** - Open and save configuration files, print options and exit the software.
- **Datalog** - Setup, Start, Stop and Close real-time log sessions.
- **Download** - Download logged readings and view log status.
- **View** - View alarm log, reading history and open previously logged data files.
- **Graph** - Change graph settings, zoom in and out, and print the current graph.
- **Tools** - Synchronize thermometer readout date/time, access logging labels, configure thermometer readout and probe settings, change LogWare options.
- **Window** - Rearrange and switch between open windows.
- **Help** - Display help topics and About dialog.

The toolbar allows quick access to the most common menu functions. The toolbar is divided into sections that correspond to related functions. Toolbar sections may be rearranged as desired by clicking on the vertical bar on the left side of the section and dragging it to the desired location.

The following toolbars are available:

File Toolbar



Figure 2 File Toolbar

New - Create a new real-time log session configuration

Open - Open a real-time log session configuration file

Save As - Save the current real-time log session configuration to a configuration file

Datalog Toolbar



Figure 3 Datalog Toolbar

Setup - Display the Log Session Settings dialog to configure a real-time log session

Start - Start a real-time log session

Stop - Stop a real-time log session

Close - Close the real-time log session window

Mute All Alarms - Mute/unmute all audible real-time log session alarms

Graph Toolbar



Figure 4 Graph Toolbar

Graph Settings - Change graph settings, add trend lines, alarm lines, titles, captions, etc.

Zoom In - Zoom in to view a section of the graph

Restore (Zoom Out) - Restore the graph to its original zoom setting

Tools Toolbar



Figure 5 Tools Toolbar

Synchronize Date/Time - Synchronize thermometer readout's date/time with computer

Logging Labels - View/change logging labels

Thermometer Settings - View/change thermometer readout settings

Help Toolbar



Figure 6
Help Toolbar

Help - Open the LogWare Help file

Status Bar

The status bar indicates the function that the software is currently performing as well as the current system date and time.



***Note:** The date and time displayed are based on the computer's clock settings and can be modified through the Windows® Control Panel.*

1.8 LogWare options

LogWare has many features and settings that can be customized by the user to allow LogWare to operate in certain ways. Many of these settings can also be protected by an administrator passcode to prevent unauthorized persons from changing these settings. For more information on these options, please refer to Section 5.6.4, Options.

2 Real-time logging

The main purpose of LogWare is to perform real-time data acquisition from Hart Scientific multi-channel digital thermometer readouts. This section describes in detail how to configure, start and stop a real-time log session. This section also describes the features available during a real-time log session.

2.1 File menu

The File menu contains some options for creating a new real-time log session and for opening and saving real-time log session configuration files.

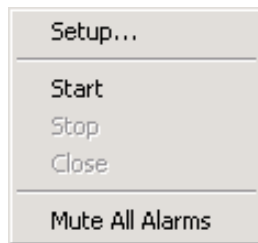


Figure 7 File Menu

2.1.1 New

The New menu option can be used to clear the current real-time log session settings and display the Log Session Settings dialog.

The  button on the toolbar corresponds to the New menu option.

When selecting the New option, a prompt is displayed to confirm that the current real-time log session settings (if any) will be reset. To abort resetting the current real-time log session settings, click the No button. Otherwise, click the Yes button to reset the current log session settings. The Log Session Settings dialog will be displayed. Please refer to Section 2, Real-Time Logging, for more information on configuring a real-time log session.

2.1.2 Open

The Open option allows a real-time log session configuration file to be opened. If the same basic configuration is frequently used to perform real-time logging, saving the log session settings to a configuration file can greatly reduce the time it takes to configure and start a real-time log session.

The  button on the toolbar corresponds to the Open menu option.



Note: When opening a configuration file, LogWare attempts to communicate with the thermometer readout to verify that the settings saved in the configuration file are compatible with the current thermometer readout configuration. Therefore, it is necessary to have the thermometer readout connected and powered on in order to successfully open a configuration file.

When selecting the Open option, the Windows® Open dialog is displayed. The \Config folder should be selected by default (or the folder that is specified for configuration files on the LogWare Options dialog). Real-time log session configuration files should have a **.rls** extension. To cancel opening a configuration file, click the Cancel button. Otherwise, select the configuration file to open and click the Open button. For more details about the real-time log session configuration file, refer to Section 7.1.1, Log Session Configuration File. Real-time log session configuration files can also be opened by clicking the Open button on the Log Session Settings dialog.

LogWare then reads the log session settings from the configuration file and verifies that the configuration of the currently connected thermometer readout is compatible with the log session settings. If there are any conflicts, or if LogWare is unable to communicate with the thermometer readout, appropriate messages are displayed.

Once the configuration file has been read in, the Log Session Settings dialog is displayed. If LogWare was unable to communicate with the thermometer readout or any other error occurred, the log session settings may be incomplete.

2.1.3 Save as

The Save As option allows the current real-time log session settings to be saved to a configuration file. If the same basic configuration is frequently used to perform real-time logging, saving the log session settings to a configuration file can greatly reduce the time it takes to configure and start a real-time log session. Before attempting to save a real-time log session configuration, use the Log Session Settings dialog to configure the log session.

The  button on the toolbar corresponds to the Save As menu option.

When selecting the Save As option, the Windows® Save As dialog is displayed. The \Config folder should be selected by default (or the folder that is specified for configuration files on the LogWare Options dialog). Real-time log session configuration files are saved with a **.rls** extension. To cancel saving a configuration file, click the Cancel button. Otherwise, enter the filename for the configuration file and click the Save button. For more details about the real-time log session configuration file, refer to Section 7.1.1, Log Session Configuration File. Real-time log session configuration files can also be saved by clicking the Save As button on the Log Session Settings dialog.

2.2 Datalog menu

The Datalog menu is used to access the Log Session Settings dialog, to start, stop and close the real-time log session and to mute audible alarms.

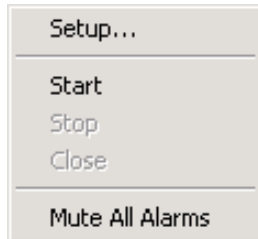


Figure 8 Datalog Menu

2.2.1 Setup


The Setup option displays the Log Session Settings dialog. This dialog is used to configure a real-time log session. If a log session has already been configured or a real-time log session configuration file has already been opened, this dialog displays the current log session settings.

The  button on the toolbar corresponds to the Setup menu option.

Please refer to Section 2.3, Log Session Settings Dialog, for more details on configuring a real-time log session.

2.2.2 Start

The Start option starts a real-time log session. This option is disabled until a log session has been completely configured.

The  button on the toolbar corresponds to the Start menu option.

Please refer to Section 2, Real-Time Logging, for more details on a real-time log session.

2.2.3 Stop

The Stop option stops the current real-time log session. This option is only enabled when a real-time log session is running. A log session may be stopped at any time by selecting this option.

The  button on the toolbar corresponds to the Stop menu option.

When selecting the Stop option, a message is displayed to confirm stopping the log session.

2.2.4 Close

The Close option closes the current real-time log session window. This option is only enabled once the real-time log session has stopped and the log session window is still open.

The  button on the toolbar corresponds to the Close menu option.

2.2.5 Mute all alarms

The Mute All Alarms option can be used to mute any audible alarms associated with the current real-time log session. When the alarms are muted, a check mark appears next to this option. Muting the alarms does not prevent alarm events from happening.

The  button on the toolbar corresponds to the Mute All Alarms menu option.

When alarms are muted, the toolbar button appears as .

2.3 Log session settings dialog

The Log Session Settings dialog is used to configure the real-time log session settings. These settings include the thermometer readout to acquire data from, communications settings, log session sample interval and start and end options, and thermometer readout channel configuration. This dialog can be displayed

by selecting the Setup option in the Datalog menu or by clicking the .

button on the toolbar. This dialog is also displayed after opening a log session configuration file.

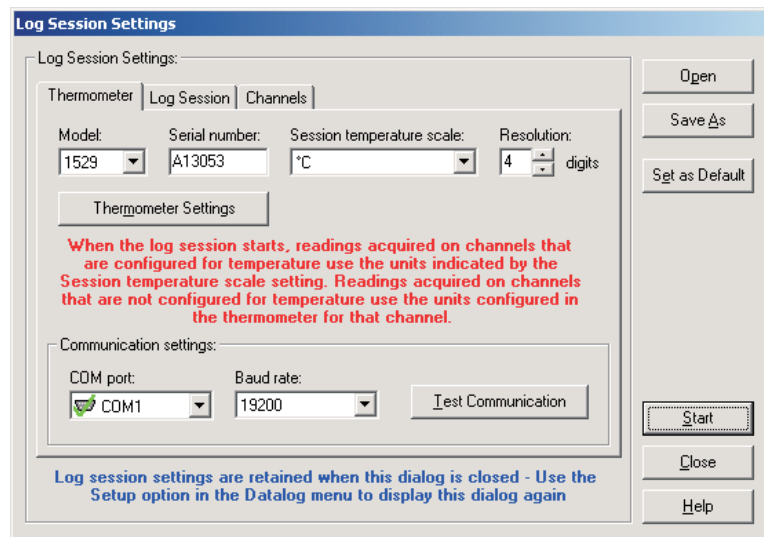


Figure 9 Log Session Settings dialog



Note: This dialog displays the current real-time log session settings. Clicking the Close button on this dialog does **not** cause these settings to be lost. To display this dialog again, select the Setup option in the Datalog menu.

This dialog has three tabs: Thermometer, Log Session and Channels. All of the information on each of these tabs must be configured before a real-time log session can be started. The following sections describe in detail each of the settings on these tabs.

Once the log session has been completely configured, the settings can be saved to a configuration file by clicking the Save As button. If these same settings may be used again at a later date, saving the settings to a configuration file greatly reduces the amount of time required to get a log session started.

When clicking the Save As button, the Windows® Save As dialog is displayed. The \Config folder should be selected by default (or the folder that is specified for configuration files on the LogWare Options dialog). Real-time log session configuration files are saved with a **.rls** extension. To cancel saving a configuration file, click the Cancel button. Otherwise, enter the filename for the configuration file and click the Save button. For more details about the real-time log session configuration file, refer to Section 7.1.1, Log Session Configuration File. Real-time log session configuration files can also be saved by selecting the Save As option in the File menu.

A log session settings configuration file that has been saved previously can be opened by clicking the Open button.



Note: *When opening a configuration file, LogWare attempts to communicate with the thermometer to verify that the settings saved in the configuration file are compatible with the current thermometer configuration. Therefore, it is necessary to have the thermometer connected and powered on in order to successfully open a configuration file.*

When clicking the Open button, the Windows® Open dialog is displayed. The \Config folder should be selected by default (or the folder that is specified for configuration files on the LogWare Options dialog). Real-time log session configuration files should have a **.rls** extension. To cancel opening a configuration file, click the Cancel button. Otherwise, select the configuration file to open and click the Open button. For more details about the real-time log session configuration file, refer to Section 7.1.1, Log Session Configuration File. Real-time log session configuration files can also be opened by selecting the Open option in the File menu.

LogWare then reads the log session settings from the configuration file and verifies that the configuration of the currently connected thermometer readout is compatible with the log session settings. If there are any conflicts, or if LogWare is unable to communicate with the thermometer readout, appropriate messages are displayed.


If LogWare was unable to communicate with the thermometer readout or any other error occurred, the log session settings may be incomplete.

LogWare has a feature which allows a log session configuration file to be defined as the default configuration file. The default configuration file is automatically loaded every time LogWare is started. To set the current configuration as the default configuration, select the Set as Default button. If the current configuration has not been saved yet, a prompt is displayed indicating that the configuration must be saved to a file before it can be set as the default. The next time LogWare is started, the default configuration file will be loaded.



Note: *Real-time log session configuration files contain information about the current thermometer readout configuration. If the thermometer readout configuration changes, the configuration file needs to be updated also.*

Once the log session settings are complete, click the Start button to start a log session. Please refer to Section 2.4, Real-Time Log Session, for more details on a real-time log session.

If the log session does not need to be started yet, click the Close button. The Log Session Settings dialog is closed but all settings are retained. To display the Log Session Settings dialog again, select the Setup option in the Datalog menu or click the  button on the toolbar.

Click the Help button to display the appropriate help topic for this dialog.

2.3.1 Thermometer tab

The Thermometer tab allows the thermometer readout that will be used to acquire data to be identified and configured.

The screenshot shows the 'Thermometer' tab of the 'Log Session Settings' dialog. At the top, there are three tabs: 'Thermometer', 'Log Session', and 'Channels'. The 'Thermometer' tab is selected. Below the tabs, there are four input fields: 'Model' (a dropdown menu showing '1529'), 'Serial number' (a text box containing 'A13053'), 'Session temperature scale' (a dropdown menu showing '°C'), and 'Resolution' (a spinner box showing '4' followed by the text 'digits'). Below these fields is a button labeled 'Thermometer Settings'. Underneath the button is a red text warning: 'When the log session starts, readings acquired on channels that are configured for temperature use the units indicated by the Session temperature scale setting. Readings acquired on channels that are not configured for temperature use the units configured in the thermometer for that channel.' Below the warning is a section titled 'Communication settings:' which contains two dropdown menus: 'COM port' (showing 'COM1') and 'Baud rate' (showing '19200'). To the right of these is a button labeled 'Test Communication'.

Figure 10 Log Session Settings - Thermometer tab

Use the Model drop-down list to select the type of thermometer readout to use to acquire data. All other controls on this tab will be disabled until a valid model is selected.

Enter the serial number of the thermometer readout to use in the Serial number box. The thermometer readout's serial number should be located on a label, generally on the back or bottom of the thermometer readout.



Note: Optionally, you may use the Thermometer Settings button to automatically determine the serial number of the thermometer readout and fill in this box.

Select the scale of the readings to be acquired from channels that are configured to read temperature using the Session temperature scale drop-down list. LogWare supports the following temperature scales:

- °C
- °F
- K (Kelvin)

When a log session starts, LogWare will automatically set the thermometer readout's system units to use the selected scale. However, if the <keep existing> option is selected, LogWare will query the thermometer readout to determine the current system units.









Note: When starting a log session, LogWare checks the conversion setting for each enabled channel. If a channel is configured to read temperature, the units that were selected in the Session temperature scale drop-down list are used. If a channel is not configured to read temperature (i.e. resistance or voltage), LogWare automatically determines the non-temperature units for that channel.

Enter the resolution (number of digits after the decimal point) that LogWare should use to format all of the measurements acquired. This setting is common to all channels from which measurements are acquired. Where applicable, LogWare will automatically set the thermometer readout's resolution to this setting when a log session starts. Where not applicable, LogWare will round more precise measurements or add 0's to less precise measurements to obtain the desired format.



Note: The resolution setting should be set prior to configuring the alarm settings for each channel since the alarm values will also use this resolution. Changing the resolution setting after configuring the alarms **does not** reformat the alarm values!!

Select the COM port on the computer to which the thermometer readout is connected using the COM port drop-down list. Only COM ports that the operating system reports as being available can be used. The following symbols are used to show the status of each COM port:

-  - Indicates a COM port that is currently available.
-  - Indicates the COM port that is currently selected by LogWare.
-  - Indicates a COM port that is currently being used by another application.
-  - Indicates an IR port (virtual COM port) that is currently available.
-  - Indicates an IR port (virtual COM port) that is currently selected by LogWare.
-  - Indicates an IR port (virtual COM port) that is currently being used by another application.



Note: For a real-time log session, LogWare requires the use of one COM port. An IR port (virtual COM port) cannot be selected as the port to use.

Select the baud rate that the computer should use to communicate over the COM port with the thermometer readout using the Baud rate drop-down list. The baud rate setting of the thermometer readout must be set to the same value as this setting. For best results, use the maximum allowable baud rate setting when possible.

After the model, COM port and baud rate settings are configured, click the Test Communication button to make sure LogWare is able to communicate with the thermometer readout. At this time, the thermometer readout must be connected to the COM port on the computer using a null modem cable. For more information, please refer to the thermometer readout's User's Guide.

When testing communications with the thermometer readout, there are three possible results:

- The thermometer readout is communicating at the correct baud rate on the selected COM port.
- The thermometer readout is communicating, but possibly at the wrong baud rate or the thermometer readout model does not match the selected model.
- The thermometer readout failed to communicate at all. Check the COM port and baud rate settings, make sure the cable is a null modem cable, and make sure the thermometer readout is plugged in and turned on.

Optional: The Thermometer Settings button may be clicked at this time to gain access to various thermometer readout settings as applicable to the selected thermometer readout. This feature is provided solely as a convenience to check the thermometer readout configuration and/or to configure each channel of the thermometer readout with appropriate conversion settings and probe coefficients. The Test Communication button should be clicked prior to clicking the Thermometer Settings button to make sure the thermometer readout is communicating successfully.



Warning: *When using the Thermometer Settings tool, make sure that the thermometer readout is not currently in any menu. Using the thermometer readout's menus while using the Thermometer Settings tool may cause the thermometer readout to fail to accept new settings.*



Note: *When starting a log session, LogWare will override many of the current thermometer readout settings as needed to acquire data from the configured channels using the specified log session settings!!*



Note: *If the serial number and session temperature scale settings have not been configured prior to clicking the Thermometer Settings button, these settings will be configured automatically after successfully viewing the thermometer readout settings dialog.*

2.3.2 Log Session tab

After the thermometer readout and communication settings are configured, the log session settings may be configured.

The screenshot shows a software window with three tabs: 'Thermometer', 'Log Session', and 'Channels'. The 'Log Session' tab is selected. The settings are as follows:

- Interval between readings:** 1 seconds
- Interval between scan sequences:** 1 minutes
- Start log session:** Three buttons: 'Start logging immediately', 'Delayed Start...', and 'Reset'
- End log session:** After specified time elapses
- After:** 24 hours

Figure 11 Log Session Settings - Log Session tab

Select the interval at which LogWare should take readings on subsequent channels using the Interval between readings settings. The range of possible settings is limited by the capabilities of the selected thermometer readout. For most Hart Scientific multi-channel digital thermometer readouts, the minimum sample interval is 2 seconds.



Note: Due to the latency inherent in the measurement process, a thermometer readout may not be able to take readings at the exact selected interval (i.e. the 1560 thermometer readout allows a minimum interval of 2 seconds, but actually takes about 2.3 seconds to make one measurement). For this reason, using the minimum interval supported by the thermometer readout may cause some duplicate readings to be taken. To avoid this problem, see the Wait until new measurement is ready (when applicable) option on the LogWare Options dialog. Please refer to Section 5.6.4, Options, for more details.

If this interval is set to a value of 0, LogWare will attempt to take simultaneous readings on all enabled channels of the thermometer readout. If this feature is used, you should use the Interval between scan sequences setting to limit how often all channels are sampled.



Note: when using this feature, the Wait until new measurement is ready option on the General tab of the Options dialog **MUST** be disabled (unchecked).

Typically, LogWare is used to take continuous measurements on all enabled channels. However, LogWare has a feature that allows one measurement to be taken on each enabled channel at a set interval. For example, LogWare could be used to monitor the gradients in a liquid bath using 10 probes at various locations in the bath. To measure gradients, readings should be taken simultaneously, or at least as close together as possible, but maybe a set of readings only needs to be taken once every minute, or once every five minutes. This can be accomplished using LogWare by enabling the Interval between scan sequences option. The Interval between scan sequences settings are then enabled.

A scan sequence is defined by LogWare as taking a single reading on each enabled channel. Select the interval to use. LogWare allows this interval to be set anywhere between 1 minute and 60 hours at 1 minute or 1 hour increments. Leaving this feature disabled causes LogWare to begin a new scan sequence immediately following completion of the previous scan sequence.

The Start log session box indicates the current setting for starting a log session. By default, this setting is configured to start a log session immediately after the Start button is clicked. LogWare may optionally be configured to start a log session unattended on any date at any time. To set the start date and time, click the Delayed Start button. The Delayed Start Date/Time dialog is displayed. Please refer to Section 2.3.4, Delayed Start Date/Time Dialog, for details on using this dialog.

Clicking the Reset button resets the start date/time setting. The log session will start immediately after the Start button is clicked.

The End log session settings allow the log session to be configured to automatically stop logging at a certain time or after a certain number of readings. By default, a log session will continue until manually stopped by the user. Select the method to use to end the log session using the drop-down list. Choices are:

- By user (default)
- After specified number of readings
- After specified time elapses

If the After specified number of readings or After specified time elapses option is selected, the end log criteria must be selected. Enter either the number of readings to acquire on each channel or the desired elapsed time setting. The range of choices is from 1 to 1,000,000 readings (10,000 default) or 1 second to 480 hours (1 hour default).



Note: *If either the After specified number of readings or After specified time elapses option is selected and the Interval between scan sequences feature is enabled, the log session will stop automatically when the end log criteria is met and the current scan sequence is complete.*

2.3.3 Channels tab

After the thermometer readout and communication settings are configured, the channels to acquire data from may be configured.

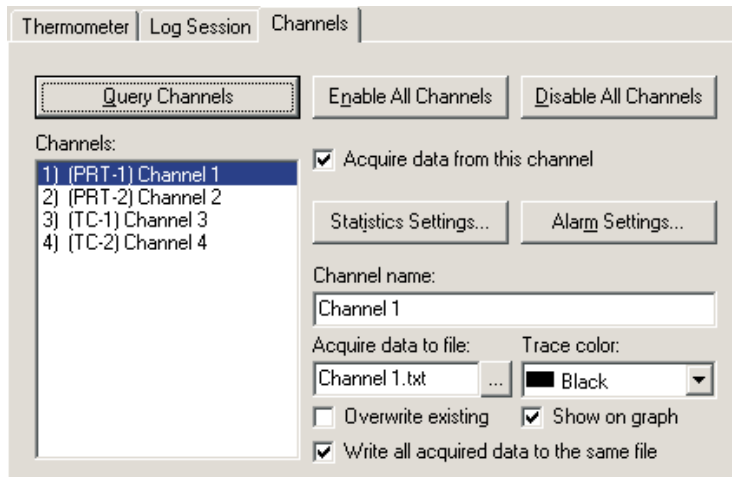


Figure 12 Log Session Settings - Channels Tab

All controls on the Channels tab (with the exception of the Query Channels button) are disabled until the channel information has been queried from the thermometer readout.

Click the Query Channels button to allow LogWare to communicate with the thermometer readout to collect as much information regarding the current channel configuration as possible.

After querying the thermometer readout for channel information, the Channels list displays one item for each channel on the thermometer readout. The channel number, module name, module channel number and currently assigned channel name are displayed in the list.

By default, all channels are enabled, meaning data should be acquired from all channels. The Enable All Channels and Disable All Channels buttons can be used to enable or disable acquiring data from all channels of the thermometer readout respectively.

To configure a channel, select the channel from the Channels list. The selected channel's settings are displayed in the other controls on this tab.

To acquire data from the selected channel, make sure the Acquire data from this channel box is checked. If this box is not checked, all other channel controls are disabled and data will not be acquired for this channel.

Click the Statistics Settings button to display the Statistics Settings dialog. This dialog is used to specify which statistics are to be maintained and displayed for the selected channel. Please refer to Section 2.3.6, Statistics Settings Dialog, for more details on the Statistics Setting dialog. By default, the statistics that are enabled are: Minimum, Maximum, Spread, Average and Standard deviation.



***Note:** Statistics are maintained for each channel independent of all other channels. As a result, each channel has its own settings which must be configured separately.*

Each channel can be configured to trip an alarm if the reading on that channel ever exceeds specific low and/or high values. To enable alarms for the selected channel, click the Alarm Settings button. The Alarm Settings dialog is displayed. This dialog is used to specify the alarm values and the actions associated with each alarm. Please refer to Section 2.3.5, Alarm Settings Dialog, for more details on the Alarm Settings dialog. By default, alarms are disabled.



***Note:** The alarm settings for each channel are independent of all other channels. As a result, each channel must be configured separately.*

Enter the name for the selected channel in the Channel name box. The channel name should be something unique that identifies the purpose, location or other information about the data being acquired on that channel. When a channel name is changed, the Channels list is updated to reflect the change. By default each channel is assigned the generic name 'Channel *n*' (where *n* is the channel number). This name appears in the Channels grid on the real-time log session window while the log session is running.

Enter the filename to which the data acquired from this channel is to be written using the Acquire data to file box. The default filename for each channel is Channel *n*.txt (where *n* is the channel number). The file is saved to the \Output folder or the folder specified for output files on the LogWare Options dialog. The location for this file can be changed by selecting the "..." (Browse) button.

Check the Overwrite existing check box to allow LogWare to automatically overwrite an existing file of the same name in the same folder. If this box is not checked and a file of the same name exists, a prompt is displayed and the option to overwrite the existing file or choose another filename is given.

If you want to write the acquired data for all enabled channels to the same output file, check the Write all acquired data to the same file option. If this option is checked, LogWare will use the filename and path of the first enabled channel for all channels. The header of the output file will then contain information for all enabled channels, and an additional column is added to the file for the channel number for each reading.



Note: This feature should typically only be used if you are using other software to sort and analyze the logged data. The output file created when using this feature can still be opened by LogWare and displayed on-screen, however all readings are displayed on the graph as if they came from the same channel (single trace).

Select the color to use for this channel's trace on the graph using the Trace color drop-down list. LogWare supports 15 different trace colors. By default, each channel is assigned one of the 15 colors.

If the trace for the selected channel is to be shown on the graph, check the Show on graph check box. Once the log session starts, each channel can be shown or hidden at any time. This setting only determines which channels are shown initially when the log session starts.

Continue configuring the other channels as needed by selecting the next channel in the list and following this process again.

2.3.4 Delayed Start Date/Time dialog

The Delayed Start Date/Time dialog allows the real-time log session to be scheduled to start at a future date and/or time.

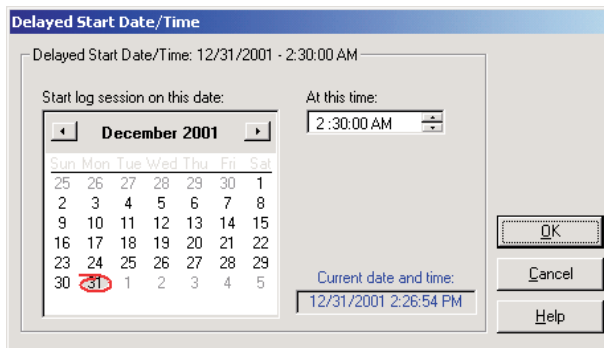


Figure 13 Delayed Start Date/Time Dialog

Select the date on which the log session should start by clicking on the appropriate date in the calendar. Today's date is indicated by the red outline circle. The currently selected date is indicated by a gray filled circle. Use the left and right arrow buttons at the top of the calendar to scroll through the months and years.

Enter the time at which the log session should start using the At this time box.



Note: When using the Delayed Start feature, make sure the Prompt to synchronize real-time log session option on the LogWare Options dialog is not checked! Otherwise the log session will not start as expected! Please refer to Section 5.6.4, Options, for more information.

Once the start date and/or time have been selected, click the OK button to close this dialog.

To cancel setting the delayed start date and/or time, click the Cancel button.

Click the Help button to display the appropriate help topic for this dialog.

2.3.5 Alarm Settings dialog

The Alarm Settings dialog is used to configure low and/or high alarms. The current low alarm or high alarm settings can be changed by selecting the Low alarm or High alarm tab respectively. Low and high alarms are independent of each other and must be configured independently.

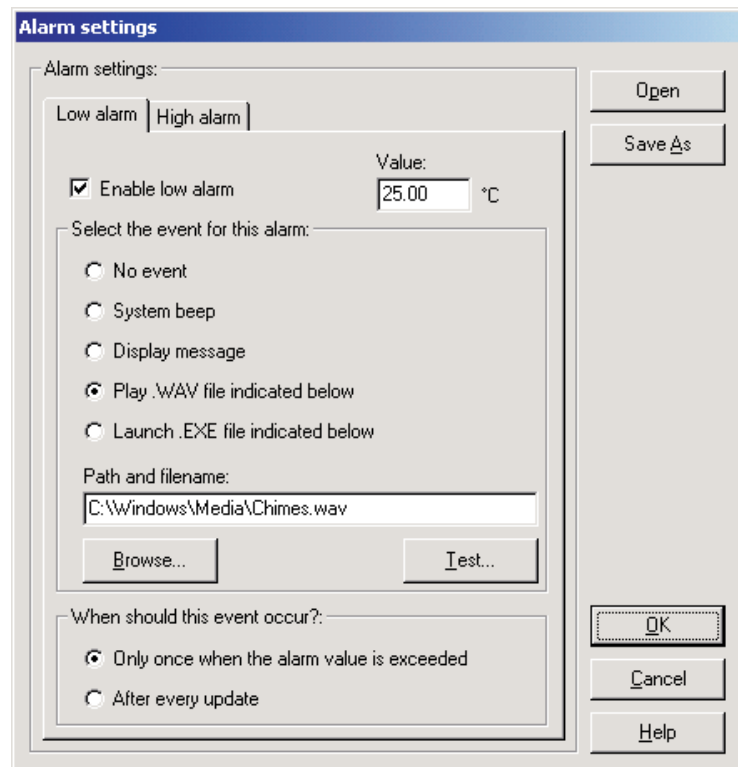


Figure 14 Alarm Settings dialog

The Low alarm tab and High alarm tab indicate the current low alarm and high alarm settings for the selected channel respectively.

To enable an alarm for the selected channel, check the Enable alarm check box on the appropriate tab. If this check box is unchecked, all other settings on the tab are disabled.

Enter the value to be used for this alarm in the Value box. The low alarm will be tripped if the reading on the selected channel meets or drops below the value on the Low alarm tab. The high alarm will be tripped if the reading on the selected channel meets or exceeds the value on the High alarm tab.

Select the event that should occur when each alarm is tripped. The choices are:

- No event
- System beep
- Display message
- Play .WAV file indicated below
- Launch .EXE file indicated below

If either the Play .WAV file indicated below or Launch .EXE file indicated below option is selected, the Path and filename box and Browse button are enabled and the appropriate file must be entered or selected. Use the Browse button to browse for the file. The Test button is enabled once a file has been entered. Select the Test button to make sure the selected file can be played or launched.



Note: *The file entered in the Path and filename box MUST exist! When the OK button is clicked, a search is performed to locate the file. If the file is not found, a warning message is displayed. If the selected file resides on a network resource, the resource must be available and properly connected when the alarm event occurs!*

LogWare ships with a few sample .WAV files that can be used. These files are located in the \Config folder.

The Launch .EXE file indicated below option can be used to launch any application such as pager software which could be configured to dial a phone number automatically and transmit a fixed message to a pager indicating that an alarm has tripped.



Note: It is not recommended that the Launch .EXE file indicated below option be used (especially in conjunction with the After every update option) until the user is familiar with how this feature works. Unless the .EXE file terminates itself automatically (or is limited to a single instance), this software will continue to launch new instances of the .EXE file every time the alarm trips, possibly causing the computer to run out of resources.

Select when the specified event for each alarm should occur. The choices are:

- Only once when the alarm value is exceeded
- After every update

If the Only once when the alarm value is exceeded option is selected, the chosen event will occur on the first reading that trips the alarm. Subsequent readings that also exceed the alarm value will not cause the event to occur again until one or more readings are taken that are back within the alarm value (alarm reset).

If the After every update option is selected, every reading that is taken that exceeds the alarm value causes the event to happen again.

The alarm settings can be saved to a configuration file. If the same alarm settings are to be used on multiple channels, this provides a method to quickly configure the alarms on subsequent channels.

When clicking the Save As button, the Windows® Save As dialog is displayed. The \Config folder should be selected by default (or the folder that is specified for configuration files on the LogWare Options dialog). Alarm configuration files are saved with a **.alm** extension. To cancel saving a configuration file, click the Cancel button. Otherwise, enter the filename for the configuration file and click the Save button. For more details about the alarm configuration file, refer to Section 7.1.2, Alarm Configuration File.

An alarm configuration file that has been saved previously can be opened by clicking the Open button.

When clicking the Open button, the Windows® Open dialog is displayed. The \Config folder should be selected by default (or the folder that is specified for configuration files on the LogWare Options dialog). Alarm configuration files should have a **.alm** extension. To cancel opening a configuration file, click the Cancel button. Otherwise, select the configuration file to open and click the Open button. For more details about the alarm configuration file, refer to Section 7.1.2, Alarm Configuration File.

Once the alarm settings are complete, click the OK button to save these settings and close this dialog.

Click the Cancel button to discard any changes you have made and close this dialog.

Click the Help button to display the appropriate help topic for this dialog.

2.3.6 Statistics Settings dialog

The Statistics Settings dialog allows specific statistical functions to be enabled or disabled for the selected channel.

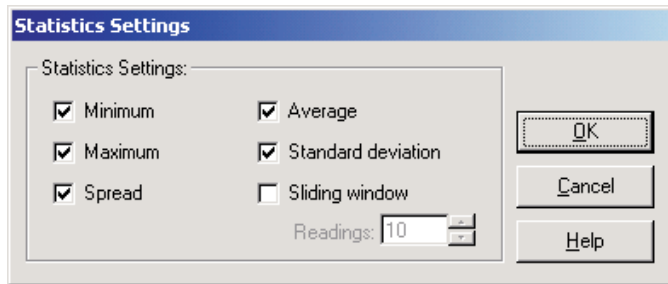


Figure 15 Statistics Settings dialog

To enable a statistical function for the selected channel, check the box next to the appropriate function.

The statistical functions supported are:

- Minimum
- Maximum
- Spread
- Average
- Standard deviation
- Sliding window

If the Sliding window function is enabled, a value must be entered into the Readings box. All other statistical values are calculated based on the most recent n measurements as indicated by the value in the Readings box. The readings setting can be set to any value between 10 and 1,000.

If the Sliding window function is disabled, the statistical values are calculated based on **all** readings since the last reset was performed.

Once the statistics settings are complete, click the OK button to save these settings and close this dialog.

Click the Cancel button to discard any changes you have made and close this dialog.

Click the Help button to display the appropriate help topic for this dialog.

2.4 Real-time log session

The Log Session window is displayed when a real-time log session starts. This window is divided into three main sections: channels grid, statistics and graph. These three sections can be resized by dragging the bars that divide the sections with the mouse.

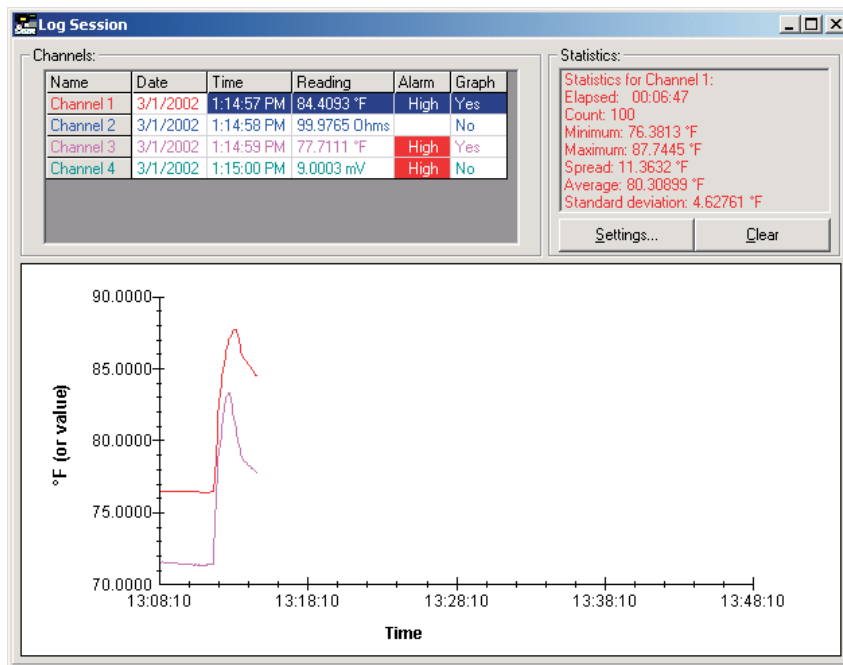


Figure 16 Real-time log session window

LogWare was designed to communicate with only one thermometer readout at a time. Therefore, only one real-time log session can be running at a time. In order to optimize data acquisition, many menu options are disabled during a real-time log session.

When a log session is started, LogWare prepares to acquire data from the enabled channels. The current conversion settings for each channel and system units are determined at this time. The scale of the readings for each channel depends on the conversion setting for each channel and the session temperature scale as set on the Log Session Settings dialog.

LogWare attempts to create the output file for each enabled channel. If LogWare is unable to create an output file, or if there is an existing file of the same name and the overwrite option was not selected, a prompt is displayed and a new path and/or filename may be entered for the output file.

If a delayed start time was selected, a dialog appears indicating the time that the log session will begin. Otherwise, a dialog appears indicating that the log session will begin momentarily while LogWare initializes the thermometer readout and finishes preparing to start the log session.




Note: *If a delayed start time was selected and the Prompt to synchronize real-time log session option is selected on the LogWare Options dialog, a prompt is displayed asking if the synchronize feature should be disabled. If the synchronize feature is not disabled, the log session will not start as expected! For more details, please refer to Section 5.6.4, Options.*

Once logging starts, one measurement is taken on each enabled channel at the interval specified on the Log Session Settings dialog. If the Interval between scan sequences feature was enabled, LogWare then waits the specified time and repeats this process. Otherwise, this process is repeated without delay.




Note: *Be aware that selecting any menu option while a real-time log session is running may cause the measurement process to halt temporarily until the menu is closed or the dialog that is displayed is closed! Also, running other applications on the computer while logging real-time data, especially those that require a lot of processing time or that access the disk frequently, may cause the measurement process to halt temporarily while the other application is busy. It is recommended that, once a real-time log session is started, the software and the computer be left alone to log the data without interruption.*

If the log session was configured to stop logging at a specified time or after a specific number of readings, LogWare continues to acquire data until the end log criteria is met. The log session can be stopped manually at any time by selecting the Stop option in the Datalog menu or by clicking the  button on the toolbar.

When a log session is stopped manually, a prompt is displayed asking if the log session window should be closed. The log session window may be closed at this time or left open. In order to start a new log session, the current log session window must be closed. Select Yes to close the log session window immediately or select No to leave the window open.

To close the real-time log session window, select the Close option in the

Datalog menu or click the  button on the toolbar. If the Show Log Session Settings dialog after closing log session option on the LogWare Options dialog is checked, the Log Session Settings dialog is automatically displayed. To prevent LogWare from displaying the Log Session Settings dialog after the log session window is closed, uncheck this option. For more details, please refer to Section 5.6.4, Options.

Once the log session has been stopped, the output file which contains all of the acquired readings can be opened by selecting the Open Previously Logged Data File option in the View menu. For more details on opening previously logged data files, please refer to Section 5.4.7, Open Previously Logged Data File.

2.4.1 Channels grid

The Channels grid is located in the upper left corner of the log session window and displays the most recent data for each enabled channel. This data includes the channel name, date, time and value and scale of the most recent reading, the current alarm state and whether or not the trace for the channel is currently displayed on the graph. The color of the data in the Channels grid matches the selected trace color for that channel. The currently selected channel is highlighted.

The first column in the Channels grid indicates the channel names as entered on the Log Session Settings dialog. The second and third columns indicate the date and time that the last reading was taken on each channel. The fourth column indicates the last reading value and scale on each channel. The fifth column indicates the current alarm state and the sixth column indicates whether the trace for each channel is currently shown on the graph. Each of these columns may be resized as desired by dragging the lines that separate each column with the mouse.

The current alarm state is determined by comparing the most recent reading on each channel to that channel's alarm settings. If the low and/or high alarms have been configured, the reading is compared to the alarm values. If the low alarm is enabled and the reading is less than or equal to the low alarm value, the Alarm box in the Channel grid flashes **Low** to indicate the low alarm has "tripped". If the high alarm is enabled and the reading is greater than or equal to the high alarm value, the Alarm box flashes **High** to indicate the high alarm has "tripped". If the reading falls between the low and high alarm values or if the alarms are not enabled, the Alarms box remains blank (alarm reset).

Double-clicking on a row in the Channels grid toggles the state of the trace on the graph. The current state for each channel is indicated in the Graph column.

Right-clicking on the Channels grid displays a popup menu. This menu consists of some of the more common options for working with channels during a

real-time log session. The menu options apply only to the currently selected channel.

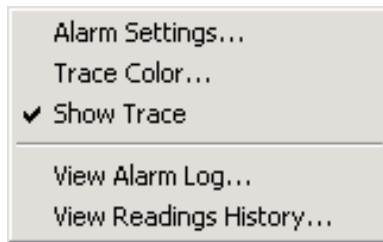


Figure 17 Channels popup menu

The alarm settings for the selected channel can be viewed by selecting the Alarm Settings option in the popup menu, or by selecting the Alarm Settings option in the View menu. If the selected channel does not have any alarms configured, this menu option is disabled. Selecting the Alarm Settings option displays the Current Alarm Settings dialog. Please refer to Section 2.4.4, Current Alarm Settings Dialog, for more information on this dialog.

The graph trace color for the selected channel can be changed while the log session is running by selecting the Trace Color option in the popup menu, or by selecting the Trace Color option in the Graph menu. Please refer to Section 2.4.7, Trace Color Dialog, for more details on using the Trace Color dialog. The color of the channel in the Channels grid, statistics list and the graph trace are all updated to the new color.

The Show Trace option in the popup menu indicates whether the trace for this channel is currently shown on the graph. A check appears next to this menu option if the trace is visible. The Graph column of the Channels grid also indicates whether the trace is currently shown on the graph. To toggle showing the trace for the selected channel on the graph, select the Show Trace option in the popup menu, or select the Show Trace option in the View menu.

The View Alarm Log option in the popup menu displays the alarm log file for this log session. Every time an alarm is tripped or reset, an entry is written to the alarm log file. This file is named "alarms.log" and is located in the default output folder. The alarm log file can also be viewed by selecting the Alarm Log option in the View menu. The alarm log file is displayed in the File Viewer window. Please refer to Section 7.2.3, Alarm Log File, for more information on the contents of the alarm log file. Please refer to Section 2.4.8, File Viewer Window, for more information on using the File Viewer window.



Note: When viewing the alarm log file while a real-time log session is running, the File Viewer window displays only the information that had been written to the file at the time the file was opened. Click the Refresh button on the File Viewer window to update the file. The alarm log file can also be printed by clicking the Print button on the File Viewer window.

The output file for each channel can be opened and viewed by selecting the View Readings History option in the popup menu or by selecting the Readings History option in the View menu. The output file contains all of the data that has been logged for the selected channel during the current log session. The output file is displayed in the File Viewer window. Please refer to Section 7.2.1, Output File, for more information on the contents of the output file. Please refer to Section 2.4.8, File Viewer Window, for more information on using the File Viewer window.



Note: If the data is being acquired using the Write all acquired data to the same file option, the View Readings History feature displays all readings for all channels!



Note: When viewing the output file while a real-time log session is running, the File Viewer window displays only the information that had been written to the file at the time the file was opened. Click the Refresh button on the File Viewer window to update the file. To see the statistics for the output file, click the Statistics button on the File Viewer window. The output file can also be printed by clicking the Print button on the File Viewer window or by selecting the Print | Readings option in the File menu.

2.4.2 Statistics list

The Statistics list is located in the upper right corner of the log session window and displays statistical information for the currently selected channel. The color of the statistics data matches the trace color for the selected channel. Each of the statistical settings can be enabled or disabled by clicking the Settings button. Only the enabled statistics are displayed in the Statistics list. Please refer to Section 2.3.6, Statistics Settings Dialog, for more information on the Statistics Settings dialog.

Clicking the Clear button resets all statistical values for the selected channel and begins calculating the statistical data again starting with the next reading acquired on the selected channel.

2.4.3 Graph

The graph is located at the bottom of the log session window. The graph gives a visual representation of the readings on all enabled channels.

The trace for each enabled channel can be toggled on or off. The Graph column in the Channels grid indicates the current state of the trace for each channel. To

show or hide the trace for the selected channel, either double-click on the channel, or right-click on the Channels grid and select the Show Trace option in the popup menu or select the Show trace option in the View menu.

The Y-axis of the graph is always labeled in the session temperature scale that was selected on the Log Session Settings dialog. The Y-axis is automatically scaled to keep the entire visible portion of all traces within the graph bounds. As traces are shown and/or hidden, the graph should automatically rescale to maximize the resolution.



Note: For readings on all channels that are configured for non-temperature units (i.e. Ohms or mV), the units of the Y-axis are assumed to be in the appropriate units for that channel.

The X-axis of the graph displays the current time according to the computer's clock. As the traces reach the right side of the graph, the traces shift to the left and the left-most part of the traces are pushed off the graph. The graph is set to display 600 readings on the X-axis at a time. This setting is fixed and cannot be changed.

Gridlines can be automatically displayed on the graph by checking the Show gridlines on graphs option of the General tab of the LogWare Options dialog. Please refer to Section 5.6.4.1, General Tab, for more details.

Right-clicking on the graph displays a popup menu. This menu is identical to the Graph menu and consists of options for working with the graph during a real-time log session. Some of the menu options apply only to the currently selected channel.

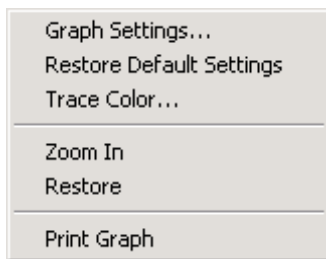


Figure 18 Graph popup menu

There are many customizable settings for the graph such as titles, a legend, background and text colors, grid lines and trend lines. All of these features can be accessed by selecting the Graph Settings option in the popup menu or in the Graph menu. The default graph settings can be restored by selecting the Restore Default Settings option in the popup menu or in the Graph menu. For more details on customizing the graph settings, please refer to Section 5.5.1, Graph Settings Dialog.

The Zoom In and Restore options are not applicable to a real-time log session and are disabled.

The graph may be printed at any time by selecting the Print Graph option in the popup menu or in the Graph menu, or by selecting the Print | Graph option in the File menu.

2.4.4 Current Alarm Settings dialog

The Current Alarm Settings dialog displays the current alarm settings.

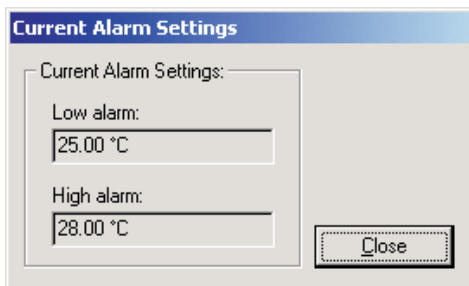


Figure 19 Current Alarm Settings dialog

The Low alarm box indicates the current low alarm value and units. If the low alarm is not enabled, the Low alarm box is disabled and no value is displayed.

The High alarm box indicates the current high alarm value and units. If the high alarm is not enabled, the High alarm box is disabled and no value is displayed.

Click the Close button to close this dialog.

2.4.5 Output file

The data that is acquired during a real-time log session is written to a file, called an output file. This file contains information about the thermometer read-out used to acquire the data, alarm settings (if applicable), and the date, time, value, scale and alarm state for each reading.

To view the contents of an output file during a real-time log session, select the Readings History option in the View menu, or right-click on the Channels grid and select the View Readings History option in the popup menu.

The output file is described in detail in Section 7.2.1, Output File.

2.4.6 Alarm log file

During a real-time log session, a log of all alarm activity is recorded in a file call the alarm log. This file contains alarm setting information for each channel

and the date, time, value, scale and alarm state for each time an alarm tripped or reset.

To view the contents of the alarm log file during a real-time log session, select the Alarm Log option in the View menu, or right-click on the Channels grid and select the View Alarm Log option in the popup menu.

The alarm log file is described in detail in Section 7.2.3, Alarm Log File.

2.4.7 Trace color dialog

The Trace Color dialog allows a new color to be selected for the graph trace.

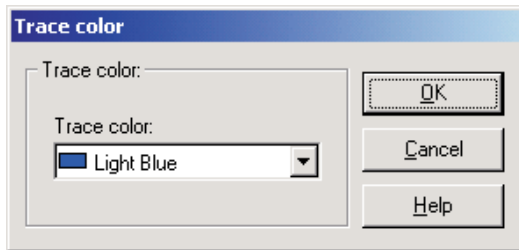


Figure 20 Trace color dialog

The current trace color is selected when the dialog is displayed. Select the new color using the Trace color drop-down list.

Click the OK button to save the settings you have selected and to close the dialog.

Click the Cancel button to discard any changes and close the dialog.

Click the Help button to display the appropriate help topic for this dialog.

2.4.8 File Viewer window

The File Viewer window displays the contents of the indicated file.

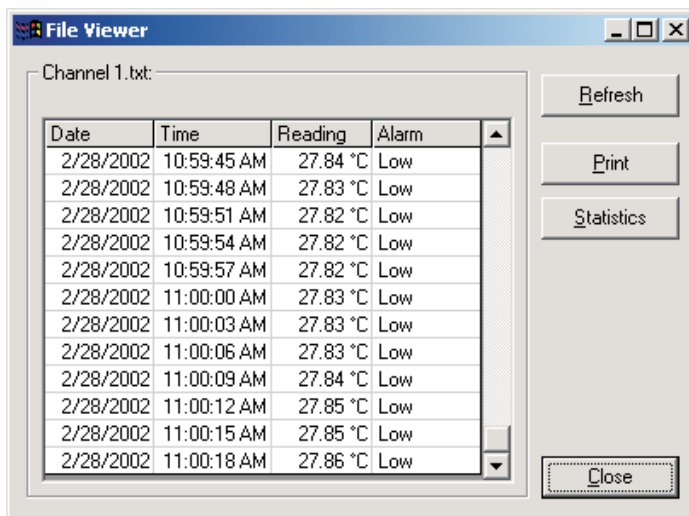


Figure 21 File Viewer window

The data grid displays the contents of the file at the time the viewer window is displayed. The data displayed in the data grid will vary depending on the contents of the file being viewed.

To update the data in the data grid, click the Refresh button. The contents of file will be read in again and the data grid is updated.

To print the contents of the indicated file to the default printer, click the Print button.

The Statistics button is shown if there is statistical data associated with the displayed data. To see the statistical data, click the Statistics button. The Statistics dialog is displayed. Please refer to Section 2.4.9, Statistics Dialog, for more details on this dialog.

The Close button closes this window.

2.4.9 Statistics dialog

The Statistics dialog displays the current statistical values. Only the values for the enabled statistics are displayed.

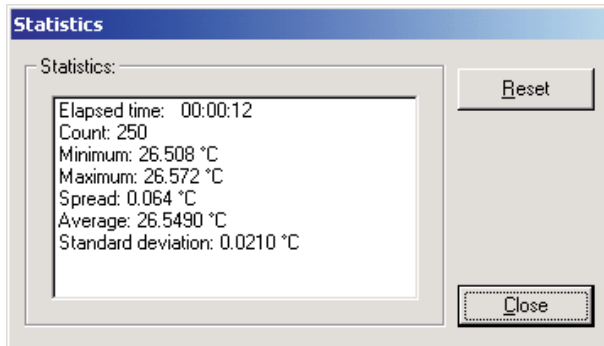


Figure 22 Statistics dialog

If the Sliding window statistical setting is enabled, the statistical values are based on the number of readings specified for the sliding window. Otherwise, the statistical values are based on all readings since the statistics were last reset.

Click the Reset button to reset the statistical values.

Click the Close button to close this dialog.

2.5 Previous log session

The Previous Log Session window allows multiple output files or download files to be opened and viewed. This window is divided into three main sections:

channels, statistics and graph. These three sections can be resized by dragging the bars that divide the sections with the mouse.

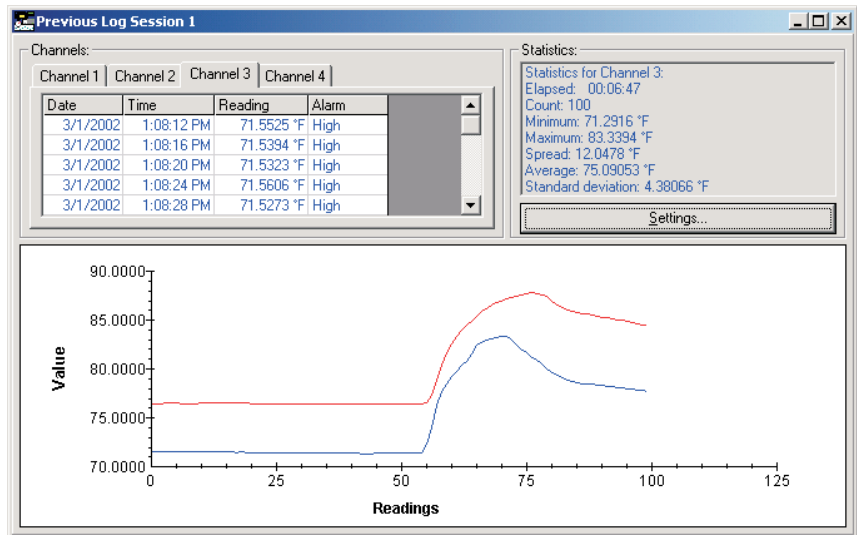


Figure 23 Previous Log Session window

A Previous Log Session window can be opened by selecting the Open Previously Logged Data File option in the View menu or by selecting one of the files in the Most Recently Used (MRU) section at the bottom of the View menu. Up to 8 Previous Log Session windows can be open at a time.

When selecting the Open Previously Logged Data File option, the Windows® Open dialog is displayed. The \Config folder should be selected by default (or the folder that is specified for configuration files on the LogWare Options dialog). Both output files and download files can be opened in the Previous Log Session window. These files should have a .txt extension. To cancel opening a file, click the Cancel button. Otherwise, select up to 10 files to open and click the Open button. If more than 10 files are selected, only the first 10 files are opened and the rest are ignored. For more details about output files, please re-

fer to Section 7.2.1, Output file. For more details about download files, please refer to Section 7.2.2, Download file.




Note: To open an output file that was created when using the Write all acquired data to the same file option (multiple data sets in the same file), you must not select more than one file on the Windows® Open dialog. If the file contains data from more than 10 channels, only the data from the first 10 channels will be displayed. Data sets from each channel are separated into individual files and may be opened in another window if desired. A message is displayed if the selected file contains more than 10 data sets and indicates the names and path where the files are located.

A progress dialog is displayed as each of the selected files are read in and analyzed. As each of the files is read in, a tab is created in the Channels section of the window and statistical data is calculated. Each channel is automatically assigned a trace color that can be changed at any time. The X-axis of the graph automatically grows to accommodate the file with the most readings. The Y-axis of the graph is automatically scaled to display all traces on the graph.



Note: The graph allows a maximum of 32,767 readings per channel. If any one file contains more readings than that, the first 32,767 readings are read in and the rest are ignored. The graph also limits the total number of data points to 128,000 for all channels. So, for 10 channels, a maximum of 12,800 readings per channel can be displayed.

To close the Previous Log Session window, click the  button in the title bar.

2.5.1 Channels

The Channels section is located in the upper left corner of the previous log session window and displays one tab for each file that was opened. Each tab is given the name of the file and the grid on the tab contains all of the data that was read in from the file. Depending on the type of file being opened, this data can include but is not necessarily limited to, the date, time, value and scale of each reading as well as the alarm state at the time that reading was taken.



Note: When opening files that contain data that was downloaded from some of the thermometers supported by LogWare, no date or time information is available. Download files never contain alarm information.

Selecting a tab in the Channels section displays the data associated with that channel. The color of the data in the grid matches the graph trace color assigned to this channel.

Right-clicking on any of the channel tabs displays a popup menu. This menu

allows channels to be added to or removed from this session, and also allows the trace color to be changed, or the trace can be hidden.

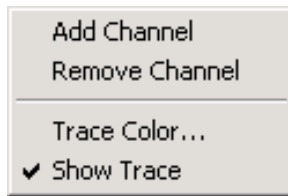


Figure 24 Channels popup menu

LogWare is capable of opening up to 10 files (channels) in the same Previous Log Session window. To add another channel to the window, select the Add Channel option in the popup menu. The Windows® Open dialog is displayed. The \Config folder should be selected by default (or the folder that is specified for configuration files on the LogWare Options dialog). To cancel opening a file, click the Cancel button. Otherwise, select up to 9 more files to open and click the Open button. Since LogWare only allows a total of 10 files to be opened, any additional files are ignored. The Add Channel option is disabled if the log session already has 10 channels open.

To remove a channel from the Previous Log Session window, select the tab for the channel and select the Remove Channel option in the popup menu. Any channel can be removed **except the first channel**. The first channel cannot be removed. When a channel is removed, the corresponding tab, grid and graph trace are also removed and the graph Y-axis is rescaled as needed.

To change the color of the trace on the graph for any channel, select the tab for the channel and select the Trace Color option in the popup menu, or by selecting the Trace Color option in the Graph menu. The Trace Color dialog is displayed. Please refer to Section 2.4.7, Trace Color Dialog, for more information on this dialog.

The Show Trace option in the popup menu indicates whether the trace for the selected channel is currently shown on the graph. A check appears next to this option if the trace is visible. The trace on the graph for any channel can be toggled on or off at any time. To toggle the trace for any channel, select the tab for the channel and select the Show Trace option in the popup menu or select the Show trace option in the View menu. If a check appears next to this option, the trace is currently shown. If no check appears, the trace is hidden.

2.5.2

Statistics list

The Statistics list is located in the upper right corner of the Previous Log Session window and displays statistical information for the currently selected channel. The statistics are calculated when the data is read in from the file. The color of the statistics data matches the trace color for the selected channel.

Each of the statistical settings can be enabled or disabled by clicking the Settings button. Only the enabled statistics are displayed in the Statistics list. Please refer to Section 2.3.6, Statistics Settings Dialog, for more information on the Statistics Settings dialog.

2.5.3 Graph

The graph is located at the bottom of the Previous Log Session window. The graph gives a visual representation of the readings on all channels.

The trace for each channel can be toggled on or off. To show or hide the trace for the selected channel, right-click on the tab in the Channels grid and select the Show Trace option in the popup menu or select the Show Trace option in the View menu.

The Y-axis of the graph is automatically scaled to display all visible traces in their entirety. The Y-axis is labeled Value by default and the resolution of the values is extracted from the data read in from the channels. The X-axis of the graph grows automatically to display the entire trace of the channel with the most readings. The X-axis is labeled Readings by default and is numbered according to the number of readings in each file.



Note: *The graph allows a maximum of 32,767 readings per channel. If any one file contains more readings than that, the first 32,767 readings are read in and the rest are ignored. The graph also limits the total number of data points to 128,000 for all channels. So, for 10 channels, a maximum of 12,800 readings per channel can be displayed.*

Gridlines can be automatically displayed on the graph by checking the show gridlines on graphs option of the General tab of the LogWare Options dialog. Please refer to Section 5.6.4.1, General Tab, for more details.

Right-clicking on the graph displays a popup menu. This menu is identical to the Graph menu and allows quick access to graph-related functions. Some of the options apply only to the currently selected channel.

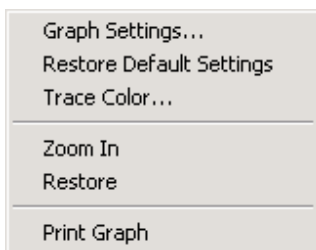


Figure 25 Graph popup menu


There are many customizable settings for the graph such as titles, a legend, background and text colors, grid lines and trend lines. All of these features can be accessed by selecting the Graph Settings option in the popup menu or in the Graph menu. The default graph settings can be restored by selecting the Restore Default Settings option in the popup menu or in the Graph menu. For more details on customizing the graph settings, please refer to Section 5.5.1, Graph Settings Dialog.

The Zoom In and Restore options are can be used to zoom in to a small section of the graph and to restore the graph to its default state respectively. Please refer to Section 2.5.4, Graph Zooming, for more details on using the graph zooming feature.

The graph may be printed at any time by selecting the Print Graph option in the popup menu or in the Graph menu, or by selecting the Print | Graph option in the File menu.

2.5.4 Graph zooming

LogWare allows the graph in the Previous Log Session window to be zoomed to view a small section of the graph in more detail. The zoom feature can be enabled by selecting the Zoom In option in the Graph menu, or by right-clicking on the graph and selecting the Zoom In option in the popup menu, or by clicking

the  button on the toolbar. When the zoom feature is enabled, the mouse pointer turns into a crosshair when positioned over the graph, the toolbar button remains depressed and a check mark appears next to the Zoom In option indicating that the graph is now in “zoom mode”.

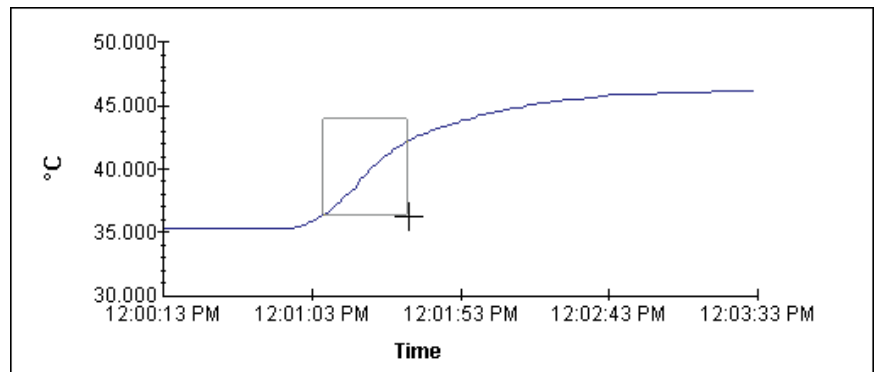



Figure 26 Graph zooming

To zoom in to a portion of the graph, use the mouse to select the portion of the graph to zoom by clicking the left mouse button and holding it down while dragging the mouse to draw a rectangular box around the desired area. Release-

ing the mouse button causes the graph to zoom in to the selected area. Repeat this process to zoom in closer and closer.

To restore the graph to its normal state, select the Restore option in the Graph menu, or right-click on the graph and select the Restore option in the popup menu, or click the  button on the toolbar. The check mark is then removed in the Zoom In option and the Zoom In toolbar button no longer appears depressed.

3 Downloading logged data

LogWare provides tools to download data that has been logged to the thermometer readout's memory or logs and to view the current status of the thermometer readout's logs or memory. These tools are accessible by selecting options in the Download menu. The following sections describe the download and log status dialogs for each of the thermometer readouts supported by LogWare.

3.1 Download menu

The Download menu contains options for downloading logged data from the thermometer readout's memory or logs and viewing the current status of the thermometer readout's logs or memory.

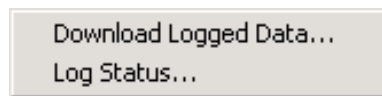


Figure 27 Download menu

3.1.1 Download logged data

The Download Logged Data option can be used to download data that has been logged to the thermometer readout's logs or memory. All thermometer readouts supported by LogWare support logging of some type. Please refer to Section 1.4.4, Firmware Requirements, for more information.

When selecting the Download Logged Data option, the Select Model dialog is displayed. LogWare needs to know the model of the thermometer readout in order to display appropriate options dialog. Please refer to Section 3.2, Select Model Dialog, for more details.

After the thermometer readout model and communication settings have been selected, the appropriate Download Options dialog for the selected thermometer readout is displayed. The Download Options dialogs are described in more detail in Section 3.3, Download Options Dialogs.

3.1.2 Log status

The Log Status option can be used to view the current status of the thermometer readout's logs or memory. This feature is only supported by some of the thermometer readouts supported by LogWare.

When selecting the Log Status option, the Select Model dialog is displayed. LogWare needs to know the model of the thermometer readout in order to display appropriate status dialog. Please refer to Section 3.2, Select Model Dialog, for more details on the Select Model dialog.

After the thermometer readout model and communication settings have been selected, the appropriate Log Status dialog for the selected thermometer readout is displayed. The Log Status dialogs are described in more detail in Section 3.4, Log Status Dialogs.

3.2 Select Model dialog

When any of the Download menu options is selected, the Select Model dialog is displayed. LogWare needs to know the model of the thermometer readout to interface with.

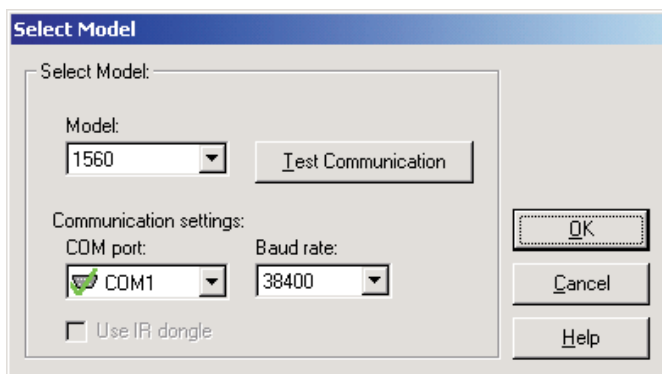








Figure 28 Select Model dialog

If this dialog has been used previously to select a thermometer readout model and communication settings, the previous settings are restored.

Use the Model drop-down list to select the type of thermometer readout being used. All other controls on this dialog are disabled until a model is selected. The list of available thermometer readouts depends on the Download menu option selected. Only thermometer readouts that support the selected feature are listed.

Select the COM port on the computer to which the thermometer readout is connected using the COM port drop-down list. Only COM ports that the operating system reports as being available can be used. The following symbols are used to show the status of each COM port:

-  - Indicates a COM port that is currently available.
-  - Indicates the COM port that is currently selected by LogWare.
-  - Indicates a COM port that is currently being used by another application.
-  - Indicates an IR port (virtual COM port) that is currently available.

-  - Indicates an IR port (virtual COM port) that is currently selected by LogWare.
-  - Indicates an IR port (virtual COM port) that is currently being used by another application.

Select the baud rate that the computer should use to communicate over the COM port with the thermometer readout using the Baud rate drop-down list. The baud rate setting of the thermometer readout must be set to the same value as this setting. For best results, use the maximum allowable baud rate setting when possible.

Check the Use IR dongle option to use the thermometer readout's IR port instead of an RS-232 cable. This option is only enabled when the selected thermometer readout has an IR port and the selected download feature allows using the IR port to communicate with the thermometer readout. For more information on using the IR dongle, please refer to Section 6.13, Using an IR Dongle with LogWare.

After the model, COM port, baud rate and/or IR settings are configured, click the Test Communication button to make sure communication with the thermometer readout is successful. At this time, the thermometer readout must be connected to the COM port on the computer using a null modem cable or, if the IR dongle is being used, the IR dongle must be connected to the selected COM port and the dongle must be aligned with the thermometer readout's IR port. For more information, please refer to the thermometer's User's Guide.

When testing communication with the thermometer readout, there are three possible results:

- The thermometer readout is communicating at the correct baud rate on the selected COM port.
- The thermometer readout is communicating, but possibly at the wrong baud rate or the thermometer readout model does not match the selected model.
- The thermometer readout failed to communicate at all. Check the COM port and baud rate settings, make sure the cable is a null modem cable, and make sure the thermometer readout is plugged in and turned on. If using the IR dongle, make sure the dongle is connected to the selected COM port and the dongle is aligned properly with the thermometer readout's IR port.

After all of the settings on this dialog are configured and communication with the thermometer readout is successful, click the OK button to accept the settings and continue.

Click the Cancel button close this dialog and abort using the selected feature.

Click the Help button to display the appropriate help topic for this dialog.

3.3 Download options dialogs

LogWare provides a tool to download data from the thermometer readout's memory or logs. This tool is accessible by selecting the Download Logged Data option in the Download menu. The following sections describe the dialogs for each of the thermometer readouts that support this feature.

3.3.1 1523 Handheld download options dialog

The Model 1523 Handheld Thermometer readout can store readings in a log, called the demand log. LogWare can download the readings from this log and write the readings into download files. For more information on download files, please refer to Section 7.2.2, Download File on page 231. When the Download Logged Data option in the Download menu is selected and the 1523 Thermometer readout is selected on the Select Model dialog, the 1523 Download Options dialog is displayed.

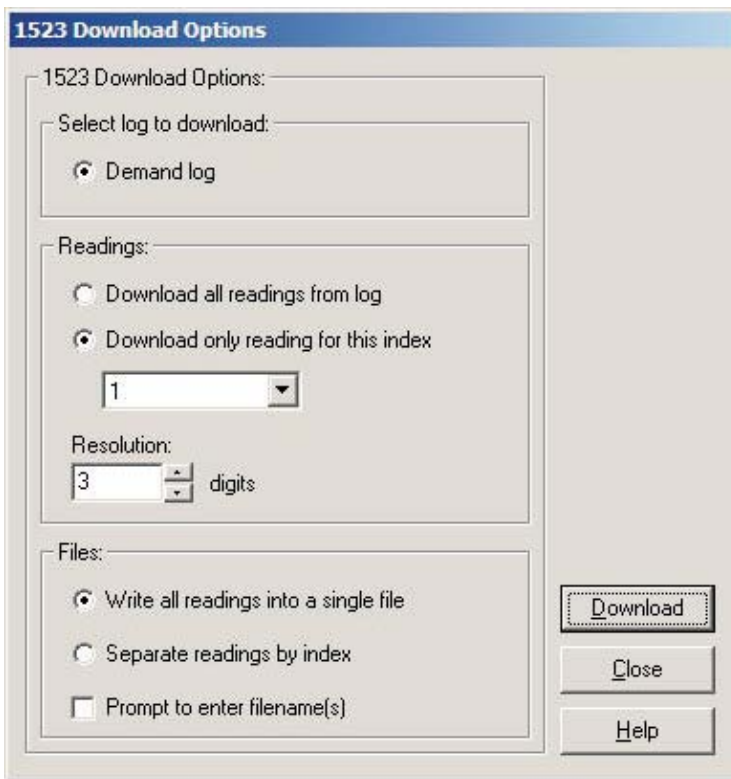


Figure 29 1523 Download Options dialog

The demand log stores up to 25 readings. For more details on the demand log, see the 1523 User's Guide.

The Readings section allows the readings to be downloaded to be specified.

Selecting the Download all readings from log option indicates that all of the readings in the selected log should be downloaded.

Selecting the Download only reading for this index option allows only the reading at the selected index (1 to 25) to be downloaded. If this option is selected, use the drop-down list to select the appropriate index of the reading to download.

Select the resolution to use to format the readings when they are written to the download file. This value can be set between 0 and 3 digits (places after decimal point).

Select the method to use when writing the downloaded readings to a file using the options in the Files section.

Selecting the Write all readings into a single file option causes all of the downloaded readings to be written to one file.

Selecting the Separate readings by index option causes one file to be created for each index for which there is a reading available. Each file will contain a single reading.

If the Prompt to enter filename(s) check box is checked, a prompt is displayed after the data is downloaded for each file that needs to be created, depending on the other selections above. If this option is not checked, the default filenames are used.

After selecting the appropriate options on this dialog, click the Download button to start the download process. For more details on the download process, please refer to Section 8, Download Process.

To close this dialog, click the Close button.

Click the Help button to display the appropriate help topic for this dialog.

3.3.2 1524 Handheld download options dialog

The Model 1524 Handheld Thermometer readout can store readings in any of two logs: demand log and auto log. *LogWare* can download the readings from these logs and write the readings into download files. For more information on download files, please refer to Section 7.2.2, Download File. When the Download Logged Data option in the Download menu is selected and the 1524

Thermometer readout is selected on the Select Model dialog, the 1524 Download Options dialog is displayed.

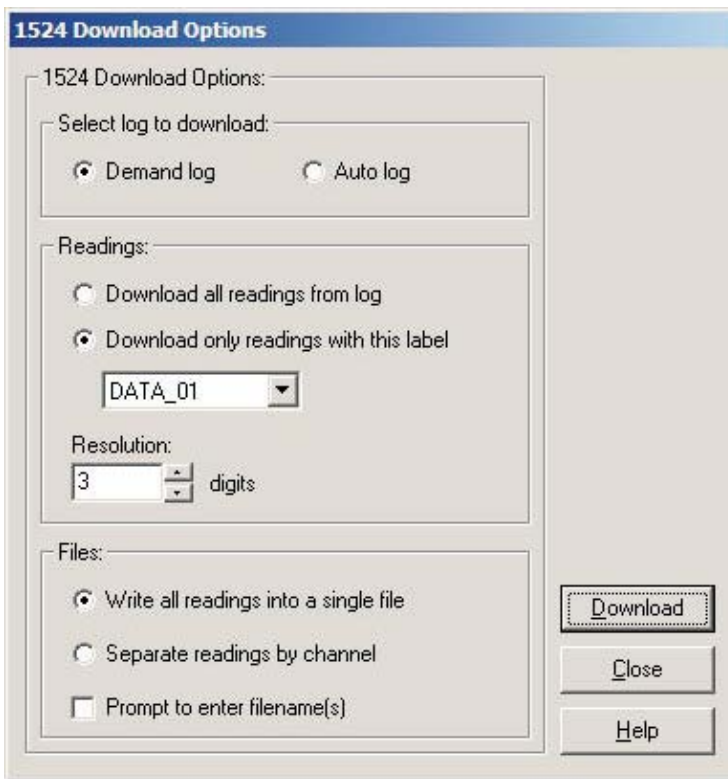


Figure 30 1524 Download Options dialog

The Select log to download section allows the log to be specified. The 1524 Handheld Thermometer readout has two logs: demand log and auto log. The demand log stores up to 25 readings while the auto log stores up to 15,000 readings. For more details on the demand and auto logs, see the 1524 User's Guide.

The Readings section allows the readings to be downloaded to be specified.

Selecting the Download all readings from log option indicates that all of the readings in the selected log should be downloaded.

Selecting the Download only readings with this label option allows only readings that are assigned the label indicated below to be downloaded from the selected log. If this option is selected, a list of current labels is read in from the thermometer readout. Use the drop-down list to select the appropriate label.

Select the resolution to use to format the readings when they are written to the download file. This value can be set between 0 and 3 digits (places after decimal point).

Select the method to use when writing the downloaded readings to a file using the options in the Files section.

Selecting the Write all readings into a single file option causes all of the downloaded readings to be written to one file.

Selecting the Separate readings by channel option causes all of the downloaded readings to be separated according to the channel that each reading was taken from and written into separate files.

If the Prompt to enter filename(s) check box is checked, a prompt is displayed after the data is downloaded for each file that needs to be created, depending on the other selections above. If this option is not checked, the default filenames are used.

After selecting the appropriate options on this dialog, click the Download button to start the download process. For more details on the download process, please refer to Section 8, Download Process.

To close this dialog, click the Close button.

Click the Help button to display the appropriate help topic for this dialog.

3.3.3 1529 Chub-E4 download options dialog

The Model 1529 Chub-E4 Thermometer readout (including models 1529, 1529-R and 1529-T) can store readings in any of two logs: demand log and auto log. LogWare can download the readings from these logs and write the readings into download files. For more information on download files, please refer to Section 7.2.2, Download File. When the Download Logged Data option in the Download menu is selected and the 1529 thermometer readout is selected on the Select Model dialog, the 1529 Chub-E4 Download Options dialog is displayed (Figure 31).

The Select log to download section allows the log to be specified. The 1529 Chub-E4 Thermometer readout has two logs: demand log and auto log. The demand log stores up to 100 readings while the auto log stores up to 8,160 readings. For more details on the demand and auto logs, see the 1529 Chub-E4 Thermometer readout User's Guide.

The Readings section allows the number of readings to be downloaded to be specified.

Selecting the Download all readings from log option indicates that all of the readings in the selected log should be downloaded.

Selecting the Download only readings with this label option allows only readings that are assigned the label indicated below to be downloaded from the selected log. If this option is selected, a list of current labels is read in from the thermometer readout. Use the drop-down list to select the appropriate label.

Enter the resolution to use to format the readings when they are written to the download file in the Resolution box. This value can be set between 0 and 4 digits (places after decimal point).

Select the method to use when writing the downloaded readings to a file using the options in the Files section.

Selecting the Write all readings into a single file option causes all of the downloaded readings to be written to one file.

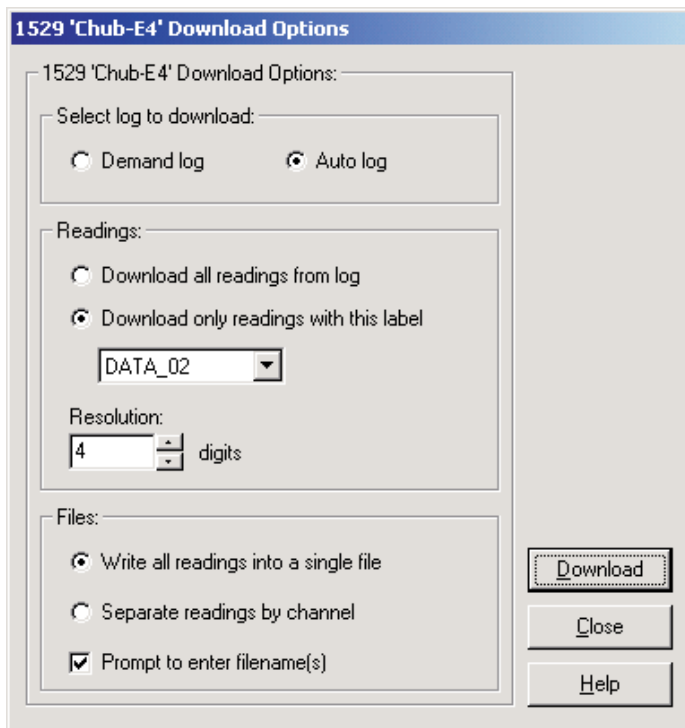


Figure 31 1529 Chub-E4 Download Options dialog

Selecting the Separate readings by channel option causes all of the downloaded readings to be separated according to the channel that each reading was taken from and written into separate files.

If the Prompt to enter filename(s) check box is checked, a prompt is displayed after the data is downloaded for each file that needs to be created, depending on the other selections above. If this option is not checked, the default filenames are used.

After selecting the appropriate options on this dialog, click the Download button to start the download process. For more details on the download process, please refer to Section 8, Download Process.

To close this dialog, click the Close button.

Click the Help button to display the appropriate help topic for this dialog.

3.3.4 1560 Black Stack download options dialog

The Model 1560 Black Stack Thermometer readout stores the values of up to the last 1,000 readings taken in the thermometer readout's memory, and is referred to by LogWare as the history log. LogWare can download the readings from the history log and write the readings into download files. For more information on download files, please refer to Section 7.2.2, Download File. When the Download Logged Data option in the Download menu is selected and the 1560 Black Stack Thermometer readout is selected on the Select Model dialog, the 1560 Black Stack Download Options dialog is displayed.

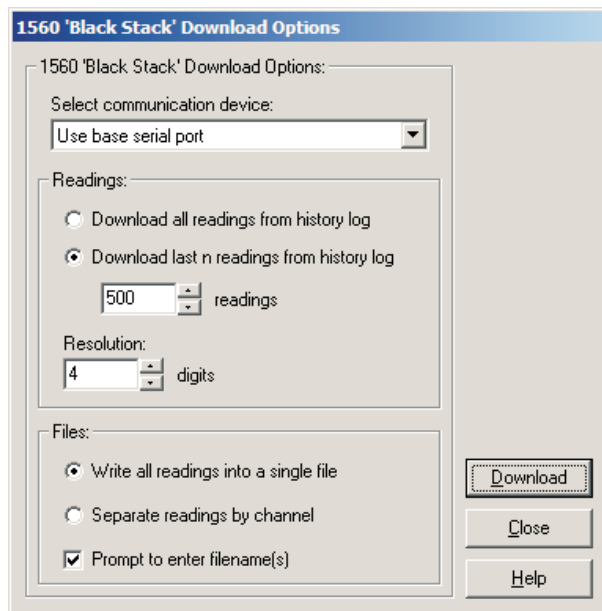


Figure 32 1560 Black Stack Download Options dialog

The Select communication device setting determines the device (serial port) to which the readings will be downloaded.

- **Use base serial port** - Select this option if the serial cable is connected to the serial port on the base unit of the 1560 (immediately behind the display).
- **Use communication module serial port** - Select this option if the serial cable is connected to the serial port on the Model 3560 Communication Module.

The Readings section allows the number of readings to be downloaded to be specified. The history log can contain up to 1,000 readings.



Note: *The actual number of readings in the history log depends on the number of channels on the thermometer readout. Typically, the history log can hold 1,000 readings as long as there are not more than 63 total input channels on the thermometer readout. For every channel over 63, the number of readings that the history log can hold decreases by approximately 16.*

Selecting the Download all readings from history log option indicates that the entire history log should be downloaded.

Selecting the Download last n readings from history log option allows a specific number of readings to be downloaded from the history log. If this option is selected, enter the number of readings to download in the readings box. This value can be set between 1 and the total number of readings in the history log.

Enter the resolution to use to format the readings when they are written to the download file in the Resolution box. This value can be set between 0 and 12 digits (places after decimal point).

Select the method to use when writing the downloaded readings to a file using the options in the Files section.

Selecting the Write all readings into a single file option causes all of the downloaded readings to be written to one file.

Selecting the Separate readings by channel option causes all of the downloaded readings to be separated according to the channel that each reading was taken from and written into separate files.

If the Prompt to enter filename(s) check box is checked, a prompt is displayed after the data is downloaded for each file that needs to be created, depending on the other selections above. If this option is not checked, the default filenames are used.



Note: *When downloading data from the 1560, the measurement process in the thermometer readout is halted until downloading is complete!*

After selecting the appropriate options on this dialog, click the Download button to start the download process. For more details on the download process, please refer to Section 8, Download Process.

To close this dialog, click the Close button.

Click the Help button to display the appropriate help topic for this dialog.

3.3.5 1575A Super-Thermometer download options dialog

The Model 1575A Super-Thermometer readout can store readings in any of two logs: history log and memory log. LogWare can download the readings from these logs and write the readings into download files. For more information on download files, please refer to Section 7.2.2, Download File. When the Download Logged Data option in the Download menu is selected and the 1575A Super-Thermometer readout is selected on the Select Model dialog, the 1575A Super-Thermometer Download Options dialog is displayed.

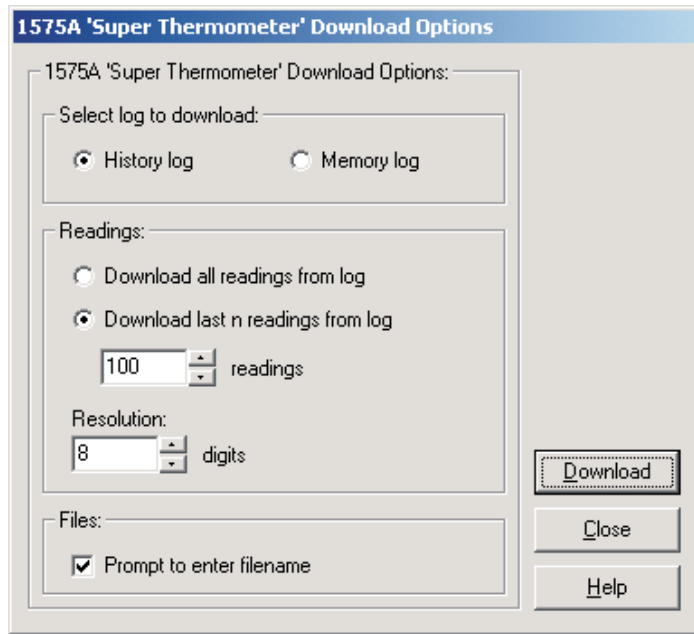


Figure 33 1575A 'Super Thermometer' Download Options dialog

The Select log to download section allows the log to be specified. The 1575A Super-Thermometer readout has two logs: history log and memory log. The history log stores the most recent 100 readings taken by the thermometer readout. The memory log can be used to store up to 100 readings manually. For more details on the history and memory logs, see the 1575A Super-Thermometer readout User's Guide.

The Readings section allows the number of readings to be downloaded to be specified.

Selecting the Download all readings from log option indicates that all of the readings in the selected log should be downloaded.

Selecting the Download last n readings from log option allows a specific number of readings to be downloaded from the selected log. If this option is se-

lected, enter the number of readings to download in the readings box. This value can be set between 1 and 100.



Note: *It is not possible for LogWare to determine exactly how many readings are currently stored in the memory log. Therefore, memory locations that do not have a value stored in them always return a value of 0 °C.*

Enter the resolution to use to format the readings when they are written to the download file in the Resolution box. This value can be set between 0 and 8 digits (places after decimal point).

Select the options regarding the file to create in the Files section.

If the Prompt to enter filename check box is checked, a prompt is displayed after the data is downloaded for the file that needs to be created. If this option is not checked, the default filename is used.

After selecting the appropriate options on this dialog, click the Download button to start the download process. For more details on the download process, please refer to Section 8, Download Process.

To close this dialog, click the Close button.

Click the Help button to display the appropriate help topic for this dialog.

3.3.6 1590 Super-Thermometer II download options

The Model 1590 Super-Thermometer II readout can store readings in any of two logs: history log and memory log. LogWare can download the readings from these logs and write the readings into download files. For more information on download files, please refer to Section 7.2.2, Download File. When the Download Logged Data option in the Download menu is selected and the 1590 Super-Thermometer II readout is selected on the Select Model dialog, the 1590 Super-Thermometer II Download Options dialog is displayed (Figure 34).

The Select log to download section allows the log to be specified. The 1590 Super-Thermometer II readout has two logs: history log and memory log. The history log stores the most recent 100 readings taken by the thermometer readout. The memory log can be used to store up to 100 readings manually. For more details on the history and memory logs, see the 1590 Super-Thermometer II readout User's Guide.

The Readings section allows the number of readings to be downloaded to be specified.

Selecting the Download all readings from log option indicates that all of the readings in the selected log should be downloaded.

Selecting the Download last n readings from log option allows a specific number of readings to be downloaded from the selected log. If this option is selected, enter the number of readings to download in the readings box. This value can be set between 1 and 100.



Note: It is not possible for LogWare to determine exactly how many readings are currently stored in the memory log. Therefore, memory locations that do not have a value stored in them always return a value of 0 °C.

Enter the resolution to use to format the readings when they are written to the download file in the Resolution box. This value can be set between 0 and 8 digits (places after decimal point).

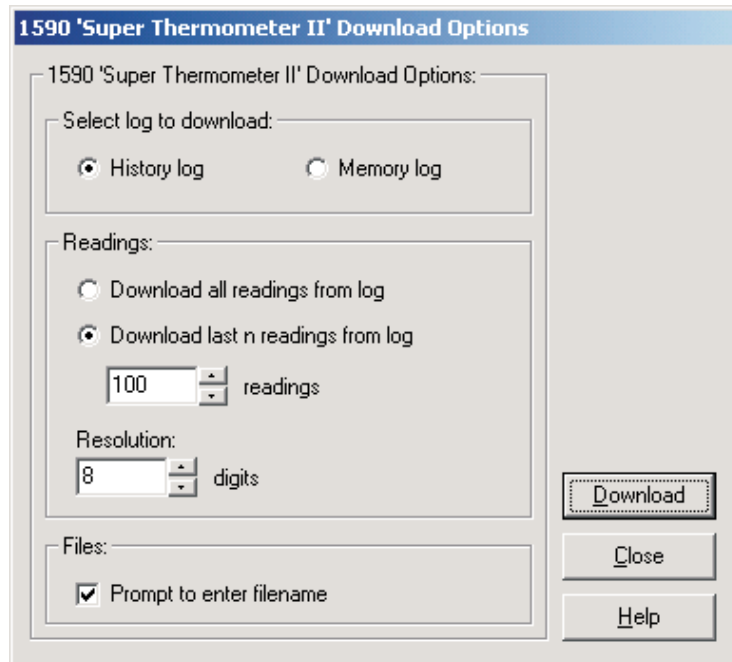


Figure 34 1590 Super-Thermometer II Download Options dialog

Select the options regarding the file to create in the Files section.

If the Prompt to enter filename check box is checked, a prompt is displayed after the data is downloaded for the file that needs to be created. If this option is not checked, the default filename is used.

After selecting the appropriate options on this dialog, click the Download button to start the download process. For more details on the download process, please refer to Section 8, Download Process.

To close this dialog, click the Close button.

Click the Help button to display the appropriate help topic for this dialog.

3.4 Log status dialogs

LogWare provides a tool to view the status of the thermometer readout's memory or logs. This tool is accessible by selecting the Log Status option in the Download menu. The only thermometer readouts supported by LogWare that have this feature are the 1529 Chub-E4 Thermometer readout and the 1560 Black Stack Thermometer readout. The following section describes the dialogs for viewing the log status for each supported thermometer readout.

3.4.1 1523 Handheld log status dialog

The Model 1523 Handheld Thermometer readout can store readings in a log, called the demand log. LogWare can display the current status of this log and provides a way to clear all readings from the log. When the Log Status option in the Download menu is selected and the 1523 Thermometer readout is selected on the Select Model dialog, the 1523 Log Status dialog is displayed.

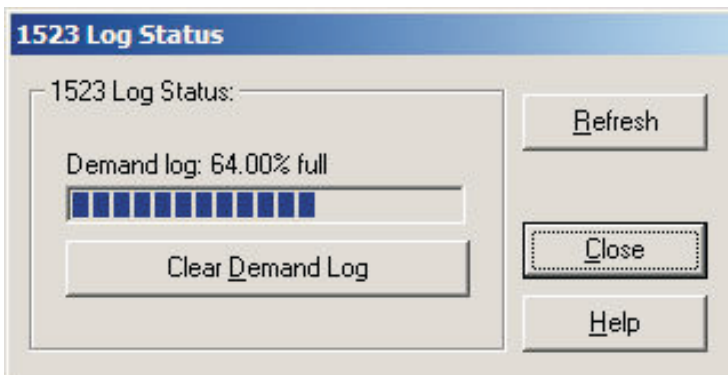


Figure 35 1523 Log Status dialog

The Demand log progress bar indicates the current state of the demand log. The demand log in the 1523 Handheld Thermometer readout holds up to 25 readings.

To clear all readings from the demand log, click the Clear Demand Log button.

To refresh the log status information on this dialog, click the Refresh button. The log status is read in from the thermometer readout again and the dialog is updated.

To close this dialog, click the Close button.

Click the Help button to display the appropriate help topic for this dialog.

3.4.2 1524 Handheld log status dialog

The Model 1524 Handheld Thermometer readout can store readings in any of two logs: demand log and auto log. LogWare can display the current status of these logs and provides a way to clear all readings from these logs. When the Log Status option in the Download menu is selected and the 1524 Thermometer readout is selected on the Select Model dialog, the 1524 Log Status dialog is displayed.

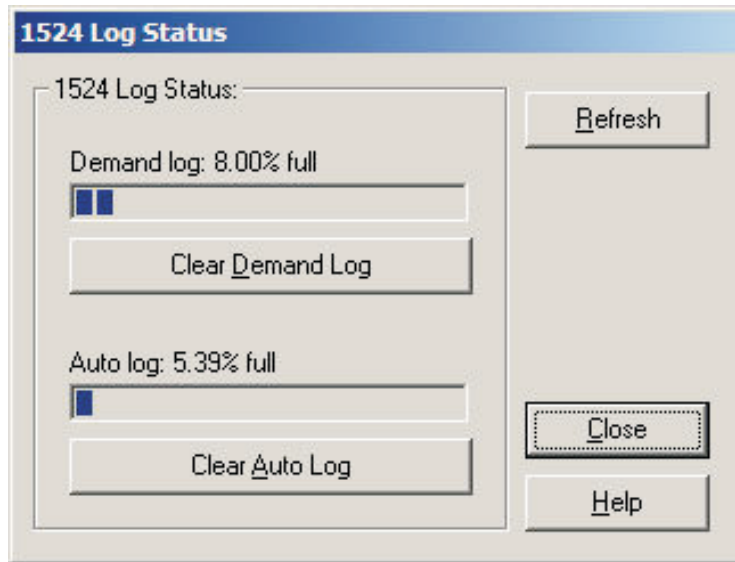


Figure 36 1524 Log Status dialog

The Demand log progress bar indicates the current state of the demand log. The demand log in the 1524 Handheld Thermometer readout holds up to 25 readings.

To clear all readings from the demand log, click the Clear Demand Log button.

The Auto log progress bar indicates the current state of the auto log. The auto log in the 1524 Handheld Thermometer readout holds up to 15,000 readings.

To clear all readings from the auto log, click the Clear Auto Log button.

To refresh the log status information on this dialog, click the Refresh button. The log status is read in from the thermometer readout again and the dialog is updated.

To close this dialog, click the Close button.

Click the Help button to display the appropriate help topic for this dialog.

3.4.3 1529 Chub-E4 log status dialog

The Model 1529 Chub-E4 Thermometer readout (including models 1529, 1529-R and 1529-T) can store readings in any of two logs: demand log and auto log. LogWare can display the current status of these logs and provides a way to clear all readings from these logs. When the Log Status option in the Download menu is selected and the 1529 Chub-E4 Thermometer readout is selected on the Select Model dialog, the 1529 Chub-E4 Log Status dialog is displayed.

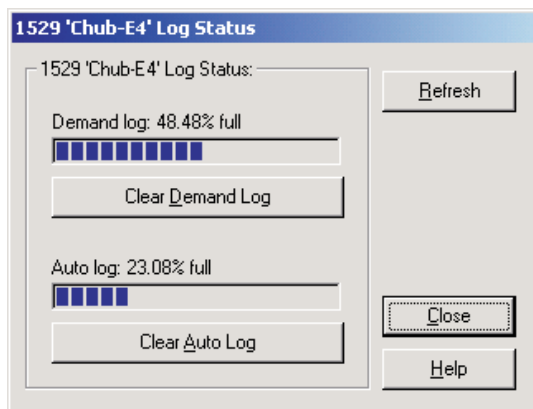


Figure 37 1529 Chub-E4 Log Status dialog

The Demand log progress bar indicates the current status of the demand log. The demand log in the 1529 Chub-E4 Thermometer readout can hold up to 100 readings.

To clear all readings from the demand log, click the Clear Demand Log button.

The Auto log progress bar indicates the current state of the auto log. The auto log in the 1529 Chub-E4 Thermometer readout can hold up to 8,160 readings.

To clear all readings from the auto log, click the Clear Auto Log button.

To refresh the log status information on this dialog, click the Refresh button. The log status is read in from the thermometer readout again and the dialog is updated.

To close this dialog, click the Close button.

Click the Help button to display the appropriate help topic for this dialog.

3.4.4 1560 Black Stack log status dialog

The Model 1560 Black Stack Thermometer readout stores the values of up to the last 1,000 readings taken in the thermometer's memory, and is referred to by LogWare as the history log. LogWare can display the current status of the

history log. When the Log Status option in the Download menu is selected and the 1560 Black Stack Thermometer readout is selected on the Select Model dialog, the 1560 Black Stack Log Status dialog is displayed.

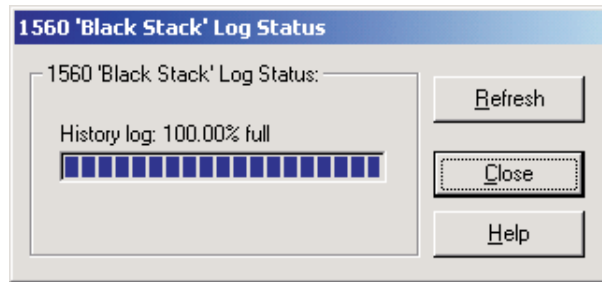


Figure 38 1560 Black Stack Log Status dialog

The history log progress bar indicates the current state of the history log. The history log in the 1560 Black Stack Thermometer readout stores the values of up to the last 1,000 readings taken. The history log of the 1560 Black Stack Thermometer readout will nearly always indicate 100% full. It is not possible for LogWare to clear the history log.



Note: The actual number of readings in the history log depends on the number of channels on the thermometer readout. Typically, the history log can hold 1,000 readings as long as there are not more than 63 total input channels on the thermometer readout. For every channel over 63, the number of readings that the history log can hold decreases by approximately 16.

To refresh the log status information on this dialog, click the Refresh button. The log status is read in from the thermometer readout again and the dialog is updated.

To close this dialog, click the Close button.

Click the Help button to display the appropriate help topic for this dialog.

4 Tools

LogWare provides tools to synchronize the thermometer readout's clock, view/modify logging labels and view/modify various thermometer readout settings including filtering, sampling, measurement, calibration, module, channel and probe settings. These tools are accessible by selecting options in the Tools menu. The following sections describe the dialogs for these tools for each of the thermometer readouts supported by LogWare.

4.1 Tools menu

The Tools menu contains options for synchronizing the thermometer readout's clock, viewing/modifying logging labels and viewing/modifying thermometer readout settings.

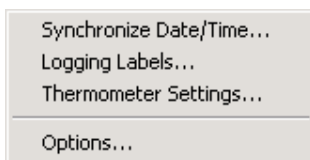


Figure 39 Tools menu

4.1.1 Synchronize date/time

The Synchronize Date/Time option can be used to synchronize the thermometer readout's clock to the computer's clock. All thermometer readouts supported by LogWare support this feature.

When selecting the Synchronize Date/Time option, the Select Model dialog is displayed. LogWare needs to know the model of the thermometer readout in order to send the proper commands to synchronize the thermometer readout's clock. Please refer to Section 4.2, Select Model Dialog, for more details on the Select Model dialog.

After the thermometer readout model and communication settings have been selected, the appropriate commands are sent to the thermometer readout to update its date and time settings (as applicable to the selected thermometer readout). After the commands have been sent, a message is displayed indicating that the thermometer readout's clock has been updated.



Note: Before synchronizing the thermometer readout's clock, make sure the computer's clock is accurate. Use the Windows® Control Panel to adjust the date and time as necessary.



Note: *The thermometer readout's display may not immediately reflect the updated time.*

4.1.2 Logging labels

The Logging Labels option can be used to view and/or modify the labels that the thermometer readout assigns to readings that are stored in the thermometer readout's log(s). This feature is only supported by the 1529 Chub-E4 Thermometer readout.

When selecting the Logging Labels option, the Select Model dialog is displayed. LogWare needs to know the model of the thermometer readout. Please refer to Section 4.2, Select Model Dialog, for more details on the Select Model dialog.

After the thermometer readout model and communication settings have been selected, the Logging Labels dialog for the selected thermometer readout is displayed. The Logging Labels dialog is described in more detail in Section 4.3, Logging Labels Dialogs.

4.1.3 Thermometer settings

The Thermometer Settings option can be used to view and/or modify many thermometer readout, module, and probe settings including filtering, sampling, measurement, calibration, and conversion settings. All thermometer readouts supported by LogWare support this feature. Please refer to Section 1.4.4, Firmware Requirements, for more information.

When selecting the Thermometer Settings option, the Select Model dialog is displayed. LogWare needs to know the model of the thermometer readout in order to display the appropriate thermometer readout settings dialogs. Please refer to Section 4.2, Select Model Dialog, for more details on the Select Model dialog.

After the thermometer readout model and communication settings have been selected, the appropriate Thermometer Settings dialog for the selected thermometer readout is displayed. The Thermometer Settings dialogs are described in more detail in Section 4.4, Thermometer Settings Dialogs.



Warning: *When using the Thermometer Settings tool, make sure that the thermometer readout is not currently in any menu. Using the thermometer readout's menus while using the Thermometer Settings tool may cause the thermometer readout to fail to accept new settings.*

4.2 Select Model dialog

When any of the Tools menu options is selected, the Select Model dialog is displayed. LogWare needs to know the model of the thermometer readout to interface with.

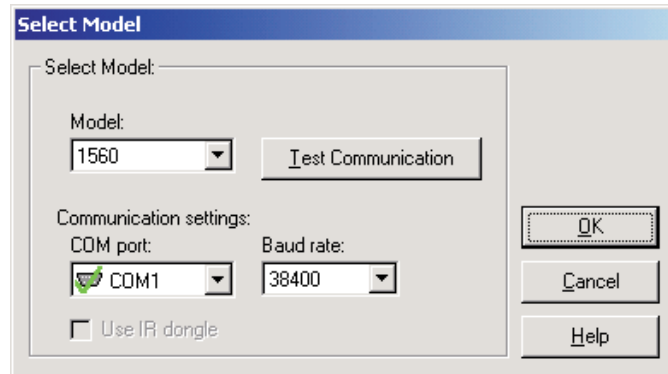







Figure 40 Select Model dialog

If this dialog has been used previously to select a thermometer readout model and communication settings, the previous settings are restored.

Use the Model drop-down list to select the type of thermometer readout being used. All other controls on this dialog are disabled until a model is selected. The list of available thermometer readouts depends on the Tools menu option selected. Only thermometer readouts that support the selected feature are listed.

Select the COM port on the computer to which the thermometer readout is connected using the COM port drop-down list. Only COM ports that the operating system reports as being available can be used. The following symbols are used to show the status of each COM port:

-  - Indicates a COM port that is currently available.
-  - Indicates the COM port that is currently selected by LogWare.
-  - Indicates a COM port that is currently being used by another application.
-  - Indicates an IR port (virtual COM port) that is currently available.
-  - Indicates an IR port (virtual COM port) that is currently selected by LogWare.

-  - Indicates an IR port (virtual COM port) that is currently being used by another application.

Select the baud rate that the computer should use to communicate over the COM port with the thermometer readout using the Baud rate drop-down list. The baud rate setting of the thermometer readout must be set to the same value as this setting. For best results, use the maximum allowable baud rate setting when possible.

Check the Use IR dongle option to use the thermometer readout's IR port instead of an RS-232 cable. This option is only enabled when the selected thermometer readout has an IR port and the selected feature allows using the IR port to communicate with the thermometer readout. For more information on using the IR dongle, please refer to Section 6.13, Using an IR Dongle with LogWare.

After the model, COM port, baud rate and/or IR settings are configured, click the Test Communication button to make sure communication with the thermometer readout is successful. At this time, the thermometer readout must be connected to the COM port on the computer using a null modem cable or, if the IR dongle is being used, the IR dongle must be connected to the selected COM port and the dongle must be aligned with the thermometer readout's IR port. For more information, please refer to the thermometer readout's User's Guide.

When testing communication with the thermometer readout, there are three possible results:

- The thermometer readout is communicating at the correct baud rate on the selected COM port.
- The thermometer readout is communicating, but possibly at the wrong baud rate or the thermometer readout model does not match the selected model.
- The thermometer readout failed to communicate at all. Check the COM port and baud rate settings, make sure the cable is a null modem cable, and make sure the thermometer readout is plugged in and turned on. If using the IR dongle, make sure the dongle is connected to the selected COM port and the dongle is aligned properly with the thermometer readout's IR port.

After all of the settings on this dialog are configured and communication with the thermometer readout is successful, click the OK button to accept the settings and continue.

Click the Cancel button close this dialog and abort using the selected feature.

Click the Help button to display the appropriate help topic for this dialog.

4.3 Logging labels dialogs

LogWare provides a tool to view and/or modify the labels that the thermometer readout assigns to readings that are stored in the thermometer readout's log(s). This tool is accessible by selecting the Logging Labels option in the Tools menu. The following sections describe the dialogs for each of the thermometer readouts that support this feature.

4.3.1 1523 Handheld logging label dialog

The Model 1523 Handheld Thermometer readout can store readings into a log, called the demand log. When readings are stored in the log, a label can be assigned to each reading to indicate the reason, purpose or to otherwise identify the reading(s). LogWare provides a tool that can be used to view and/or modify the current logging label. When the Logging Labels option in the Tools menu is selected and the 1523 Thermometer readout is selected on the Select Model dialog, the 1523 Logging Label dialog is displayed.



Figure 41 1523 Logging Label dialog

When the 1523 Logging Label dialog is displayed, the current label name is read in from the thermometer readout.

The 1523 Handheld Thermometer readout has a single logging label for the demand log. The Label name box indicates the current name for this label. Enter the new name for the demand log label.

After the label has been modified as needed, click the OK button to write the new label name to the thermometer readout. A progress dialog is displayed while the new label name is being written to the thermometer readout.

If the label was not modified, or to close this dialog without writing the new label name to the thermometer readout, click the Close button.

Click the Help button to display the appropriate help topic for this dialog.

4.3.2 1524 Handheld logging labels dialog

The Model 1524 Handheld Thermometer readout can store readings into logs. When readings are stored in a log, a label can be assigned to each reading or set of readings to indicate the reason, purpose or to otherwise identify the reading(s). LogWare provides a tool that can be used to view and/or modify the current logging label settings. When the Logging Labels option in the Tools menu is selected and the 1524 Thermometer readout is selected on the Select Model dialog, the 1524 Logging Labels dialog is displayed.

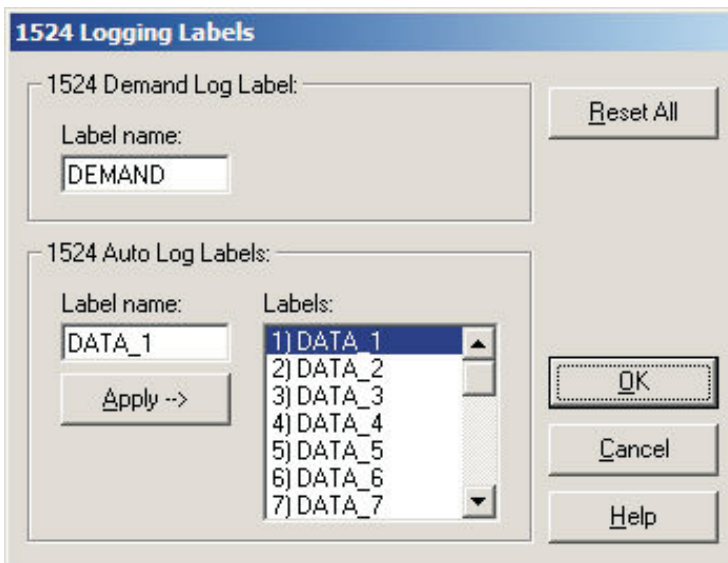


Figure 42 1524 Logging Labels dialog

When the 1524 Logging Labels dialog is displayed, all of the current label names are read in from the thermometer readout. The label used for the demand log is displayed in the 1524 Demand Log Label section. The labels used for the auto log are displayed in the 1524 Auto Log Labels section.

The 1524 Handheld Thermometer readout has a single logging label for the demand log. The Label name box indicates the current name for this label. Enter the new name for the demand log label.

The Labels list in the 1524 Auto Log Labels section indicates the current names of all auto log labels. The 1524 Handheld Thermometer readout has 25 logging labels for the auto log.

To modify any of one of the labels, click on the appropriate label in the Labels list. The Label name box indicates the name of the label that is currently selected. Enter the new name for the selected label and press the Enter key or click the Apply —> button. The Labels list is updated to reflect the change. If

the Enter key was pressed, the next label in the list will automatically be selected and the name of the next label can be entered in the Label name box.

Repeat this process for all labels that need to be modified.

To reset all labels to their default names, click the Reset All button. The Labels list is updated to show the default names for all labels.

The 1524 Handheld Thermometer readout has a single logging label for the demand log. The Demand Log Label box indicates the current name for this label. Enter the new name for the demand log label.

After all labels have been modified as needed, click the OK button to write the new label names to the thermometer readout. A progress dialog is displayed while the new label names are being written to the thermometer readout.

If no labels were modified, or to close this dialog without writing the new label names to the thermometer readout, click the Close button.

Click the Help button to display the appropriate help topic for this dialog.

4.3.3 1529 Chub-E4 logging labels

The Model 1529 Chub-E4 Thermometer readout (including models 1529, 1529-R and 1529-T) can store readings into logs. When readings are stored in a log, a label can be assigned to each reading or set of readings to indicate the reason, purpose or to otherwise identify the reading(s). LogWare provides a tool that can be used to view and/or modify the current logging label settings. When the Logging Labels option in the Tools menu is selected and the 1529 Chub-E4 Thermometer readout is selected on the Select Model dialog, the 1529 Chub-E4 Logging Labels dialog is displayed.

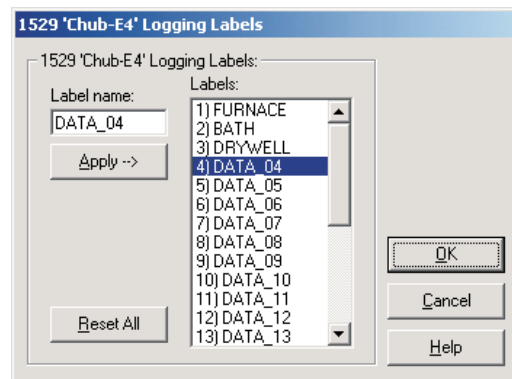


Figure 43 1529 Chub-E4 Logging Labels dialog

When the Logging Labels dialog is displayed, the current logging labels are read in from the thermometer readout. The Labels list indicates the current logging labels. The 1529 Chub-E4 Thermometer readout has 25 logging labels.

To modify any of one of the labels, click on the appropriate label in the Labels list. The Label name box indicates the name of the label that is currently selected. Enter the new name for the selected label and press the Enter key or click the Apply --> button. The Labels list is updated to reflect the change. If the Enter key was pressed, the next label in the list will automatically be selected and the name of the next label can be entered in the Label name box.

Repeat this process for all labels that need to be modified.

To reset all labels to their default names, click the Reset All button. The Labels list is updated to show the default names for all labels.

After all labels have been modified as needed, click the OK button to write the new label names to the thermometer readout. A progress dialog is displayed while the new label names are being written to the thermometer readout.

If no labels were modified, or to close this dialog without writing the new label names to the thermometer readout, click the Close button.

Click the Help button to display the appropriate help topic for this dialog.

4.4 Thermometer settings dialogs

LogWare provides tools to view and/or modify various thermometer readout settings including filtering, sampling, measurement, calibration, module, channel, and probe settings. These tools are accessible by selecting the Thermometer Settings option in the Tools menu. The following sections describe the dialogs for each of the thermometer readouts that support this feature.

4.4.1 1523/1524 Handheld Settings

The 1523/1524 Settings dialog displays information about the thermometer readout including model number, serial number, identification information and

the total number of channels. It also allows many thermometer readout settings to be viewed and modified.

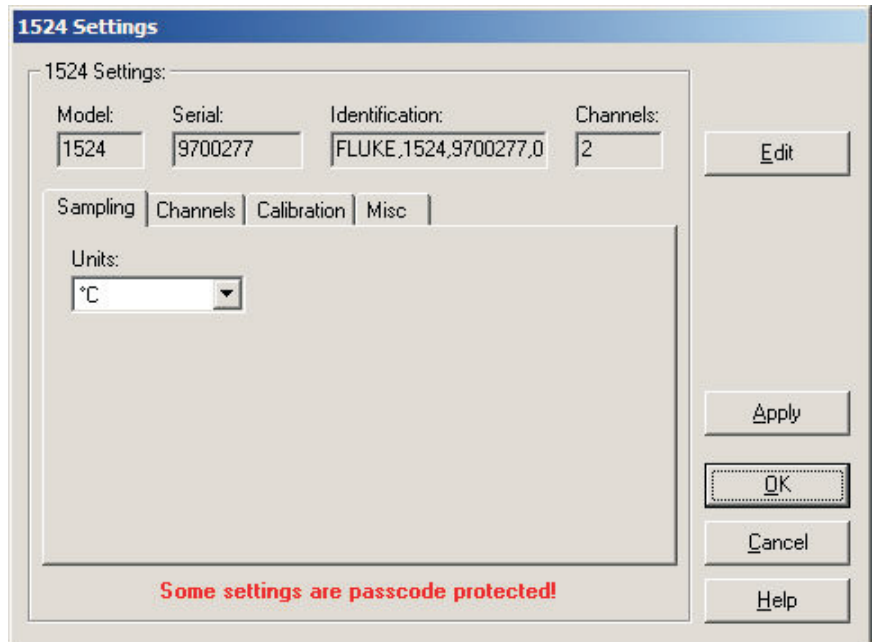


Figure 44 1523/1524 Settings dialog

The 1523/1524 Handheld Thermometer readout must be connected to an available COM port and the communication settings must be properly configured so that this information can be read in from the thermometer readout. If communication with the thermometer readout is not successful, an error message is displayed and this dialog is blank.

The Model box indicates the model number of the thermometer readout. The Serial box indicates the serial number assigned to this thermometer readout. The Identification box indicates the manufacturer name, model number, serial number and firmware version installed in this thermometer readout. The Channels box indicates the total number of input channels on this thermometer readout.

The sampling settings can be viewed and modified by selecting the Sampling tab. The Channels tab displays a list of channels and allows access to channel and probe settings dialogs. The thermometer readout's calibration settings can be viewed and modified by selecting the Calibration tab. Other settings can be viewed and modified by selecting the Misc tab.

Some of the settings on the 1523/1524 Settings dialog may be passcode protected. If they are, a message appears at the bottom of the dialog and the Edit button is displayed. The controls for the passcode protected settings are dis-

abled. To edit the protected settings, click the Edit button and enter the current passcode for the thermometer readout. If the correct passcode is entered, the controls are enabled and the values may be changed. The controls remain enabled until this dialog is closed or until the current changes are applied by clicking the Apply button.

After modifying any of the settings on this dialog, click the Apply or OK button to write the new settings to the thermometer readout. If the Apply button is clicked, this dialog remains open. If the OK button is clicked, this dialog closes after writing the settings.

LogWare has a feature that causes all of the thermometer readout settings to be read back in from the thermometer readout after the Apply button is clicked and the new settings have been written to the thermometer readout. This feature can be useful to verify that the desired settings were properly applied to the thermometer readout. To enable or disable this feature, see Section 5.6.4, Options.

To close this dialog without writing the settings to the thermometer readout, click the Cancel button.

Click the Help button to open the appropriate help topic for this dialog.

4.4.1.1 Sampling tab

Select the Sampling tab to view and/or modify sampling settings.



Figure 45 1523/1524 Settings – Sampling tab

Select the units to use when displaying readings in temperature using the Units drop-down list. Temperature readings can be displayed in either °C or °F. This setting applies to all channels that are configured to display readings in temperature.

Select the Channels tab to view a list of channels and to access the channel and probe settings dialogs to view and/or modify channel and probe specific settings.

4.4.1.2 Channels tab

Select the channel to view or modify from the Channels list. This list indicates the channel number, channel type and the serial number of the probe on each channel.

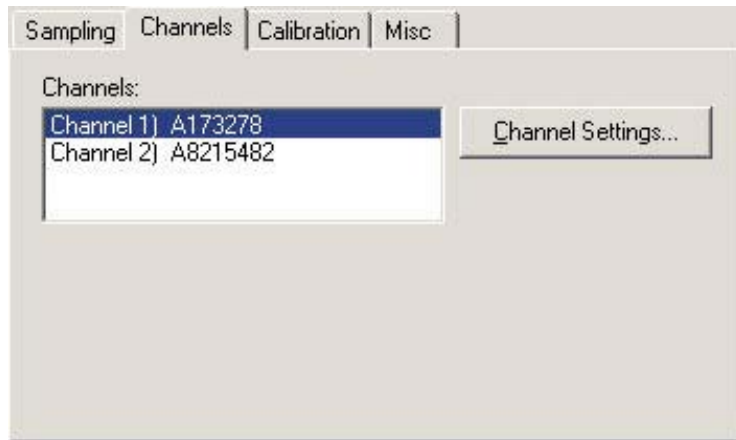


Figure 46 1523/1524 Settings – Channels tab

Click the Channel Settings button to display the appropriate channel settings dialog. When this button is clicked, the settings for the selected channel and probe are read in from the thermometer readout.

4.4.1.3 Calibration tab

Select the Calibration tab to view and/or modify the thermometer readout's calibration settings.

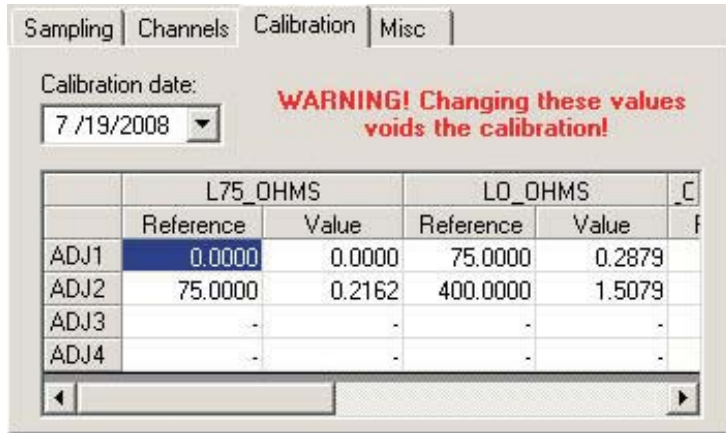


Figure 47 1523/1524 Settings – Calibration tab



CAUTION!! DO NOT CHANGE THESE VALUES UNLESS THIS READOUT HAS BEEN RECALIBRATED OR UNLESS THESE VALUES DO NOT MATCH THE VALUES LOCATED ON THE LATEST REPORT OF CALIBRATION! CHANGING THESE VALUES VOIDS THE CURRENT CALIBRATION!

Select the date this readout was last calibrated using the drop-down calendar or by entering the date. This date should be located on the Report of Calibration.

Enter the reference and offset values for each calibration range in the grid. The calibration ranges are indicated in the first row of the grid. Each calibration range can have up to 4 adjustment pairs consisting of a reference and an offset value. Reference and offset values for unused adjustment pairs must be left blank. These values can be found on the Report of Calibration for the readout only if this readout received an accredited calibration.

4.4.1.4 Misc tab

Select the Misc tab to view and/or modify other miscellaneous thermometer readout settings.

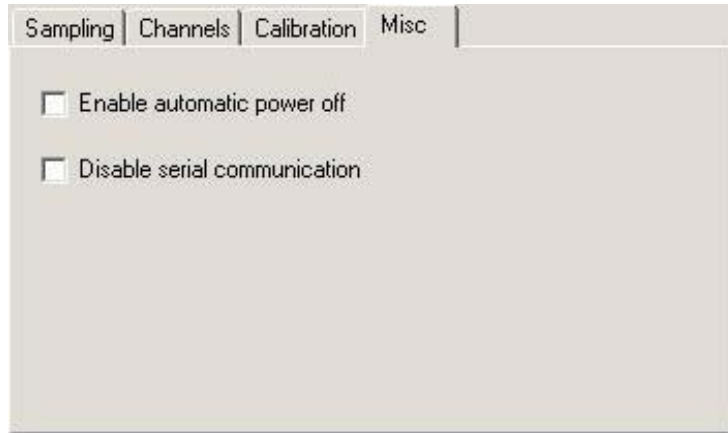


Figure 48 1523/1524 Settings – Misc tab

The Enable automatic power off setting determines whether the thermometer readout will automatically shut itself off after a specific amount of time. By checking this option, the time period is set to 30 minutes.

The Disable serial communication setting determines whether the serial port of the thermometer readout is enabled. Disabling the serial port increases battery life. If the serial port is disabled, it can only be re-enabled using the thermometer readout's SETUP menu.

4.4.1.5 1523/1524 PRT/thermistor channel settings dialog

The 1523/1524 PRT/Thermistor Channel Settings dialog displays information about the selected channel.

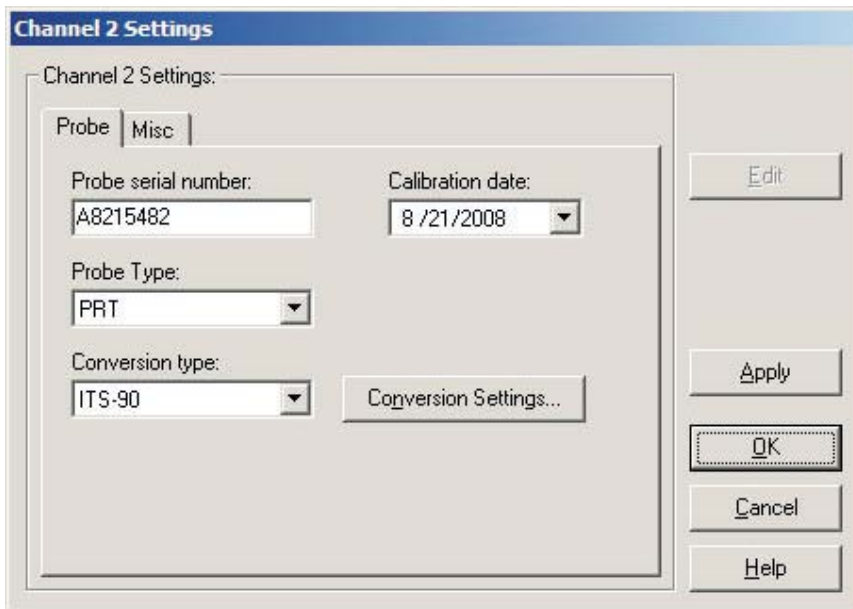


Figure 49 1523/1524 PRT/Thermistor Channel Settings dialog

Some or all of the settings on this dialog may be passcode protected. If so, a message appears at the bottom of the Probe tab and the Edit button is displayed. The protected controls are disabled. To edit the protected settings, click the Edit button and enter the current passcode for the thermometer readout. If the correct passcode is entered, the protected controls are enabled and the values may be changed. The controls remain enabled until this dialog is closed or until the current changes are applied by clicking the Apply button.

After modifying any of the settings on this dialog, click the Apply or OK button to write the new settings to the thermometer readout. If the Apply button is clicked, this dialog remains open. If the OK button is clicked, this dialog closes after writing the settings.

LogWare has a feature that causes all of the channel settings to be read back in from the thermometer readout after the Apply button is clicked and the new settings have been written to the thermometer readout. This feature can be useful to verify that the desired settings were properly applied to the thermometer readout. To enable or disable this feature, see Section 5.6.4, Options.

To close this dialog without writing the channel settings to the thermometer readout, click the Cancel button.

Click the Help button to open the appropriate help topic for this dialog.

4.4.1.5.1 Probe tab

Select the Probe tab to view and/or modify probe settings.

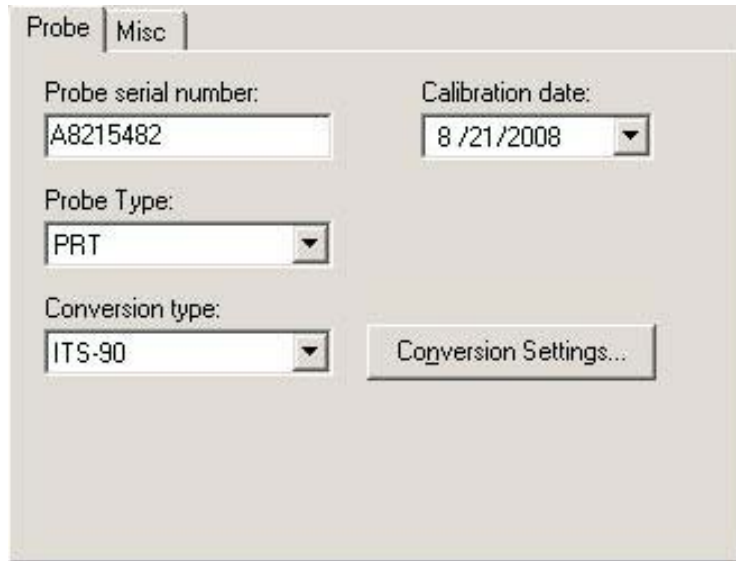


Figure 50 1523/1524 PRT/Thermistor Channel Settings – Probe tab

Enter the serial number of the probe attached to this channel in the probe serial number box. The serial number can be any combination of numbers and/or letters up to 8 characters in length. The underscore “_” character can also be used in serial numbers. Select the date this probe was last calibrated using the drop-down calendar or by entering the date. This date should be located on the Report of Calibration for the probe.

Select the appropriate probe type for this probe using the Probe Type drop-down list. The choices are PRT or Thermistor.

Select the conversion type to use for the probe on this channel. The conversion types available in the drop-down list are dependent on the Probe Type selection.

For PRT probe types, the available conversion types are:

- ITS-90
- ITS-90 Subrange 5
- PT100
- Callendar-Van Dusen
- Ohms

For Thermistor probe types, the available conversion types are:

- R(T)
- Ohms

Refer to the Report of Calibration for the probe for information on the conversion type that the probe was calibrated for.

If the selected conversion type requires coefficients or other settings, the Conversion Settings button is enabled. Click the Conversion Settings button to display the dialog for the selected conversion type. Otherwise, the Conversion Settings button is disabled.



Note: The conversion settings are not written to the thermometer readout until the Apply or OK button is clicked on this dialog!

4.4.1.5.2

Misc tab

Select the Misc tab to view and/or modify other probe settings.

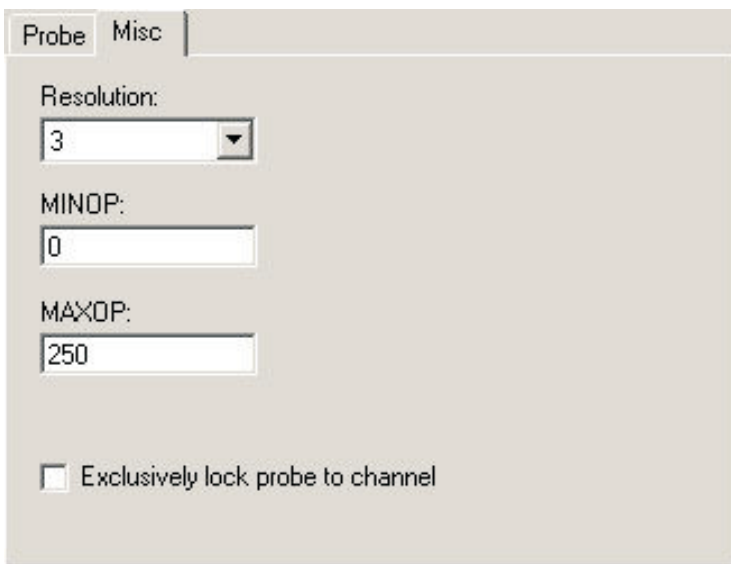


Figure 51 1523/1524 PRT/Thermistor Channel Settings – Misc tab

Select the resolution for the readings for this probe using the Resolution drop-down list. The resolution can be set to any value between 0 and 3, or Auto.

Enter the minimum operating temperature for this probe in the MINOP box. If the temperature measured by the thermometer readout is less than this setting, the display indicates an overload condition (OL).

Enter the maximum operating temperature for this probe in the MAXOP box. If the temperature measured by the thermometer readout is greater than this setting, the display indicates an overload condition (OL).

To exclusively lock the connected probe to this channel, check the Exclusively lock probe to channel checkbox. Locking a probe to a channel prevents any other probe from being used on this channel, and also prevents this probe from being used on any other channel, including channels of other thermometer readouts.

4.4.1.6 1523/1524 thermocouple channel settings dialog

The Channel Settings dialog for thermocouple probes displays information about the selected channel.

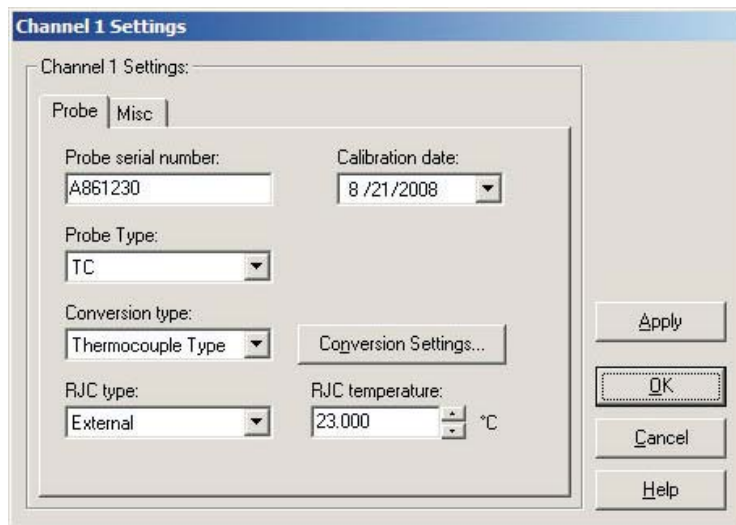


Figure 52 1523/1524 Thermocouple Channel Settings dialog

Some or all of the settings on this dialog may be passcode protected. If so, a message appears at the bottom of the Probe tab and the Edit button is displayed. The protected controls are disabled. To edit the protected settings, click the Edit button and enter the current passcode for the thermometer readout. If the correct passcode is entered, the protected controls are enabled and the values may be changed. The controls remain enabled until this dialog is closed or until the current changes are applied by clicking the Apply button.

After modifying any of the settings on this dialog, click the Apply or OK button to write the new settings to the thermometer readout. If the Apply button is

clicked, this dialog remains open. If the OK button is clicked, this dialog closes after writing the settings.

LogWare has a feature that causes all of the channel settings to be read back in from the thermometer readout after the Apply button is clicked and the new settings have been written to the thermometer readout. This feature can be useful to verify that the desired settings were properly applied to the thermometer readout. To enable or disable this feature, see Section 5.6.4, Options.

To close this dialog without writing the channel settings to the thermometer readout, click the Cancel button.

Click the Help button to open the appropriate help topic for this dialog.

4.4.1.6.1 Probe tab

Select the Probe tab to view and/or modify probe settings.

Figure 53 1523/1524 Thermocouple Channel Settings – Probe tab

Enter the serial number of the probe attached to this channel in the probe serial number box. The serial number can be any combination of numbers and/or letters up to 8 characters in length. The underscore “_” character can also be used in serial numbers. Select the date this probe was last calibrated using the drop-down calendar or by entering the date. This date should be located on the Report of Calibration for the probe.

Select the appropriate probe type for this probe using the Probe Type drop-down list. Currently, the only choice is TC.

Select the conversion type to use for the probe on this channel. For TC probe types, the available conversion types are:

- Thermocouple Type
- Volts

Refer to the Report of Calibration for the probe for information on the conversion type that the probe was calibrated for.

If the selected conversion type requires coefficients or other settings, the Conversion Settings button is enabled. Click the Conversion Settings button to display the dialog for the selected conversion type. Otherwise, the Conversion Settings button is disabled.



***Note:** The conversion settings are not written to the thermometer readout until the Apply or OK button is clicked on this dialog!*

Select the type of reference junction compensation to use on this channel using the RJC type drop-down list. Available choices are Internal and External. If External is chosen as the RJC type, the external RJC temperature must be entered in the RJC temperature box.


Enter the external RJC temperature value in the RJC temperature box. This box is disabled if the RJC type is set to Internal.



***Note:** The RJC temperature value must always be in °C!*

4.4.1.6.2 Misc tab

Select the Misc tab to view and/or modify other probe settings.

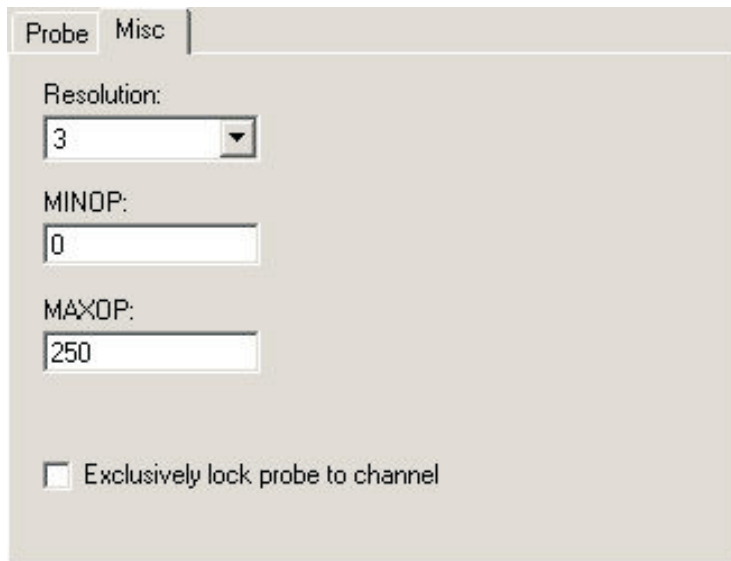

The image shows a software dialog box titled "1523/1524 Thermocouple Channel Settings" with two tabs: "Probe" and "Misc". The "Misc" tab is active. It contains the following controls:

- Resolution:** A drop-down menu with the value "3" selected.
- MINOP:** A text input field containing the value "0".
- MAXOP:** A text input field containing the value "250".
- Exclusively lock probe to channel:** An unchecked checkbox.

Figure 54 1523/1524 Thermocouple Channel Settings – Misc tab

Select the resolution for the readings for this probe using the Resolution drop-down list. The resolution can be set to any value between 0 and 3, or Auto.

Enter the minimum operating temperature for this probe in the MINOP box. If the temperature measured by the thermometer readout is less than this setting, the display indicates an overload condition (OL).

Enter the maximum operating temperature for this probe in the MAXOP box. If the temperature measured by the thermometer readout is greater than this setting, the display indicates an overload condition (OL).

To exclusively lock the connected probe to this channel, check the Exclusively lock probe to channel checkbox. Locking a probe to a channel prevents any other probe from being used on this channel, and also prevents this probe from being used on any other channel, including channels of other thermometer readouts.

4.4.2 1529 Chub-E4 settings dialog

The 1529 Chub-E4 Settings dialog displays information about the thermometer readout including model number, serial number, identification information, and

the total number of channels. It also allows many thermometer readout settings to be viewed and/or modified.

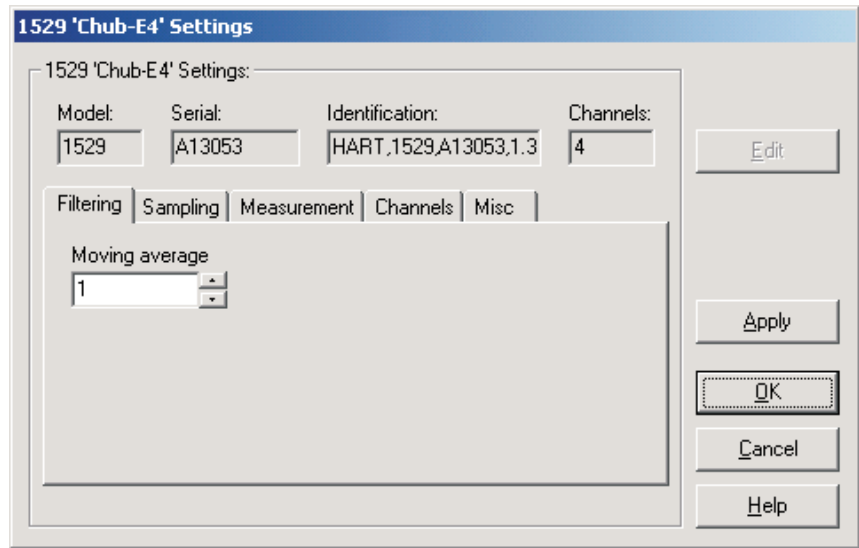


Figure 55 1529 Chub-E4 Settings dialog

The 1529 Chub-E4 must be connected to an available COM port or IR dongle and the communication settings must be properly configured so that this information can be read in from the thermometer readout. If communication with the thermometer readout is not successful, an error message is displayed and this dialog is blank. A progress dialog is displayed while the current thermometer readout settings are being read in.

The Model box indicates the model number of the thermometer readout. The Serial box indicates the serial number assigned to this thermometer readout. The Identification box indicates the manufacturer name, model number, serial number, and firmware version installed in this thermometer readout. The Channels box indicates the total number of input channels on this thermometer readout.

The filtering settings can be viewed and modified by selecting the Filtering tab. The sampling settings can be viewed and modified by selecting the Sampling tab. The measurement settings can be viewed and modified by selecting the Measurement tab. The Channels tab displays a list of channels and allows access to channel and probe settings dialogs. The settings on each of these tabs are described in detail in the following sections. For more details on these settings, please refer to the thermometer readout's User's Guide.

Some of the settings on the 1529 Chub-E4 Settings dialog may be passcode protected. If they are, a message appears at the bottom of the dialog and the Edit button is displayed. The controls for the passcode protected settings are

disabled. To edit the protected settings, click the Edit button and enter the current passcode for the thermometer readout. If the correct passcode is entered, the controls are enabled and the values may be changed. The controls remain enabled until this dialog is closed or until the current changes are applied by clicking the Apply button.

After modifying any of the settings on this dialog, click the Apply or OK button to write the new settings to the thermometer readout. If the Apply button is clicked, this dialog remains open. If the OK button is clicked, this dialog closes after writing the settings.

LogWare has a feature that causes all of the thermometer readout settings to be read back in from the thermometer readout after the Apply button is clicked and the new settings have been written to the thermometer readout. This feature can be useful to verify that the desired settings were properly applied to the thermometer readout. To enable or disable this feature, see Section 5.6.4, Options.

To close this dialog without writing the settings to the thermometer readout, click the Cancel button.

Click the Help button to open the appropriate help topic for this dialog.

4.4.2.1 Filtering tab

Select the Filtering tab to view and/or modify the filter settings.

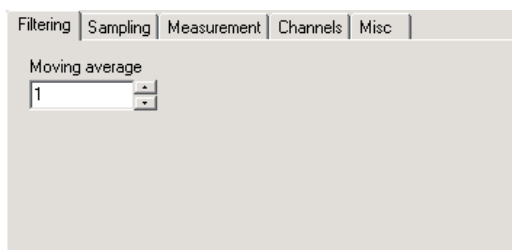


Figure 56 1529 Chub-E4 Settings - Filtering tab

The Moving average setting is sometimes also called a sliding window filter. Enter the number of readings to average (size of window) in the box provided. Setting this option to any value other than 1 causes the thermometer readout to display the average of the last n readings on a specific channel as the current reading for that channel. If the moving average is set to 1, the most recent measurement on each channel is displayed as the current reading. The moving average can be set to any value between 1 and 10.

4.4.2.2 Sampling tab

Select the Sampling tab to view and/or modify sampling settings.

Figure 57 1529 Chub-E4 Settings - Sampling tab

Select the units to use when displaying readings in temperature using the Units drop-down list. Temperature readings can be displayed in either °C, °F or K. This setting applies to all channels that are configured to display readings in temperature.

Select the resolution for the readings using the Resolution drop-down list. The resolution can be set to any value between 0 and 4, or Auto.

Select the time interval between subsequent measurements using the Sample interval drop-down list. The sample interval can be set to any one of the following settings:

- 0.1, 0.2 or 0.5 seconds
- 1, 2, 5, 10 or 30 seconds
- 1, 2, 5, 10 or 30 minutes
- 1 hour



Note: Setting the sample interval to 0.1 or 0.2 seconds causes the 1529 to enter a special “fast” sampling mode. Please refer to the 1529 Chub-E4 User’s Guide for more information on using sample interval settings of less than 0.5 seconds.

Select the decimal format character using the Decimal format drop-down list. The 1529 Chub-E4 Thermometer readout supports using either the period (.) or comma (,) as the decimal separator.

Select the date format using the Date format drop-down list. The 1529 Chub-E4 Thermometer readout supports displaying the date in one of four formats:

- MM-DD-YY
- MM-DD-YYYY
- DD/MM/YY

- DD/MM/YYYY

Select the time format using the Time format drop-down list. The 1529 Chub-E4 Thermometer readout supports displaying the time in either 12-hour or 24-hour format.

Checking the Stamp all measurements with date and time option causes the 1529 Chub-E4 Thermometer readout to always send date and time information with the measurement when querying measurements using an RS-232, IR, or IEEE interface.

4.4.2.3 Measurement tab

Select the Measurement tab to view and/or modify measurement related settings.

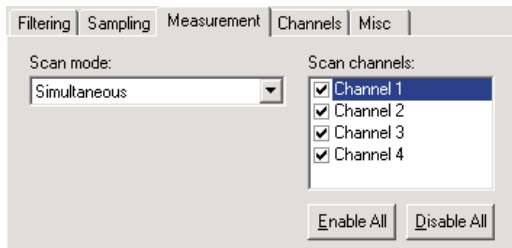


Figure 58 1529 Chub-E4 Settings - Measurement Tab

Select the scan mode for the 1529 Chub-E4 Thermometer readout using the Scan mode drop-down list. The scan mode can be set to one of the following options:

- Scan
- Simultaneous

Select the channels that the thermometer readout should measure. Each channel can be enabled or disabled by checking or unchecking the box next to the channel number independent of all other channels.

To enable all channels, click the Enable All button. This causes the box next to every channel to be checked.

To disable all channels, click the Disable All button. This causes the box next to every channel to be unchecked.

4.4.2.4 Channels Tab

Select the Channels tab to view a list of channels and to access the channel and

probe settings dialogs to view and/or modify channel and probe specific settings.

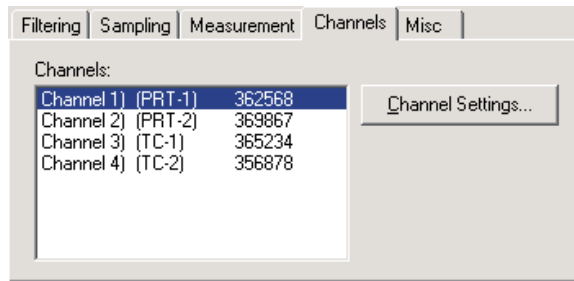


Figure 60 1529 Chub-E4 Settings - Channels tab

Select the channel to view and/or modify from the Channels list. This list indicates the channel number, channel type and the serial number of the probe on each channel.

Click the Channel Settings button to display the appropriate Channel Settings dialog. When this button is clicked, the settings for the selected channel and probe are read in from the thermometer readout. The Channel Settings dialogs are described in detail in Section 4.4.2.6, 1529 Chub-E4 PRT/Thermistor Channel Settings Dialog and Section 4.4.2.7, 1529 Chub-E4 Thermocouple Channel Settings Dialog.

4.4.2.5 Misc tab

Select the Misc tab to view and/or modify other miscellaneous thermometer readout settings.

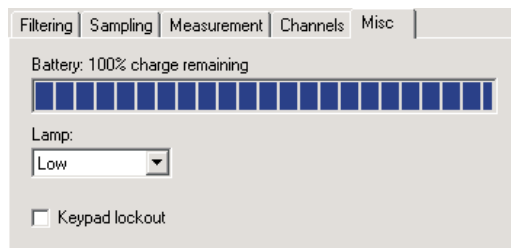


Figure 59 1529 Chub-E4 Settings - Misc tab

The Battery indicator displays the remaining charge available in the battery.

Select the setting for the lamp on the thermometer readout's display using the Lamp drop-down list. The lamp can be set to one of the following settings:

- Off

- Low
- Medium
- High

To prevent the buttons on the front panel of the thermometer readout from being used to view and/or modify thermometer readout settings, check the Keypad lockout box. Leaving this box unchecked allows the buttons on the front panel to be used to view and/or modify thermometer readout settings.

4.4.2.6 1529 Chub-E4 PRT/thermistor channel settings dialog

The 1529 Chub-E4 PRT/Thermistor Channel Settings dialog displays information about the selected PRT/Thermistor channel.

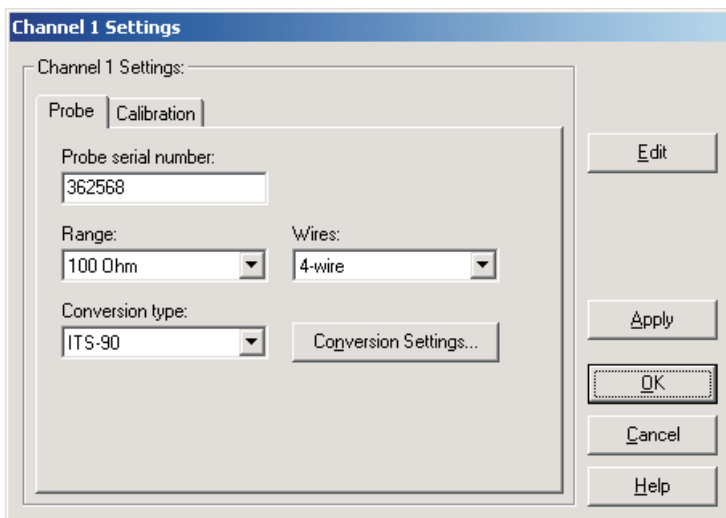


Figure 61 1529 Chub-E4 PRT/thermistor channel settings dialog

Some or all of the settings on this dialog may be passcode protected. If so, a message appears at the bottom of the Probe tab and/or Calibration tab and the Edit button is displayed. The protected controls are disabled. To edit the protected settings, click the Edit button and enter the current passcode for the thermometer readout. If the correct passcode is entered, the protected controls are enabled and the values may be changed. The controls remain enabled until this dialog is closed or until the current changes are applied by clicking the Apply button.

After modifying any of the settings on this dialog, click the Apply or OK button to write the new settings to the thermometer readout. If the Apply button is

clicked, this dialog remains open. If the OK button is clicked, this dialog closes after writing the settings.

LogWare has a feature that causes all of the channel settings to be read back in from the thermometer readout after the Apply button is clicked and the new settings have been written to the thermometer readout. This feature can be useful to verify that the desired settings were properly applied to the thermometer readout. To enable or disable this feature, see Section 5.6.4, Options.

To close this dialog without writing the channel settings to the thermometer readout, click the Cancel button.

Click the Help button to open the appropriate help topic for this dialog.

4.4.2.6.1 Probe tab

Select the Probe tab to view and/or modify probe settings.

Figure 62 1529 Chub-E4 PRT/thermistor channel settings - Probe tab

Enter the serial number of the probe attached to this channel in the Probe serial number box. The serial number can be any combination of numbers and/or letters up to 8 characters in length. The underscore “_” character can also be used in serial numbers.

Select the resistance range for this channel using the Range drop-down list. The choices for the resistance range are 100 Ohm or 10K Ohm.

Select the wires for this channel using the Wires drop-down list. The choices for the wires setting are 2-wire, 3-wire or 4-wire.

Select the conversion type to use for the probe on this channel. The Model 1529 Chub-E4 Thermometer readout supports the following conversion types:

- ITS-90

- ITS-SR5 (ITS-90 subrange 5 only)
- CVD (Callendar-Van Dusen)
- PT100
- R(T)
- T(R)
- YSI-400
- Ohms

Refer to the Report of Calibration for the probe for information on the conversion type that the probe was calibrated for.

If the selected conversion type requires coefficients or other settings, the Conversion Settings button is enabled. Click the Conversion Settings button to display the dialog for the selected conversion type. Otherwise, the Conversion Settings button is disabled. Please refer to Section 9, Conversion Settings Dialogs, for a detailed description of the Conversion Settings dialogs.



Note: *The conversion settings are not written to the thermometer readout until the Apply or OK button is clicked on this dialog!*

4.4.2.6.2 Calibration tab

Select the Calibration tab to view and/or modify channel calibration information.

Figure 63 1529 Chub-E4 PRT/thermistor channel settings - Calibration tab



Caution: Do not change these values unless this channel has been recalibrated or unless these values do not match the values located on the latest Report of Calibration for this channel! Changing these values voids the current calibration!

Select the date this channel was last calibrated using the drop-down calendar or by entering the date. This date should be located on the Report of Calibration for this channel.

Select the date this channel is due for recalibration using the drop-down calendar or by entering the date.

Enter the values for the calibration constants in the boxes provided. The calibration constants required by a PRT/Thermistor channel are: C0, C100, C400, C0K, C10K and C100K. The values of these constants should be located on the Report of Calibration for this channel.

4.4.2.7 1529 Chub-E4 thermocouple channel settings

The 1529 Chub-E4 Thermocouple Channel Settings dialog displays information about the selected thermocouple channel.

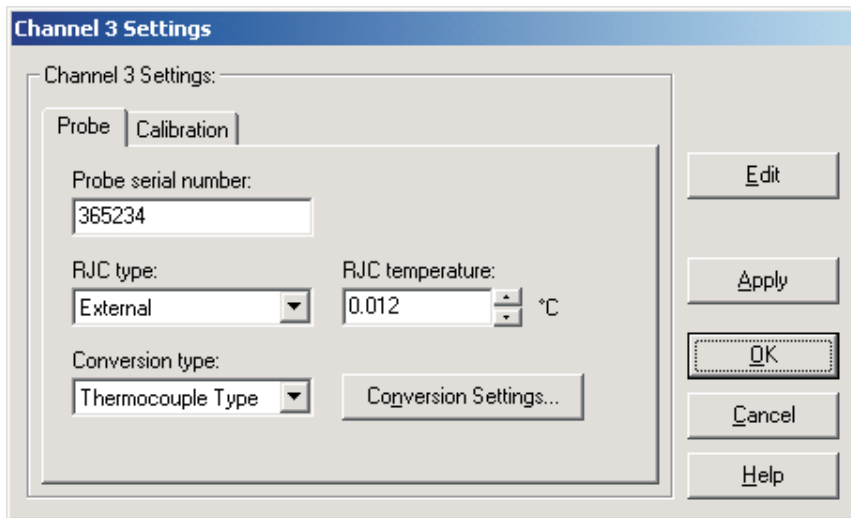


Figure 64 1529 Chub-E4 Thermocouple channel settings dialog

Some or all of the settings on this dialog may be passcode protected. If so, a message appears at the bottom of the Probe tab and/or Calibration tab and the Edit button is displayed. The protected controls are disabled. To edit the protected settings, click the Edit button and enter the current passcode for the thermometer readout. If the correct passcode is entered, the protected controls are enabled and the values may be changed. The controls remain enabled until this dialog is closed or until the current changes are applied by clicking the Apply button.

After modifying any of the settings on this dialog, click the Apply or OK button to write the new settings to the thermometer readout. If the Apply button is clicked, this dialog remains open. If the OK button is clicked, this dialog closes after writing the settings.

LogWare has a feature that causes all of the channel settings to be read back in from the thermometer readout after the Apply button is clicked and the new settings have been written to the thermometer readout. This feature can be useful to verify that the desired settings were properly applied to the thermometer readout. To enable or disable this feature, see Section 5.6.4, Options.

To close this dialog without writing the channel settings to the thermometer readout, click the Cancel button.

Click the Help button to open the appropriate help topic for this dialog.

4.4.2.7.1 Probe tab

Select the Probe tab to view and/or modify probe settings.

Figure 65 1529 Chub-E4 thermocouple channel settings - Probe tab

Enter the serial number of the probe attached to this channel in the Probe serial number box. The serial number can be any combination of numbers and/or letters up to 8 characters in length. The underscore “_” character can also be used in serial numbers.

Select the type of reference junction compensation to use on this channel using the RJC type drop-down list. Available choices are Internal and External. If External is chosen as the RJC type, the external RJC temperature must be entered in the RJC temperature box.

Enter the external RJC temperature value in the RJC temperature box. This box is disabled if the RJC type is set to Internal.



Note: The RJC temperature value must always be in °C!

Select the conversion type to use for the probe on this channel. The Model 1529 Chub-E4 Thermometer readout supports the following conversion types:

- Polynomial
- Thermocouple Type (B, E, J, K, N, R, S, T or AuPt)
- Volts

Refer to the Report of Calibration for the probe for information on the conversion type that the probe was calibrated for.

If the selected conversion type requires coefficients or other settings, the Conversion Settings button is enabled. Click the Conversion Settings button to display the dialog for the selected conversion type. Otherwise, the Conversion

Settings button is disabled. Please refer to Section 9, Conversion Settings Dialogs, for a detailed description of the Conversion Settings dialogs.



Note: *The conversion settings are not written to the thermometer readout until the Apply or OK button is clicked on this dialog!*

4.4.2.7.2 Calibration tab

Select the Calibration tab to view and/or modify channel calibration information.

Figure 66 1529 Chub-E4 thermocouple channel settings - Probe tab



Caution: *Do not change these values unless this channel has been recalibrated or unless these values do not match the values located on the latest Report of Calibration for this channel! Changing these values voids the current calibration!*

Select the date this channel was last calibrated using the drop-down calendar or by entering the date. This date should be located on the Report of Calibration for this channel.

Select the date this channel is due for recalibration using the drop-down calendar or by entering the date.

Enter the values for the calibration constants in the boxes provided. The calibration constants required by a Thermocouple channel are: C0, C100 and CRJ. The values of these constants should be located on the Report of Calibration for this channel.

4.4.3 1560 Black Stack settings dialog

The 1560 Black Stack Settings dialog displays information about the thermometer readout including model number, serial number, identification information,

and the total number of channels. It also allows many thermometer readout settings to be viewed and/or modified.

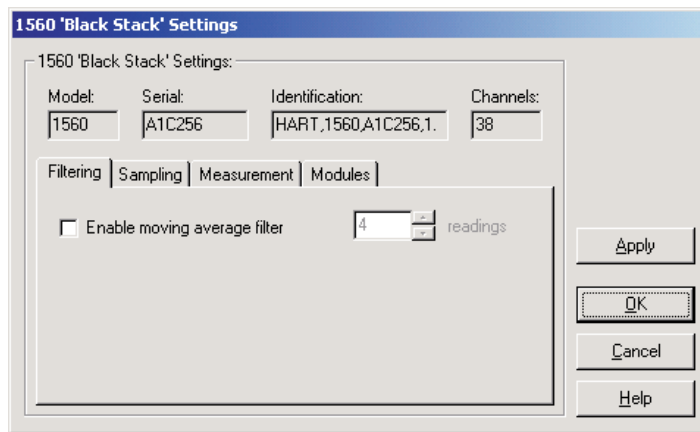


Figure 67 1560 Black Stack Settings dialog

The 1560 Black Stack must be connected to an available COM port and the communication settings must be properly configured so that this information can be read in from the thermometer readout. If communication with the thermometer readout is not successful, an error message is displayed and this dialog is blank. A progress dialog is displayed while the current thermometer readout settings are being read in.

The Model box indicates the model number of the thermometer readout. The Serial box indicates the serial number assigned to this thermometer readout. The Identification box indicates the manufacturer name, model number, serial number, and firmware version installed in this thermometer readout. The Channels box indicates the total number of input channels on this thermometer readout.

The filtering settings can be viewed and/or modified by selecting the Filtering tab. The sampling settings can be viewed and/or modified by selecting the Sampling tab. The measurement settings can be viewed and/or modified by selecting the Measurement tab. The Modules tab displays a list of modules currently attached to this thermometer readout and allows access to module settings dialogs. The settings on each of these tabs are described in detail in the following sections. For more details on these settings, please refer to the thermometer readout's User's Guide.

After modifying any of the settings on this dialog, click the Apply or OK button to write the new settings to the thermometer readout. If the Apply button is clicked, this dialog remains open. If the OK button is clicked, this dialog closes after writing the settings.

LogWare has a feature that causes all of the thermometer readout settings to be read back in from the thermometer readout after the Apply button is clicked and the new settings have been written to the thermometer readout. This feature can be useful to verify that the desired settings were properly applied to the thermometer readout. To enable or disable this feature, see Section 5.6.4, Options.

To close this dialog without writing the settings to the thermometer readout, click the Cancel button.

Click the Help button to open the appropriate help topic for this dialog.

4.4.3.1 Filtering tab

Select the Filtering tab to view and/or modify the filter settings.

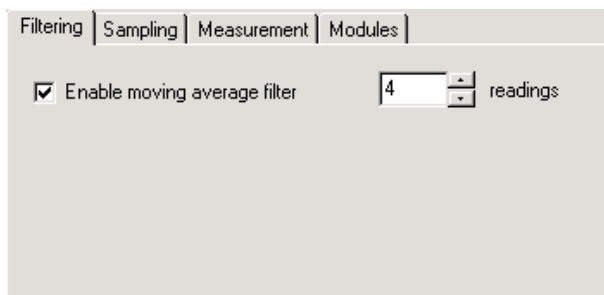


Figure 68 1560 Black Stack settings - Filtering tab

Check the Enable moving average filter box to enable the moving average filter. This filter is sometimes also called a sliding window filter. When this box is checked, the number of readings to average (size of window) must be entered in the box provided. Enabling this option causes the thermometer readout to display the average of the last n readings on a specific channel as the current reading for that channel. If this feature is disabled, the most recent measurement on each channel is displayed as the current reading. The moving average filter can be set to any value between 1 and 10.

4.4.3.2 Sampling tab

Select the Sampling tab to view and/or modify sampling settings.

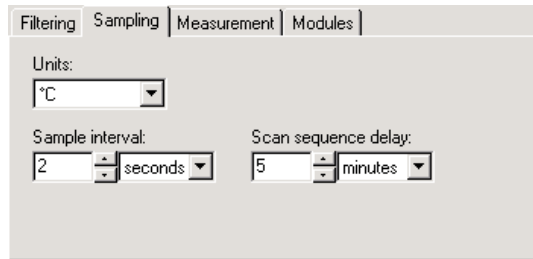


Figure 69 1560 Black Stack Settings – Sampling tab

Select the units to use when displaying readings in temperature using the Units drop-down list. Temperature readings can be displayed in either °C, °F or K. This setting applies to all channels that are configured to display readings in temperature.

Select the time interval between subsequent measurements using the Sample interval controls. The sample interval can be set to any value from 0 to 32,767 seconds (0 to 546 minutes or 0 to 9 hours).



Note: Setting the sample interval to 0 seconds or 1 second is equivalent to setting it to 2 seconds because the 1560 Black Stack cannot sample at a rate faster than once every 2 seconds. This is the default behavior of the 1560 Black Stack Thermometer readout.

Select the time interval between subsequent scans through a scan sequence using the Scan sequence delay controls. The scan sequence delay can be set to any value from 0 to 10,000 seconds (0 to 166 minutes or 0 to 2 hours).



Note: Setting the scan sequence delay to an interval of less than the setting of the sample interval multiplied by the number of channels in the scan sequence effectively disables the scan sequence delay feature.

4.4.3.3 Measurement tab

Select the Measurement tab to view and/or modify measurement related settings.

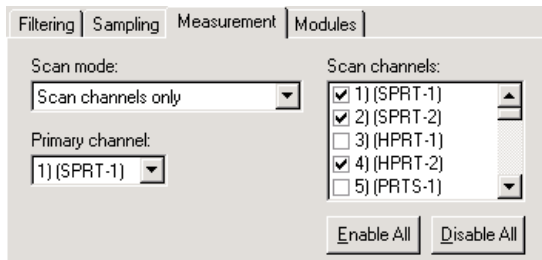


Figure 70 1560 Black Stack settings - Measurement tab

Select the scan mode for the 1560 Black Stack Thermometer readout using the Scan mode drop-down list. The scan mode can be set to one of the following options:

- Primary channel only
- Alternate scan/primary channels
- Scan channels only

If the Primary channel only option is selected, the thermometer readout will be set to measure only the channel selected in the Primary channel drop-down list below.

If the Alternate scan/primary channels option is selected, the thermometer readout will be set to alternate measurements on the primary channel and each of the enabled scan channels.

If the Scan channels only option is selected, the thermometer readout will be set to scan through each of the enabled channels in the Scan channels list.

Select the channel to set as the primary channel using the Primary channel drop-down list. This list indicates the channel number, module name and module input number for all channels on the thermometer readout. The primary channel setting is only enabled if the scan mode setting is set to the Primary channel only or the Alternate scan/primary channels option.

Select the channels to include in the thermometer readout's scan sequence. The Scan channels list indicates the channel number, module name and module input number for all channels on the thermometer readout. Each channel can be enabled or disabled by checking or unchecking the box next to the channel number independent of all other channels. The Scan channels list is only en-

abled if the scan mode setting is set to either the Alternate scan/primary channels or the Scan channels only option.

To include all channels in the scan sequence, click the Enable All button. This causes the box next to every channel to be checked.

To exclude all channels from the scan sequence, click the Disable All button. This causes the box next to every channel to be unchecked.

4.4.3.4 Modules tab

Select the Modules tab to view a list of modules currently attached to the 1560 Black Stack and to access the Module Settings dialogs to view and/or modify module specific settings.

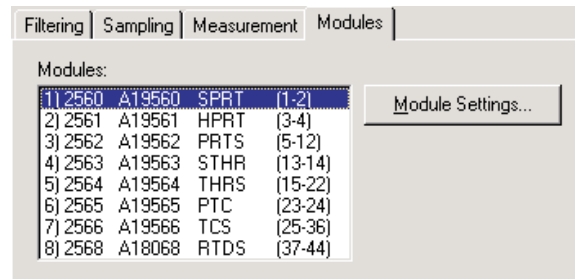


Figure 71 1560 Black Stack settings - Modules tab

Select the module to view and/or modify from the Modules list. This list indicates the module positions, model numbers, serial numbers, module names and the range of channels on each module.

Click the Module Settings button to display the appropriate Module Settings dialog. When this button is clicked, the settings for the selected module are read in from the thermometer readout. The Module Settings dialogs are described in detail in the following sections.

4.4.3.5 2560 SPRT Module Settings Dialog

The 2560 Module Settings dialog displays information about the 2560 module including model number, serial number, identification information and the

number of channels on this module. It also allows many module specific settings to be viewed and/or modified.

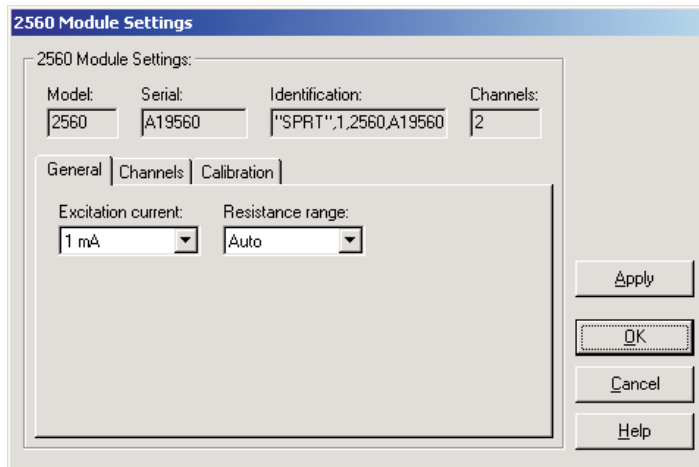


Figure 72 2560 SPRT Module Settings dialog

The Model box indicates the model number of this module. The Serial box indicates the serial number assigned to this module. The Identification box indicates the module name, number of devices, model number, serial number and firmware version installed in this module. The Channels box indicates the number of input channels on this module.

General module settings can be viewed and modified by selecting the General tab. The Channels tab displays a list of the channels on this module and allows access to channel settings dialogs. The module calibration information can be viewed and modified by selecting the Calibration tab.

After modifying any of the settings on this dialog, click the Apply or OK button to write the new module settings to the thermometer readout. If the Apply button is clicked, this dialog remains open. If the OK button is clicked, this dialog closes after writing the settings.

LogWare has a feature that causes all of the module settings to be read back in from the thermometer readout after the Apply button is clicked and the new settings have been written to the thermometer readout. This feature can be useful to verify that the desired settings were properly applied to the thermometer readout. To enable or disable this feature, see Section 5.6.4, Options.

To close this dialog without writing the module settings to the thermometer readout, click the Cancel button.

Click the Help button to open the appropriate help topic for this dialog.

4.4.3.5.1 General tab

Select the General tab to view and/or modify general module settings.

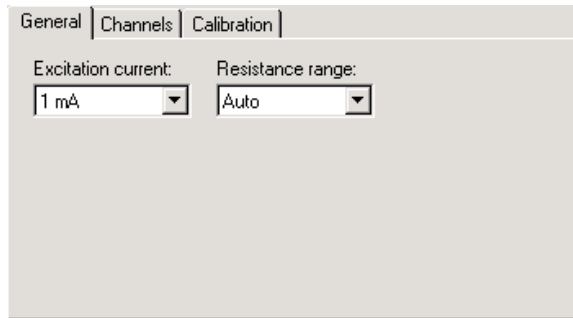


Figure 73 2560 SPRT Module Settings - General tab

Select the excitation current for this module using the Excitation current drop-down list. The 2560 module can use either a 1 mA or a 1.414 mA excitation current. This setting applies to all channels on this module.

Select the resistance range for this module using the Resistance range drop-down list. The 2560 module allows the selection of either Auto, High or Low. This setting applies to all channels on this module.

4.4.3.5.2 Channels tab

Select the Channels tab to view a list of channels on this module and to access the channel settings dialogs to view and/or modify channel specific settings.

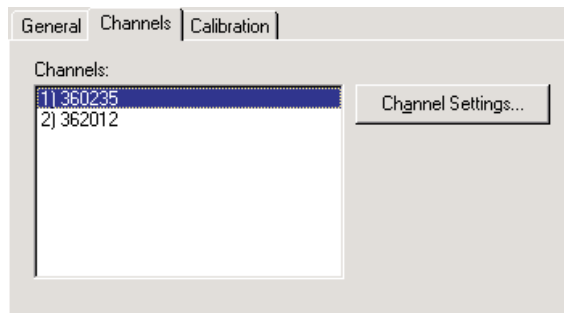


Figure 74 2560 SPRT Module Settings - Channels tab

Select the channel to view and/or modify from the Channels list. This list indicates the channel number (with respect to this module) and the serial number of the probe on each channel.

Click the Channel Settings button to display the appropriate Channel Settings dialog. When this button is clicked, the settings for the selected channel are read in from the thermometer readout. The Channel Settings dialog is described in detail in Section 4.4.3.14, 1560 Black Stack PRT/Thermistor Channel Settings Dialog.

4.4.3.5.3 Calibration tab

Select the Calibration tab to view and/or modify module calibration information.

Figure 75 2560 SPRT Module Settings - Calibration tab

The settings on the Calibration tab may be passcode protected. If they are, a message appears at the bottom of the Calibration tab and the Edit button is disabled. The other controls on this tab are disabled. To edit the calibration parameters, click the Edit button and enter the current passcode for the thermometer readout. If the correct passcode is entered, the calibration parameter controls are enabled and the values may be changed. The controls remain enabled until this dialog is closed or until the current changes are applied by clicking the Apply button.



Caution: Do not change these values unless this module has been recalibrated or unless these values do not match the values located on the latest Report of Calibration for this module! Changing these values voids the current calibration!

Select the date this module was last calibrated using the drop-down calendar or by entering the date. This date should be located on the Report of Calibration for this module.

Enter the values for the calibration constants in the boxes provided. The calibration constants required by the 2560 module are: 0 ADJ, 100 ADJ and 400 ADJ. The values of these constants should be located on the Report of Calibration for this module.

4.4.3.6 2561 HPRT Module Settings dialog

The 2561 Module Settings dialog displays information about the 2561 module including model number, serial number, identification information and the number of channels on this module. It also allows many module specific settings to be viewed and/or modified.

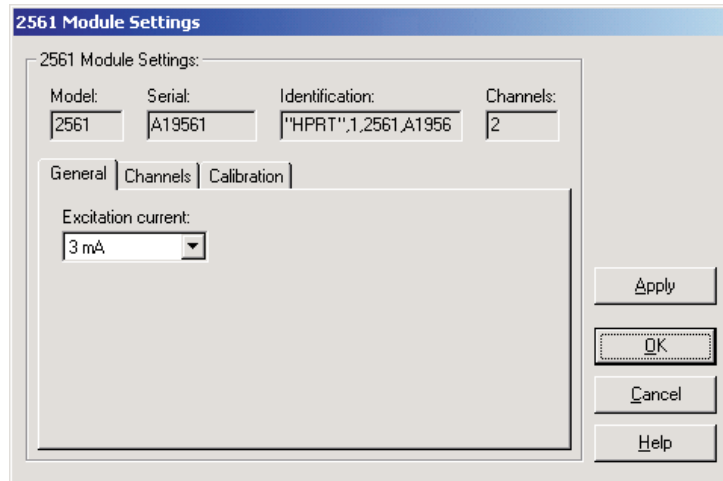


Figure 76 2561 HPRT Module Settings dialog

The Model box indicates the model number of this module. The Serial box indicates the serial number assigned to this module. The Identification box indicates the module name, number of devices, model number, serial number and firmware version installed in this module. The Channels box indicates the number of input channels on this module.

General module settings can be viewed and modified by selecting the General tab. The Channels tab displays a list of the channels on this module and allows access to channel settings dialogs. The module calibration information can be viewed and modified by selecting the Calibration tab.

After modifying any of the settings on this dialog, click the Apply or OK button to write the new module settings to the thermometer readout. If the Apply button is clicked, this dialog remains open. If the OK button is clicked, this dialog closes after writing the settings.

LogWare has a feature that causes all of the module settings to be read back in from the thermometer readout after the Apply button is clicked and the new settings have been written to the thermometer readout. This feature can be useful to verify that the desired settings were properly applied to the thermometer readout. To enable or disable this feature, see Section 5.6.4, Options.

To close this dialog without writing the module settings to the thermometer readout, click the Cancel button.

Click the Help button to open the appropriate help topic for this dialog.

4.4.3.6.1 General tab

Select the General tab to view and/or modify general module settings.

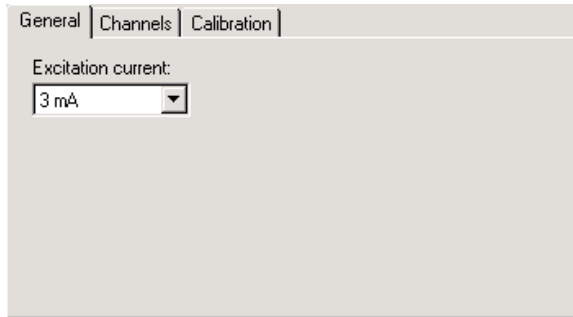


Figure 77 2561 HPRT Module Settings - General tab

Select the excitation current for this module using the Excitation current drop-down list. The 2561 module can use either a 3 mA or a 5 mA excitation current. This setting applies to all channels on this module.

4.4.3.6.2 Channels tab

Select the Channels tab to view a list of channels on this module and to access the channel settings dialogs to view and/or modify channel specific settings.

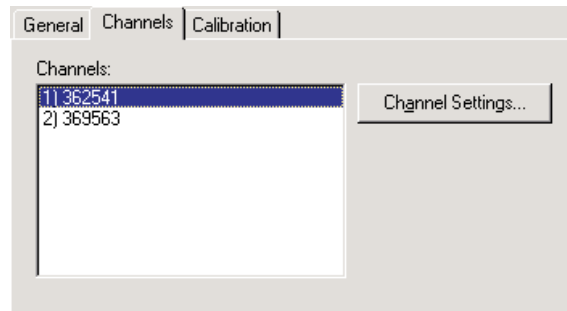


Figure 78 2561 HPRT Module Settings - Channels tab

Select the channel to view and/or modify from the Channels list. This list indicates the channel number (with respect to this module) and the serial number of the probe on each channel.

Click the Channel Settings button to display the appropriate Channel Settings dialog. When this button is clicked, the settings for the selected channel are read in from the thermometer readout. The Channel Settings dialog is described in detail in Section 4.4.3.14, 1560 Black Stack PRT/Thermistor Channel Settings Dialog.

4.4.3.6.3 Calibration tab

Select the Calibration tab to view and/or modify module calibration information.

Figure 79 2561 HPRT Module Settings - Calibration tab

The settings on the Calibration tab may be passcode protected. If they are, a message appears at the bottom of the Calibration tab and the Edit button is displayed. The other controls on this tab are disabled. To edit the calibration parameters, click the Edit button and enter the current passcode for the thermometer readout. If the correct passcode is entered, the calibration parameter controls are enabled and the values may be changed. The controls remain enabled until this dialog is closed or until the current changes are applied by clicking the Apply button.



Caution: Do not change these values unless this module has been recalibrated or unless these values do not match the values located on the latest Report of Calibration for this module! Changing these values voids the current calibration!

Select the date this module was last calibrated using the drop-down calendar or by entering the date. This date should be located on the Report of Calibration for this module.

Enter the values for the calibration constants in the boxes provided. The calibration constants required by the 2561 module are: 0 ADJ and 10 ADJ. The

values of these constants should be located on the Report of Calibration for this module.

4.4.3.7 2562 PRTS Module Settings dialog

The 2562 Module Settings dialog displays information about the 2562 module including model number, serial number, identification information and the number of channels on this module. It also allows many module specific settings to be viewed and/or modified.

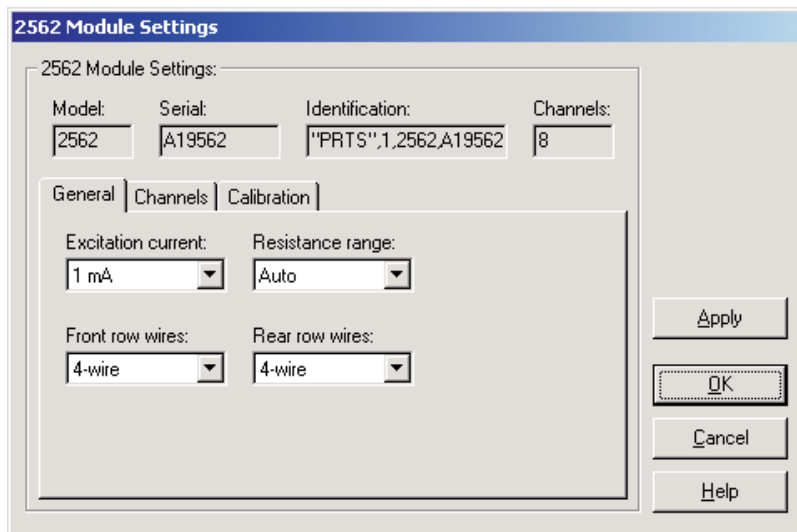


Figure 80 2562 'PRTS' Module Settings dialog

The Model box indicates the model number of this module. The Serial box indicates the serial number assigned to this module. The Identification box indicates the module name, number of devices, model number, serial number, and firmware version installed in this module. The Channels box indicates the number of input channels on this module.

General module settings can be viewed and modified by selecting the General tab. The Channels tab displays a list of the channels on this module and allows access to channel settings dialogs. The module calibration information can be viewed and modified by selecting the Calibration tab.

After modifying any of the settings on this dialog, click the Apply or OK button to write the new module settings to the thermometer readout. If the Apply button is clicked, this dialog remains open. If the OK button is clicked, this dialog closes after writing the settings.

LogWare has a feature that causes all of the module settings to be read back in from the thermometer readout after the Apply button is clicked and the new set-


tings have been written to the thermometer readout. This feature can be useful to verify that the desired settings were properly applied to the thermometer readout. To enable or disable this feature, see Section 5.6.4, Options.

To close this dialog without writing the module settings to the thermometer readout, click the Cancel button.

Click the Help button to open the appropriate help topic for this dialog.

4.4.3.7.1 General tab

Select the General tab to view and/or modify general module settings.

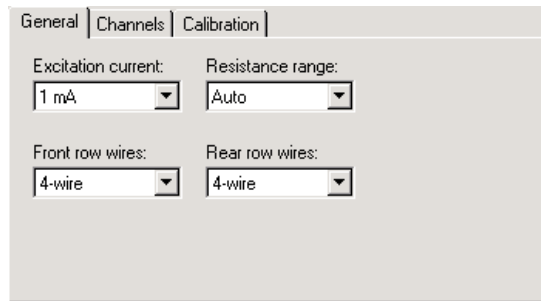

The screenshot shows a dialog box titled "2562 PRTS Module Settings" with three tabs: "General", "Channels", and "Calibration". The "General" tab is active. It contains four settings, each with a drop-down menu:

- Excitation current: 1 mA
- Resistance range: Auto
- Front row wires: 4-wire
- Rear row wires: 4-wire

Figure 81 2562 PRTS Module Settings - General tab

Select the excitation current for this module using the Excitation current drop-down list. The 2562 module can use either a 1 mA or a 1.414 mA excitation current. This setting applies to all channels on this module.

Select the resistance range for this module using the Resistance range drop-down list. The 2562 module allows the selection of either Auto, High or Low. This setting applies to all channels on this module.

Select the front row wires option for this module using the Front row wires drop-down list. The 2562 module allows the selection of either 4-wire or 3-wire. This setting applies only to the four channels on the front row of this module.

Select the rear row wires option for this module using the Rear row wires drop-down list. The 2562 module allows the selection of either 4-wire or 3-wire. This setting applies only to the four channels on the rear row of this module.

4.4.3.7.2 Channels tab

Select the Channels tab to view a list of channels on this module and to access the channel settings dialogs to view and/or modify channel specific settings.

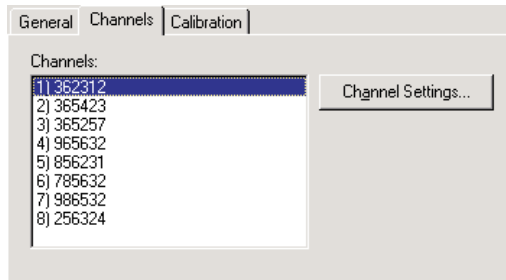


Figure 82 2562 'PRTS' Module Settings - Channels tab

Select the channel to view and/or modify from the Channels list. This list indicates the channel number (with respect to this module) and the serial number of the probe on each channel.

Click the Channel Settings button to display the appropriate Channel Settings dialog. When this button is clicked, the settings for the selected channel are read in from the thermometer readout. The Channel Settings dialog is described in detail in Section 4.4.3.14, 1560 Black Stack PRT/Thermistor Channel Settings Dialog.

4.4.3.7.3 Calibration tab

Select the Calibration tab to view and/or modify module calibration information.

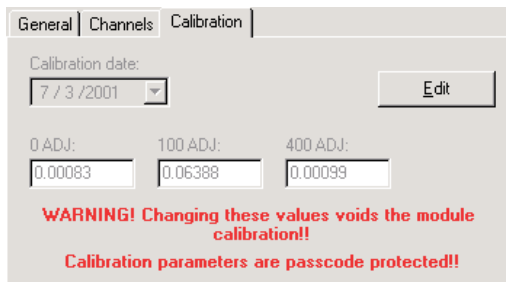


Figure 83 2562 'PRTS' Module Settings - Calibration tab

The settings on the Calibration tab may be passcode protected. If they are, a message appears at the bottom of the Calibration tab and the Edit button is dis-

played. The other controls on this tab are disabled. To edit the calibration parameters, click the Edit button and enter the current passcode for the thermometer readout. If the correct passcode is entered, the calibration parameter controls are enabled and the values may be changed. The controls remain enabled until this dialog is closed or until the current changes are applied by clicking the Apply button.



Caution: Do not change these values unless this module has been recalibrated or unless these values do not match the values located on the latest Report of Calibration for this module! Changing these values voids the current calibration!

Select the date this module was last calibrated using the drop-down calendar or by entering the date. This date should be located on the Report of Calibration for this module.

Enter the values for the calibration constants in the boxes provided. The calibration constants required by the 2562 module are: 0 ADJ, 100 ADJ and 400 ADJ. The values of these constants should be located on the Report of Calibration for this module.

4.4.3.8 2563 STHR Module Settings dialog

The 2563 Module Settings dialog displays information about the 2563 module including model number, serial number, identification information and the number of channels on this module. It also allows many module specific settings to be viewed and/or modified.

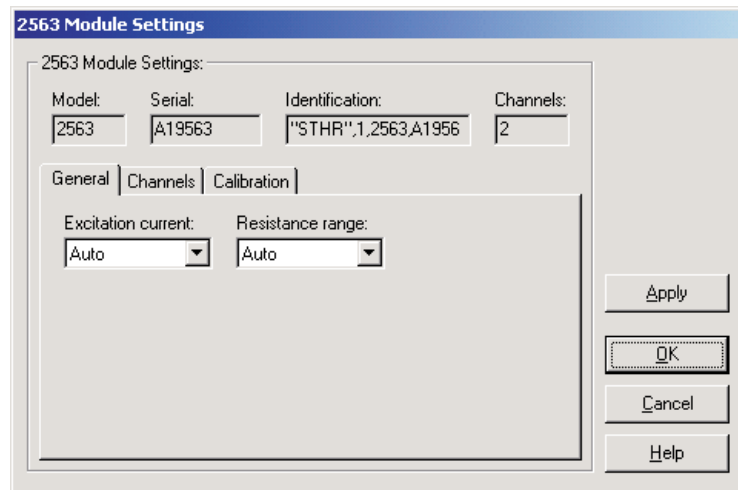


Figure 84 2563 STHR Module Settings dialog

The Model box indicates the model number of this module. The Serial box indicates the serial number assigned to this module. The Identification box indicates the module name, number of devices, model number, serial number and firmware version installed in this module. The Channels box indicates the number of input channels on this module.

General module settings can be viewed and modified by selecting the General tab. The Channels tab displays a list of the channels on this module and allows access to channel settings dialogs. The module calibration information can be viewed and modified by selecting the Calibration tab.

After modifying any of the settings on this dialog, click the Apply or OK button to write the new module settings to the thermometer readout. If the Apply button is clicked, this dialog remains open. If the OK button is clicked, this dialog closes after writing the settings.

LogWare has a feature that causes all of the module settings to be read back in from the thermometer readout after the Apply button is clicked and the new settings have been written to the thermometer readout. This feature can be useful to verify that the desired settings were properly applied to the thermometer readout. To enable or disable this feature, see Section 5.6.4, Options.

To close this dialog without writing the module settings to the thermometer readout, click the Cancel button.

Click the Help button to open the appropriate help topic for this dialog.

4.4.3.8.1 General tab

Select the General tab to view and/or modify general module settings.

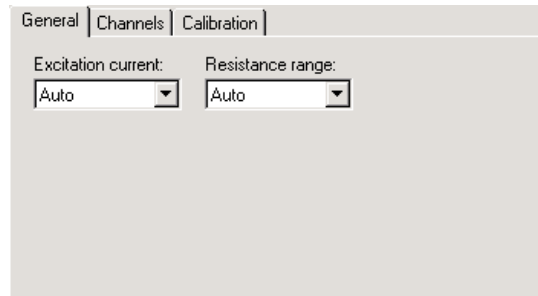


Figure 85 2563 STHR Module Settings - General tab

Select the excitation current for this module using the Excitation current drop-down list. The 2563 module can use either a 2 μA or a 10 μA excitation current, or Auto. This setting applies to all channels on this module.

Select the resistance range for this module using the Resistance range drop-down list. The 2563 module allows the selection of either Auto, High or Low. This setting applies to all channels on this module.

4.4.3.8.2 Channels tab

Select the Channels tab to view a list of channels on this module and to access the channel settings dialogs to view and/or modify channel specific settings.

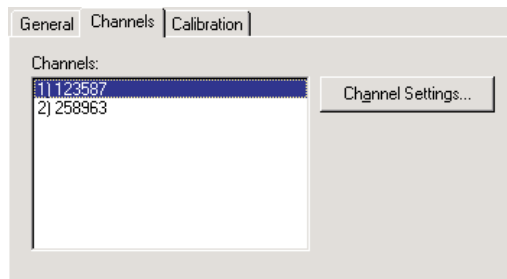


Figure 86 2563 STHR Module Settings - Channels tab

Select the channel to view and/or modify from the Channels list. This list indicates the channel number (with respect to this module) and the serial number of the probe on each channel.

Click the Channel Settings button to display the appropriate channel settings dialog. When this button is clicked, the settings for the selected channel are read in from the thermometer readout. The Channel Settings dialog is described in detail in Section 4.4.3.14, 1560 Black Stack PRT/Thermistor Channel Settings Dialog.

4.4.3.8.3 Calibration tab

Select the Calibration tab to view and/or modify module calibration information.

Figure 87 2563 STHR Module Settings - Calibration tab

The settings on the Calibration tab may be passcode protected. If they are, a message appears at the bottom of the Calibration tab and the Edit button is disabled. The other controls on this tab are disabled. To edit the calibration parameters, click the Edit button and enter the current passcode for the thermometer readout. If the correct passcode is entered, the calibration parameter controls are enabled and the values may be changed. The controls remain enabled until this dialog is closed or until the current changes are applied by clicking the Apply button.



Caution: Do not change these values unless this module has been recalibrated or unless these values do not match the values located on the latest Report of Calibration for this module! Changing these values voids the current calibration!

Select the date this module was last calibrated using the drop-down calendar or by entering the date. This date should be located on the Report of Calibration for this module.

Enter the values for the calibration constants in the boxes provided. The calibration constants required by the 2563 module are: 0 ADJ, 10K ADJ and 100K ADJ. The values of these constants should be located on the Report of Calibration for this module.

4.4.3.9 2564 THRS Module Settings dialog

The 2564 Module Settings dialog displays information about the 2564 module including model number, serial number, identification information and the number of channels on this module. It also allows many module specific settings to be viewed and/or modified.

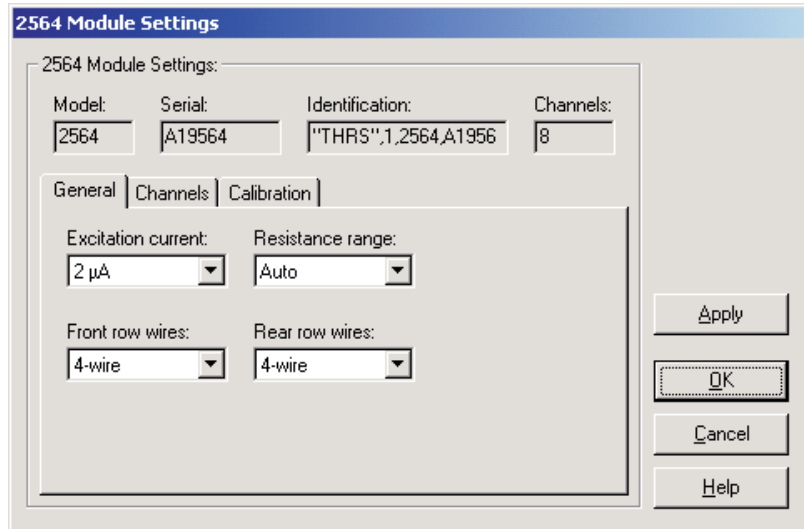


Figure 88 2564 THRS Module Settings dialog

The Model box indicates the model number of this module. The Serial box indicates the serial number assigned to this module. The Identification box indicates the module name, number of devices, model number, serial number and firmware version installed in this module. The Channels box indicates the number of input channels on this module.

General module settings can be viewed and modified by selecting the General tab. The Channels tab displays a list of the channels on this module and allows access to channel settings dialogs. The module calibration information can be viewed and modified by selecting the Calibration tab.

After modifying any of the settings on this dialog, click the Apply or OK button to write the new module settings to the thermometer readout. If the Apply button is clicked, this dialog remains open. If the OK button is clicked, this dialog closes after writing the settings.

LogWare has a feature that causes all of the module settings to be read back in from the thermometer readout after the Apply button is clicked and the new settings have been written to the thermometer readout. This feature can be useful to verify that the desired settings were properly applied to the thermometer readout. To enable or disable this feature, see Section 5.6.4, Options.

To close this dialog without writing the module settings to the thermometer readout, click the Cancel button.

Click the Help button to open the appropriate help topic for this dialog.

4.4.3.9.1 General tab

Select the General tab to view and/or modify general module settings.

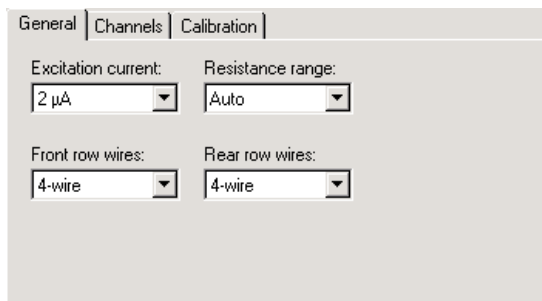


Figure 89 2564 THRS Module Settings - General tab

Select the excitation current for this module using the Excitation current drop-down list. The 2564 module can use either a 2 µA or a 10 µA excitation current, or Auto. This setting applies to all channels on this module.

Select the resistance range for this module using the Resistance range drop-down list. The 2564 module allows the selection of either Auto, High or Low. This setting applies to all channels on this module.

Select the front row wires option for this module using the Front row wires drop-down list. The 2564 module allows the selection of either 4-wire or 3-wire. This setting applies only to the four channels on the front row of this module.

Select the rear row wires option for this module using the Rear row wires drop-down list. The 2564 module allows the selection of either 4-wire or 3-wire. This setting applies only to the four channels on the rear row of this module.

4.4.3.9.2 Channels tab

Select the Channels tab to view a list of channels on this module and to access the channel settings dialogs to view and/or modify channel specific settings.

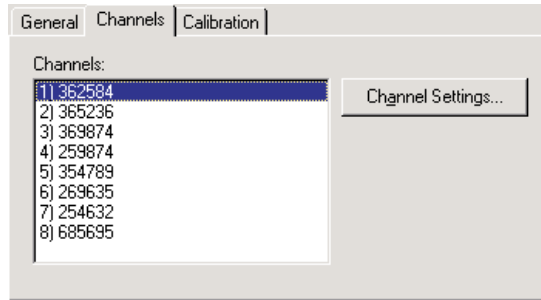


Figure 90 2564 THRS Module Settings - Channels tab

Select the channel to view and/or modify from the Channels list. This list indicates the channel number (with respect to this module) and the serial number of the probe on each channel.

Click the Channel Settings button to display the appropriate channel settings dialog. When this button is clicked, the settings for the selected channel are read in from the thermometer readout. The Channel Settings dialog is described in detail in Section 4.4.3.14, 1560 Black Stack PRT/Thermistor Channel Settings Dialog.

4.4.3.9.3 Calibration tab

Select the Calibration tab to view and/or modify module calibration information.

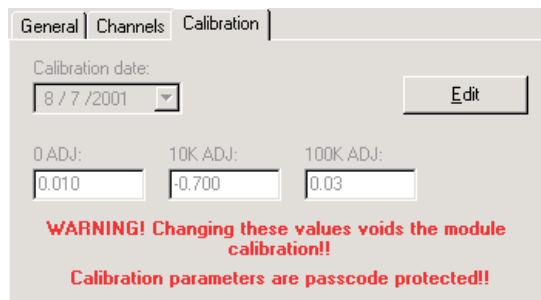


Figure 91 2564 THRS Module Settings - Calibration tab

The settings on the Calibration tab may be passcode protected. If they are, a message appears at the bottom of the Calibration tab and the Edit button is displayed. The other controls on this tab are disabled. To edit the calibration parameters, click the Edit button and enter the current passcode for the thermometer readout. If the correct passcode is entered, the calibration parameter controls are enabled and the values may be changed. The controls remain enabled until this dialog is closed or until the current changes are applied by clicking the Apply button.



Caution: *Do not change these values unless this module has been recalibrated or unless these values do not match the values located on the latest Report of Calibration for this module! Changing these values voids the current calibration!*

Select the date this module was last calibrated using the drop-down calendar or by entering the date. This date should be located on the Report of Calibration for this module.

Enter the values for the calibration constants in the boxes provided. The calibration constants required by the 2564 module are: 0 ADJ, 10K ADJ and 100K ADJ. The values of these constants should be located on the Report of Calibration for this module.

4.4.3.10 2565 PTC Module Settings dialog

The 2565 Module Settings dialog displays information about the 2565 module including model number, serial number, identification information and the

number of channels on this module. It also allows many module specific settings to be viewed and/or modified.

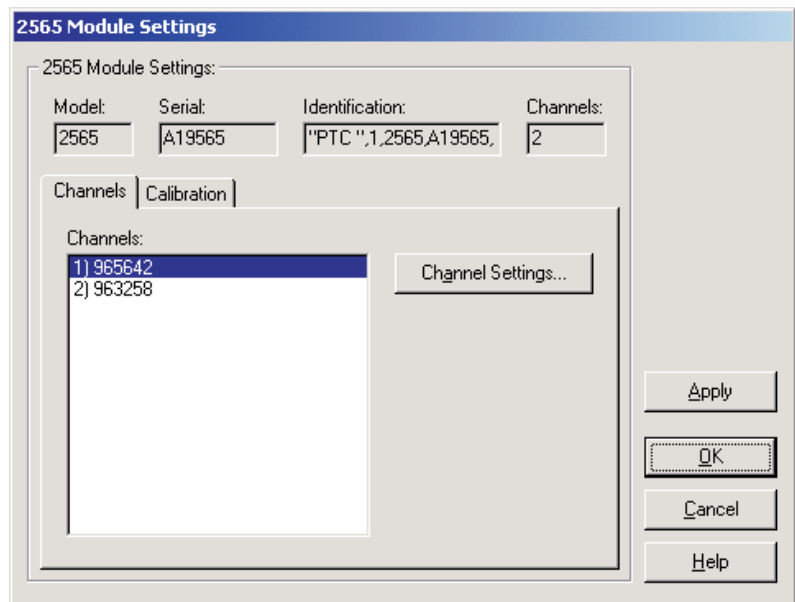


Figure 92 2565 PTC Module Settings dialog

The Model box indicates the model number of this module. The Serial box indicates the serial number assigned to this module. The Identification box indicates the module name, number of devices, model number, serial number and firmware version installed in this module. The Channels box indicates the number of input channels on this module.

The Channels tab displays a list of the channels on this module and allows access to channel settings dialogs. The module calibration information can be viewed and modified by selecting the Calibration tab.

After modifying any of the settings on this dialog, click the Apply or OK button to write the new module settings to the thermometer readout. If the Apply button is clicked, this dialog remains open. If the OK button is clicked, this dialog closes after writing the settings.

LogWare has a feature that causes all of the module settings to be read back in from the thermometer readout after the Apply button is clicked and the new settings have been written to the thermometer readout. This feature can be useful to verify that the desired settings were properly applied to the thermometer readout. To enable or disable this feature, see Section 5.6.4, Options.

To close this dialog without writing the module settings to the thermometer readout, click the Cancel button.

Click the Help button to open the appropriate help topic for this dialog.

4.4.3.10.1 Channels tab

Select the Channels tab to view a list of channels on this module and to access the channel settings dialogs to view and/or modify channel specific settings.

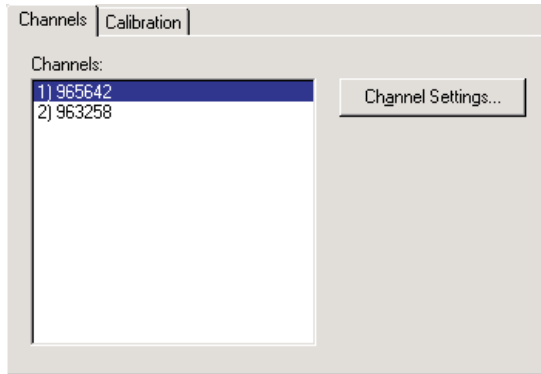


Figure 93 2565 PTC Module Settings - Channels tab

Select the channel to view and/or modify from the Channels list. This list indicates the channel number (with respect to this module) and the serial number of the probe on each channel.

Click the Channel Settings button to display the appropriate channel settings dialog. When this button is clicked, the settings for the selected channel are read in from the thermometer readout. The Channel Settings dialog is described in detail in Section 4.4.3.15, 1560 Black Stack Thermocouple Channel Settings Dialog.

4.4.3.10.2 Calibration tab

Select the Calibration tab to view and/or modify module calibration information.

Figure 94 2565 PTC Module Settings - Calibration tab

The settings on the Calibration tab may be passcode protected. If they are, a message appears at the bottom of the Calibration tab and the Edit button is displayed. The other controls on this tab are disabled. To edit the calibration parameters, click the Edit button and enter the current passcode for the thermometer readout. If the correct passcode is entered, the calibration parameter controls are enabled and the values may be changed. The controls remain enabled until this dialog is closed or until the current changes are applied by clicking the Apply button.



Caution: Do not change these values unless this module has been recalibrated or unless these values do not match the values located on the latest Report of Calibration for this module! Changing these values voids the current calibration!

The AMP GA constant should always be set to 1 and cannot be changed.

Select the date this module was last calibrated using the drop-down calendar or by entering the date. This date should be located on the Report of Calibration for this module.

Enter the values for the calibration constants in the boxes provided. The calibration constants required by the 2565 module are: OS 1, GA 1 and CJ 1 (for channel 1) and OS 2, GA 2 and CJ 2 (for channel 2). The values of these constants should be located on the Report of Calibration for this module.

4.4.3.11 2566 TCS Module Settings dialog

The 2566 Module Settings dialog displays information about the 2566 module including model number, serial number, identification information and the number of channels on this module. It also allows many module specific settings to be viewed and/or modified.

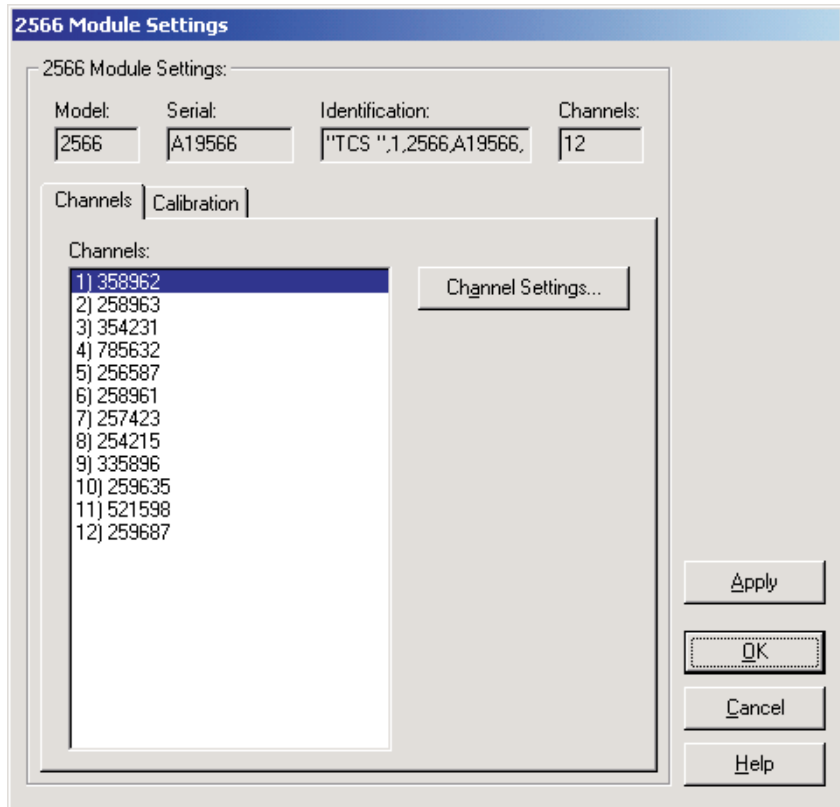


Figure 95 2566 TCS Module Settings dialog

The Model box indicates the model number of this module. The Serial box indicates the serial number assigned to this module. The Identification box indicates the module name, number of devices, model number, serial number and firmware version installed in this module. The Channels box indicates the number of input channels on this module.

The Channels tab displays a list of the channels on this module and allows access to channel settings dialogs. The module calibration information can be viewed and/or modified by selecting the Calibration tab.

After modifying any of the settings on this dialog, click the Apply or OK button to write the new module settings to the thermometer readout. If the Apply

button is clicked, this dialog remains open. If the OK button is clicked, this dialog closes after writing the settings.

LogWare has a feature that causes all of the module settings to be read back in from the thermometer readout after the Apply button is clicked and the new settings have been written to the thermometer readout. This feature can be useful to verify that the desired settings were properly applied to the thermometer readout. To enable or disable this feature, see Section 5.6.4, Options.

To close this dialog without writing the module settings to the thermometer readout, click the Cancel button.

Click the Help button to open the appropriate help topic for this dialog.

4.4.3.11.1 Channels tab

Select the Channels tab to view a list of channels on this module and to access the channel settings dialogs to view and/or modify channel specific settings.

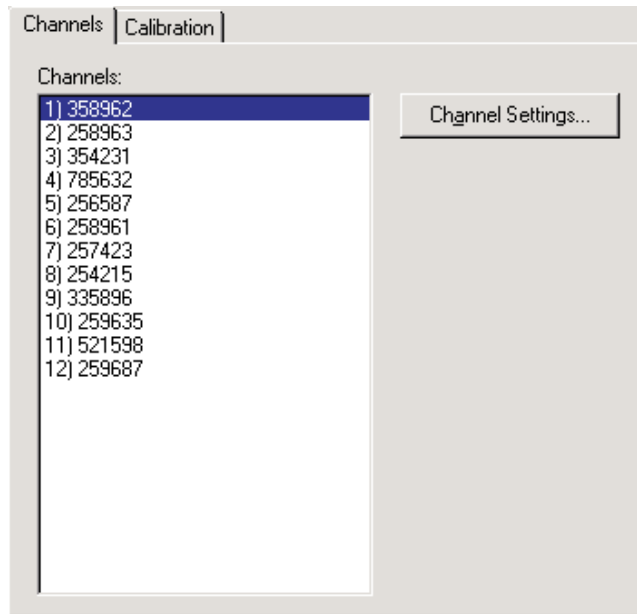


Figure 96 2566 TCS Module Settings - Channels tab

Select the channel to view and/or modify from the Channels list. This list indicates the channel number (with respect to this module) and the serial number of the probe on each channel.

Click the Channel Settings button to display the appropriate channel settings dialog. When this button is clicked, the settings for the selected channel are read

in from the thermometer readout. The Channel Settings dialog is described in detail in Section 4.4.3.15, 1560 Black Stack Thermocouple Channel Settings Dialog.

4.4.3.11.2 Calibration tab

Select the Calibration tab to view and/or modify module calibration information.

Figure 97 2566 TCS Module Settings - Calibration tab

The settings on the Calibration tab may be passcode protected. If they are, a message appears at the bottom of the Calibration tab and the Edit button is disabled. The other controls on this tab are disabled. To edit the calibration parameters, click the Edit button and enter the current passcode for the thermometer readout. If the correct passcode is entered, the calibration parameter controls are enabled and the values may be changed. The controls remain enabled until this dialog is closed or until the current changes are applied by clicking the Apply button.



Caution: Do not change these values unless this module has been recalibrated or unless these values do not match the values located on the latest Report of Calibration for this module! Changing these values voids the current calibration!

The AMP GA constant should always be set to 1 and cannot be changed.

Select the date this module was last calibrated using the drop-down calendar or by entering the date. This date should be located on the Report of Calibration for this module.

The calibration constants for this module are divided onto two tabs: Inputs 1 to 6 and Inputs 7 to 12.

4.4.3.11.2.1 Inputs 1 to 6 tab

Select the Inputs 1 to 6 tab to view and/or modify the calibration constants for inputs (channels) 1 to 6 (front row) of this module.

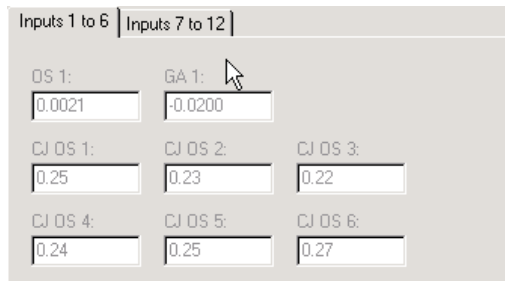


Figure 98 2566 TCS Module Settings - Inputs 1 to 6 tab

Enter the values for the calibration constants for inputs 1 to 6 in the boxes provided. The calibration constants required by the 2566 module are: OS 1, GA 1, CJ OS 1, CJ OS 2, CJ OS 3, CJ OS 4, CJ OS 5 and CJ OS 6. The values of these constants should be located on the Report of Calibration for this module.

4.4.3.11.2 Inputs 7 to 12 tab

Select the Inputs 7 to 12 tab to view and/or modify the calibration constants for inputs (channels) 7 to 12 (rear row) of this module.

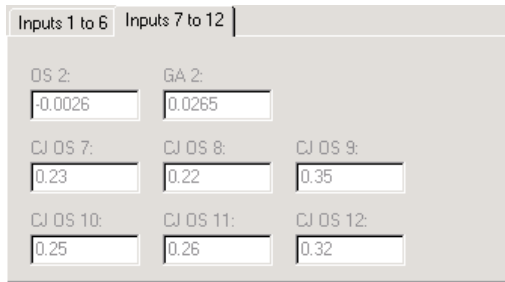


Figure 99 2566 TCS Module Settings - Inputs 7 to 12 tab

Enter the values for the calibration constants for inputs 7 to 12 in the boxes provided. The calibration constants required by the 2566 module are: OS 2, GA 2, CJ OS 7, CJ OS 8, CJ OS 9, CJ OS 10, CJ OS 11 and CJ OS 12. The values of these constants should be located on the Report of Calibration for this module.

4.4.3.12 2567 RTD Module Settings dialog

The 2567 Module Settings dialog displays information about the 2567 module including model number, serial number, identification information and the



number of channels on this module. It also allows many module specific settings to be viewed and/or modified.

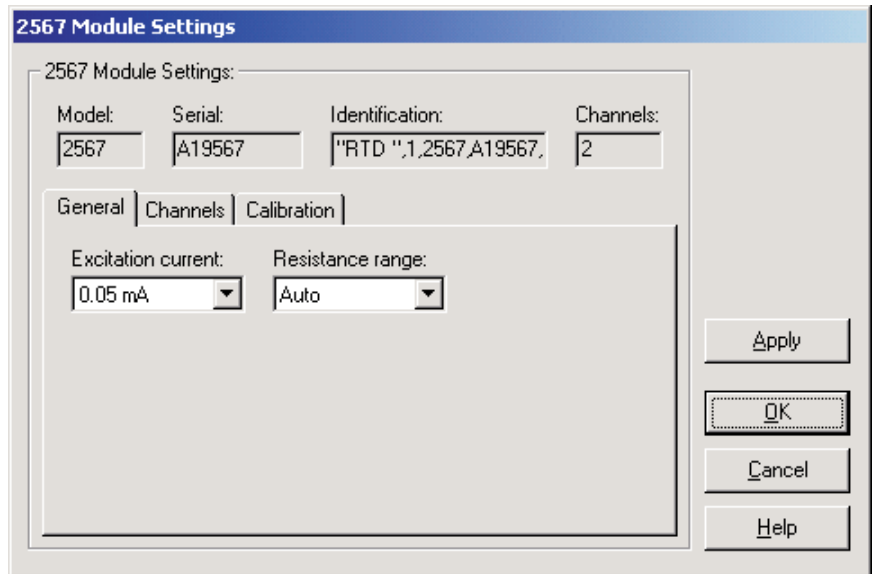


Figure 100 2567 RTD Module Settings dialog

The Model box indicates the model number of this module. The Serial box indicates the serial number assigned to this module. The Identification box indicates the module name, number of devices, model number, serial number and firmware version installed in this module. The Channels box indicates the number of input channels on this module.

General module settings can be viewed and/or modified by selecting the General tab. The Channels tab displays a list of the channels on this module and allows access to channel settings dialogs. The module calibration information can be viewed and/or modified by selecting the Calibration tab.

After modifying any of the settings on this dialog, click the Apply or OK button to write the new module settings to the thermometer readout. If the Apply button is clicked, this dialog remains open. If the OK button is clicked, this dialog closes after writing the settings.

LogWare has a feature that causes all of the module settings to be read back in from the thermometer readout after the Apply button is clicked and the new settings have been written to the thermometer readout. This feature can be useful to verify that the desired settings were properly applied to the thermometer readout. To enable or disable this feature, see Section 5.6.4, Options.

To close this dialog without writing the module settings to the thermometer readout, click the Cancel button.

Click the Help button to open the appropriate help topic for this dialog.

4.4.3.12.1 General tab

Select the General tab to view and/or modify general module settings.

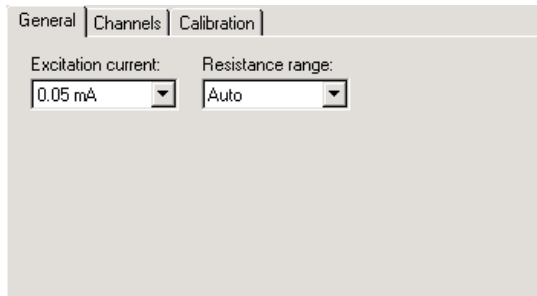


Figure 101 2567 RTD Module Settings - General tab

Select the excitation current for this module using the Excitation current drop-down list. The 2567 module can use either a 0.05 mA or a 0.1 mA excitation current. This setting applies to all channels on this module.

Select the resistance range for this module using the Resistance range drop-down list. The 2567 module allows the selection of either Auto, High or Low. This setting applies to all channels on this module.

4.4.3.12.2 Channels tab

Select the Channels tab to view a list of channels on this module and to access the channel settings dialogs to view and/or modify channel specific settings.

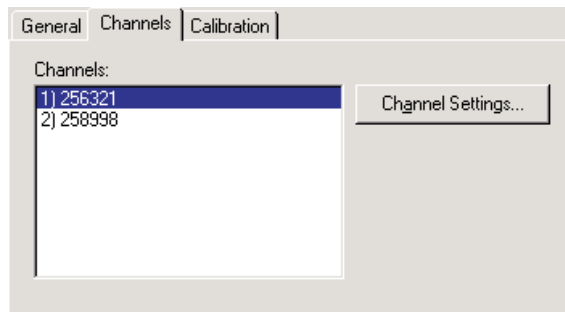


Figure 102 2567 RTD Module Settings - Channels tab

Select the channel to view and/or modify from the Channels list. This list indicates the channel number (with respect to this module) and the serial number of the probe on each channel.

Click the Channel Settings button to display the appropriate channel settings dialog. When this button is clicked, the settings for the selected channel are read in from the thermometer readout. The Channel Settings dialog is described in detail in Section 4.4.3.14, 1560 Black Stack PRT/Thermistor Channel Settings Dialog.

4.4.3.12.3 Calibration tab

Select the Calibration tab to view and/or modify module calibration information.

Figure 103 2567 RTD Module Settings - Calibration tab

The settings on the Calibration tab may be passcode protected. If they are, a message appears at the bottom of the Calibration tab and the Edit button is displayed. The other controls on this tab are disabled. To edit the calibration parameters, click the Edit button and enter the current passcode for the thermometer readout. If the correct passcode is entered, the calibration parameter controls are enabled and the values may be changed. The controls remain enabled until this dialog is closed or until the current changes are applied by clicking the Apply button.



Caution: Do not change these values unless this module has been recalibrated or unless these values do not match the values located on the latest Report of Calibration for this module! Changing these values voids the current calibration!

Select the date this module was last calibrated using the drop-down calendar or by entering the date. This date should be located on the Report of Calibration for this module.

Enter the values for the calibration constants in the boxes provided. The calibration constants required by the 2567 module are: 0 ADJ, 1K ADJ and 4K

ADJ. The values of these constants should be located on the Report of Calibration for this module.

4.4.3.13 2568 RTDS Module Settings

The 2568 Module Settings dialog displays information about the 2568 module including model number, serial number, identification information and the number of channels on this module. It also allows many module specific settings to be viewed and/or modified.

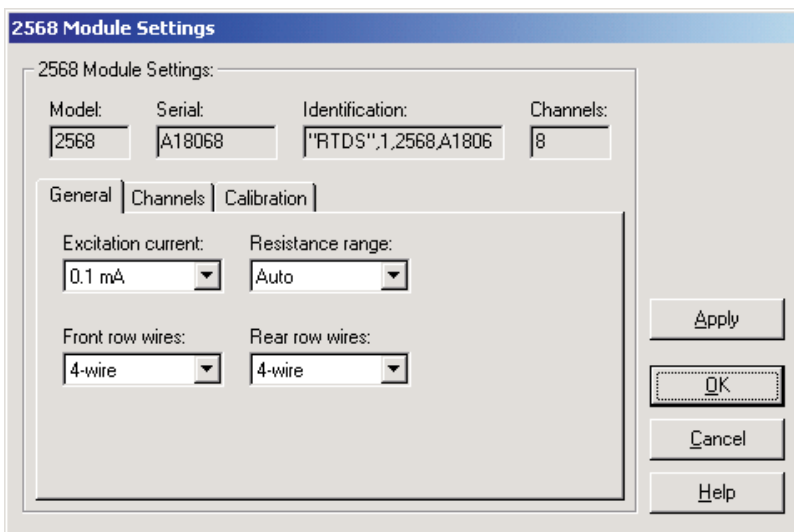


Figure 104 2568 RTDS Module Settings dialog

The Model box indicates the model number of this module. The Serial box indicates the serial number assigned to this module. The Identification box indicates the module name, number of devices, model number, serial number, and firmware version installed in this module. The Channels box indicates the number of input channels on this module.

General module settings can be viewed and/or modified by selecting the General tab. The Channels tab displays a list of the channels on this module and allows access to channel settings dialogs. The module calibration information can be viewed and/or modified by selecting the Calibration tab.

After modifying any of the settings on this dialog, click the Apply or OK button to write the new module settings to the thermometer readout. If the Apply button is clicked, this dialog remains open. If the OK button is clicked, this dialog closes after writing the settings.

LogWare has a feature that causes all of the module settings to be read back in from the thermometer readout after the Apply button is clicked and the new settings have been written to the thermometer readout. This feature can be useful

to verify that the desired settings were properly applied to the thermometer readout. To enable or disable this feature, see Section 5.6.4, Options.

To close this dialog without writing the module settings to the thermometer readout, click the Cancel button.

Click the Help button to open the appropriate help topic for this dialog.

4.4.3.13.1 General tab

Select the General tab to view and/or modify general module settings.

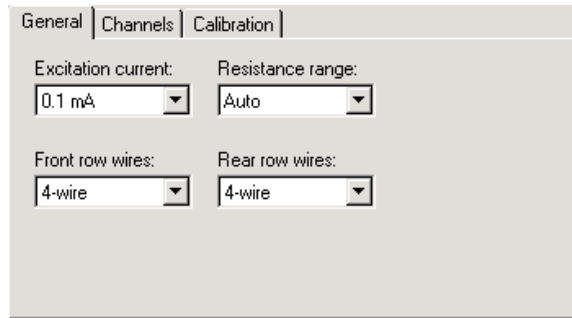


Figure 105 2568 RTDS Module Settings - General tab

Select the excitation current for this module using the Excitation current drop-down list. The 2568 module can use either a 0.05 mA or a 0.1 mA excitation current. This setting applies to all channels on this module.

Select the resistance range for this module using the Resistance range drop-down list. The 2568 module allows the selection of either Auto, High or Low. This setting applies to all channels on this module.

Select the front row wires option for this module using the Front row wires drop-down list. The 2568 module allows the selection of either 4-wire or 3-wire. This setting applies only to the four channels on the front row of this module.

Select the rear row wires option for this module using the Rear row wires drop-down list. The 2568 module allows the selection of either 4-wire or 3-wire. This setting applies only to the four channels on the rear row of this module.

4.4.3.13.2 Channels tab

Select the Channels tab to view a list of channels on this module and to access the channel settings dialogs to view and/or modify channel specific settings.

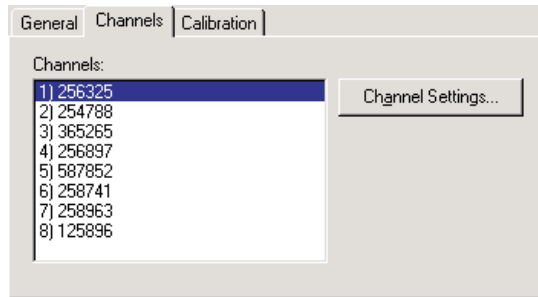


Figure 106 2568 RTDS Module Settings - Channels tab

Select the channel to view and/or modify from the Channels list. This list indicates the channel number (with respect to this module) and the serial number of the probe on each channel.

Click the Channel Settings button to display the appropriate channel settings dialog. When this button is clicked, the settings for the selected channel are read in from the thermometer readout. The Channel Settings dialog is described in detail in Section 4.4.3.14, 1560 Black Stack PRT/Thermistor Channel Settings Dialog.

4.4.3.13.3 Calibration tab

Select the Calibration tab to view and/or modify module calibration information.

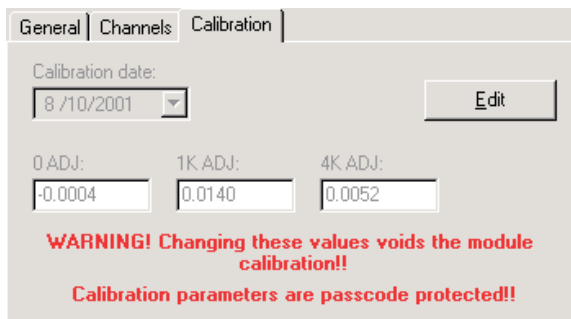


Figure 107 2568 RTDS Module Settings - Calibration tab

The settings on the Calibration tab may be passcode protected. If they are, a message appears at the bottom of the Calibration tab and the Edit button is displayed. The other controls on this tab are disabled. To edit the calibration parameters, click the Edit button and enter the current passcode for the thermometer readout. If the correct passcode is entered, the calibration parameter controls are enabled and the values may be changed. The controls remain enabled until this dialog is closed or until the current changes are applied by clicking the Apply button.



Caution: Do not change these values unless this module has been recalibrated or unless these values do not match the values located on the latest Report of Calibration for this module! Changing these values voids the current calibration!

Select the date this module was last calibrated using the drop-down calendar or by entering the date. This date should be located on the Report of Calibration for this module.

Enter the values for the calibration constants in the boxes provided. The calibration constants required by the 2568 module are: 0 ADJ, 1K ADJ and 4K ADJ. The values of these constants should be located on the Report of Calibration for this module.

4.4.3.14 1560 Black Stack PRT/Thermistor Channel Settings dialog

The 1560 Black Stack PRT/Thermistor Channel Settings dialog displays information about the selected channel.

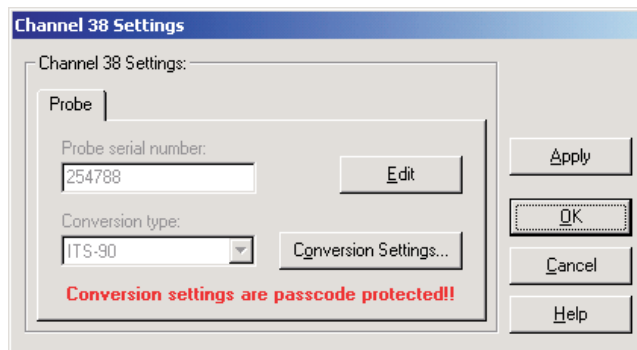


Figure 108 1560 Black Stack PRT/Thermistor Channel Settings dialog

After modifying any of the settings on this dialog, click the Apply or OK button to write the new settings to the thermometer readout. If the Apply button is clicked, this dialog remains open. If the OK button is clicked, this dialog closes after writing the settings.

LogWare has a feature that causes all of the channel settings to be read back in from the thermometer readout after the Apply button is clicked and the new settings have been written to the thermometer readout. This feature can be useful to verify that the desired settings were properly applied to the thermometer readout. To enable or disable this feature, see Section 5.6.4, Options.

To close this dialog without writing the channel settings to the thermometer readout, click the Cancel button.

Click the Help button to open the appropriate help topic for this dialog.

4.4.3.14.1 Probe tab

Select the Probe tab to view and/or modify probe settings.

The screenshot shows a dialog box titled 'Probe'. It contains two main sections. The first section is labeled 'Probe serial number:' and has a text input field containing '254788' and an 'Edit' button to its right. The second section is labeled 'Conversion type:' and has a dropdown menu showing 'ITS-90' and a 'Conversion Settings...' button to its right. At the bottom of the dialog, there is a red text message: 'Conversion settings are passcode protected!'.

Figure 109 1560 Black Stack PRT/Thermistor Channel Settings - Probe tab

The settings on the Probe tab may be passcode protected. If they are, a message appears at the bottom of the Probe tab and the Edit button is displayed. The other controls on this tab are disabled except the Conversion Settings button. If the selected conversion type does not have any settings, this button should also be disabled. To edit the probe settings, click the Edit button and enter the current passcode for the thermometer readout. If the correct passcode is entered, the probe setting controls are enabled and the values may be changed. The controls remain enabled until this dialog is closed or until the current changes are applied by clicking the Apply button.

Enter the serial number of the probe attached to this channel in the Probe serial number box. The serial number can be any combination of numbers and/or letters up to 8 characters in length. The dash “-” character can also be used in serial numbers.

Select the conversion type to use for the probe on this channel. Available conversion types will vary depending on the module that this channel belongs to. Refer to the Report of Calibration for the probe attached to this channel for information on the conversion type that the probe was calibrated for.

If the selected conversion type requires coefficients or other settings, the Conversion Settings button is enabled. Click the Conversion Settings button to display the dialog for the selected conversion type. Otherwise, the Conversion

Settings button is disabled. Please refer to Section 9, Conversion Settings Dialogs, for a detailed description of the Conversion Settings dialogs.



Note: The conversion settings are not written to the thermometer readout until the Apply or OK button is clicked on this dialog!

4.4.3.15 1560 Black Stack Thermocouple Channel Settings dialog

The 1560 Black Stack Thermocouple Channel Settings dialog displays information about the selected channel.

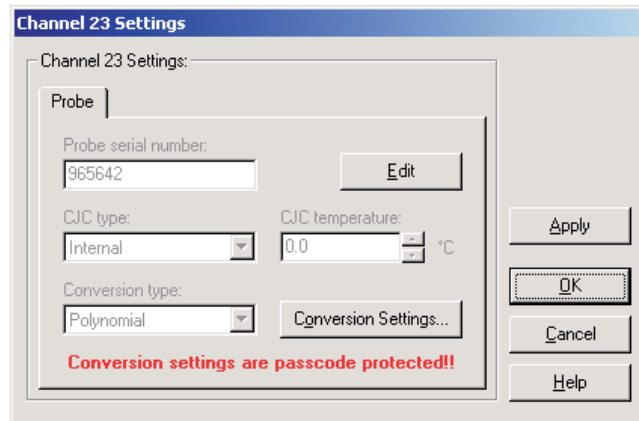


Figure 110 1560 Black Stack Thermocouple Channel Settings dialog

After modifying any of the settings on this dialog, click the Apply or OK button to write the new settings to the thermometer readout. If the Apply button is clicked, this dialog remains open. If the OK button is clicked, this dialog closes after writing the settings.

LogWare has a feature that causes all of the channel settings to be read back in from the thermometer readout after the Apply button is clicked and the new settings have been written to the thermometer readout. This feature can be useful to verify that the desired settings were properly applied to the thermometer readout. To enable or disable this feature, see Section 5.6.4, Options.

To close this dialog without writing the channel settings to the thermometer readout, click the Cancel button.

Click the Help button to open the appropriate help topic for this dialog.

4.4.3.15.1 Probe tab

Select the Probe tab to view and/or modify probe settings.

Figure 111 1560 Black Stack PRT/Thermistor Channel Settings - Probe tab

The settings on the Probe tab may be passcode protected. If they are, a message appears at the bottom of the Probe tab and the Edit button is displayed. The other controls on this tab are disabled except the Conversion Settings button. If the selected conversion type does not have any settings, this button should also be disabled. To edit the probe settings, click the Edit button and enter the current passcode for the thermometer readout. If the correct passcode is entered, the probe setting controls are enabled and the values may be changed. The controls remain enabled until this dialog is closed or until the current changes are applied by clicking the Apply button.

Enter the serial number of the probe attached to this channel in the probe serial number box. The serial number can be any combination of numbers and/or letters up to 8 characters in length. The dash “-” character can also be used in serial numbers.

Select the type of cold junction compensation to use on the selected channel using the CJC type drop-down list. Available choices are Internal and External. If External is chosen as the CJC type, the external CJC temperature must be entered in the CJC temperature box.

Enter the external CJC temperature value in the CJC temperature box. This box is disabled if the CJC type is set to Internal.



Note: The CJC temperature value must always be in °C!

Select the conversion type to use for the probe on this channel. Available conversion types will vary depending on the module that this channel belongs to. Refer to the Report of Calibration for the probe attached to this channel for information on the conversion type that the probe was calibrated for.

If the selected conversion type requires coefficients or other settings, the Conversion Settings button is enabled. Click the Conversion Settings button to display the dialog for the selected conversion type. Otherwise, the Conversion Settings button is disabled. Please refer to Section 9, Conversion Settings Dialogs, for a detailed description of the Conversion Settings dialogs.



Note: The conversion settings are not written to the thermometer readout until the Apply or OK button is clicked on this dialog!

4.4.4 1575A Super-Thermometer Settings

The 1575A Super-Thermometer Settings dialog displays information about the thermometer readout including model number, identification information and the total number of channels. It also allows many thermometer readout settings to be viewed and/or modified.

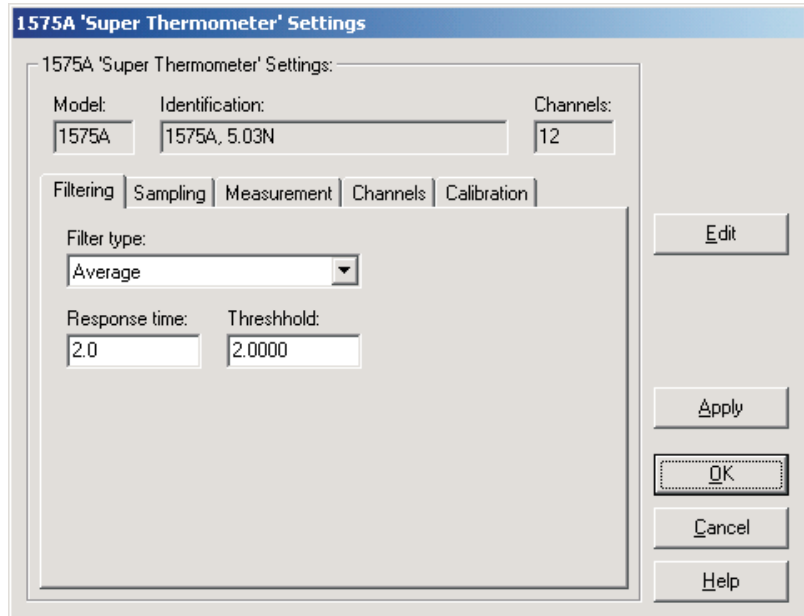


Figure 112 1575A Super-Thermometer Settings dialog

The 1575A Super-Thermometer readout must be connected to an available COM port and the communication settings must be properly configured so that this information can be read in from the thermometer readout. If communication with the thermometer readout is not successful, an error message is displayed and this dialog is blank. A progress dialog is displayed while the current thermometer readout settings are being read in.

The Model box indicates the model number of the thermometer readout. The Identification box indicates the model number and firmware version installed in this thermometer readout. The Channels box indicates the total number of input channels on this thermometer readout.

The filtering settings can be viewed and modified by selecting the Filtering tab. The sampling settings can be viewed and/or modified by selecting the Sampling tab. The measurement settings can be viewed and/or modified by selecting the Measurement tab. The Channels tab displays a list of channels and allows access to channel settings dialogs. The thermometer readout calibration information can be viewed and/or modified by selecting the Calibration tab. The settings on each of these tabs are described in detail in the following sections. For more details on these settings, please refer to the thermometer readout's User's Guide.

If any of the settings on the 1575A Super-Thermometer Settings dialog are passcode protected, a message appears at the bottom of the appropriate tab and the Edit button is displayed. The controls for the passcode protected settings are disabled. To edit the protected settings, click the Edit button and enter the current passcode for the thermometer readout. If the correct passcode is entered, the controls are enabled and the values may be changed. The controls remain enabled until this dialog is closed or until the current changes are applied by clicking the Apply button.

After modifying any of the settings on this dialog, click the Apply or OK button to write the new settings to the thermometer readout. If the Apply button is clicked, this dialog remains open. If the OK button is clicked, this dialog closes after writing the settings.


LogWare has a feature that causes all of the thermometer readout settings to be read back in from the thermometer readout after the Apply button is clicked and the new settings have been written to the thermometer readout. This feature can be useful to verify that the desired settings were properly applied to the thermometer readout. To enable or disable this feature, see Section 5.6.4, Options.

To close this dialog without writing the settings to the thermometer readout, click the Cancel button.

Click the Help button to open the appropriate help topic for this dialog.

4.4.4.1 Filtering tab

Select the Filtering tab to view and/or modify the filter settings.

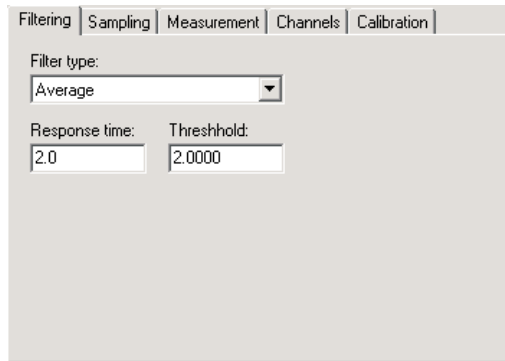

The screenshot shows a software dialog box with five tabs: Filtering, Sampling, Measurement, Channels, and Calibration. The 'Filtering' tab is selected. Inside the dialog, there is a 'Filter type:' label followed by a dropdown menu currently showing 'Average'. Below this are two text input fields: 'Response time:' containing the value '2.0' and 'Threshold:' containing the value '2.0000'.

Figure 113 1575A Super-Thermometer Settings - Filtering tab

Use the Filter type drop-down list to select the type of filtering to use. The filter type can be set to one of the following options:

- None
- Average
- Exponential

If the filter type is set to None, the Response time and Threshold settings are not applicable and are disabled. If the filter type is set to Average or Exponential, the Response time and Threshold settings are enabled.

Enter the response time for the filter in the Response time box. The response time can be set to any value between 2 and 200 seconds.

Enter the threshold to use for the selected filter in the Threshold box. The threshold can be set to any value between .0001 and 1,000,000.

4.4.4.2 Sampling tab

Select the Sampling tab to view and/or modify sampling settings.

The screenshot shows the 'Sampling' tab of the 1575A Super-Thermometer settings. It includes the following controls:

- Units:** A drop-down menu set to °C.
- Sample action:** A drop-down menu set to Run.
- Sample n:** A numeric input field set to 1.
- Sample timing:** Three numeric input fields: Sample interval (2), Integration time (2), and Conversion time (2).
- Resolutions:** Three numeric input fields: Temperature (8), Resistance (8), and Ratio (8).

A red warning message at the bottom of the tab reads: "Some sampling settings are passcode protected!!"

Figure 114 1575A Super-Thermometer Settings - Sampling tab

Select the units using the Units drop-down list. Readings can be displayed in either °C, °F, K, Ohms, Ratio or W.



Note: This setting may override the selected conversion type for each channel. Please refer to the thermometer readout's User's Guide for more information on setting the system units.

Select the sample action from the Sample action drop-down list. The sample action can be set to one of the following options:

- Run
- Stop
- Sample n

If the sample action is set to Run or Stop, the Sample n setting is not applicable and is disabled. If the filter type is set to Sample n , the Sample n setting is enabled.

If the sample action is set to Sample n , enter the number of samples in the Sample n box. The Sample n setting can be set to any value between 1 and 32,000.

The Sample timing controls allow the thermometer readout's sample timing settings to be viewed and/or modified. The sample timing settings include the Sample interval, Integration time and Conversion time.

Enter the sample interval value in the Sample interval box provided or use the up/down buttons. The sample interval can be set to any value between 2 and 99,999 seconds.



Note: *The sample interval must be greater than or equal to the integration time, which in turn must be greater than or equal to the conversion time. If the sample interval is set to a value that is less than the current integration time, the integration time and conversion time settings are automatically set to match the sample interval setting!*

Enter the integration time value in the Integration time box provided or use the up/down buttons. The integration time can be set to any value between 2 and 9,999 seconds.



Note: *The integration time must be greater than or equal to the conversion time. If the integration time is set to a value that is less than the current conversion time, the conversion time setting is automatically set to match the integration time setting!*

Enter the conversion time value in the Conversion time box provided or use the up/down buttons. The conversion time can be set to any value between 2 and 10 seconds.

The Resolutions controls allow the thermometer readout's resolution settings to be viewed and/or modified. The resolution settings include the Temperature, Resistance and Ratio settings.

Enter the resolution for temperature readings in the Temperature box provided or use the up/down buttons. The temperature resolution can be set to any value between 0 and 8 digits (places after decimal point).

Enter the resolution for resistance readings in the Resistance box provided or use the up/down buttons. The resistance resolution can be set to any value between 0 and 8 digits (places after decimal point).

Enter the resolution for ratio readings in the Ratio box provided or use the up/down buttons. The ratio resolution can be set to any value between 0 and 8 digits (places after decimal point).

4.4.4.3 Measurement tab

Select the Measurement tab to view and/or modify measurement settings.

The screenshot displays the Measurement tab of the 1575A Super-Thermometer settings. The interface includes the following elements:

- Channel mode:** A drop-down menu set to "Scanner sequence".
- Repetitions:** A numeric input field set to "1" with up/down arrows.
- Scanner channel:** A drop-down menu set to "1".
- Scanner sequence mode:** A drop-down menu set to "User-defined scanner sequence".
- Seq. channels:** A numeric input field set to "10" with up/down arrows.
- User-defined scanner sequence:**
 - Steps:** A numeric input field set to "100" with up/down arrows.
 - Scanner sequence:** A list box containing:
 - 1) Input 2
 - 2) Scanner 1
 - 3) Scanner 2
 - 4) Scanner 3
 - 5) Scanner 4
- Step channel:** A drop-down menu set to "Input 2".

Figure 115 1575A Super-Thermometer Settings - Measurement tab

Use the Channel mode drop-down list to select the channels to take measurements from. The channel mode can be set to one of the following options:

- Alternate input 1/input 2
- Input 1
- Input 2
- Scanner channel **(only if Model 2575 Mighty Mux Multiplexer is connected)**
- Scanner sequence **(only if Model 2575 Mighty Mux Multiplexer is connected)**

If the channel mode is set to Alternate input 1/input 2, Input 1 or Input 2, the scanner channel and scanner sequence mode settings are not applicable and are disabled. If the channel mode is set to Scanner channel, the Scanner channel setting is enabled. If the channel mode is set to Scanner sequence, the scanner sequence mode setting is enabled.

Enter the repetitions setting in the Repetitions box provided or use the up/down buttons. The repetitions can be set to any value between 1 and 32,000 readings.

If the channel mode is set to Scanner channel, select the scanner channel to activate using the Scanner channel drop-down list.

If the channel mode is set to Scanner sequence, select the scanner sequence mode using the Scanner sequence mode drop-down list. The scanner sequence mode can be set to one of the following options:

- Scanner channels only
- Input 2 and scanner channels
- Alternate input 2/scanner channels
- Alternate scanner channel 1/scanner channels
- User-defined scanner sequence

If the scanner sequence mode is set to Scanner channels only, Input 2 and scanner channels, Alternate input 2/scanner channels or Alternate scanner channel 1/scanner channels, the Seq. channels setting must be set and is enabled. If the scanner sequence mode is set to User-defined scanner sequence, the user-defined scanner sequence settings must be set and are enabled.

If the scanner sequence mode is set to Scanner channels only, Input 2 and scanner channels, Alternate input 2/scanner channels or Alternate scanner channel 1/scanner channels, enter the number of scanner channels to include in the scanner sequence in the Seq. channels box provided or use the up/down buttons.

If the scanner sequence mode is set to User-defined scanner sequence, the user-defined scanner sequence can be configured using the Steps, Scanner sequence and Step channel controls.

Enter the number of steps in the user-defined scanner sequence in the Steps box provided or use the up/down buttons. The number of steps in a user-defined scanner sequence can be set to any value between 1 and 100.

The Scanner sequence list indicates the current channel assignments for each step in the user-defined scanner sequence. To change the current channel assigned to a step, select the step in the list. The current channel assignment appears in the Step channel control.

Select the channel assignment for the currently selected step using the Step Channel drop-down list.

4.4.4.4 Channels tab

Select the Channels tab to view a list of channels on the 1575A Super-Ther-

thermometer readout and to access the channel settings dialogs to view and/or modify channel specific settings.

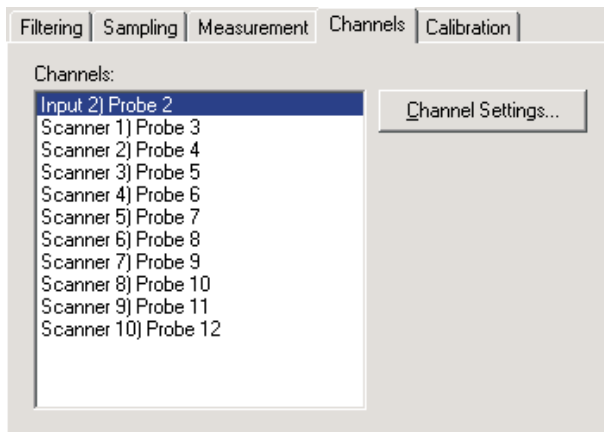


Figure 116 1575A Super-Thermometer Settings - Channels tab

Select the channel to view and/or modify from the Channels list. This list indicates the channel number and the current probe that is assigned to each channel.

Click the Channel Settings button to display the appropriate channel settings dialog. When this button is clicked, the settings for the selected channel are read in from the thermometer readout. The Channel Settings dialog is described in detail in Section 4.4.4.6, 1575A Super-Thermometer/2575 Mighty Mux Channel Settings Dialog.

4.4.4.5 Calibration tab

Select the Calibration tab to view and/or modify thermometer readout calibration information.

Figure 117 1575A Super-Thermometer Settings - Calibration tab



Caution: Do not change these values unless this thermometer readout has been recalibrated or unless these values do not match the values located on the latest Report of Calibration for this thermometer readout! Changing these values voids the current calibration!

Enter the value of the external reference resistor in the External resistor box provided. If no external reference resistor is being used, the value in this box is ignored by the thermometer readout.

Enter the values for the internal reference resistors in the boxes provided. The Model 1575A Super-Thermometer readout has four internal reference resistors: 1 Ohm, 10 Ohm, 100 Ohm and 10K Ohm. The values of these reference resistors as well as the ADC constant value should be located on the Report of Calibration for this thermometer readout.

Select the state of the internal reference resistor oven using the Oven drop-down list. The oven can either be turned On or Off. Typically, the oven should be set On.

4.4.4.6 1575A Super Thermometer/2575 Mighty Mux Channel Settings

The 1575A 'Super Thermomete' /2575 Mighty Mux Channel Settings dialog displays information about the selected PRT/Thermistor channel.

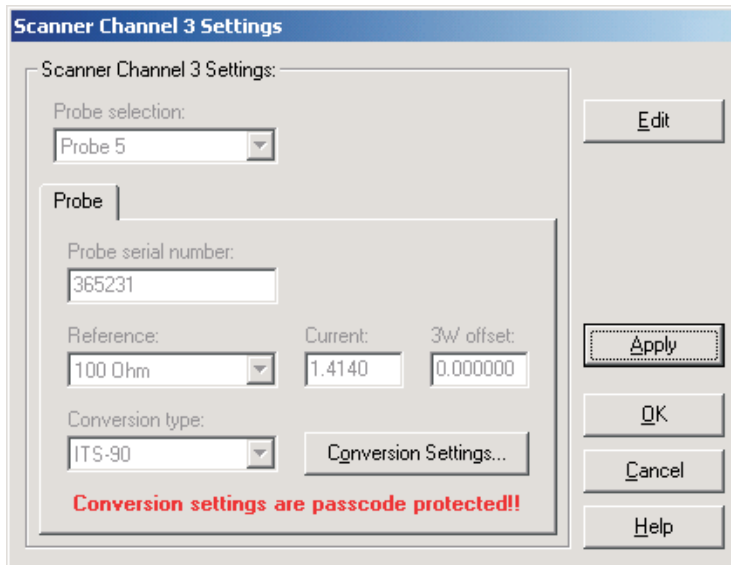


Figure 118 1575A Super Thermometer/2575 Mighty Mux Channel Settings dialog

Some or all of the settings on this dialog may be passcode protected. If so, a message appears at the bottom of the Probe tab and/or Calibration tab and the Edit button is displayed. The protected controls are disabled. To edit the protected settings, click the Edit button and enter the current passcode for the thermometer readout. If the correct passcode is entered, the protected controls are enabled and the values may be changed. The controls remain enabled until this dialog is closed or until the current changes are applied by clicking the Apply button.

Select the probe to be assigned to this channel using the Probe selection drop-down list. The Model 1575A Super-Thermometer readout can hold information for up to 16 probes simultaneously. Any one of these probes can then be assigned to any channel.

When a new probe is selected, the current settings for that probe are read in from the thermometer readout. A progress dialog is displayed while the probe settings are being read in.

After modifying any of the settings on this dialog, click the Apply or OK button to write the new settings to the thermometer readout. If the Apply button is clicked, this dialog remains open. If the OK button is clicked, this dialog closes after writing the settings.

LogWare has a feature that causes all of the channel settings to be read back in from the thermometer readout after the Apply button is clicked and the new settings have been written to the thermometer readout. This feature can be useful to verify that the desired settings were properly applied to the thermometer readout. To enable or disable this feature, see Section 5.6.4, Options.

To close this dialog without writing the channel settings to the thermometer readout, click the Cancel button.

Click the Help button to open the appropriate help topic for this dialog.

4.4.4.6.1 Probe tab

Select the Probe tab to view and/or modify probe settings.

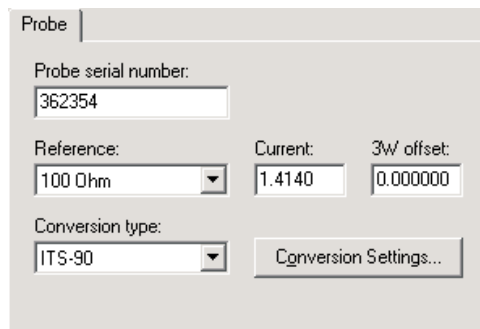


Figure 119 1575A Super-Thermometer/2575 Mighty Mux Channel Settings - Probe tab

Enter the serial number of the selected probe in the Probe serial number box. The serial number can be any combination of numbers and/or letters up to 8 characters in length. The dash “-” character can also be used in serial numbers.

Select the reference resistor for the selected probe using the Reference drop-down list. The choices for the reference resistor are:

- External resistor
- 1 Ohm
- 10 Ohm
- 100 Ohm
- 10K Ohm



Note: The External resistor option is not available for Input 2 because the external reference resistor must be connected to Input 2!

Enter the excitation current for the selected probe in the Current box provided. The excitation current can be set to any value between 0.0001 and 21.0 mA.

Enter the 3-wire offset value for the selected probe in the 3W offset box provided. The 3-wire offset can be set to any value between 0 and 1,000,000 ohms.

Select the conversion type to use for the selected probe. The Model 1575A Super-Thermometer readout supports the following conversion types:

- ITS-90
- IPTS-68
- CVD (Callendar-Van Dusen)
- W(T90)
- R(T)
- T(R)
- Polynomial
- Ohms
- Ratio

Refer to the Report of Calibration for the selected probe for information on the conversion type that the probe was calibrated for.

If the selected conversion type requires coefficients or other settings, the Conversion Settings button is enabled. Otherwise, the Conversion Settings button is disabled. Click the Conversion Settings button to display the dialog for the selected conversion type. Please refer to Section 9, Conversion Settings Dialogs, for a detailed description of the Conversion Settings dialogs.



***Note:** The conversion settings are not written to the thermometer readout until the Apply or OK button is clicked on this dialog!*

4.4.5 1590 Super-Thermometer II Settings dialog

The 1590 Super-Thermometer II Settings dialog displays information about the thermometer readout including model number, identification information and

the total number of channels. It also allows many thermometer readout settings to be viewed and/or modified.

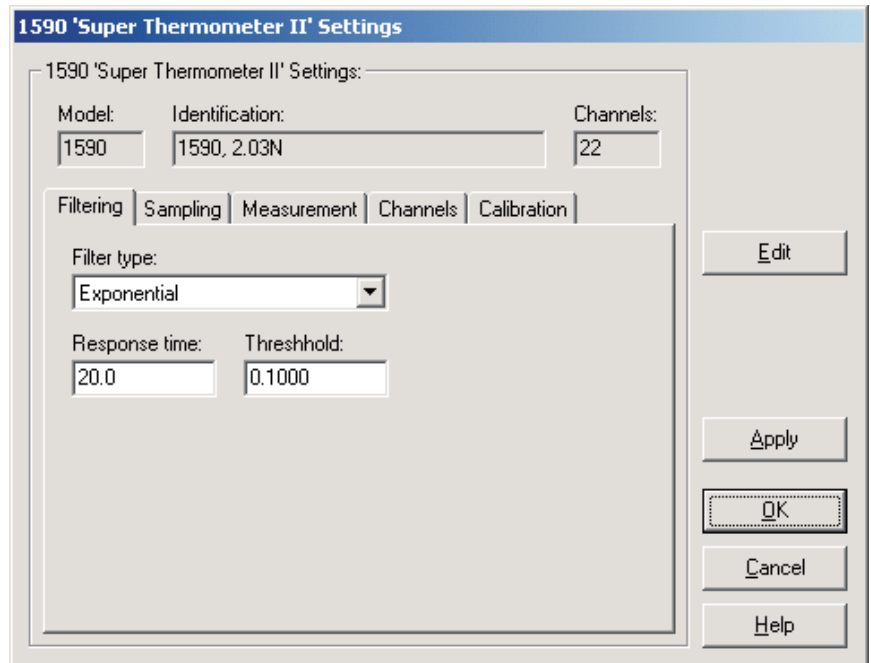


Figure 120 1590 Super-Thermometer II Settings dialog

The 1590 Super-Thermometer II readout must be connected to an available COM port and the communication settings must be properly configured so that this information can be read in from the thermometer readout. If communication with the thermometer readout is not successful, an error message is displayed and this dialog is blank. A progress dialog is displayed while the current thermometer readout settings are being read in.

The Model box indicates the model number of the thermometer readout. The Identification box indicates the model number and firmware version installed in this thermometer readout. The Channels box indicates the total number of input channels on this thermometer readout.

The filtering settings can be viewed and/or modified by selecting the Filtering tab. The sampling settings can be viewed and/or modified by selecting the Sampling tab. The measurement settings can be viewed and/or modified by selecting the Measurement tab. The Channels tab displays a list of channels and allows access to channel settings dialogs. The thermometer readout calibration information can be viewed and/or modified by selecting the Calibration tab. The settings on each of these tabs are described in detail in the following sec-

tions. For more details on these settings, please refer to the thermometer readout's User's Guide.

If any of the settings on the 1590 Super-Thermometer II Settings dialog are passcode protected, a message appears at the bottom of the appropriate tab and the Edit button is displayed. The controls for the passcode protected settings are disabled. To edit the protected settings, click the Edit button and enter the current passcode for the thermometer readout. If the correct passcode is entered, the controls are enabled and the values may be changed. The controls remain enabled until this dialog is closed or until the current changes are applied by clicking the Apply button.

After modifying any of the settings on this dialog, click the Apply or OK button to write the new settings to the thermometer readout. If the Apply button is clicked, this dialog remains open. If the OK button is clicked, this dialog closes after writing the settings.

LogWare has a feature that causes all of the thermometer readout settings to be read back in from the thermometer readout after the Apply button is clicked and the new settings have been written to the thermometer readout. This feature can be useful to verify that the desired settings were properly applied to the thermometer readout. To enable or disable this feature, see Section 5.6.4, Options.

To close this dialog without writing the settings to the thermometer readout, click the Cancel button.

Click the Help button to open the appropriate help topic for this dialog.

4.4.5.1 Filtering tab

Select the Filtering tab to view and/or modify the filter settings.

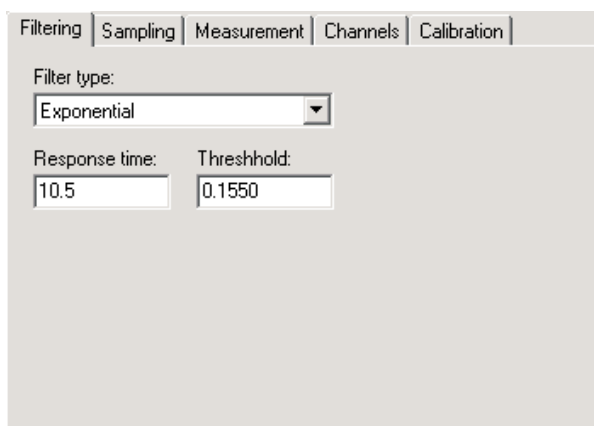


Figure 121 1590 Super-Thermometer II Settings - Filtering tab

Use the Filter type drop-down list to select the type of filtering to use. The filter type can be set to one of the following options:

- None
- Average
- Exponential

If the filter type is set to None, the Response time and Threshold settings are not applicable and are disabled. If the filter type is set to Average or Exponential, the Response time and Threshold settings are enabled.

Enter the response time for the filter in the Response time box. The response time can be set to any value between 2 and 200 seconds.

Enter the threshold to use for the selected filter in the Threshold box. The threshold can be set to any value between .0001 and 1,000,000.

4.4.5.2 Sampling tab

Select the Sampling tab to view and/or modify sampling settings.

Figure 122 1590 Super-Thermometer II Settings - Sampling tab

Select the units using the Units drop-down list. Readings can be displayed in either °C, °F, K, Ohms, Ratio or W.



Note: This setting may override the selected conversion type for each channel. Please refer to the thermometer readout's User's Guide for more information on setting the system units.

Select the sample action from the Sample action drop-down list. The sample action can be set to one of the following options:

- Run
- Stop
- Sample n

If the sample action is set to Run or Stop, the Sample n setting is not applicable and is disabled. If the filter type is set to Sample n , the Sample n setting is enabled.

If the sample action is set to Sample n , enter the number of samples in the Sample n box. The Sample n setting can be set to any value between 1 and 32,000.

The Sample timing controls allow the thermometer readout's sample timing settings to be viewed and/or modified. The sample timing settings include the Sample interval, Integration time and Conversion time.

Enter the sample interval value in the Sample interval box provided or use the up/down buttons. The sample interval can be set to any value between 2 and 99,999 seconds.



***Note:** The sample interval must be greater than or equal to the integration time, which in turn must be greater than or equal to the conversion time. If the sample interval is set to a value that is less than the current integration time, the integration time and conversion time settings are automatically set to match the sample interval setting!*

Enter the integration time value in the Integration time box provided or use the up/down buttons. The integration time can be set to any value between 2 and 9,999 seconds.



***Note:** The integration time must be greater than or equal to the conversion time. If the integration time is set to a value that is less than the current conversion time, the conversion time setting is automatically set to match the integration time setting!*

Enter the conversion time value in the Conversion time box provided or use the up/down buttons. The conversion time can be set to any value between 2 and 10 seconds.

The Resolutions controls allow the thermometer readout's resolution settings to be viewed and/or modified. The resolution settings include the Temperature, Resistance and Ratio settings.

Enter the resolution for temperature readings in the Temperature box provided or use the up/down buttons. The temperature resolution can be set to any value between 0 and 8 digits (places after decimal point).

Enter the resolution for resistance readings in the Resistance box provided or use the up/down buttons. The resistance resolution can be set to any value between 0 and 8 digits (places after decimal point).

Enter the resolution for ratio readings in the Ratio box provided or use the up/down buttons. The ratio resolution can be set to any value between 0 and 8 digits (places after decimal point).

4.4.5.3 Measurement tab

Select the Measurement tab to view and/or modify measurement settings.

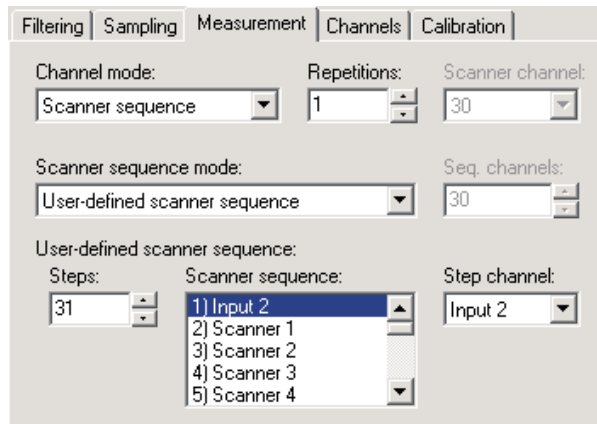


Figure 123 1590 Super-Thermometer II Settings - Measurement tab

Use the Channel mode drop-down list to select the channels to take measurements from. The channel mode can be set to one of the following options:

- Alternate input 1/input 2
- Input 1
- Input 2
- Scanner channel (only if Model 2590 ‘Mighty Mux II’ Multiplexer(s) is/are connected)
- Scanner sequence (only if Model 2590 ‘Mighty Mux II’ Multiplexer(s) is/are connected)

If the channel mode is set to Alternate Input 1/Input 2, Input 1 or Input 2, the scanner channel and scanner sequence mode settings are not applicable and are disabled. If the channel mode is set to Scanner channel, the Scanner channel setting is enabled. If the channel mode is set to Scanner sequence, the scanner sequence mode setting is enabled.

Enter the repetitions setting in the Repetitions box provided or use the up/down buttons. The repetitions can be set to any value between 1 and 32,000 readings.

If the channel mode is set to Scanner channel, select the scanner channel to activate using the Scanner channel drop-down list.

If the channel mode is set to Scanner sequence, select the scanner sequence mode using the Scanner sequence mode drop-down list. The scanner sequence mode can be set to on of the following options:

- Scanner channels only
- Input 2 and scanner channels
- Alternate input 2/scanner channels
- Alternate scanner channel 1/scanner channels
- User-defined scanner sequence

If the scanner sequence mode is set to Scanner channels only, Input 2 and scanner channels, Alternate input 2/scanner channels or Alternate scanner channel 1/scanner channels, the Seq. channels setting must be set and is enabled. If the scanner sequence mode is set to User-defined scanner sequence, the user-defined scanner sequence settings must be set and are enabled.

If the scanner sequence mode is set to Scanner channels only, Input 2 and scanner channels, Alternate input 2/scanner channels or Alternate scanner channel 1/scanner channels, enter the number of scanner channels to include in the scanner sequence in the Seq. channels box provided or use the up/down buttons.

If the scanner sequence mode is set to User-defined scanner sequence, the user-defined scanner sequence can be configured using the Steps, Scanner sequence, and Step channel controls.

Enter the number of steps in the user-defined scanner sequence in the Steps box provided or use the up/down buttons. The number of steps in a user-defined scanner sequence can be set to any value between 1 and 100.

The Scanner sequence list indicates the current channel assignments for each step in the user-defined scanner sequence. To change the current channel assigned to a step, select the step in the list. The current channel assignment appears in the Step channel control.

Select the channel assignment for the currently selected step using the Step Channel drop-down list.

4.4.5.4 Channels tab

Select the Channels tab to view a list of channels on the 1590 Super-Thermom-

eter II readout and to access the channel settings dialogs to view and/or modify channel specific settings.

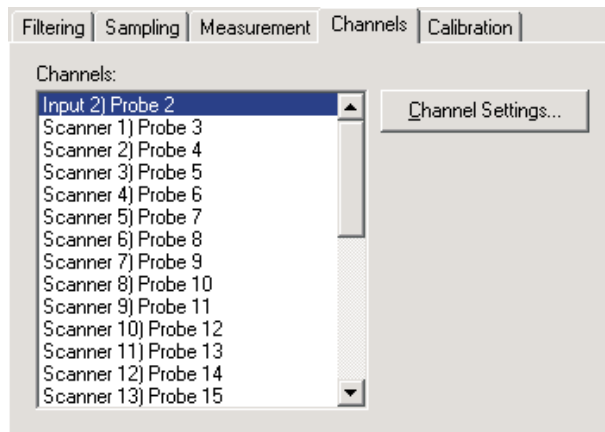


Figure 124 1590 Super-Thermometer II Settings - Channels tab

Select the channel to view and/or modify from the Channels list. This list indicates the channel number and the current probe that is assigned to each channel.

Click the Channel Settings button to display the appropriate channel settings dialog. When this button is clicked, the settings for the selected channel are read in from the thermometer readout. The Channel Settings dialog is described in detail in Section 4.4.5.6, 1590 Super-Thermometer II/2590 'Mighty Mux II' Channel Settings Dialog.

4.4.5.5 Calibration tab

Select the Calibration tab to view and/or modify thermometer readout calibration information.

The screenshot shows the Calibration tab with the following settings:

- External resistor: 100.00121231300
- Internal resistors:
 - 1 Ohm: 0.9999544564600
 - 10 Ohm: 10.001212346465
 - 100 Ohm: 100.00000484600
 - 10K Ohm: 1.000245647E+04
- ADC: -0.37
- Oven: On

WARNING! Changing these values voids the thermometer calibration!!
Calibration parameters are passcode protected!!

Figure 125 1590 Super-Thermometer II Settings - Calibration tab



Caution: Do not change these values unless this thermometer readout has been recalibrated or unless these values do not match the values located on the latest Report of Calibration for this thermometer readout! Changing these values voids the current calibration!

Enter the value of the external reference resistor in the External resistor box provided. If no external reference resistor is being used, the value in this box is ignored by the thermometer readout.

Enter the values for the internal reference resistors in the boxes provided. The Model 1590 Super-Thermometer II readout has four internal reference resistors: 1 Ohm, 10 Ohm, 100 Ohm and 10K Ohm. The values of these reference resistors as well as the ADC constant value should be located on the Report of Calibration for this thermometer readout.

Select the state of the internal reference resistor oven using the Oven drop-down list. The oven can either be turned On or Off. Typically, the oven should be set On.

4.4.5.6 1590 Super-Thermometer II/2590 'Mighty Mux II' Channel Settings dialog

The 1590 Super-Thermometer II/2590 'Mighty Mux II' Channel Settings dialog displays information about the selected PRT/Thermistor channel.

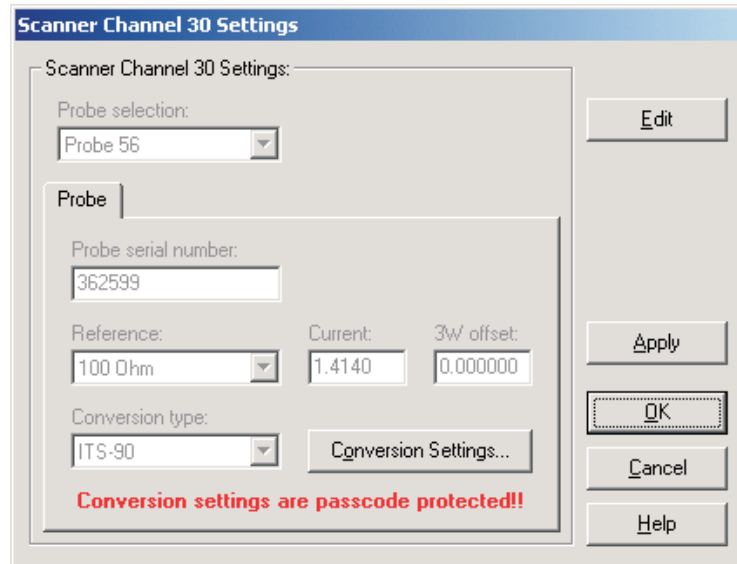


Figure 126 1590 Super-Thermometer II/2590 'Mighty Mux II' Channel Settings dialog

Some or all of the settings on this dialog may be passcode protected. If so, a message appears at the bottom of the Probe tab and/or Calibration tab and the Edit button is displayed. The protected controls are disabled. To edit the protected settings, click the Edit button and enter the current passcode for the thermometer readout. If the correct passcode is entered, the protected controls are enabled and the values may be changed. The controls remain enabled until this dialog is closed or until the current changes are applied by clicking the Apply button.

Select the probe to be assigned to this channel using the Probe selection drop-down list. The Model 1590 Super-Thermometer II readout can hold information for up to 56 probes simultaneously. Any one of these probes can then be assigned to any channel.

When a new probe is selected, the current settings for that probe are read in from the thermometer readout. A progress dialog is displayed while the probe settings are being read in.

After modifying any of the settings on this dialog, click the Apply or OK button to write the new settings to the thermometer readout. If the Apply button is

clicked, this dialog remains open. If the OK button is clicked, this dialog closes after writing the settings.


LogWare has a feature that causes all of the channel settings to be read back in from the thermometer readout after the Apply button is clicked and the new settings have been written to the thermometer readout. This feature can be useful to verify that the desired settings were properly applied to the thermometer readout. To enable or disable this feature, see Section 5.6.4, Options.

To close this dialog without writing the channel settings to the thermometer readout, click the Cancel button.

Click the Help button to open the appropriate help topic for this dialog.

4.4.5.6.1 Probe tab

Select the Probe tab to view and/or modify probe settings.

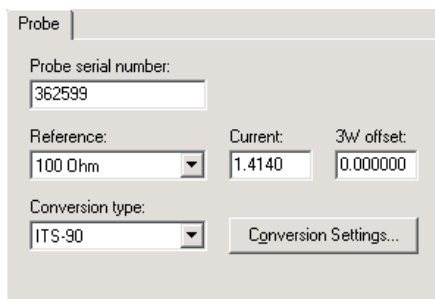

The screenshot shows a dialog box titled "Probe" with the following fields and controls:

- Probe serial number: 362599
- Reference: 100 Ohm (dropdown menu)
- Current: 1.4140
- 3W offset: 0.000000
- Conversion type: ITS-90 (dropdown menu)
- Conversion Settings... (button)

Figure 127 1590 Super-Thermometer II/2590 'Mighty Mux II' Channel Settings - Probe tab

Enter the serial number of the selected probe in the Probe serial number box. The serial number can be any combination of numbers and/or letters up to 8 characters in length. The dash “-” character can also be used in serial numbers.

Select the reference resistor for the selected probe using the Reference drop-down list. The choices for the reference resistor are:

- External resistor
- 1 Ohm
- 10 Ohm
- 100 Ohm
- 10K Ohm



Note: The External resistor option is not available for Input 2 because the external reference resistor must be connected to Input 2!

Enter the excitation current for the selected probe in the Current box provided. The excitation current can be set to any value between 0.0001 and 21.0 mA.

Enter the 3-wire offset value for the selected probe in the 3W offset box provided. The 3-wire offset can be set to any value between 0 and 1,000,000 ohms.

Select the conversion type to use for the selected probe. The Model 1590 Super-Thermometer II readout supports the following conversion types:

- ITS-90
- IPTS-68
- CVD (Callendar-Van Dusen)
- W(T90)
- R(T)
- T(R)
- Polynomial
- Ohms
- Ratio

Refer to the Report of Calibration for the selected probe for information on the conversion type that the probe was calibrated for.

If the selected conversion type requires coefficients or other settings, the Conversion Settings button is enabled. Otherwise, the Conversion Settings button is disabled. Click the Conversion Settings button to display the dialog for the selected conversion type. Please refer to Section 9, Conversion Settings Dialogs, for a detailed description of the Conversion Settings dialogs.



Note: The conversion settings are not written to the thermometer readout until the Apply or OK button is clicked on this dialog!

4.4.6 1594A/1595A Hart Super-Thermometer Settings dialog

The 1594A/1595A Hart Super-Thermometer Settings dialog displays information about the thermometer readout including model number, identification in-

formation and the total number of channels. It also allows many thermometer readout settings to be viewed and modified.

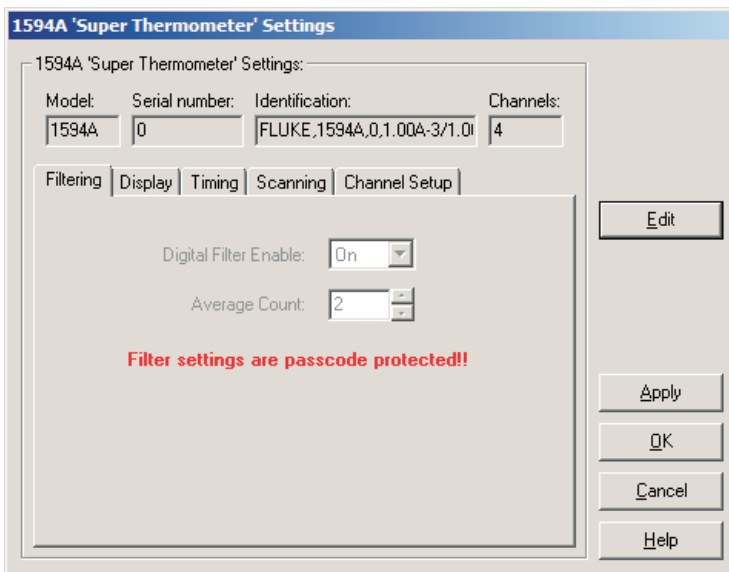


Figure 128 1594A/1595A Hart Super-Thermometer settings

The 1594A/1595A Hart Super-Thermometer readout must be connected to an available COM port and the communication settings must be properly configured so that this information can be read in from the thermometer readout. If communication with the thermometer readout is not successful, an error message is displayed and this dialog is blank. A progress dialog is displayed while the current thermometer readout settings are being read in.

The Model box indicates the model number of the thermometer readout. The Serial number box indicates the serial number assigned to this thermometer readout. The Identification box indicates the model number, serial number and firmware version installed in this thermometer readout. The Channels box indicates the total number of input channels on this thermometer readout.

The filtering settings can be viewed and modified by selecting the Filtering tab. The display settings can be viewed and modified by selecting the Display tab. The timing settings can be viewed and modified by selecting the Timing tab. The scan settings can be viewed and modified by selecting the Scanning tab. The Channel Setup tab displays a list of channels and allows access to probe settings and channel settings dialogs. The settings on each of these tabs are described in detail in the following sections. For more detail on these settings, please refer to the thermometer readout's User's Guide.

If any of the settings on the 1594A/1595A Hart Super-Thermometer Settings dialog are passcode protected, a message appears at the bottom of the appropriate tab and the Edit button is displayed. The controls for the passcode protected settings are disabled. To edit the protected settings, click the Edit button and enter the current passcode for the thermometer readout. If the correct passcode is entered, the controls are enabled and the values may be changed.

After modifying any of the settings on this dialog, click the Apply or OK button to write the new settings to the thermometer readout. If the Apply buttons clicked, this dialog remains open. If the OK button is clicked, this dialog closes after writing the settings.

LogWare has a feature that causes all of the thermometer readout settings to be read back in from the thermometer readout after the Apply button is clicked and the new settings have been written to the thermometer readout. This feature can be useful to verify that the desired settings were properly applied to the thermometer readout. To enable or disable this feature, see Section 5.6.4, Options.

To close this dialog without writing the settings to the thermometer readout, click the Cancel button.

Click the Help button to open the appropriate help topic for this dialog.

4.4.6.1 Filtering tab

Select the Filtering tab to view and/or modify the filter settings.

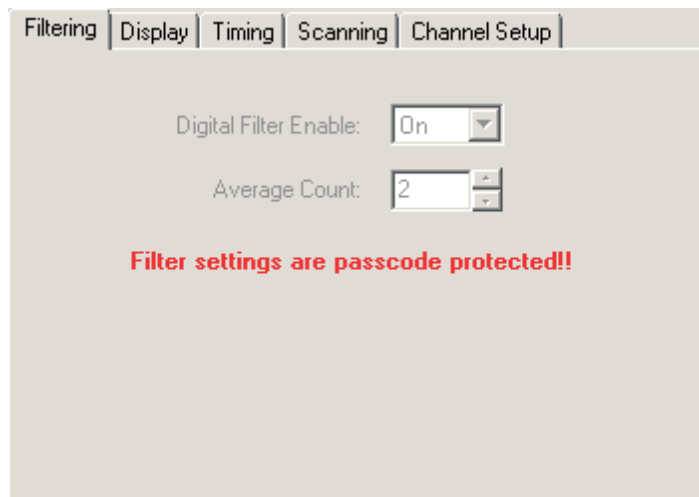


Figure 129 1594A/1595A Hart Super-Thermometer Filtering tab

Use the Digital Filter Enable drop-down list to select whether the digital filter is enabled (On) or disabled (Off). If the setting is Off, the Average Count setting is not applicable and is disabled.

The Average Count set the number of Display Interval measurements that are averaged by the digital filter. Possible values are between 2 and 100. If the Digital Filter Enable is Off, this settings is not applicable and is disabled.

4.4.6.2 Display tab

Select the Display tab to view and/or modify display settings.

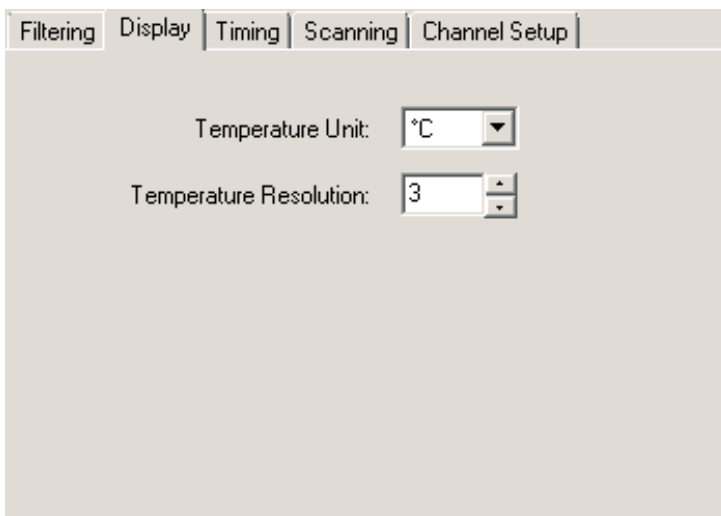


Figure 130 1594A/1595A Hart Super-Thermometer Display tab

Select the units using the Units drop-down list. Readings can be displayed in either °C, °F or K.



Note: This setting may be overridden depending on the selected probe conversion type and channel settings for each channel. Please refer to the thermometer readout's User's Guide for more information on setting the system units.

Enter the resolution for temperature readings in the Temperature Resolution box provided or click the up/down buttons. The temperature resolution can be set to any value between 0 and 6 digits (places after decimal point).

4.4.6.3 4.4.6.3 Timing tab

Select the Timing tab to view and/or modify timing settings.

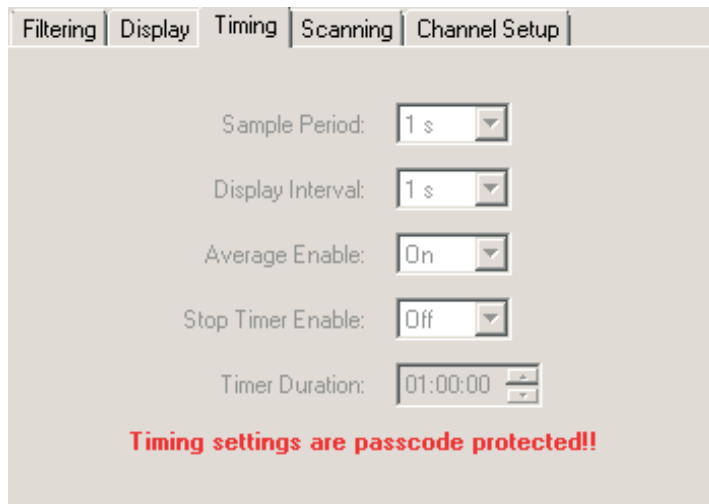


Figure 131 1594A/1595A Hart Super-Thermometer Timing tab

Use the Sample Period drop-down list to select the length of time required to obtain an individual measurement. Possible values are 1 s, 2 s, 5 s and 10 s.



Note: Changing this setting may affect the Display Interval setting, as Sample Period cannot be greater than Display Interval.

Use the Display Interval drop-down list to select how often the instrument's display, statistics fields, memory and communications ports are updated. Possible values are 1 s, 2 s, 10 s, 30 s and 60 s.



Note: Changing this setting may affect the Sample Period setting, as Sample Period cannot be greater than Display Interval.

Use the Average Enable drop-down list to select whether averaging is enabled (On) or disabled (Off). If the setting is On, all measurements acquired at the Sample Period rate are averaged prior to updating the instrument's display. If the setting is Off, the latest measurement acquired at the Sample Period rate is used to update the instrument's display.

Use the Stop Timer Enable drop-down list to select whether the stop timer is enabled (On) or disabled (Off). If the setting is On, a timer begins counting down when a measurement is started and stops the measurement automatically

when the timer elapses. If the setting is Off, the Timer Duration is not applicable and is disabled.

The Timer Duration determines the amount of time a measurement runs. If the Stop Timer Enable is Off, this setting is not applicable and is disabled. The timer duration can be set to any time between 00:00:01 and 23:59:59 with the time format set as hh:mm:ss.

4.4.6.4 Scanning tab

Select the Scanning tab to view and/or modify scan settings.

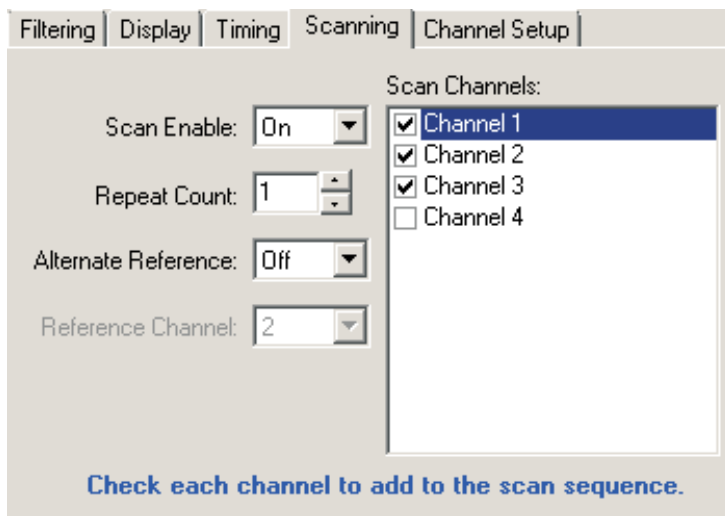


Figure 132 1594A/1595A Hart Super-Thermometer Scanning tab

Use the Scan Enable drop-down list to select whether scanning is enabled (On) or disabled (Off). If Scan Enable is Off, Repeat Count, Alternate Reference and Reference Channel are not applicable and are disabled and only one channel can be selected in the Scan Channels list.

The Repeat Count determines how many times a channel is measured before the instrument continues to the next channel to be measured. If Scan Enable is Off, this setting is not applicable and is disabled.

Use the Alternate Reference drop-down list to select whether the reference channel is automatically measured after each individual channel is measured. If this setting is Off, the Reference Channel is not applicable and is disabled. If Scan Enable is Off, this setting is not applicable and is disabled.

When the Alternate Reference is enabled (On), this setting determines which channel shall be used for the reference. If Scan Enable or Alternate Reference is Off, this setting is not applicable and is disabled.

The Scan Channels list shows all available channels in 1594A/1595A Hart Super-Thermometer readout. If Scan Enable is Off, a check next to a channel indicates the measurement channel. Click the checkbox next to a channel to change the measurement channel. If Scan Enable is On, a check next to a channel indicates the channel is part of the instrument's scan sequence. Click the checkbox next to a channel to change the scan sequence.

4.4.6.5 Channel Setup tab

Select the Channel Setup tab to view a list of channels on the 1594A/1595A Hart Super-Thermometer readout and to access the channel and probe settings for each channel.

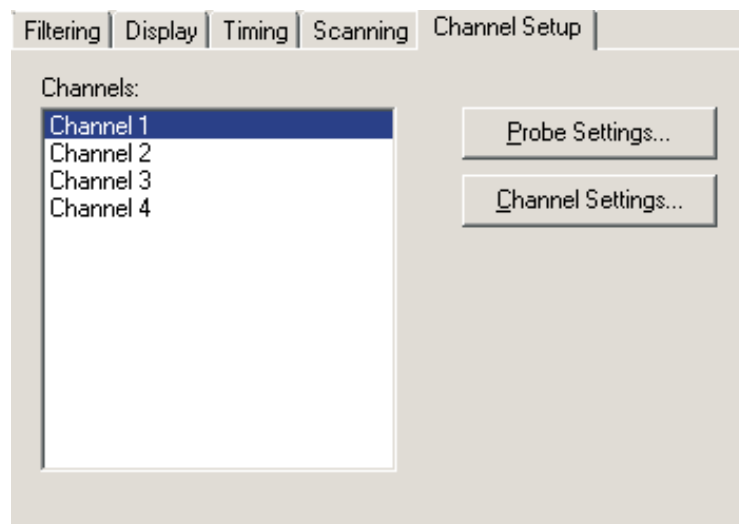


Figure 133 1594A/1595A Hart Super-Thermometer Channel setup tab

The Channels list shows all available channels in 1594A/1595A Hart Super-Thermometer readout. Select a channel to view or modify the channel or probe settings.

Click the Probe Settings button to display the appropriate probe settings for the selected channel. When this button is clicked, the probe settings are read in from the thermometer readout.

Click the Channel Settings button to display the appropriate channel settings for the selected channel. When this button is clicked, the channel settings are read in from the thermometer readout.

4.4.6.6 1594A/1595A Hart Super-Thermometer Probe Settings dialog

The 1594A/1595A Hart Super-Thermometer probe settings dialog displays probe information pertaining to the current channel selection.

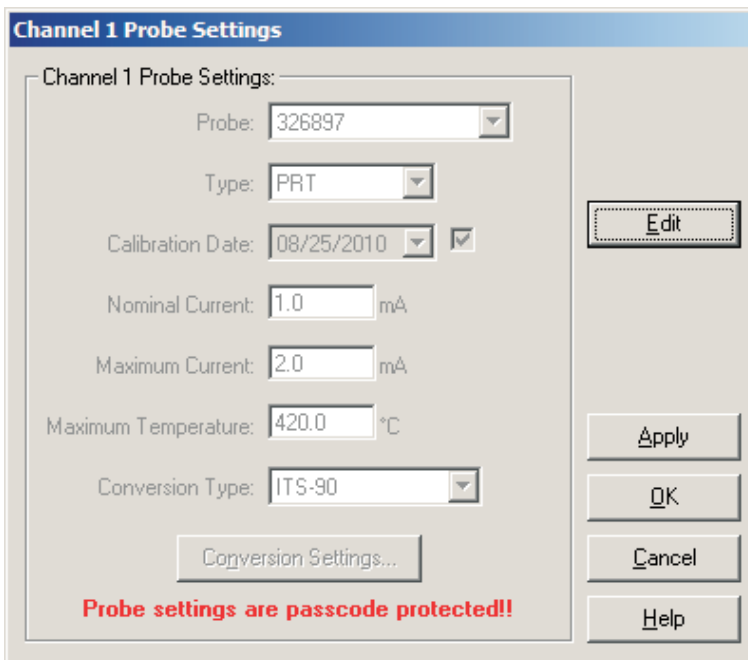


Figure 134 1594A/1595A Hart Super-Thermometer Probe settings dialog

The settings on the Probe Settings dialog may be passcode protected. If so, a message appears at the bottom of the dialog and the Edit button is displayed. The controls on this dialog are disabled when the settings are protected. To edit the probe settings, click the Edit button and enter the current passcode for the thermometer readout. If the correct passcode is entered, the controls are enabled and the values may be changed.

Select the probe to be assigned to this channel using the Probe selection drop-down list. The Model 1594A/1595A Hart Super-Thermometer readout can hold information for up to 100 probes. Any one of these probes can then be assigned to any channel.

When a new probe is selected, the current settings for that probe are read in from the thermometer readout. A progress dialog is displayed while the probe settings are being read in.

After modifying any of the settings on this dialog, click the Apply or OK button to write the new settings to the thermometer readout. If the Apply button is

clicked, this dialog remains open. If the OK button is clicked, this dialog closes after writing the settings.

To close this dialog without writing the channel settings to the thermometer readout, click the Cancel button.

Click the Help button to open the appropriate help topic for this dialog.

Use the Type drop-down list to select the probe type for the selected probe. Possible probe types are PRT, Thermistor or Resistor.

The Calibration Date may be used to record the date the probe was calibrated. The calibration date is an optional parameter. The checkbox next to the calibration date field indicates whether the probe's calibration date is saved in the instrument. If the checkbox is checked, the calibration date is saved in the instrument. If the checkbox is unchecked, the probe's calibration date is set to zeros in the instrument.

The Nominal Current is the usual sensing current that should be used to measure the resistance of the probe. Possible values depend on the probe type selected.

- PRT probe can be set to any value between 0.01 and 20.0 mA
- Thermistor probe can be set to any value between 0.001 and 0.1 mA
- Resistor probe can be set to any value between 0.001 and 20 mA.

The Maximum Current sets a limit for the probe's sensing current. Possible values depend on the probe type selected.

- PRT probe can be set to any value between 0.01 and 20.0 mA
- Thermistor probe can be set to any value between 0.001 and 0.1 mA
- Resistor probe can be set to any value between 0.001 and 20 mA.

The Maximum Temperature sets the temperature limit for the probe. Possible values depend on the probe type selected.

- PRT probe can be set to any value between 25.0 and 2000.0 °C
- Thermistor probe can be set to any value between 25.0 and 1000.0 °C
- Maximum temperature is not available for a resistor probe.

Select the conversion type to use for the selected probe. The Model 1594A/1595A Hart Super-Thermometer readout supports the following conversion types:

- ITS-90
- CVD-ALPHA
- CVD-ABC
- PT-100
- R(T)
- T(R)

- Polynomial
- Ohms

Refer to the Report of Calibration for the selected probe for information on the conversion type that the probe was calibrated for.

If the selected conversion type requires coefficients or other settings, the Conversion Settings button is enabled. Click the Conversion Settings button to display the dialog for the selected conversion type. Otherwise, the Conversion Settings button is disabled.



Note: The conversion settings are not written to the thermometer readout until the Apply or OK button is clicked on this dialog!

4.4.6.7 1594A/1595A Hart Super-Thermometer Channel Settings dialog

The 1594A/1595A Hart Super-Thermometer channel settings dialog displays information pertaining to the current channel selection.

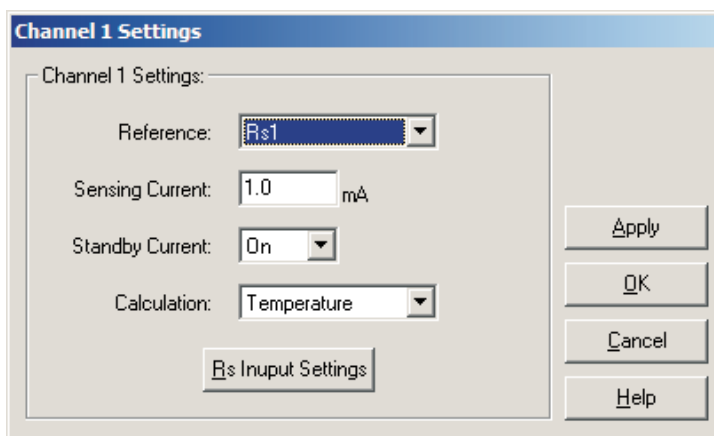


Figure 135 1594A/1595A Hart Super-Thermometer Channel settings

After modifying any of the settings on this dialog, click the Apply or OK button to write the new settings to the thermometer readout. If the Apply button is clicked, this dialog remains open. If the OK button is clicked, this dialog closes after writing the settings.

To close this dialog without writing the channel settings to the thermometer readout, click the Cancel button.

Click the Help button to open the appropriate help topic for this dialog.

Use the Reference drop-down list to select the reference resistor for the selected channel. Possible internal reference resistor values are 1 ohm, 10 ohm,

25 ohm, 100 ohm and 10,000 ohm. In addition, an external reference resistor connected to Rs1, Rs2, CH2 or CH4 can be selected.

Sensing Current

Sensing Current the sensing current used with the channel and probe. Possible values depend on the reference resistor selected.

- 1 ohm and 10 ohm internal reference can have the sensing current set to any value between 0.0001 and 20.0 mA
- 25 ohm internal reference can have the sensing current set to any value between 0.0001 and 10.0 mA
- 100 ohm internal reference can have the sensing current set to any value between 0.0001 and 5.0 mA
- 10,000 ohm internal reference can have the sensing current set to any value between 0.0001 and 0.1 mA
- The external reference can have the sensing current set to any value between 0.0001 and 20.0 mA

The sensing current is also limited by the probe's or reference resistor's maximum current value.

Use the Standby Current drop-down list to select whether the standby current is enabled (On) or disabled (Off) for the selected channel.

Use the Calculation drop-down list to determine the type of measurement calculation that is performed for the selected channel. Possible options are Ratio, Resistance and Temperature.

The Temperature option is not available if any of the following are true:

- The probe assignment for the selected channel is Undefined
- The probe conversion type is set to None
- The probe type is set to Resistor
- The reference resistor is set to Variable Resistance

The Resistance option is not available if the reference resistor is set to Variable Resistance.

If the selected reference is an external resistor, the Rs Input Settings button is enabled. Click the Rs Input Settings button to display the dialog for the reference channel. Otherwise, the Rs Input Settings button is disabled.

4.4.6.8 1594A/1595A Hart Super-Thermometer Rs Settings dialog

The 1594A/1595A Hart Super-Thermometer Rs Settings dialog displays reference resistor information pertaining to the current channel selection.

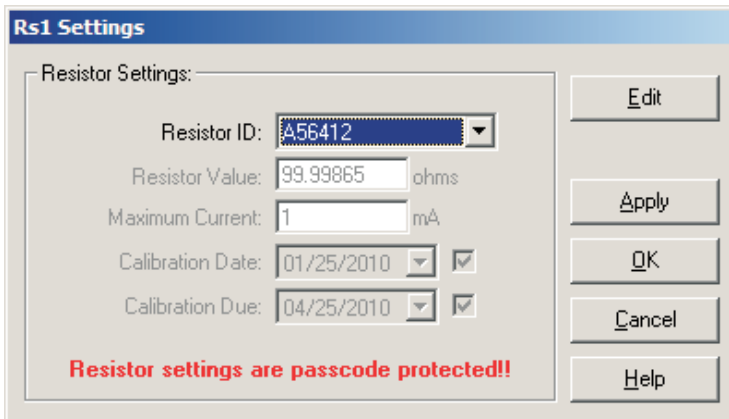


Figure 136 1594A/1595A Hart Super-Thermometer Rs1 settings

The settings on the Rs Settings dialog may be passcode protected. If they are, a message appears at the bottom of the dialog and the Edit button is displayed. The controls on this dialog are disabled, except the Resistor ID. To edit the resistor settings, click the Edit button and enter the current passcode for the thermometer readout. If the correct passcode is entered, the controls are enabled and the values may be changed.

Select the resistor to be assigned to this Rs channel using the Resistor ID selection drop-down list. The Model 1594A/1595A Hart Super-Thermometer readout can hold information for up to 50 resistors. Any one of these resistors can then be assigned to any Rs channel.

When a different resistor is selected, the current settings for that resistor are read in from the thermometer readout. A progress dialog is displayed while the probe settings are being read in.

After modifying any of the settings on this dialog, click the Apply or OK button to write the new settings to the thermometer readout. If the Apply button is clicked, this dialog remains open. If the OK button is clicked, this dialog closes after writing the settings.

To close this dialog without writing the channel settings to the thermometer readout, click the Cancel button.

Click the Help button to open the appropriate help topic for this dialog.

Use the Type drop-down list to select the probe type for the selected probe. Possible probe types are PRT, Thermistor or Resistor.

The Resistor Value is the calibrated resistance of the resistor.

The Maximum Current sets a limit for the resistor's sensing current. Possible values are any value between 0.01 and 20.0 mA.

The Calibration Date may be used to record the date the resistor was calibrated. The calibration date is an optional parameter. The checkbox next to the calibration date field indicates whether the resistor's calibration date is saved in the instrument. If the checkbox is checked, the resistor's calibration date is saved in the instrument. If the checkbox is unchecked, the resistor's calibration date is set to zeros in the instrument.

The Calibration Due date may be used to record the date the resistor is due for recalibration. The calibration due date is an optional parameter. The checkbox next to this field indicates whether the resistor's calibration due date is saved in the instrument. If the checkbox is checked, the resistor's calibration due date is saved in the instrument. If the checkbox is unchecked, the resistor's calibration due date is set to zeros in the instrument.

5 Menus

LogWare's features and tools are all accessible through menu options. This section describes each of LogWare's menus.

5.1 File menu

The File menu provides options for creating a new real-time configuration, opening and saving real-time configuration files, printing and exiting LogWare.

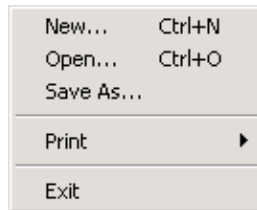



Figure 137 File menu

5.1.1 New

The New option can be used to clear the current real-time log session settings and display the Log Session Settings dialog.


The  button on the toolbar corresponds to the New option.

When selecting the New option, a prompt is displayed to confirm that the current real-time log session settings (if any) will be reset. To abort resetting the current real-time log session settings, click the No button. Otherwise, click the Yes button to reset the current log session settings. The Log Session Settings dialog will be displayed. Please refer to Section 2.3, Log Session Settings Dialog, for more information on configuring a real-time log session.

The New option and toolbar button are disabled while a real-time log session window is open.

5.1.2 Open

The Open option allows a real-time log session configuration file to be opened. If the same basic configuration is frequently used to perform real-time logging, saving the log session settings to a configuration file can greatly reduce the time it takes to configure and start a real-time log session.

The  button on the toolbar corresponds to the Open option.



Note: When opening a configuration file, LogWare attempts to communicate with the thermometer readout to verify that the settings saved in the configuration file are compatible with the current thermometer readout configuration. Therefore, it is necessary to have the thermometer readout connected and powered on in order to successfully open a configuration file.

When selecting the Open option, the Windows® Open dialog is displayed. The \Config folder should be selected by default (or the folder that is specified for configuration files on the LogWare Options dialog). Real-time log session configuration files should have a **.rls** extension. To cancel opening a configuration file, click the Cancel button. Otherwise, select the configuration file to open and click the Open button. For more details about the real-time log session configuration file, refer to Section 7.1.1, Log Session Configuration File. Real-time log session configuration files can also be opened by selecting the Open button on the Log Session Settings dialog.


LogWare then reads the log session settings from the configuration file and verifies that the configuration of the currently connected thermometer readout is compatible with the log session settings. If there are any conflicts, or if LogWare is unable to communicate with the thermometer readout, appropriate messages are displayed.

Once the configuration file has been read in, the Log Session Settings dialog is displayed. If LogWare was unable to communicate with the thermometer readout or any other error occurred, the log session settings may be incomplete.

The Open option and toolbar button are disabled while a real-time log session window is open.

5.1.3 Save as

The Save As option allows the current real-time log session settings to be saved to a configuration file. If the same basic configuration is frequently used to perform real-time logging, saving the log session settings to a configuration file can greatly reduce the time it takes to configure and start a real-time log session. Before attempting to save a real-time log session configuration, use the Log Session Settings dialog to configure the log session.

The  button on the toolbar corresponds to the Save As option.

When selecting the Save As option, the Windows® Save As dialog is displayed. The \Config folder should be selected by default (or the folder that is specified for configuration files on the LogWare Options dialog). Real-time log session configuration files are saved with a **.rls** extension. To cancel saving a configuration file, click the Cancel button. Otherwise, enter the filename for the configuration file and click the Save button. For more details about the real-time log session configuration file, refer to Section 7.1.1, Log Session Configuration File. Real-time log session configuration files can also be saved by selecting the Save As button on the Log Session Settings dialog.

The Save As option and toolbar button are disabled while a real-time log session window is open.

5.1.4 Print

The Print option allows the current graph or log file to be printed. The Print option has 2 submenu options: Readings and Graph. The submenu options are disabled if there is no real-time or previous log session window open.

5.1.4.1 Readings

The Readings option allows the output file for the currently selected channel to be printed. This option applies to both real-time and previous log sessions.

5.1.4.2 Graph

The Graph option allows the graph in the currently selected window to be printed. This option applies to both real-time and previous log sessions.

The graph can also be printed by selecting the Print Graph option in the Graph menu.

5.1.5 Exit

The Exit option closes the LogWare Software.

The Exit option is disabled while a real-time log session window is open.

5.2 Datalog menu

The Datalog menu provides options for working with configuring, starting, stopping and closing a real-time log session.

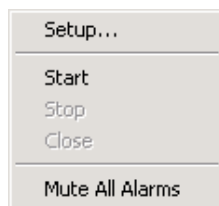



Figure 138 Datalog menu

5.2.1 Setup

The Setup option displays the Log Session Settings dialog. This dialog is used to configure a real-time log session. If a log session has already been configured or a real-time log session configuration file has already been opened, this dialog displays the current log session settings.


The  button on the toolbar corresponds to the Setup option.

Please refer to Section 2.3, Log Session Settings Dialog, for more details on configuring a real-time log session.

This option is disabled while a real-time log session window is open.

5.2.2 Start

The Start option starts a real-time log session. This option is disabled until a log session has been completely configured.

The  button on the toolbar corresponds to the Start option.

Please refer to Section 2.4, Real-Time Log Session, for more details on a real-time log session.

5.2.3 Stop


The Stop option stops the current real-time log session. This option is only enabled when a real-time log session is running. A log session may be stopped at any time by selecting this option.

The  button on the toolbar corresponds to the Stop option.

When selecting the Stop option, a message is displayed to confirm stopping the log session.

5.2.4 Close


The Close option closes the current real-time log session window. This option is only enabled once the real-time log session has stopped and the log session window is still open.

The  button on the toolbar corresponds to the Close option.

5.2.5 Mute all alarms

The Mute All Alarms option can be used to mute any audible alarms associated with the current real-time log session. When the alarms are muted, a check mark appears next to this option. Muting the alarms does not prevent alarm events from happening.

The  button on the toolbar corresponds to the Mute All Alarms option.

When alarms are muted, the toolbar button appears as  .

5.3 Download menu

The Download menu contains options for downloading logged data from the thermometer readout's memory or logs and viewing the current status of the thermometer readout's logs or memory.

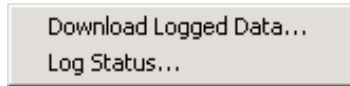


Figure 139 Download Menu

5.3.1 Download logged data

The Download Logged Data option can be used to download data that has been logged to the thermometer readout's logs or memory. All thermometer readouts supported by LogWare support logging of some type. Please refer to Section 1.4.4, Firmware Requirements, for more information.

When selecting the Download Logged Data option, the Select Model dialog is displayed. LogWare needs to know the model of the thermometer readout in order to display appropriate options dialog. Please refer to Section 3.2, Select Model Dialog, for more details on the Select Model dialog.

After the thermometer readout model and communication settings have been selected, the appropriate Download Options dialog for the selected thermometer readout is displayed. The Download Options dialogs are described in more detail in Section 3.3, Download Options Dialogs.

This option is disabled while a real-time log session is running.

5.3.2 Log status

The Log Status option can be used to view the current status of the thermometer readout's logs or memory. This feature is only supported by some of the thermometer readouts supported by LogWare.

When selecting the Log Status option, the Select Model dialog is displayed. LogWare needs to know the model of the thermometer readout in order to display appropriate status dialog. Please refer to Section 3.2, Select Model Dialog, for more details on the Select Model dialog.

After the thermometer readout model and communication settings have been selected, the appropriate Download Options dialog for the selected thermometer readout is displayed. The Log Status dialogs are described in more detail in Section 3.4, Log Status Dialogs.

This option is disabled while a real-time log session is running.

5.4 View menu

The View menu provides options for showing or hiding the toolbar and status bar, viewing current alarm settings, showing or hiding graph traces, viewing output and alarm log files, and for opening download files or output files from previous log sessions.

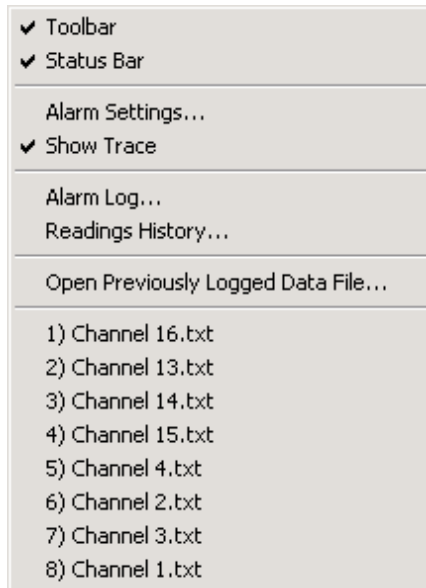


Figure 140 View Menu

5.4.1 Toolbar

The Toolbar option toggles the display of the toolbar at the top of the main window. A check mark appears next to this option if the toolbars are currently displayed.

5.4.2 Status bar

The Status bar option toggles the display of the status bar at the bottom of the main window. A check mark appears next to this option if the status bar is currently displayed.

5.4.3 Alarm settings

The Alarm Settings option displays the current alarm settings for the selected channel. This option only applies to channels in a real-time log session.

If no alarms were configured for the selected channel, this option is disabled.

5.4.4 Show trace

The Show Trace option toggles displaying the trace on the graph for the selected channel. This option applies to both real-time and previous log session windows. A check mark appears next to this option if the trace for the selected channel is currently displayed.

This option is only enabled when a real-time or previous log session window is displayed.

5.4.5 Alarm log

The Alarm Log option displays the alarm log for the current real-time log session. Please refer to Section 7.2.3, Alarm Log File, for more details on alarm log files.

This option is only enabled while a real-time log session window is open.

5.4.6 Readings history

The Readings History option displays the output file for the selected channel of a real-time log session. Please refer to Section 7.2.1, Output File, for more details on output files.

This option is only enabled while a real-time log session window is open.

5.4.7 Open previously logged data file

The Open Previously Logged Data File option allows file(s) containing logged or downloaded data to be opened. Please refer to Section 2.5, Previous Log Session, for more information on opening previous log session files.

When selecting the Open Previously Logged Data File option, the Windows® Open dialog is displayed. The \Config folder should be selected by default (or the folder that is specified for configuration files on the LogWare Options dialog). Both output files and download files can be opened in the Previous Log Session window. These files should have a **.txt** extension. To cancel opening a file, click the Cancel button. Otherwise, select up to 10 files to open and click the Open button. If more than 10 files are selected, only the first 10 files are opened and the rest are ignored. For more details about output files, please refer to Section 7.2.1, Output File. For more details about download files, please refer to Section 7.2.2, Download File.

5.4.8 MRU list

The View menu MRU list indicates up to the last 8 output or download files that were opened in a Previous Log Session window. Click on an item to open that file in a new Previous Log Session window.

5.5 Graph menu

The Graph menu provides options for customizing graph settings, changing the graph trace colors, accessing graph zooming features and printing the graph.

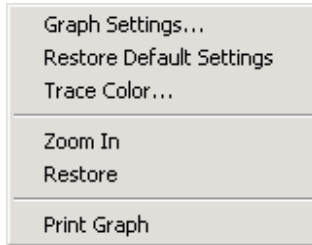



Figure 141 Graph Menu

5.5.1 Graph settings dialog

The Graph Settings option displays the Graph Settings dialog.

The  button on the toolbar corresponds to the Graph Settings option. The Graph Settings dialog allows many of the graph settings to be changed including titles, axis settings, grids, colors, trend lines and legends. The Graph Settings option applies to the current log session window.

The Graph Settings option and toolbar button are only enabled if a real-time or previous log session window is open.

The following sections describe in detail each of the customizable graph settings.

After modifying any of the settings on this dialog, click the Apply Now or OK button to apply the new settings to the graph. If the Apply Now button is clicked, this dialog remains open. If the OK button is clicked, this dialog closes.

To close this dialog without saving new settings, click the Cancel button.

Click the Help button to open the appropriate help topic for this dialog.

5.5.1.1 Titles tab

The Titles tab can be used to enter the titles to appear on the top, bottom, left and right sides of the graph.

The screenshot shows the 'Titles' tab of the Graph Settings dialog. It features a tabbed interface with 'Titles' selected. The 'Titles' section includes four text input fields: 'Graph Title' (empty), 'Bottom Title' (containing 'Time'), 'Left Title' (containing '°C (or value)'), and 'Right Title' (empty). Below the 'Left Title' field are three radio buttons: 'Horizontal', 'Up' (selected), and 'Down'. Below the 'Right Title' field are three radio buttons: 'Horizontal' (selected), 'Up', and 'Down'.

Figure 142 Graph Settings - Titles tab

When you enter text for a title, the Graph control adjusts the rest of the graphing window to provide space—either redrawing the graph and associated objects at a smaller size or decreasing the space between objects. When you clear the text box for a title, you disable it and provide more space for the rest of the graph.

If you enter a title that's too long to appear in a single line, the Graph control automatically word-wraps it. If a title doesn't display at all, it's because the Graph control can't make the font small enough to fit all the text in the space provided. Increase the size of the graphing window to make the title appear.



Note: To restore the titles to their default settings, select the *Restore Default Settings* option in the Graph menu.

Enter the top title for the graph in the Graph Title box. The graph title can contain up to 80 characters. This title appears centered at the top of the graph. By default, the graph title is blank.

Enter the bottom title for the graph in the Bottom Title box. The bottom title can contain up to 80 characters. This title appears centered at the bottom of the

graph and is generally used to denote the scale of the X-axis. By default, this title is set to "Time".

Enter the title for the left Y-axis of the graph in the Left Title box. The left title can contain up to 80 characters. This title appears centered vertically on the left side of the graph and is generally used to denote the scale of the Y-axis. The title can be displayed horizontally or vertically running up or down. By default, this title is set to the Readings Scale with the Up orientation option selected.

Enter the title for the right Y-axis of the graph in the Right Title box. The right title can contain up to 80 characters. This title appears centered vertically on the right side of the graph and is generally used to denote the scale of a secondary Y-axis. The title can be displayed horizontally or vertically running up or down. By default, the right title is blank.



Note: *LogWare only supports on Y-axis, on the left side of the graph. However, a title can still be placed on the right side of the graph.*

5.5.1.2 Axis tab

The Axis tab can be used to customize the graph's X- and Y-axis settings, colors, ticks and gridlines.

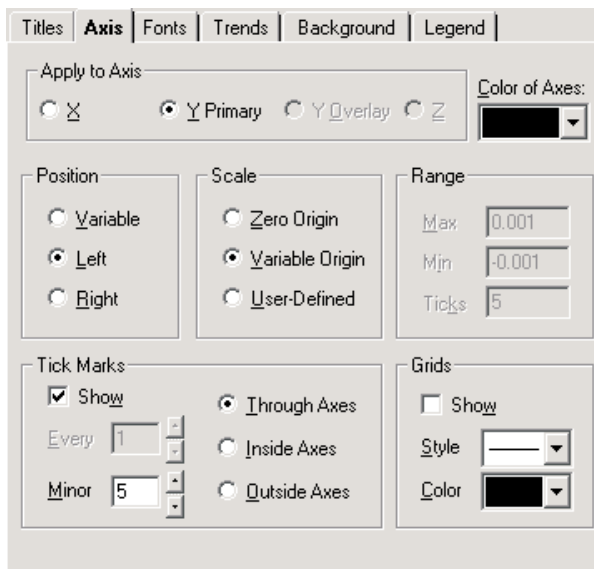


Figure 143 Graph Settings - Axis tab



Note: To restore the axes to their default settings, select the Restore Default Settings option in the Graph menu.

Select the axis for which changes are to be made using the Apply to Axis option. The other controls on this tab change to reflect the current settings of the selected axis. By default, the Y Overlay and Z axes options do not apply to the graph and are not enabled.

Select the color for the X-and Y-axes by selecting a color bar from the Color of Axes drop-down list. This setting applies to both axes and is therefore independent of the current Apply to Axis setting. By default, this option is set to black.

The Position options allow the X-axis to be positioned on either the top or bottom of the graph and the Y-axis to be positioned on either the left or right side of the graph. Selecting the Variable option allows the graph itself to determine whether to position the axes on the top, bottom, left or right side.

By default, the X-axis is set to Bottom and the Y-axis is set to Left.

The Scale options allow the user to determine the minimum and maximum values for an axis. Each option is described below:

- **Zero Origin** - This option draws the selected axis including the origin (zero) and extending far enough in the positive and negative directions to include all of the graph's values
- **Variable Origin** - Select this option to have the axis "zoom in" on the range of the graph's actual values, whether or not that range includes zero.
- **User-Defined** - This option allows the user to set values for the minimum, maximum, and number of ticks for the axis. When this option is selected, the settings in the Range group are enabled

By default, the X-axis and Y-axis are both set to Variable Origin.



*Note: Changing the Scale settings (especially for the X-axis) while acquiring real-time data could possibly cause graphing problems to occur. Therefore, it is recommended that these settings **NOT** be changed by the user when acquiring real-time data.*

The Range options are only enabled if the User-Defined option is selected in the Scale options for the selected axis. Enter the Maximum and Minimum values for the selected axis. The selected axis is rescaled to the settings entered.

The Ticks option allows the user to determine how many major tick marks should be displayed between the minimum and maximum values on the selected axis. Each major tick mark is labeled with the appropriate value. The tick marks themselves are only displayed if the Show check box in the Tick Marks controls is checked.

By default, these settings are not used by the software because of the default Scale settings.

The Tick Marks options allow the user to specify whether or not to show tick marks on the selected axis. To show tick marks, check the Show check box. If this check box is not checked, the tick marks are not shown on the selected axis. Valid settings are between 1 and 1,000.

The Every option is only enabled for the X-axis. This option determines how often major tick marks appear on the X-axis. When acquiring real-time data, the graph displays the most recent 600 readings acquired. The Every option is set to 150 meaning that a major tick mark appears every 150 readings, or $600 \div 150 = 4$ major tick marks appear on the X-axis. When analyzing previously acquired data, the Every option is calculated by the software depending on the number of readings being displayed. The X-axis must end with a tick mark. If you set an Every value that doesn't include the last value on the axis, the graph will extend the X-axis so that it ends on a tick mark.

The Minor option allows the user to determine how many minor tick marks appear between each major tick mark. By default, the X-axis and Y-axis are set to display 5 minor tick marks between major tick marks. Valid settings are between 0 and 100.

The Through Axes, Inside Axes and Outside Axes options allow the user to specify how the tick marks should be drawn on the axis. By default, both the X-axis and Y-axis are set to Through Axes.

The Grids options allow the user to specify whether or not to display grid lines on the graph. To show the grid lines for the selected axis, check the Show check box. Select the list Style and Color for the grid lines. Both axes are set to the same style and color, but each axis can be set to show or hide grid lines independent of the other.

By default, grid lines are not shown on the graph.

5.5.1.3 Fonts tab

The Fonts tab can be used to change the typefaces and sizes of fonts used on the graph.

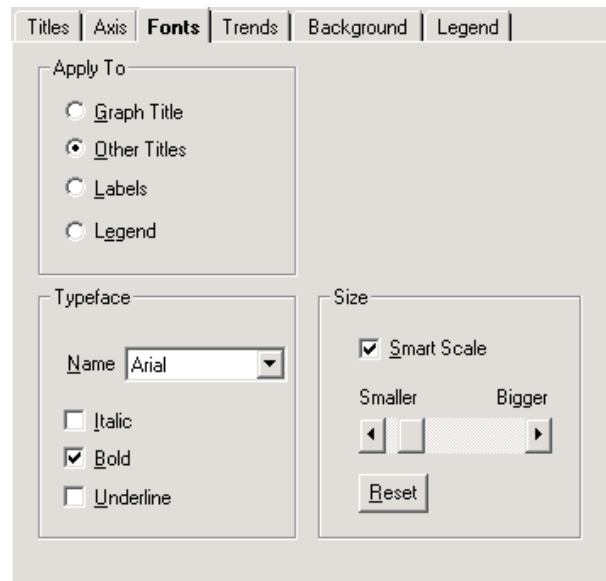


Figure 144 Graph Settings - Fonts tab



Note: To restore the fonts to their default settings, select the Restore Default Settings option in the Graph menu.

Select the title, label or legend option to which the font settings are to be applied using the Apply To option. The Typeface and Size controls display the current settings for the selected option.

Select the font name from the Typeface drop-down list and choose the styles to apply to the selected title, label or legend.

Select the font size to apply to the selected title, label, or legend by clicking on the scroll bar or dragging the scroll bar's thumb. Selecting the Smart Scale option allows the graph to automatically reduce the size of the selected font if the dimensions of the graph become too small to display the title, label or legend in the selected font size. Click the Reset button to restore the size to its default setting.

5.5.1.4 Trends tab

The Trends tab can be used to superimpose trend lines (Min, Max, Average, etc.), curve fit and limit (alarm) lines on the graph.

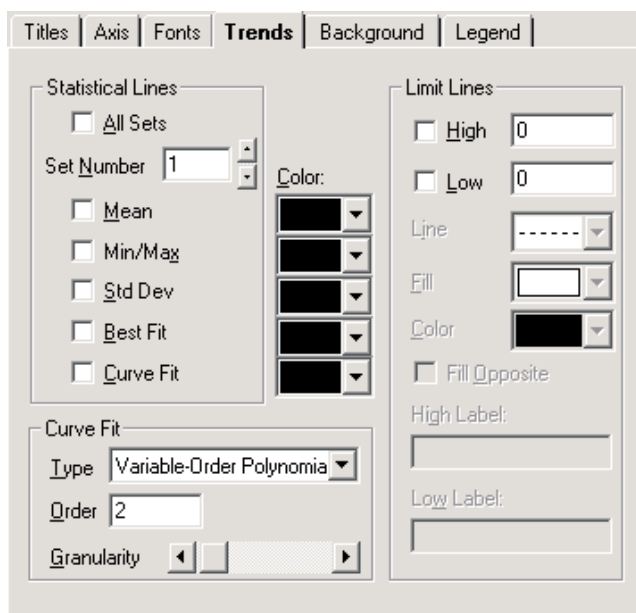


Figure 145 Graph Settings - Trends tab



Note: To restore the trend settings to their default settings, select the Restore Default Settings option in the Graph menu.

The Statistical Lines options allow the user to specify which statistical lines should be displayed on the graph.

Check the All Sets option to apply the statistical lines settings to all of the sets (traces) on the graph. When this check box is checked, the Set Number controls are disabled.

The Set Number box allows the user to select which set (trace) to which to apply the statistical lines settings. This setting can only be changed if the All Sets check box is not checked.



Note: *This software only has one set of data. When acquiring real-time data, this set is the data being acquired. When analyzing previously acquired data, this set is the data which was previously acquired.*

Check the Mean check box to enable a mean line, which is drawn horizontally through the average value of all data points in the selected set.

Check the Min/Max check box to enable min and max lines, which are drawn horizontally through the lowest and highest data points in the selected set.

Check the Std Dev check box to enable a pair of standard-deviation lines, which are drawn horizontally through the standard deviation from the mean (in both the positive and negative directions) for the selected set.

Check the Best Fit check box to enable a best-fit line, which is a straight line indicating the trend of data points (a first-order polynomial curve) for the selected set.

Check the Curve Fit check box if you want to fit a curve through your data points for the selected set. You can choose the type, order and granularity of the curve in the Curve Fit options below. These settings will apply to all curve fit lines drawn on the graph.

By default, no trend lines are enabled.

Select the Curve Fit options for the curve fit statistical line. The following options are available:

Curve type	Description/formula
Variable-Order Polynomial	Polynomial curve of variable order
Logarithmic	$y = a + b * \log(x)$
Exponential 1	$y = a * \exp(b * x)$
Exponential 2	$y = a * x * \exp(-b * x)$
Power	$y = a * (x ^ b)$
Inverse 1	$y = a + b / x$
Inverse 2	$y = a / (b + x)$
Inverse 3	$y = 1 / (a + b * x)$
Inverse 4	$y = x / (a * x + b)$
Inverse 5	$y = 1 / (a + b * x) ^ 2$
Spline	Spline fit through all points
Moving Average Mid	Moving average plotted at midpoint of averaged group
Moving Average End	Moving average plotted at end point of averaged group

The Order option applies only to three curve types:

- For variable-order polynomial curves, Order is the order of the polynomial used in curve fitting. A setting of 1 produces a straight line (the same as a best-fit line); a setting one less than the number of points produces a curve that passes through every point.
- For moving-average mid and moving-average end curves, Order is the range of data points over which moving averages are averaged, beginning with the first point.

The Granularity scroll bar sets the granularity of all curve types except moving-average. The granularity is the number of “steps,” or straight line segments, making up the curve. Higher values create smoother curves, but require more drawing time. The default Granularity setting is 50 curve steps, which generally creates a smooth-looking curve at a high drawing speed. Each click to the left decreases the number of steps by 2 (to a minimum of 10), and each click to the right increases the number of steps by 2 (to a maximum of 1000). With spline curves, you generally need higher granularities than normal—up to 10 times the number of points in the graph.

Select the colors for each of the statistical lines to display on the graph using the Color drop-down lists. The selected color for each statistical line applies to that statistical line for all sets of data.

Select the limit (alarm) lines options to display on the graph using the Limit Lines options. To display the high alarm line, check the High check box. To display the low alarm line, check the Low check box.



Note: Changing the low and high limit line settings on this dialog **DOES NOT** change the low and high alarm settings entered on the Alarm Settings dialog and therefore does not affect the alarm events options. These settings are only used to display the alarm lines on the graph. Only change these settings if absolutely necessary.

Select the line style for the limit (alarm) lines using the Line drop-down list. A fill pattern can also be selected using the Fill drop-down list. Select the color of the limit (alarm) lines and the fill pattern using the Color drop-down list.

Select the Fill Opposite check box to fill the area between the High and Low limit (alarm) lines with the selected pattern. If this option is not checked, the areas above the High limit value and below the Low limit value are filled in with the selected pattern.

Enter the text to appear as a label on the High and Low limit (alarm) lines in the High Label and Low Label boxes. By default, these labels are set to “High Alarm” and “Low Alarm” respectively.



Note: Unlike statistical lines, the graph can only display one set of limit (alarm) lines for all sets of data (traces).

5.5.1.5 Background tab

The Background tab can be used to set the background colors of titles, the legend and the graph itself.

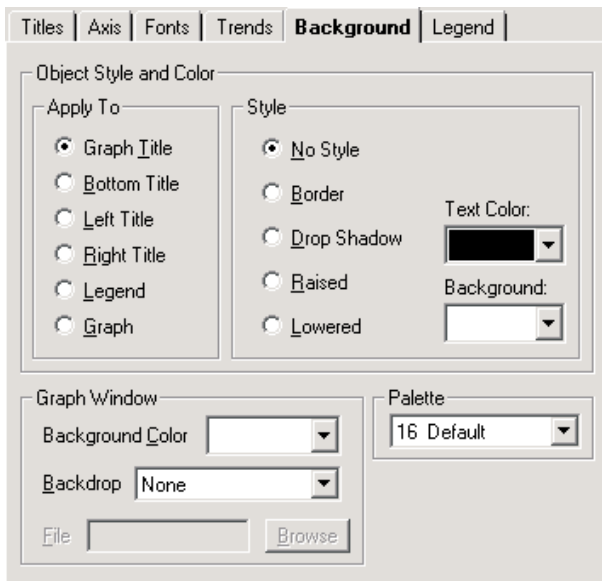


Figure 146 Graph Settings - Background tab



Note: To restore the background settings to their default settings, select the Restore Default Settings option in the Graph menu.

Select the title, legend, or graph option to which the background settings are to be applied using the Apply To options. The other controls on this tab change to show the current settings for the selected item.

Select the background style options to apply to the selected title, legend or graph using the Style options. Select the text and background color for the selected item using the drop-down lists provided.

Select the background options for the graph window using the Graph Window options. Use the Background Color drop-down list to select the background color for the plotting area of the graph.

The Backdrop option can be used to display a graphic (.BMP or .WMF) file such as a company logo or texture pattern behind the graph. If any other option except "None" is selected, the File and Browse button controls are enabled. Enter the name of the file or browse for the filename using the Browse button.

Select the color palette to use using the Palette drop-down list. Using this drop-down list, choose a 16-color or 128-color palette for the screen. All of the Color drop-down lists for the graph settings use this palette to determine the available choices.

- 16-color palettes consist of 16 differentiated colors—default, pastel, or grayscale
- 128-color palettes always consist of the 16 colors from the default palette, followed by 96 colors that vary according to the palette you select. The graph reserves 16 additional colors for special graphic needs. These colors cannot be selected

By default, the “16 Default” palette is selected.



Note: Selecting any other palette will invalidate the Trace Color option on the Log Session Settings dialog. The Palette option should NOT be changed.

5.5.1.6 Legend tab

The Legend tab can be used to add a legend to the graph.

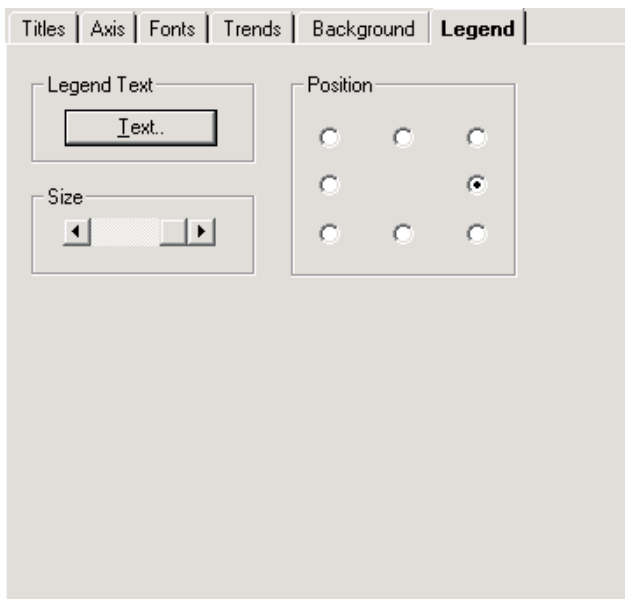


Figure 147 Graph Settings - Legend tab



Note: To restore the legend settings to their default settings, select the Restore Default Settings option in the Graph menu.

Click the Text button to enter the labels for each set (trace) of data on the graph. By entering text into the boxes provided, the legend will automatically appear on the graph. To hide the legend, remove the text from all of the boxes.

By default, all of the legend text boxes are empty causing the legend to be hidden.

Select the position with respect to the graph to place the legend using the Position options.

Select the size of the legend using the Size options. To change the font or style that the legend uses, see the Fonts tab. To set the legend colors, see the Background tab.

5.5.2 Restore default settings

The Restore Default Settings option restores the graph to its default settings and applies to the current log session window.

The Restore Default Settings option and toolbar button are only enabled if a real-time or previous log session window is open.


5.5.3 Trace color

The Trace Color option allows the trace color for the selected channel to be changed. Please refer to Section 2.4, Real-Time Log Session, for more details on changing the trace color for real-time log sessions and Section 2.5, Previous Log Session, for more details on changing the trace color for previous log sessions.

The Trace Color option is only enabled if a real-time or previous log session window is open.

5.5.4 Zoom in

The Zoom In option allows the graph to be zoomed in.

The  button on the toolbar corresponds to the Zoom In option. Please refer to Section 2.5.4, Graph Zooming, for more details on using this feature.

The Zoom In option and toolbar button are only enabled if a previous log session window is open.

5.5.5 Restore

The Restore option restores the graph to its default zoom state.

The  button on the toolbar corresponds to the Restore option. Please refer to Section 2.5.4, Graph Zooming, for more details on using this feature.

The Restore option and toolbar button are only enabled if a previous log session window is open.

5.5.6 Print graph

The Print Graph option prints the current graph on the selected printer.

The graph can also be printed by selecting the Print | Graph option in the File menu.

5.6 Tools menu

The Tools menu provides options for synchronizing the thermometer readout's clock, viewing and/or modifying logging labels, viewing and/or modifying thermometer readout settings and configuring LogWare's options.

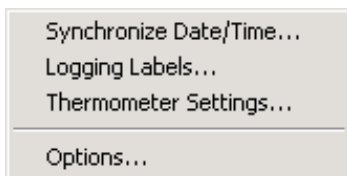



Figure 148 Tools Menu

5.6.1 Synchronize date/time

The Synchronize Date/Time option provides a utility to synchronize the thermometer readout's clock to the computer's clock.

The  button on the toolbar corresponds to the Synchronize Date/Time option. Please refer to Section 4.1.1, Synchronize Date/Time, for a description of this feature.

The Synchronize Date/Time option and toolbar button are disabled while a real-time log session is running.

5.6.2 Logging labels


The Logging Labels option provides a utility to view and/or modify logging labels.

The  button on the toolbar corresponds to the Logging Labels option. Please refer to Section 4.3, Logging Labels Dialogs, for a description of this feature.

The Logging Labels option and toolbar button are disabled while a real-time log session is running.

5.6.3 Thermometer settings

The Thermometer Settings option provides a utility to access the thermometer readout settings dialogs.

The  button on the toolbar corresponds to the Thermometer Settings option. Please refer to Section 4.4, Thermometer Settings Dialogs, for a description of these dialogs.

The Thermometer Settings option and toolbar button are disabled while a real-time log session is running.

5.6.4 Options

The Options option displays the LogWare Options dialog to view and/or modify LogWare's default settings.

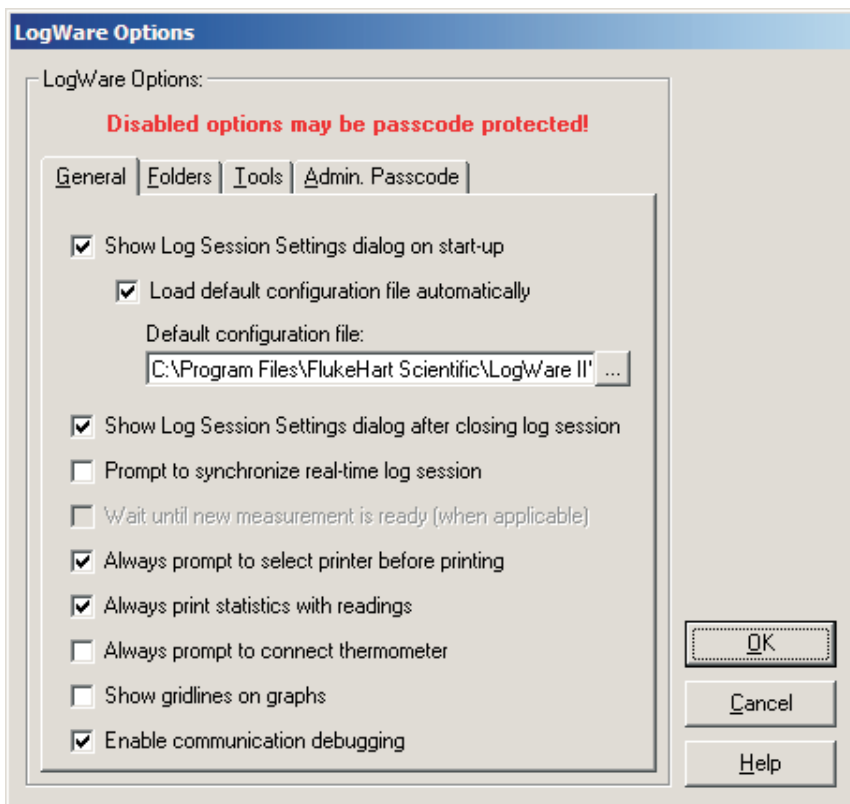


Figure 149 LogWare Options Dialog

The Options option is disabled while a real-time log session is running.

The following sections describe in detail each of the LogWare Options settings.

After modifying any of the settings on this dialog, click the OK button to save the new settings and close the dialog.

To close this dialog without saving new settings, click the Cancel button.

Click the Help button to open the appropriate help topic for this dialog.

5.6.4.1 General tab

The General tab displays options pertaining to the behavior of the LogWare software in general.

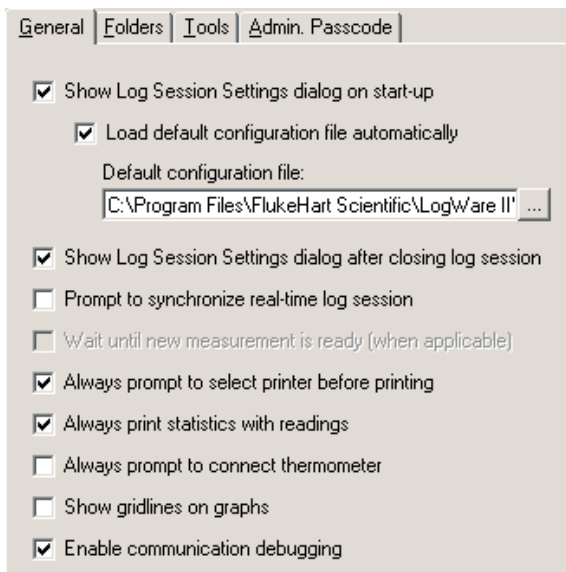


Figure 150 LogWare Options - General tab

The Show Log Session Settings dialog on start-up option causes LogWare to automatically display the Log Session Settings dialog every time LogWare starts up. If checked, the Log Session Settings dialog is displayed immediately after LogWare starts up. If not checked, LogWare does not display the Log Session Settings dialog on start-up and the other options on this tab are disabled.

The Load default configuration file automatically option allows LogWare to automatically load the specified configuration file every time LogWare starts up. This option is useful if the same hardware configuration is used repeatedly to perform real-time log sessions. When this option is checked, the configuration file to load must be specified.

The Default configuration file box indicates the configuration file that will be loaded automatically every time LogWare starts up. The configuration file will

only be loaded if the Show Log Session Settings dialog on start-up and Load default configuration file automatically options are both checked. Use the "...” (Browse) button to locate the configuration file to be used as the default configuration file.

The Show Log Session Settings dialog after closing log session option determines whether the Log Session Settings dialog is displayed after the real-time log session window is closed. All real-time log session settings are still active in LogWare after a log session is closed. To access the Log Session Settings dialog, select the Setup option in the Datalog menu. If this option is not checked, this dialog is not displayed when the log session window is closed.

The Prompt to synchronize real-time log session option determines whether or not a prompt is displayed when the real-time log session is initialized and ready to start acquiring data. This prompt can be used to facilitate synchronizing LogWare with another process or application. If this option is not checked, LogWare begins acquiring data immediately after initialization of the session is complete.



Note: *When using the delayed start feature of LogWare to start a log session unattended, make sure this setting is not checked! Otherwise, the log session may not start as expected!*

The Wait until new measurement is ready (when applicable) option determines whether LogWare should always wait until the thermometer readout indicates a new measurement is ready on the current channel before querying the measurement during a real-time log session. If this option is not checked, LogWare always queries the last measurement taken on the current channel.

Not all thermometers supported by LogWare support waiting until a new measurement is ready. For these thermometer readouts, LogWare always queries the last measurement taken on the current channel.



Note: *This option is protected by the administrator passcode. To change this setting, the correct administrator passcode must be entered. For more information, see Section 5.6.4.4, Admin. Passcode Tab.*

The Always prompt to select printer before printing option determines whether the Print Setup dialog is displayed when printing. The Print Setup dialog allows the printer to be selected as well as the paper size, source and orientation settings (as supported by the selected printer driver). If this option is not checked, LogWare always prints to the default printer using that printer's default settings.

The Always print statistics with readings option determines whether LogWare appends statistical data to the end of the readings when printing an output file. If this option is checked, the statistical data is printed.



***Note:** LogWare allows the readings history to be printed with or without statistics while a real-time log session is running. However, when printing the readings history during a real-time log session, be aware that the statistical values may not exactly match the printed readings, since the statistics may get updated after the readings have already been printed. For best results, only print the readings AFTER the real-time log session has stopped.*

The Always prompt to connect thermometer option determines whether LogWare displays a prompt before attempting to communicate with the thermometer readout. These prompts are intended to be a reminder to the user to connect the thermometer readout so that the software can communicate with it. If the thermometer readout is frequently disconnected from the computer, this option may be left checked. If the thermometer readout is always connected to the computer, these prompts may become bothersome. In this case, uncheck this option so that the prompts are not displayed.

The Show gridlines on graphs option determines whether gridlines are enabled on the real-time log session and previous log session graphs by default. The gridlines can be enabled or disabled for a specific graph at any time using the Grids controls on the Axis tab of the Graph Settings dialog. Please refer to Section 5.5.1.2, Axis Tab, for more details.

The Enable communication debugging option is used to enable debugging features in LogWare. This option should remain unchecked unless otherwise directed by a Customer Service representative.

5.6.4.2 Folders tab

The Folders tab displays options pertaining to the default folders to which *LogWare* saves various types of files.

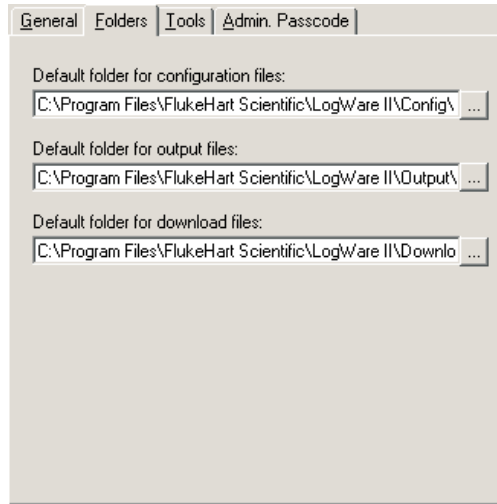


Figure 151 *LogWare Options - Folders tab*

The Default folder for configuration files option allows the default folder to which configuration files are saved to be specified. *LogWare* allows many types of configuration files to be saved including real-time log session (**.rls**), alarm (**.alm**) and conversion types (**.i90**, **.i68**, **.cvd**, **.ply**, **.hpt**, **.w90**, **.thr**, **.tcp**, **.tct**, **.tcx**). Use the “...” (Browse) button to select an existing path.

The Default folder for output files option allows the default folder to which output files are saved to be specified. Output files are the files created by a real-time log session and typically have the extension **.txt**. Use the “...” (Browse) button to select an existing path.

The Default folder for download files option allows the default folder to which download files are saved to be specified. Download files are the files created when data is downloaded from the memory or log(s) of a thermometer readout and typically have the extension **.txt**. Use the “...” (Browse) button to select an existing path.

5.6.4.3 Tools tab

The Tools tab displays options pertaining to the Tools menu options.

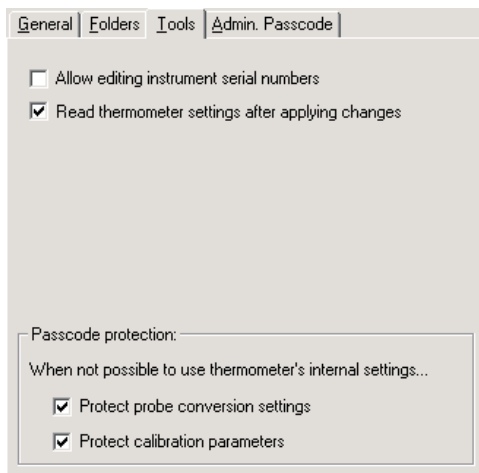


Figure 152 LogWare Options - Tools tab

The Allow editing instrument serial numbers option determines whether LogWare will allow the serial number of the thermometer readouts to be changed using the Thermometer Settings option in the Tools menu. If this option is checked, the Serial number boxes on the thermometer settings dialogs are enabled (where applicable) and the serial number of the thermometer readout can be changed.



Note: Many Hart Scientific thermometer readouts save the thermometer readout's serial number in its memory. However, many thermometer readouts still have their serial number set to 0. This option can be used to allow the thermometer readout's serial number to be set. Once the serial number is set, this option should always be left off to prevent the serial number from being changed to an incorrect setting.



Note: This option is protected by the administrator passcode. To change this setting, the correct administrator passcode must be entered. For more information, see Section 5.6.4.4, Admin. Passcode Tab.

The Read thermometer settings after applying changes option determines whether LogWare will query the thermometer readout for current settings after new settings have been applied. This option is used on the thermometer settings dialogs which are accessible from the Thermometer Settings option.

The Protect probe conversion settings option determines whether probe conversion settings (including coefficients) should be passcode protected if LogWare is unable to determine the protection settings of the thermometer readout or if the thermometer readout does not support passcode protection. The thermometer readout's protection settings are always used where possible.



Note: This option is protected by the administrator passcode. To change this setting, the correct administrator passcode must be entered. For more information, see Section 5.6.4.4, Admin. Passcode Tab.

The Protect thermometer calibration parameters option determines whether thermometer readout calibration parameters should be passcode protected if LogWare is unable to determine the protection settings of the thermometer readout or if the thermometer readout does not support passcode protection. The thermometer readout's protection settings are always used where possible.



Note: This option is protected by the administrator passcode. To change this setting, the correct administrator passcode must be entered. For more information, see Section 5.6.4.4, Admin. Passcode Tab.

5.6.4.4 Admin. Passcode tab

The Admin. Passcode tab allows options pertaining to the administrator passcode to be selected.

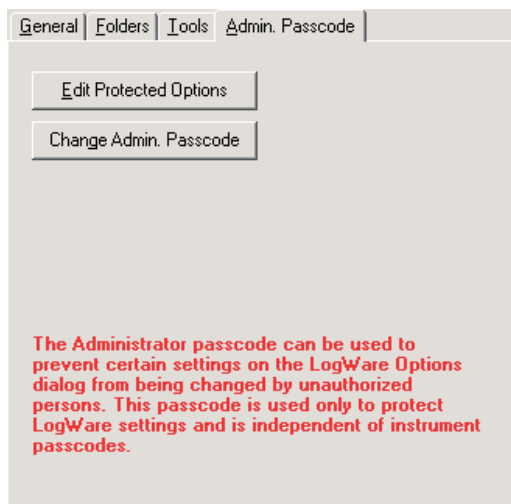


Figure 153 LogWare Options - Admin. Passcode tab

The Edit Protected Options button allows the user to change the settings of the passcode protected options. This button is disabled if an administrator passcode

has not been setup or if the correct passcode has been entered since the LogWare Options dialog was displayed.

Clicking the Edit Protected Options button displays the Enter Passcode dialog. The current administrator passcode must be entered to enable changing the protected options.

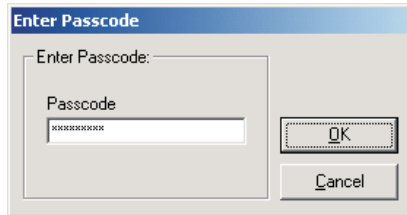


Figure 154 Enter Passcode dialog

Enter the current administrator passcode and click the OK button to continue. If the passcode entered does not match the current administrator passcode, a message is displayed and access is denied. Otherwise, if the correct passcode is entered, the Edit Protected Options button becomes disabled and all passcode protected options on the LogWare Options dialog become enabled. These options remain enabled until the LogWare Options dialog is closed or until the user attempts to change the administrator passcode.

The Change Admin. Passcode button allows the administrator passcode to be setup or changed.

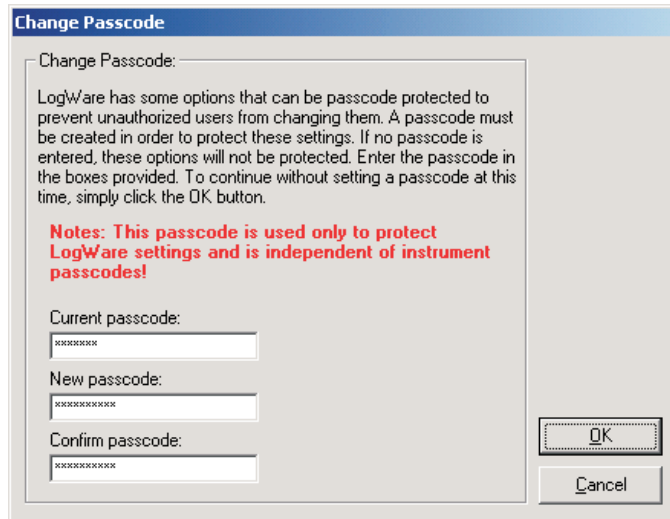


Figure 155 Change Passcode dialog



Note: The administrator passcode in LogWare is completely independent of the thermometer readout passcodes. The administrator passcode is used by LogWare to protect certain LogWare options and default settings on the LogWare Options dialog only! Thermometer readout passcodes are specific to the thermometer readout being used and are required to change certain settings in the thermometer readout itself.

If an administrator passcode has already been setup, the current passcode is required to gain access to the Change Passcode dialog and clicking the Change Admin. Passcode button displays the Enter Passcode dialog first. The current administrator passcode must be entered to enable changing the administrator passcode. Enter the current administrator passcode and click the OK button to continue. If the passcode entered does not match the current administrator passcode, a message is displayed and access is denied. Otherwise, if the correct passcode is entered, the Change Passcode dialog is displayed.

To change the current administrator passcode, enter the new passcode in the New passcode box. Then confirm the new passcode by entering the passcode again in the Confirm passcode box. Click the OK button to save the new passcode. To disable passcode protection, simply leave the New passcode and Confirm passcode boxes blank and click the OK button. To cancel changing the current passcode, click the Cancel button.

5.7 Window menu

The Window menu provides options for arranging the open log session windows and for switching between open log session windows.

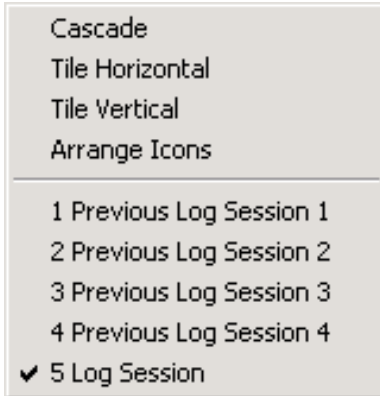


Figure 156 Window menu

5.7.1 Cascade

The Cascade option cascades all open windows.

5.7.2 Tile horizontal

The Tile Horizontal option tiles all open windows horizontally.

5.7.3 Tile vertical

The Tile Vertical option tiles all open windows vertically.

5.7.4 Arrange icons

The Arrange Icons option arranges the icons for all minimized windows along the bottom of LogWare's main display.

5.7.5 Window list

This is a list of all currently open windows. To switch to a window, click on the appropriate item in the window list. A check mark appears next to the currently selected window.

5.8 Help menu

The Help menu provides options for opening the LogWare's online help and for displaying information about LogWare.

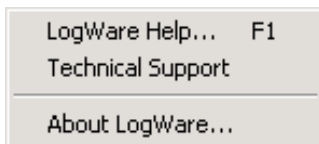


Figure 157 Help menu

5.8.1 LogWare help

The LogWare Help option opens the LogWare help file.

5.8.2 Technical support

The Technical Support option opens the LogWare help file to the Technical Support topic. For technical support information, please refer to Section 10.1, Technical Support.

5.8.3 About LogWare

The About LogWare option displays the About LogWare dialog.

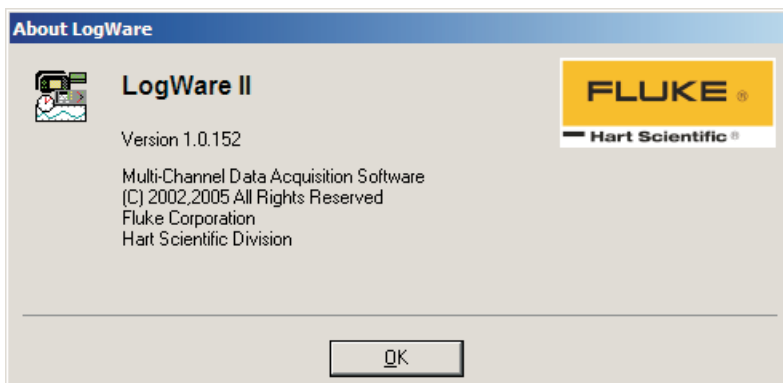


Figure 158 About LogWare dialog

The About LogWare dialog displays specific information about the currently installed version of LogWare. The software's name, version, build, and copyright information are displayed.


6 How to

The following sections describe in detail how to perform certain tasks using *LogWare*.

6.1 Setting up and starting a real-time log session

The main purpose of *LogWare* is to perform real-time data acquisition. In order to do this, *LogWare* must be configured to acquire data on the channels of the thermometer readout to be used. *LogWare* needs to know what data to acquire, when to acquire it and when to stop.

To configure a real-time log session:

1. Select the New option in the File menu.
2. This will clear any current log session settings and display the Log Session Settings dialog.
3. Configure the Thermometer, Log Session and Channels settings according to the needs of this session.
4. Once configured, either click the Start button on the Log Session Settings dialog or click the Close button, then select the Start option in the Datalog menu or click the  button on the toolbar. This will start the real-time log session and the Log Session window is displayed.
5. If a delayed start time was selected, the real-time log session will begin at the selected time. Otherwise, the log session will start immediately.

Please refer to Section 2, Real-Time Logging, for more information.

6.2 Starting a real-time log session unattended (delayed start)

The main purpose of *LogWare* is to perform real-time data acquisition. Many times, the technician is present when a log session starts. However, sometimes it is necessary or desirable to start a real-time log session at off hours when the technician is out.

By default, a real-time log session starts immediately when the Start button on the Log Session Settings dialog is clicked or when the Start option in the Datalog menu is selected. However, *LogWare* has a feature called the Delayed Start which allows a real-time log session to start at a predefined date and time.

To set the delayed start date and time:

1. Select the Setup option in the Datalog menu.
2. The Log Session Settings dialog is displayed.

3. Select the Log Session tab.
4. Click the Delayed Start button.
5. The Delayed Start Date/Time dialog is displayed.
6. Choose the date and time that the real-time log session should start. Since the computer's clock will be used to determine when the log session starts, make sure the computer's clock is accurate. Click the OK button to accept the selected date and time.



Note: To cause the real-time log session to start immediately when clicking the Start button after having set a delayed start date and time, click the Reset button.

7. Click the Start button on the Log Session Settings dialog to start the log session.
8. LogWare displays a message indicating the date and time the real-time log session will start.
9. To cancel the delayed start, click the Cancel button at any time.

After a real-time log session ends, the delayed start time is reset so that the next real-time log session will start immediately when the Start button on the Log Session Settings dialog is clicked or when the Start option in the Datalog menu is selected. To set a new delayed start time for the next real-time log session, follow the steps above.



Note: Every effort has been made to make the real-time log session take its first reading at the selected date and time. However, LogWare may or may not be able to do so depending on the processor speed of the computer, amount of available memory, free hard disk space, other running applications, etc.

Please refer to Section 2, Real-Time Logging, for more information.

6.3 Stopping a real-time log session automatically

The main purpose of LogWare is to perform real-time data acquisition. Many times, the technician is present when a log session needs to be stopped. However, sometimes it is necessary or desirable to stop a real-time log session when the technician is busy with another task, or at off hours when the technician is out.

The End Log feature of LogWare provides three methods for stopping a real-time log session:

- By user
- After specified number of readings

- After specified time elapses

The first method, By user, is the default setting.

To stop the real-time log session automatically:

1. Select the Setup option in the Datalog menu
2. The Log Session Settings dialog is displayed.
3. Select the Log Session tab.
4. Select the method to use to stop the real-time log session using the End log drop-down list:
 - By user – Log session stops when the Stop option in the Datalog menu is selected.
 - After specified number of readings – Log session stops automatically when number of readings specified have been acquired on each channel.
 - After specified time elapses – Log session stops automatically after the specified time has elapsed.
5. Specify the End Log criteria:
 - If the After specified number of readings option was selected, enter the number of readings to be acquired on each channel in the After box.
 - If the After specified time elapses option was selected, enter the amount of time to elapse in the After box.
6. Click the Start button on the Log Session Settings dialog to start the log session.
7. When the selected end log criteria are met, logging will stop automatically.

Please refer to Section 2, Real-Time Logging, for more information.

6.4 Opening a configuration file automatically on start-up

LogWare is designed to be flexible in the thermometer readouts that it can use and the features it supports. However, most users of LogWare will probably always use the same thermometer readout, scale and communication settings to do all or most of their data acquisition.

For this reason, LogWare allows the real-time log session settings to be saved to a configuration file. Opening the configuration file restores the settings, therefore eliminating the need to manually configure LogWare every time a new log session is started.

Configuration files save the following information about a real-time log session:

- Thermometer readout model, serial number, session temperature scale and resolution settings
- Communications COM port and baud rate settings
- Interval between readings and interval between scan sequences
- Start log session and End log session settings
- Channel configurations including statistical and alarm settings

In addition to allowing configuration files to be saved and opened, LogWare optionally allows one configuration file to be opened automatically every time LogWare runs. This configuration file can be specified in one of two ways: using the LogWare Options dialog or using the Log Session Settings dialog.

To specify the default configuration file using the LogWare Options dialog:

1. Select the Options option in the Tools menu.
2. The LogWare Options dialog is displayed.
3. Select the General tab.
4. Check the Show Log Session Settings dialog on start-up check box.
5. Check the Open default configuration file automatically check box.
6. Click the “...” (Browse) button to select the configuration file to open or type the path and filename in the Default configuration file box.
7. Click the OK button to save the changes.



Note: *The real-time log session settings must already be saved to a configuration file in order to accomplish these steps.*

To specify the default configuration file using the Log Session Settings dialog:

1. Select the Setup option in the Datalog menu.
2. The Log Session Settings dialog is displayed.
3. Fill in all of the information on this dialog.
4. Click the Set as Default button. A prompt is displayed indicating that the current settings must be saved to a configuration file. Click the Yes button to save the current settings to a configuration file.
5. The Windows® Save As dialog is displayed to save the current settings to a configuration file. Enter the filename for this configuration file and click the OK button to save the settings to the file.

6. A prompt is then displayed indicating that this configuration file will be made the default configuration file. Click the Yes button to set this configuration file as the default. A message is displayed indicating the current configuration is now set as the default.
7. Click the Close button to close the Log Session Settings dialog.

Test this feature to make sure it works properly by closing LogWare and running it again. The thermometer readout must be connected and powered on when opening a configuration file so that LogWare can verify that the information in the configuration file matches the thermometer readout configuration. The Log Session Settings dialog should be displayed with all of the configuration information filled in.

To disable opening the configuration file on start-up, uncheck the Open default configuration file automatically check box on the LogWare Options dialog.

Please refer to Section 5.6.4, Options, for more information.

6.5 Changing thermometer and probe settings

LogWare provides a feature which allows many of the thermometer readout and probe settings to be read in, viewed, modified and written back out to the thermometer readout. These settings vary from thermometer readout to thermometer readout but include:

- Filtering (averaging)
- Resolution
- Calibration and recalibration dates
- Calibration constants
- Probe serial number
- Probe wires and excitation current
- Probe calibration and recalibration dates
- Probe conversion type and calibration coefficients



Note: *In order to use this feature of LogWare, the thermometer readout must have a recent firmware version. If your thermometer readout does not have a recent firmware version, it may be upgraded. For more details, contact Hart Scientific Technical Support. Please refer to Section 1.4.4, Firmware Requirements, for more details.*

To view, modify and write thermometer and probe settings:

1. Select the Thermometer Settings option in the Tools menu.
2. The Select Model dialog is displayed. Select the thermometer readout model. Also, select the COM port to which the thermometer readout is connected and the baud rate to use. Make sure LogWare can communi-

cate with the thermometer readout by clicking the Test Communication button. Click the OK button when communication is successful.

3. The thermometer readout settings are then read in and displayed in the thermometer settings dialog. The thermometer settings dialog is specific to the selected thermometer readout.
4. If applicable, to view the settings for a specific module, select the module from the Modules list and click the Module Settings button.
5. To view the settings for a specific channel, select the channel from the Channels list and click the Channel Settings button.
6. The settings can be modified at this time. Some settings may be passcode protected. Passcode protected settings can be viewed, but can only be modified after entering the correct passcode.
 - a. Click the Edit button.
 - b. The Enter Passcode dialog is displayed. Enter the passcode for the connected thermometer readout and click the OK button.
 - c. If the correct passcode was entered, the passcode protected settings are enabled and may be modified.
7. To write the changes that have been made to the thermometer readout, click the OK button on any and/or all dialogs.
8. To close the dialogs without writing the new settings to the thermometer readout, click the Cancel button on any of the dialogs.



***Note:** The passcode for the thermometer readout is completely independent of the administrator passcode in LogWare. The administrator passcode is used by LogWare to protect certain LogWare options and default settings on the LogWare Options dialog only! Thermometer passcodes are specific to the thermometer readout being used and are required to change certain settings in the thermometer readout itself. Please refer to Section 4.4, Thermometer Settings Dialogs, for more information.*

6.6 Importing logged data into a spreadsheet

The data logged by LogWare is saved into ASCII text files referred to as output files. These files can be opened and graphed by LogWare by selecting the Open Previously Logged Data File option in the View menu.

However, these files may also be imported into spreadsheet software such as Microsoft® Excel for graphing or analysis.

To import an output file into Excel:



Note: This example uses Microsoft® Excel XP. These steps may or may not apply if using a different version of Excel or another spreadsheet software package.

1. In Excel, select the Open option in the File menu to display the Open dialog.
2. In the Files of type box at the bottom of the open dialog, select the Text Files (*.prn;*.txt;*.csv) option.
3. Navigate to the folder where the output file is located. Select the output file to open and click the Open button.
4. The Text Import Wizard – Step 1 of 3 dialog appears.

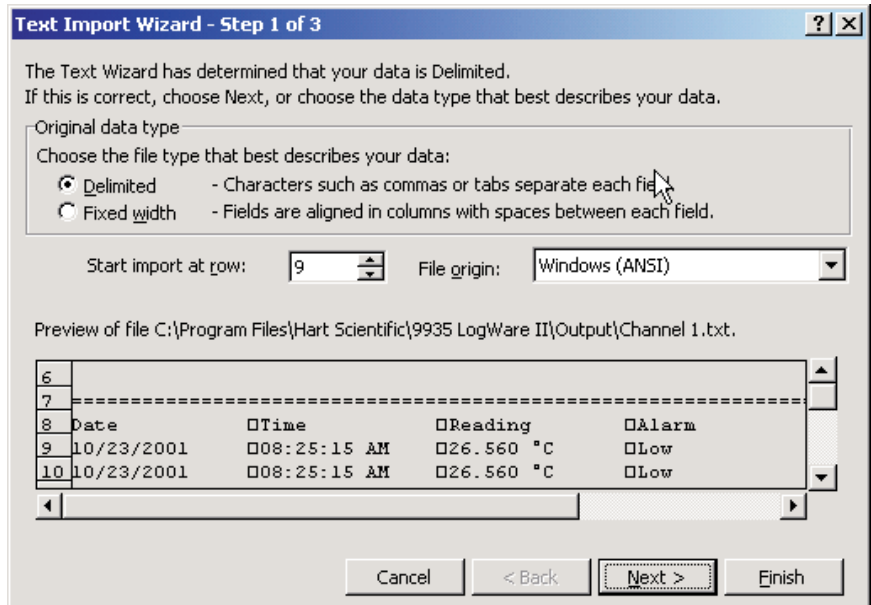


Figure 159 Excel XP Text Import Wizard – Step 1 of 3

5. Make sure the Original data type option is set to Delimited. If the thermometer and probe information is not needed, set the Start import at row option to the first row containing the logged readings (row 9 in this example). Set the File Origin option to Windows (ANSI) and click the Next button.

- The Text Import Wizard – Step 2 of 3 dialog appears.

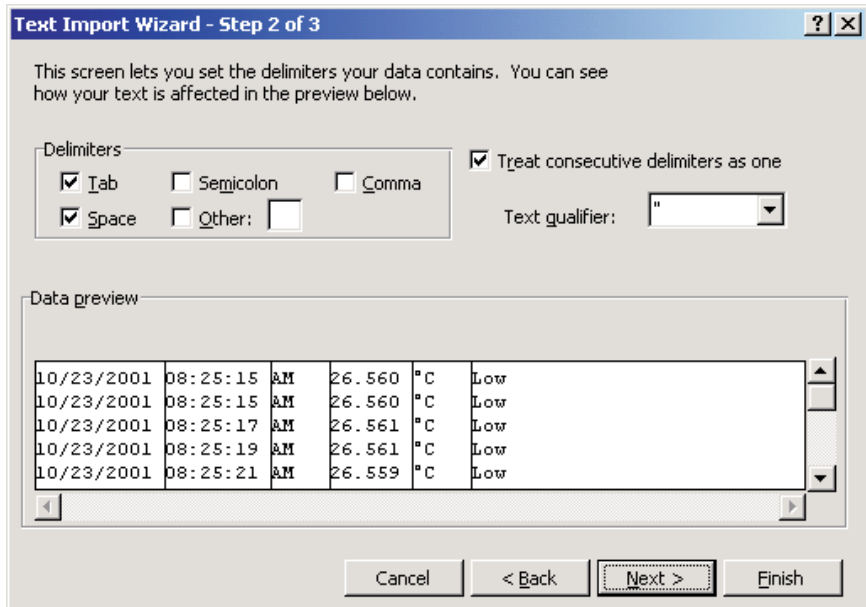


Figure 161 Excel XP Text Import Wizard – Step 2 of 3

- In the Delimiters box, select the Tab and Space options. Make sure the Treat consecutive delimiters as one check box is checked and click the Finish button.
- The output file is displayed in the spreadsheet. Resize the columns so the contents of each cell are visible.
- Select the Date, Time, and Reading columns by clicking on the gray column headers while holding down the control key labeled A, B and D, as shown below:

	A	B	C	D	E	F
1	10/23/2001	8:25:15 AM	26.56	°C	Low	
2	10/23/2001	8:25:15 AM	26.56	°C	Low	
3	10/23/2001	8:25:17 AM	26.561	°C	Low	
4	10/23/2001	8:25:19 AM	26.561	°C	Low	
5	10/23/2001	8:25:21 AM	26.559	°C	Low	
6	10/23/2001	8:25:23 AM	26.559	°C	Low	
7	10/23/2001	8:25:25 AM	26.561	°C	Low	
8	10/23/2001	8:25:27 AM	26.572	°C	Low	
9	10/23/2001	8:25:29 AM	26.572	°C	Low	
10	10/23/2001	8:25:31 AM	26.509	°C	Low	
11	10/23/2001	8:25:35 AM	26.508	°C	Low	

Figure 160 Excel XP data

10. In the Insert menu of Excel, select the Chart option. The Chart Wizard – Step 1 of 4 – Chart Type dialog is displayed.

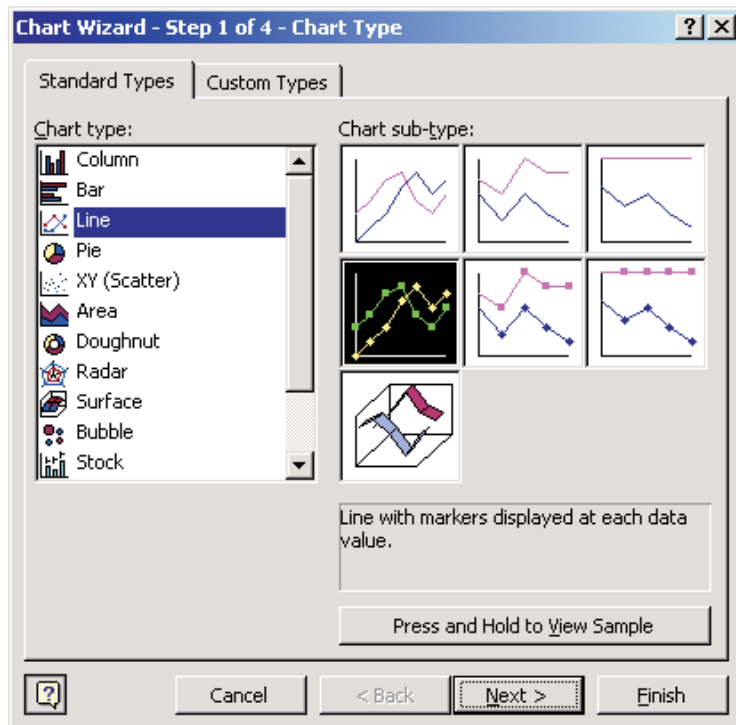


Figure 162 Excel XP Chart Wizard – Step 1 of 4 – Chart Type

11. In the Chart type list, select Line and click the Finish button.

12. The graph should be displayed as shown below:

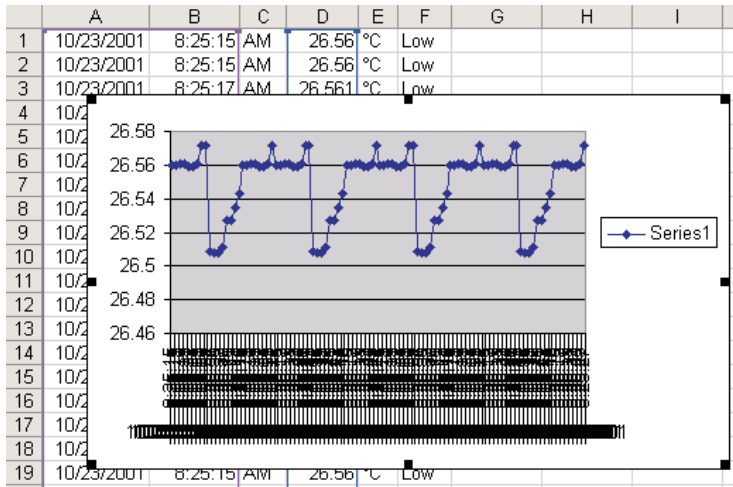


Figure 163 Excel XP graph

13. Format and size the graph as desired.

6.7 Synchronizing the thermometer readout's clock to the computer

The thermometer readouts supported by LogWare have an internal clock. LogWare provides a utility to synchronize the thermometer readout's clock with the clock on the computer.



Note: LogWare itself does not use the clock on the thermometer readout at all. All dates and times associated with readings acquired by LogWare use the computer's clock.

Before synchronizing the thermometer readout's clock, make sure the computer's clock is accurate. Use the Windows® Control Panel to adjust the date and time as necessary.

To synchronize the thermometer readout's clock with the computer's clock:

1. Select the Synchronize Date/Time option in the Tools menu or click the



button on the toolbar.

2. The Select Model dialog is displayed. Select the thermometer readout model. Also, select the COM port to which the thermometer readout is connected and the baud rate to use. Make sure LogWare can communicate with the thermometer readout by clicking the Test Communication button. Click the OK button when communication is successful.
3. LogWare attempts to set the thermometer readout's clock to match the clock on the computer.

Please refer to Section 4.1.1, Synchronize Date/Time, for more information.

6.8 Protecting calibration parameters and probe conversion settings

Some of the thermometer readouts supported by LogWare have built-in passcode protection for calibration parameters and/or probe conversion settings (including probe characterization coefficients). These settings are typically protected because altering their values changes the calibration and/or readout accuracy of the thermometer readout. Some thermometer readouts, however, lack this protection.

LogWare includes a feature that allows these settings to be passcode protected (for thermometer readouts that do not inherently include this protection) using an administrator passcode to prevent unauthorized users from changing the current values.

To protect calibration parameters and probe conversion settings:

1. Select the Options option in the Tools menu.
2. The LogWare Options dialog is displayed.
3. Select the Tools tab. The Tools tab contains options that pertain to the thermometer settings dialogs.
4. In the Passcode Protection section there are two options:
 - Protect probe conversion settings
 - Protect calibration parameters



Note: *If these settings are disabled, an administrator passcode has been setup. The administrator passcode is required to change these settings. For more information on gaining access to the options protected by the administrator passcode, see Section 6.9, Setting up an administrator passcode.*



Note: The administrator passcode in LogWare is completely independent of the thermometer readout passcodes. The administrator passcode is used by LogWare to protect certain LogWare options and default settings on the LogWare Options dialog only! Thermometer readout passcodes are specific to the thermometer readout being used and are required to change certain settings in the thermometer readout itself.

5. To passcode protect probe conversion settings, check the Protect probe conversion settings box.
6. To passcode protect thermometer readout calibration parameters, check the Protect calibration parameters check box.



Note: These options only apply to thermometer readouts that do not have built-in passcode protection for these features. For thermometer readouts that do have built-in passcode protection, the thermometer readouts settings override these settings.

Please refer to Section 5.6.4, Options, for more information.

6.9 Setting up an administrator passcode

LogWare allows the user to setup an administrator passcode to protect some of the features and options of the LogWare software. Some of these features include:

- Ability to set thermometer readout and/or module serial numbers
- Ability to specify passcode protection settings
- Real-time measurement process settings

The administrator passcode can be setup or changed at any time by following the process below.



Note: The administrator passcode in LogWare is completely independent of the thermometer readout passcodes. The administrator passcode is used by LogWare to protect certain LogWare options and default settings on the LogWare Options dialog only! Thermometer readout passcodes are specific to the thermometer readout being used and are required to change certain settings in the thermometer readout itself.

To setup or change the administrator passcode:

1. Select the Options option in the Tools menu.
2. The LogWare Options dialog is displayed.
3. Select the Admin. Passcode tab.

4. Click the Change Admin. Passcode button to setup or change the administrator passcode.
5. If a passcode has already been setup, the current passcode is required and must be entered. The Enter Passcode dialog is displayed. Enter the current administrator passcode in the Passcode box and click OK to continue.
6. The Change Passcode dialog is displayed. Enter the new administrator passcode in the New passcode box. Enter the passcode again in the Confirm passcode box. Click the OK button to accept the new passcode or click Cancel to abort setting the new passcode.



Note: DO NOT FORGET THE ADMINISTRATOR PASSCODE! Write it down and keep it in a safe place. Only reveal the passcode to users who are authorized to change the protected settings!

Once an administrator passcode has been configured, certain features and settings are disabled and can only be changed after the administrator passcode has been entered.

To gain access to passcode protected settings:

1. Select the Options option in the Tools menu.
2. The LogWare Options dialog is displayed.
3. Select the Admin. Passcode tab.
4. Click the Edit Protected Options button to enter the administrator passcode.
5. The Enter Passcode dialog is displayed. Enter the current administrator passcode in the Passcode box and click the OK button to continue.
6. If the entered passcode is correct, the features and settings on the LogWare Options dialog are enabled and may be changed. These settings are only enabled until this dialog is closed.

The administrator passcode can also be removed from LogWare, which permits anyone to change any of LogWare's options.

To remove the administrator passcode:

1. Select the Options option in the Tools menu.
2. The LogWare Options dialog is displayed.
3. Select the Admin. Passcode tab.
4. Click the Change Admin. Passcode button to setup or change the administrator passcode.

5. The Enter Passcode dialog is displayed. Enter the current administrator passcode in the Passcode box and click the OK button to continue.
6. The Change Passcode dialog is displayed. Leave the New passcode box Confirm passcode boxes empty. Click the OK button. A message is displayed warning that passcode protection is about to be disabled. Confirm disabling passcode protection by clicking the Yes button.

Please refer to Section 5.6.4, Options, for more information.

6.10 Viewing remaining memory for logging data

Some thermometer readouts that are supported by LogWare allow readings to be logged to the memory of the thermometer readout. LogWare provides a feature that indicates how much of the thermometer readout's memory is already used.

To view the status of the thermometer readout's log(s):

1. Select the Log Status option in the Download menu.
2. The Select Model dialog is displayed. Select the thermometer readout model. Also, select the COM port to which the thermometer readout is connected and the baud rate to use. Make sure LogWare can communicate with the thermometer readout by clicking the Test Communication button. Click the OK button when communication is successful.
3. The Log Status dialog for the selected thermometer readout is displayed indicating how much of the thermometer readout's memory is already used. This dialog is specific to the selected thermometer readout.
4. Click the Close button to close the Log Status dialog when finished.

Refer to Section 3.4, Log Status Dialogs, for more information.

6.11 Downloading logged data from the memory of a thermometer readout

The thermometer readouts that are supported by LogWare allow readings to be logged to the memory of the thermometer readout. LogWare provides a feature to download the logged readings from the memory of the thermometer readout and save them to a text file (or files) on the computer.

To download readings from the memory of a thermometer readout to a text file:

1. Select the Download Logged Data option in the Download menu.
2. The Select Model dialog is displayed. Select the thermometer readout model. Also, select the COM port to which the thermometer readout is connected and the baud rate to use. Make sure LogWare can communi-

cate with the thermometer readout by clicking the Test Communication button. Click the OK button when communication is successful.

3. The Download Options dialog is displayed. This dialog is specific to the selected thermometer readout.
4. Select the appropriate download options and click the Download button to download the readings.
5. Click the Close button to close the Download Options dialog when finished.

Please refer to Section 3, Downloading Logged Data, for more information.

6.12 Saving probe coefficients to a configuration file

LogWare provides a feature that allows probe coefficients to be saved to a configuration file.

To save probe coefficients to a configuration file:

1. Select the Thermometer Settings option in the Tools menu.
2. The Select Model dialog is displayed. Select the thermometer readout model. Also, select the COM port where the thermometer readout is connected and the baud rate to use. Make sure LogWare can communicate with the thermometer readout by clicking the Test Communication button. Click the OK button when communication is successful.
3. The thermometer readout settings are then read in and displayed in the Thermometer Settings dialog. This dialog is specific to the selected thermometer readout.
4. If applicable, to view the settings for a specific module, select the module from the Modules list and click the Module Settings button.
5. To view the settings for a specific channel, select the channel from the Channels list and click the Channel Settings button.
6. To view the coefficients for the probe on the selected channel, select the Conversion Settings button.
7. The conversion settings dialog for the selected conversion type is displayed.
8. Click the Save As button on the conversion settings dialog to save the coefficients for the selected channel to a configuration file.
9. The Windows® Save As dialog is displayed. Enter a filename and click the OK button to save the coefficients.

Please refer to Section 9, Conversion Settings Dialogs, for more information.

6.13 Using an IR dongle with LogWare

Some of the features of LogWare allow an IR dongle to be used to communicate with thermometer readouts that have an IR port instead of using an RS-232 cable.



Note: Hart Scientific thermometer readouts with an IR port use a proprietary protocol for communication. Because of this proprietary protocol, LogWare is not able to use IR ports that are built into a computer. Instead, LogWare requires the use of an IR dongle which can be connected to a standard RS-232 port on the computer. No special software or drivers are required to use the IR dongle. All necessary communication protocols and drivers are built into the LogWare software. To purchase an IR dongle, contact Hart Scientific Technical Support.

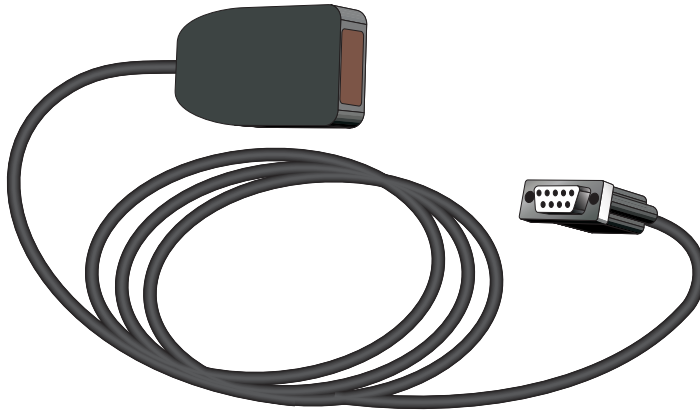


Figure 164 IR dongle

LogWare provides support for communicating over an IR port with these thermometer readouts. However, the support is limited to the following features:

- Downloading logged data from the thermometer readout
- Checking the thermometer readout's log status
- Writing the thermometer readout's logging labels
- Synchronizing the thermometer readout's date and time to the computer's clock
- Accessing the thermometer readout's settings

LogWare does not support using the IR port for performing real-time log sessions.


To use an IR dongle with LogWare:

1. Locate a 9-pin or 25-pin COM (RS-232) port on the computer. Connect the IR dongle to the COM port. If the COM port has a 25-pin connector, use a 9-25 pin adapter to connect the IR dongle.
2. Select the appropriate menu option for the feature of LogWare to use.
3. When selecting one of these menu options, the Select Model dialog appears.
4. Select the model of the thermometer readout to use and the COM port to which the IR dongle is connected.
5. If the selected feature and the selected thermometer readout model support using the IR dongle, the Use IR dongle check box is enabled. Check the Use IR dongle check box.
6. Before continuing, make sure the thermometer readout's IR port and the IR dongle are in close proximity and the path between them is unobstructed. For more details on placement of the IR dongle, refer to the thermometer readout's User's Guide.
7. Click the Test Communication button on the Select Model dialog. If LogWare succeeds in communicating with the thermometer readout using the IR dongle, a success message is displayed. Otherwise, a failure message is displayed. In case of failure, check the following:
 - ◆ The dongle is connected to the selected COM port.
 - ◆ The thermometer readout is powered up and operating.
 - ◆ The dongle is within the required range of the thermometer readout's IR port.
 - ◆ The path between the thermometer readout's IR port and the IR dongle is not obstructed.
 - ◆ The thermometer readout's IR baud rate matches the baud rate setting on the Select Model dialog.
 - ◆ The thermometer readout's IR mode is set properly to send/receive commands (see thermometer readout User's Guide).
 - ◆ Shut the thermometer readout's power off, then back on again.

6.14 Changing logging label names

Some of the thermometer readouts supported by LogWare have labels that can be assigned to data that is logged in the thermometer readout's memory. LogWare provides a utility to view and/or change the names of these logging labels.

To view and/or modify the thermometer readout's logging labels:

1. Select the Logging Labels option in the Tools menu or click the  button on the toolbar.
2. The Select Model dialog is displayed. Select the thermometer readout model. Also, select the COM port where the thermometer readout is connected and the baud rate to use. Make sure LogWare can communicate with the thermometer readout by clicking the Test Communication button. Click the OK button when communication is successful.
3. LogWare displays the Logging Labels dialog and attempts to read in the current logging label settings. This dialog is specific to the selected thermometer readout.
4. Modify the labels as desired on the Logging Labels dialog.
5. To write the new logging label settings to the thermometer readout, click the OK button.

Please refer to Section 4.3, Logging Labels Dialogs, for more information.

6.15 Unlocking a demo version of LogWare

When LogWare is purchased, a serial number is supplied on a sticker typically located on the CD pouch on the inside back cover of the LogWare User's Guide. When LogWare is installed, a dialog appears and asks for the serial number. If a valid serial number is entered, LogWare is installed as a "full working" version. If an invalid serial number is entered, LogWare is installed as a demo version.

The demo version of LogWare never "expires". The demo version is identical to the "full working" version with the following exceptions:

- Performs real-time log session, but does not write acquired data to file(s).
- Downloads data from thermometer readout log(s), but does not write downloaded data to file(s).
- Allows thermometer readout settings to be viewed, but does not write new settings to thermometer readout.
- Allows logging labels to be viewed, but does not write new logging labels to thermometer readout.
- Some example data files are installed to provide examples of the files that demo version does not write to.

The demo version of LogWare may be freely distributed and installed and is available for download on Hart Scientific's website for no charge.

The demo version may be "unlocked" at any time after obtaining a valid serial number from Hart Scientific Technical Support without having to reinstall the software.

To “unlock” a demo version of LogWare:



Note: To successfully unlock a demo version of LogWare, you must be logged into Windows as an Administrator.

1. Select the Unlock option in the Help menu. The Unlock option only appears in the Help menu for the demo version of LogWare. The Unlock Code dialog appears.

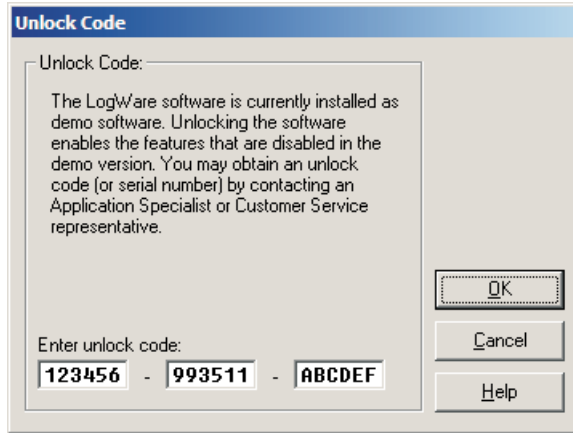


Figure 165 Unlock Code dialog

2. Enter the serial number provided in the Enter unlock code boxes at the bottom of the dialog. The serial number must have the format: XXXXXX - 9935XX - XXXXXX.
3. Click the OK button to accept the unlock code. If the serial number entered is a valid serial number, a message is displayed indicating LogWare has been unlocked. LogWare MUST be closed and restarted at this point!
4. If the serial number entered is invalid, LogWare will NOT be unlocked.
5. Click the Cancel button to close this dialog without unlocking the demo version.

6.16 Setting up LogWare to read multiple scales

LogWare was designed primarily to acquire temperature readings from digital thermometer readouts. However, LogWare is capable of acquiring temperature, resistance, and voltage readings simultaneously. This can be accomplished by configuring each channel on the thermometer readout to read in the desired units prior to starting a real-time log session.

To configure a thermometer readout to read different units on each channel:

1. Select the Setup option in the Datalog menu to display the Log Session Settings dialog.
2. Select the Thermometer tab.
3. Select the thermometer readout model using the Model drop-down list.
4. Select the COM port, and baud rate using the drop-down lists in the Communication settings section.
5. Click the Thermometer Settings button to display the Thermometer Settings dialog for the selected thermometer readout.
6. On the Thermometer Settings dialog, select the desired system units using the Units drop-down list. Click the Apply button to write the new system units setting to the thermometer readout.



Note: If a non-temperature setting is selected for the system units (i.e. Ohms, Volts, Ratio, or W), this setting will override any conversion type setting for all channels that would otherwise cause reading to be displayed in temperature!! Therefore, it is recommended that the system units be set to a temperature setting (i.e. °C, °F, K, etc.) and the conversion type be used to configure channels that are to read non-temperature values.

7. On the Thermometer Settings dialog, navigate to the first channel to configure.
 - a. For thermometer readouts that support modules, select the Modules tab, select the module from the Modules list that the desired channel is on and click the Module Settings button. The Module Settings dialog is displayed. Continue with step b. below.
 - b. Select the Channels tab, select the desired channel from the Channels list and click the Channel Settings button. The Channel Settings dialog is displayed.
8. To configure the selected channel to read temperature, set the Conversion type to the desired setting and enter the coefficients to use (as applicable) to convert the readings on that channel to temperature by clicking the Conversion Settings button.
9. To configure the selected channel to read resistance, voltage, ratio or W, set the Conversion type appropriately. Refer to the thermometer readout's User's Guide for more details on the conversion types supported by the thermometer readout.
10. Click the OK button on the Channel Settings dialog to write the new settings to the thermometer readout.
11. Repeat steps 3 to 6 for the other channels that need to be configured.

12. Click the OK button on the Thermometer Settings dialog when finished.
13. Note that the Serial number and Session temperature scale boxes are filled in automatically when returning to the Log Session Settings dialog.
14. Set the Session temperature scale to <keep existing> using the drop-down list.
15. Finish setting up the remainder of the settings on the Log Session Settings dialog for this log session.
16. When the log session starts, *LogWare* will query the thermometer read-out to determine the system temperature units (Session temperature scale) and the conversion setting of each enabled channel. The system temperature units are then automatically applied to all channels that are configured to read temperature. All other channels use the appropriate scale depending on their conversion type setting.

Please refer to Section 2, Real-Time Logging, for more information.

7 Files

This section describes all of the files that LogWare creates and/or uses. The first section describes LogWare's configuration files and the second section describes LogWare's data files.

7.1 Configuration files

Configuration files are intended to facilitate the configuration process. By default, all configuration files are saved to the \Config folder (or the folder that is specified for configuration files on the LogWare Options dialog). LogWare supports many types of configuration files:

- Real-Time Log Session configuration files
- Alarm configuration files
- Conversion configuration files



Warning: Do not attempt to manually modify the contents of any configuration file! Files that have been modified may not be able to be read by LogWare!

7.1.1 Log session configuration file

The log session configuration file is used to save the current real-time log session settings.



Warning: Do not attempt to manually modify the contents of any configuration file! Files that have been modified may not be able to be read by LogWare!

A log session configuration file contains ALL configuration information required to start a real-time log session. This information includes the thermometer information, communication settings, sample interval settings, start and end log session settings, channel, statistics and alarm settings.

Log session configuration files must always have an extension of **.rls** and can be saved and opened from any of the following locations:

- Open and Save As buttons on the Log Session Settings dialog
- Open and Save As options in the File menu
- Open and Save As buttons on the toolbar



Note: Opening a log session configuration file overwrites ALL current log session settings!

For more information on opening and saving real-time log session files, see Section 2, Real-Time Logging.

7.1.2 Alarm configuration file

The alarm configuration file is used to save the current alarm settings of a single channel.



Warning: Do not attempt to manually modify the contents of any configuration file! Files that have been modified may not be able to be read by LogWare!

An alarm configuration file contains configuration information for low and high alarms on a single channel. This information includes the alarm state (enabled or disabled), alarm values, events, and other options.

Alarm configuration files must always have an extension of **.alm** and can be saved and opened from the Alarm Settings dialog.



Note: Opening an alarm configuration file overwrites the current alarm settings!

7.1.3 Conversion configuration files

Conversion configuration files are useful for saving probe coefficients to a file. Conversion configuration files contain information concerning a single conversion type including coefficients and any other information specific to that conversion type.

Conversion configuration files have an extension that matches the type of conversion as described below:

Conversion Type	File Extension
Callendar-Van Dusen	.cvd
High Temperature PRT	.hpt
IPTS-68	.i68
ITS-90	.i90
W(T90)	.w90
Polynomial	.ply
Thermistor	.thr
Thermocouple Polynomial	.tcp
Thermocouple Table	.tct
Thermocouple Type	.tcx



Note: Opening a configuration file overwrites the current conversion settings.

Conversion configuration files can be opened and/or saved from the appropriate conversion settings dialog. For more information on conversion settings dialogs, please refer to Section 9, Conversion Settings Dialogs.

7.2 Data files

Data files are created by LogWare either during a real-time log session or when data is downloaded from a thermometer readout. LogWare creates the following types of data files:

- **Output files** – Created during real-time log session and contains acquired data for a single channel
- **Alarm Log file** – Created during real-time log session and contains information about alarms
- **Download files** – Created when data is downloaded from a thermometer readout



Warning: Do not attempt to manually modify the contents of any data file! Files that have been modified may not be able to be read by LogWare!

7.2.1 Output file

LogWare saves the readings acquired during a real-time log session to an output file. Output files also contain information about the thermometer readout, probe, and alarm settings that were configured for the log session.

During a real-time log session, the acquired readings are displayed in the Channels grid and graphed on the real-time graph. The output file can be viewed during a real-time log session by selecting the Readings History option in the View menu or by right-clicking on the Channels grid on the real-time Log Session window and selecting the View Readings History option in the popup menu.

After a real-time log session has been stopped and/or closed, the output file can be opened by selecting the Open Previously Logged Data File option in the View menu. All of the readings that were acquired are displayed in a spreadsheet-style grid and on the graph. Output files can also be opened in other analysis software packages such as spreadsheets and text editors.

The following is an example of an output file created by LogWare:

```
Channel 1.txt

Thermometer Readout Model: 1529Thermometer Serial: A13053Thermometer Readout Channel: 1
Channel name: Channel 1

Low Alarm: 79.0000 F
High Alarm: 80.0000 F
=====
Date      Time      Reading  Alarm
03/29/2010 3:21:48 PM 73.8339 F Low
03/29/2010 3:21:49 PM 73.8285 F Low
03/29/2010 3:21:50 PM 73.8312 F Low
03/29/2010 3:21:51 PM 73.8254 F Low
03/29/2010 3:21:52 PM 73.8279 F Low
03/29/2010 3:21:53 PM 73.8250 F Low
03/29/2010 3:21:54 PM 73.8236 F Low
03/29/2010 3:21:55 PM 73.8237 F Low
03/29/2010 3:21:56 PM 73.8237 F Low
03/29/2010 3:21:57 PM 73.8208 F Low
03/29/2010 3:21:58 PM 73.8214 F Low
03/29/2010 3:21:59 PM 73.8183 F Low
03/29/2010 3:22:00 PM 73.8176 F Low
03/29/2010 3:22:01 PM 73.8158 F Low
03/29/2010 3:22:02 PM 73.8192 F Low
03/29/2010 3:22:03 PM 73.8130 F Low
03/29/2010 3:22:04 PM 73.8155 F Low
03/29/2010 3:22:05 PM 73.8181 F Low
03/29/2010 3:22:06 PM 73.8150 F Low
03/29/2010 3:22:07 PM 73.8166 F Low
03/29/2010 3:22:08 PM 73.8190 F Low
03/29/2010 3:22:09 PM 73.8179 F Low
03/29/2010 3:22:10 PM 73.8191 F Low
03/29/2010 3:22:11 PM 73.8208 F Low
03/29/2010 3:22:12 PM 73.8203 F Low
```

Figure 166 Output file example

The output file contains a header indicating the filename, thermometer readout and session temperature scale information, the channel name and alarm settings (if any).

The data is formatted into columns and consist of the date, time, reading value and alarm state.



Note: If the output file was created using the Write all acquired data to the same file option, the header of the output file will then contain information for all enabled channels, and an additional column is added to the file for the channel number for each reading. This feature should typically only be used if you are using other software to sort and analyze the logged data. The output file created when using this feature can still be opened by LogWare and displayed on-screen, however all readings are displayed on the graph as if they came from the same channel (single trace).

Optionally, the statistical data for the readings may be appended to the end of the file when the output file is printed. Statistics are only appended if this feature is enabled on the LogWare Options dialog. Please refer to Section 5.6.4, Options, for more details.

7.2.2 Download file

LogWare includes a feature that allows readings that are stored in the memory of thermometer readouts that support this feature to be downloaded and written into text files, called download files.

To download readings from the memory of a thermometer readout, see the Download Logged Data option in the Download menu.

Download files can be opened by this software by selecting the Open Previously Logged Data Files option in the View menu. They can also be opened in other analysis software packages such as spreadsheets and text editors.

The format of a download file is very similar to and follows the same general format as an output file. The actual content of a download file, however, will vary slightly depending on the thermometer readout that the data was downloaded from. See the example below:

```
download1.txt
Thermometer Model: 1529
Channel: 1
-----
Date           Time           Reading         Label
11/02/2002    10:26:28       20.3548 °C      DATA_04
11/02/2002    10:26:29       20.3529 °C      DATA_04
11/02/2002    10:26:30       20.3539 °C      DATA_04
11/02/2002    10:26:31       20.3536 °C      DATA_04
11/02/2002    10:26:32       20.3531 °C      DATA_04
11/02/2002    10:26:33       20.3534 °C      DATA_04
11/02/2002    10:26:34       20.3525 °C      DATA_04
11/02/2002    10:26:35       20.3536 °C      DATA_04
```

Figure 167 Download file example

The download file contains a header indicating the filename, thermometer readout model and channel (optional). The data is formatted into columns and consist of the date, time, reading value, label and/or channel information.

7.2.3 Alarm Log File

When a real-time log session is started, LogWare creates a file called “alarms.log”. This file is located in the \Output folder by default. The purpose of this file is to maintain a history of readings that caused the alarms to trip or reset.

The alarm log file can be viewed during a real-time log session by selecting the Alarm Log menu option in the View menu or by right-clicking on the Channels grid on the real-time Log Session window and selecting the View Alarm Log option in the popup menu.

The following is an example of an alarm log file created by LogWare:

```
alarms.log
Alarm Settings:
Channel: 1      Channel 1      Low Alarm:      28.00 °C      High Alarm:      30.00 °C
Channel: 2      Channel 2
Channel: 3      Channel 3
=====
Date           Time           Name           Alarm           Reading
2/28/2002     10:49:38 AM   Channel 1     Low            25.59 °C
2/28/2002     10:49:59 AM   Channel 1     Reset         28.44 °C
2/28/2002     10:50:14 AM   Channel 1     High          30.04 °C
2/28/2002     10:50:59 AM   Channel 1     Reset         29.99 °C
```

Figure 168 Alarm log file example

The alarm log file contains a header indicating the filename and alarm settings for all enabled channels.

The data is formatted into columns and consist of the date, time, channel name, alarm state and reading value.

8 Download process

The following sections describe in detail the download process for each of the thermometer readouts supported by *LogWare*.

8.1 1523/1524 Handheld and 1529 Chub-E4 download process

The download process for the 1523/1524 Handheld and 1529 Chub-E4 Thermometer readouts consists of 3 main steps:

1. Download the data from the thermometer readout
2. Parse the data
3. Write the data to download file(s)

When the download process starts, a progress dialog appears to indicate the download progress. To abort downloading the data from the thermometer readout, click the Cancel button on the progress dialog.

If the Prompt to enter filename(s) option was selected, the Windows® Save As dialog is displayed to get the name of the file(s) that the data should be saved to once all of the data has been downloaded. The \Download folder should be selected by default (or the folder that is specified for download files on the *LogWare* Options dialog).

- If the Write all readings into a single file option was selected, all readings will be written to the specified file. By default, the filename is “download.txt”.
- If the Separate readings by channel option was selected, a prompt appears for every channel. By default, the filenames are “download n .txt” where n is the channel number.

Accept the default filename(s) or enter new filename(s) and click the Save button to continue. If a file with the selected filename already exists, confirm overwriting the existing file or enter a new filename. Clicking the Cancel button on the Windows® Save As dialog aborts the download process.

If the Prompt to enter filename(s) option was not selected, the data is saved to file(s) with the default filename(s).

Once all of the downloaded data has been saved into the specified file(s), the names of all of the files are displayed on the Downloaded Files dialog.

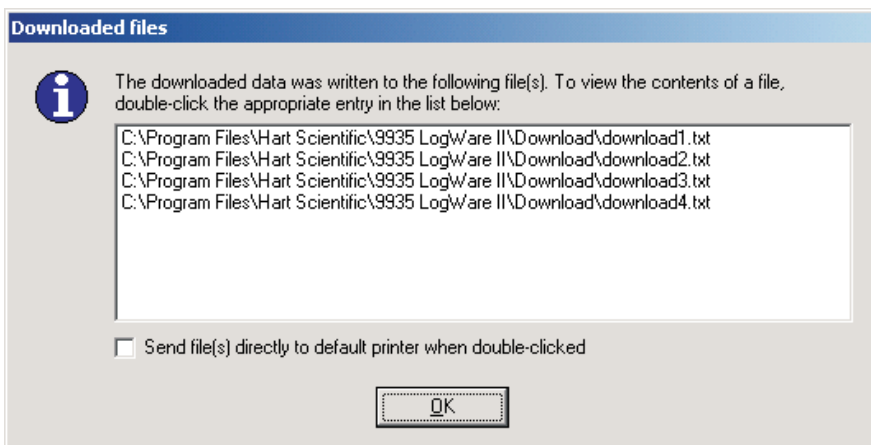


Figure 169 Downloaded Files dialog

To open a file that downloaded data was written to, double-click the appropriate entry in the list. The file will be opened in the File Viewer window. For more details on the File Viewer window, please refer to Section 2.4.8, File Viewer Window.

To print a file, check the Send file(s) directly to default printer when double-clicked option. Double-clicking an entry in the list will then cause the file to be sent to the current default printer.

Click the OK button to close this dialog.

8.2 1560 Black Stack download process

The download process for the 1560 Black Stack Thermometer readout consists of 3 main steps:

1. Download the data from the thermometer readout
2. Parse the data
3. Write the data to download file(s)

When the download process starts, a progress dialog appears to indicate the download progress. To abort downloading the data from the thermometer readout, click the Cancel button on the progress dialog.



Note: When downloading data from the 1560, the measurement process in the thermometer readout is halted until downloading is complete!

If the Prompt to enter filename(s) option was selected, the Windows® Save As dialog is displayed to get the name of the file(s) that the data should be saved to once the data has been downloaded. The \Download folder should be selected by default (or the folder that is specified for download files on the LogWare Options dialog).

- If the Write all readings into a single file option was selected, all readings will be written to the specified file. By default, the filename is “download.txt”.
- If the Separate readings by channel option was selected, a prompt appears for every channel. By default, the filenames are “download*n*.txt” where *n* is the channel number.

Accept the default filename(s) or enter new filename(s) and click the Save button to continue. If a file with the selected filename already exists, confirm overwriting the existing file or enter a new filename. Clicking the Cancel button on the Windows® Save As dialog aborts the download process.

If the Prompt to enter filename(s) option was not selected, the data is saved to file(s) with the default filename(s).

Once all of the downloaded data has been saved into the specified file(s), the names of all of the files are displayed on the Downloaded Files dialog (see Figure 169 on page 233).

To open a file that downloaded data was written to, double-click the appropriate entry in the list. The file will be opened in the File Viewer window. For more details on the File Viewer window, please refer to Section 2.4.8, File Viewer Window.

To print a file, check the Send file(s) directly to default printer when double-clicked option. Double-clicking an entry in the list will then cause the file to be sent to the current default printer.

Click the OK button to close this dialog.

8.3 1575A Super-Thermometer download process

The download process for the 1575A Super-Thermometer readout consists of 3 main steps:

1. Download the data from the thermometer readout
2. Parse the data
3. Write the data to download file

When the download process starts, a progress dialog appears to indicate the download progress. To abort downloading the data from the thermometer readout, click the Cancel button on the progress dialog.

If the Prompt to enter filename option was selected, the Windows® Save As dialog is displayed to get the name of the file that the data should be saved to once the data has been downloaded. The \Download folder should be selected by default (or the folder that is specified for download files on the LogWare Options dialog). Accept the default filename or enter new filename and click the Save button to continue. If a file with the selected filename already exists, confirm overwriting the existing file or enter a new filename. Clicking the Cancel button on the Windows® Save As dialog aborts the download process.

If the Prompt to enter filename option was not selected, the data is saved to a file with the default filename 'download.txt'.

Once all of the downloaded data has been saved into the specified file, the filename is displayed on the Downloaded Files dialog (see Figure 169 on page 233).

To open the file that downloaded data was written to, double-click the entry in the list. The file will be opened in the File Viewer window. For more details on the File Viewer window, please refer to Section 2.4.8, File Viewer Window.

To print a file, check the Send file(s) directly to default printer when double-clicked option. Double-clicking an entry in the list will then cause the file to be sent to the current default printer.

Click the OK button to close this dialog.

8.4 1590 Super-Thermometer II download process

The download process for the 1590 Super-Thermometer II readout consists of 3 main steps:

1. Download the data from the thermometer readout
2. Parse the data
3. Write the data to download file

When the download process starts, a progress dialog appears to indicate the download progress. To abort downloading the data from the thermometer readout, click the Cancel button on the progress dialog.

If the Prompt to enter filename option was selected, the Windows® Save As dialog is displayed to get the name of the file that the data should be saved to once the data has been downloaded. The \Download folder should be selected by default (or the folder that is specified for download files on the LogWare Options dialog). Accept the default filename or enter new filename and click the Save button to continue. If a file with the selected filename already exists, confirm overwriting the existing file or enter a new filename. Clicking the Cancel button on the Windows® Save As dialog aborts the download process.

If the Prompt to enter filename option was not selected, the data is saved to a file with the default filename 'download.txt'.

Once all of the downloaded data has been saved into the specified file, the file-name is displayed on the Downloaded Files dialog (see Figure 169 on page 233).

To open the file that downloaded data was written to, double-click the entry in the list. The file will be opened in the File Viewer window. For more details on the File Viewer window, please refer to Section 2.4.8, File Viewer Window.

To print a file, check the Send file(s) directly to default printer when double-clicked option. Double-clicking an entry in the list will then cause the file to be sent to the current default printer.

Click the OK button to close this dialog.

9 Conversion settings dialogs

The following sections describe in detail the conversion settings dialogs supported by LogWare. Each of these dialogs may appear slightly different depending on the features and settings supported by the thermometer readout being used.



Note: Some conversion type settings do not have an associated dialog (i.e. Ohms, Volts, Ratio, W, etc.).

9.1 ITS-90 Settings dialog

The ITS-90 Settings dialog allows the RTPW, Low Range, and High Range coefficients to be entered.

Figure 170 ITS-90 Settings dialog

If the ITS-90 coefficients were read in from the thermometer readout, the boxes are filled with the current coefficient values.

Enter the RTPW coefficient in from the Report of Calibration in the RTPW box. This box must contain a valid value.

Select the Low Range tab to select the low range and enter low range coefficients. Select the High Range tab to select the high range and enter high range coefficients.

Select the Open button to display the Windows® Open dialog to open a previously saved configuration file which contains ITS-90 coefficients. Select the Save As button to display the Windows® Save As dialog to save the current ITS-90 coefficients to a configuration file.


Click the OK button to accept the current coefficients and close this dialog.

Click the Cancel button to discard any changes you have made and close this dialog.

Click the Help button to display the appropriate help topic for this dialog.

9.1.1 Low Range tab

Select the Low Range tab to select the low range and enter low range coefficients. If the thermometer readout being used does not support ITS-90 low ranges, this tab is not displayed.

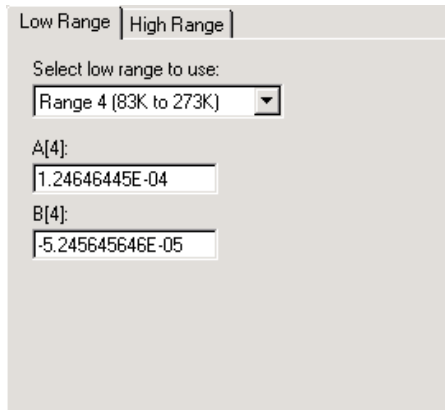

Low Range | High Range

Select low range to use:
Range 4 (83K to 273K)

A[4]:
1.24646445E-04

B[4]:
-5.245645646E-05

C[4]:

Figure 171 ITS-90 Settings – Low Range tab

Select the low range to use from the Select low range to use drop-down list. Low range choices include Range 1 to Range 5 but may be limited depending on the thermometer readout being used.

Enter the low range coefficients from the Report of Calibration in the boxes provided. The boxes displayed depend on the currently selected low range. If a coefficient is not used, set its value to 0.0.

9.1.2 High Range tab

Select the High Range tab to select the high range and enter high range coeffi-

icients. If the thermometer readout being used does not support ITS-90 high ranges, this tab is not displayed.

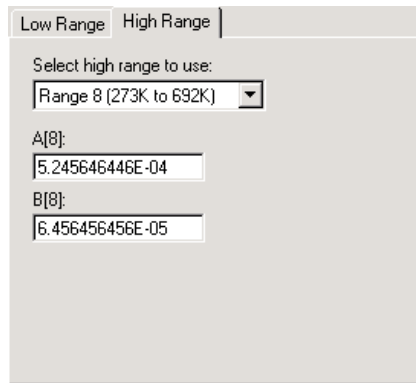


Figure 172 ITS-90 Settings – High Range tab

Select the high range to use from the Select high range to use drop-down list. High range choices include Range 6 to Range 11 but may be limited depending on the thermometer readout being used.

Enter the high range coefficients from the Report of Calibration in the boxes provided. The boxes displayed depend on the currently selected high range. If a coefficient is not used, set its value to 0.0.

9.2 IPTS-68 settings dialog

The IPTS-68 Settings dialog allows the Scale, R0, ALPHA, DELTA, A4 and C4 coefficients to be entered.

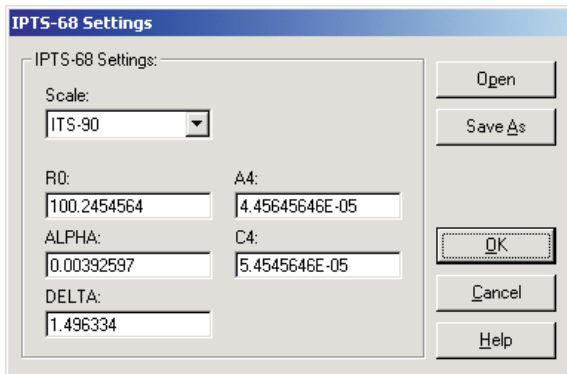


Figure 173 IPTS-68 Settings dialog

If the IPTS-68 coefficients were read in from the thermometer readout, the boxes are filled with the current coefficient values.

Select the scale to use from the Scale drop-down list. Scale choices include IPTS-68 and ITS-90. If the thermometer readout being used does not support the Scale setting, this drop-down list is not displayed.

Enter the R0, ALPHA, DELTA, A4, and C4 coefficients from the Report of Calibration in the boxes provided. All boxes must contain valid values.

Select the Open button to display the Windows® Open dialog to open a previously saved configuration file which contains IPTS-68 coefficients. Select the Save As button to display the Windows® Save As dialog to save the current IPTS-68 coefficients to a configuration file.

Click the OK button to accept the current coefficients and close this dialog.

Click the Cancel button to discard any changes you have made and close this dialog.

Click the Help button to display the appropriate help topic for this dialog.

9.3 Callendar-Van Dusen Settings dialog

The Callendar-Van Dusen Settings dialog allows the R0, ALPHA, DELTA, and BETA coefficients to be entered.

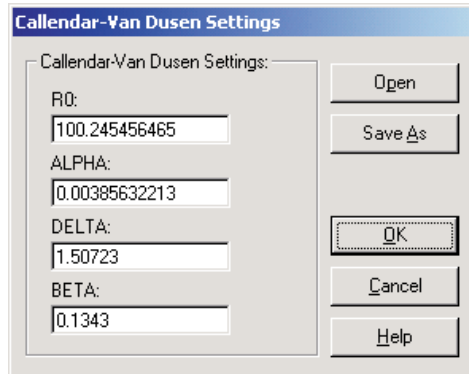


Figure 174 Callendar-Van Dusen Settings dialog

If the Callendar-Van Dusen coefficients were read in from the thermometer readout, the boxes are filled with the current coefficient values.

Enter the R0, ALPHA, DELTA, and BETA coefficients from the Report of Calibration. All boxes must contain valid values.

Select the Open button to display the Windows® Open dialog to open a previously saved configuration file which contains Callendar-Van Dusen coefficients. Select the Save As button to display the Windows® Save As dialog to save the current Callendar-Van Dusen coefficients to a configuration file.

Click the OK button to accept the current coefficients and close this dialog.

Click the Cancel button to discard any changes you have made and close this dialog.

Click the Help button to display the appropriate help topic for this dialog.

9.4 PT100 Settings dialog

The PT100 Settings dialog allows the R0 coefficient to be entered.

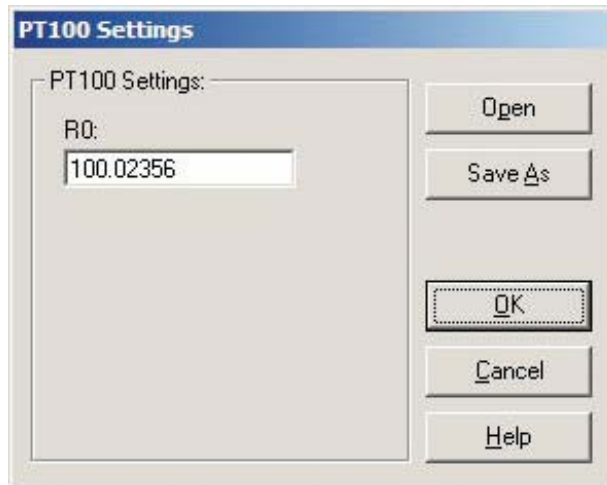


Figure 175 PT100 Settings dialog

If the PT100 coefficient was read in from the thermometer, the box is filled with the current coefficient value.

Enter the R0 coefficient from the Report of Calibration. The box must contain a valid value.



Note: Not all thermometer readouts that support the PT100 conversion type allow the R0 coefficient to be entered. This option is only available if the selected thermometer readout supports it.

Click the Open button to display the Windows® Open dialog to open a previously saved configuration file which contains a PT100 coefficient. Click the Save As button to display the Windows® Save As dialog to save the current PT100 coefficient to a configuration file.

Click the OK button to accept the current coefficients and close this dialog.

Click the Cancel button to discard any changes you have made and close this dialog.

Click the Help button to display the appropriate help topic for this dialog.

9.5 High-Temperature PRT Settings dialog

The High Temperature PRT Settings dialog allows the RTPW, A and B coefficients to be entered.

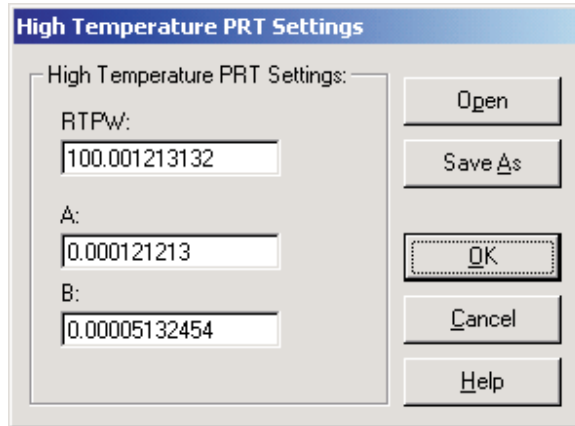


Figure 176 High-Temp PRT Settings dialog

If the High Temperature PRT coefficients were read in from the thermometer readout, the boxes are filled with the current coefficient values.

Enter the RTPW coefficient from the Report of Calibration in the RTPW box. This box must contain a valid value.

Enter the values for the A and B coefficients from the Report of Calibration in the boxes provided.

Select the Open button to display the Windows® Open dialog to open a previously saved configuration file which contains High Temperature PRT coefficients. Select the Save As button to display the Windows® Save As dialog to save the current High Temperature PRT coefficients to a configuration file.

Click the OK button to accept the current coefficients and close this dialog.

Click the Cancel button to discard any changes you have made and close this dialog.

Click the Help button to display the appropriate help topic for this dialog.

9.6 Polynomial Settings dialog

The Polynomial Settings dialog allows the A0, A1, A2,... A10 coefficients to be entered.

Label	Value
A0:	1.121313132E02
A1:	5.454644656E04
A2:	-2.45646546E05
A3:	5.465446646E07
A4:	8.456464600E09
A5:	
A6:	
A7:	
A8:	
A9:	
A10:	

Figure 177 Polynomial Settings dialog

If the Polynomial coefficients were read in from the thermometer readout, the boxes are filled with the current coefficient values.

Enter the values for all polynomial coefficients from the Report of Calibration in the boxes provided. If any coefficients are not used, leave the boxes blank or set the values to 0.0. The actual number of boxes displayed depends on the thermometer readout being used.

Select the Open button to display the Windows® Open dialog to open a previously saved configuration file which contains Polynomial coefficients. Select the Save As button to display the Windows® Save As dialog to save the current Polynomial coefficients to a configuration file.

Click the OK button to accept the current coefficients and close this dialog.

Click the Cancel button to discard any changes you have made and close this dialog.

Click the Help button to display the appropriate help topic for this dialog.

9.7 W(T90) Settings dialog

The W(T90) Settings dialog allows the RTPW coefficient to be entered.

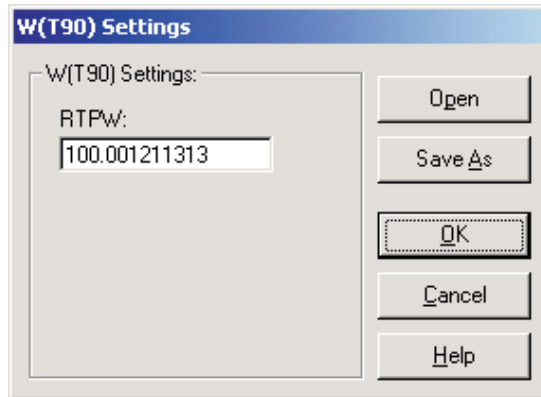


Figure 178 W(T90) Settings dialog

If the W(T90) coefficients were read in from the thermometer readout, the boxes are filled with the current coefficient values.

Enter the RTPW coefficient from the Report of Calibration in the RTPW box. This box must contain a valid value.

Select the Open button to display the Windows® Open dialog to open a previously saved configuration file which contains W(T90) coefficients. Select the Save As button to display the Windows® Save As dialog to save the current W(T90) coefficients to a configuration file.

Click the OK button to accept the current coefficients and close this dialog.

Click the Cancel button to discard any changes you have made and close this dialog.

Click the Help button to display the appropriate help topic for this dialog.

9.8 Thermistor Settings dialog

The Thermistor Settings dialog allows the Thermistor T(R) or Thermistor R(T) coefficients to be entered.

If the Thermistor coefficients were read in from the thermometer readout, the boxes are filled with the current coefficient values.

Select the Open button to display the Windows® Open dialog to open a previously saved configuration file which contains Thermistor coefficients. Select the Save As button to display the Windows® Save As dialog to save the current Thermistor coefficients to a configuration file.

Click the OK button to accept the current coefficients and close this dialog.

Click the Cancel button to discard any changes you have made and close this dialog.

Click the Help button to display the appropriate help topic for this dialog.

9.8.1 Thermistor T(R) Settings dialog

If the selected conversion type is T(R), the Thermistor T(R) Settings dialog is displayed.

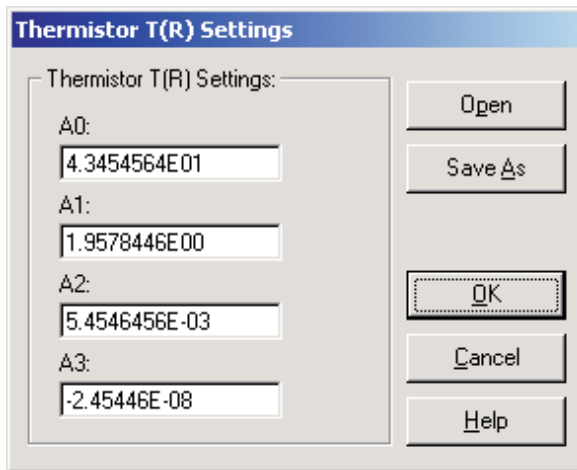


Figure 179 Thermistor T(R) Settings dialog

Enter the values for all Thermistor T(R) coefficients from the Report of Calibration in the boxes provided. If any coefficients are not used, leave the boxes blank or set the values to 0.0.

9.8.2 Thermistor R(T) Settings dialog

If the selected conversion type is R(T), the Thermistor R(T) Settings dialog is displayed.

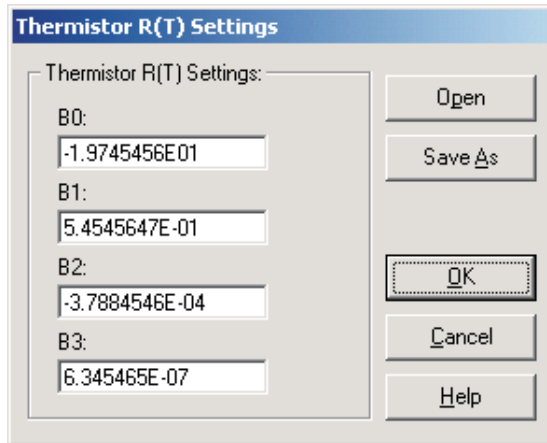


Figure 180 Thermistor R(T) Settings dialog

Enter the values for all Thermistor R(T) coefficients from the Report of Calibration in the boxes provided. If any coefficients are not used, leave the boxes blank or set the values to 0.0.

9.9 Thermocouple Polynomial Settings dialog

The Thermocouple Polynomial Settings dialog allows the mV (25C), C0, C1, C2,... C15 coefficients to be entered.

Figure 181 Thermocouple Polynomial Settings dialog

If the Thermocouple Polynomial coefficients were read in from the thermometer readout, the boxes are filled with the current coefficient values.

Enter the voltage of the thermocouple probe (in millivolts) at 25°C from the Report of Calibration in the mV (25C) box. If this value is not applicable to the thermometer readout, this box is disabled.

Enter the values for all thermocouple polynomial coefficients from the Report of Calibration in the boxes provided. If any coefficients are not used, leave the boxes blank or set the values to 0.0. The actual number of boxes displayed depends on the thermometer readout being used.

Select the Open button to display the Windows® Open dialog to open a previously saved configuration file which contains Thermocouple Polynomial coeffi-

icients. Select the Save As button to display the Windows® Save As dialog to save the current Thermocouple Polynomial coefficients to a configuration file.

Click the OK button to accept the current coefficients and close this dialog.

Click the Cancel button to discard any changes you have made and close this dialog.

Click the Help button to display the appropriate help topic for this dialog.

9.10 Thermocouple Table Settings dialog

The Thermocouple Table Settings dialog allows the Number of points and Temperature and Voltage values to be entered.

	Temperature		Voltage	
1:	-200.0	°C	-3.9	mV
2:	-100.0	°C	-1.8	mV
3:	0.0	°C	0.0	mV
4:	100.0	°C	2.7	mV
5:	200.0	°C	5.4	mV
6:	300.0	°C	8.5	mV
7:	400.0	°C	12.1	mV
8:	500.0	°C	14.5	mV
9:	600.0	°C	17.2	mV
10:	700.0	°C	20.0	mV

Figure 182 Thermocouple Table Settings dialog

If the Thermocouple Table coefficients were read in from the thermometer readout, the boxes are filled with the current coefficient values.

Select the number of points to use for the Temperature/Voltage table using the Number of points drop-down list. The choices are from 1 to 10 but may be limited depending on the thermometer readout being used.

Up to ten Temperature/Voltage pairs can be entered. Enter the temperature and voltage values at each temperature in the boxes provided. The units for the temperature and voltage values are indicated next to each box.

Select the Open button to display the Windows® Open dialog to open a previously saved configuration file which contains Thermocouple Table coefficients. Select the Save As button to display the Windows® Save As dialog to save the current Thermocouple Table coefficients to a configuration file.

Click the OK button to accept the current coefficients and close this dialog.

Click the Cancel button to discard any changes you have made and close this dialog.

Click the Help button to display the appropriate help topic for this dialog.

9.11 Thermocouple Type Settings dialog

The Thermocouple Type Settings dialog allows the Type, Temperature and Offset values, and DC1 & DC2 coefficients to be entered.

The screenshot shows the 'Thermocouple Type Settings' dialog box. It contains the following elements:

- Title Bar:** Thermocouple Type Settings
- Type:** A dropdown menu currently set to 'S'.
- Temperature and Offset:** A table with three rows and two columns.

	Temperature	Offset
1:	100.0 °C	0.0 °C
2:	200.0 °C	0.0 °C
3:	300.0 °C	0.0 °C
- DC1 and DC2:** Two input fields. DC1 contains '1.4544646E-04' and DC2 contains '-5.4568786E-05'.
- Buttons:** 'Open', 'Save As', 'OK', 'Cancel', and 'Help' are located on the right side of the dialog.

Figure 183 Thermocouple Polynomial Settings dialog

If the Thermocouple Type coefficients were read in from the thermometer readout, the boxes are filled with the current coefficient values and settings.

The Type drop-down list allows the standard thermocouple type to be selected. Thermocouple types include: B, E, J, K, N, R, S, T and AuPt. The choices in

the list may be limited depending on the thermocouple types supported by the thermometer readout being used.

The Temperature and Offset boxes provide a method to use an offset table with this thermocouple type. Up to three Temperature/Offset pairs can be entered. If no offset table is to be used, make sure the values in the Offset boxes are all set to 0.0.



Note: *If any one of the temperature values is set to 0°C, that point is ignored because the voltage on a thermocouple at 0°C is always assumed to be 0mV.*

If the selected thermocouple type is type S or type R, the DC1 and DC2 boxes are displayed. The DC1 and DC2 values are delta (Δ) values for the standard C1 and C2 polynomial coefficients. Enter the values for DC1 and DC2. These values should be located on the Report of Calibration.

Select the Open button to display the Windows® Open dialog to open a previously saved configuration file which contains Thermocouple Type coefficients. Select the Save As button to display the Windows® Save As dialog to save the current Thermocouple Type coefficients to a configuration file.

Click the OK button to accept the current coefficients and close this dialog.

Click the Cancel button to discard any changes you have made and close this dialog.

Click the Help button to display the appropriate help topic for this dialog.

9.12 Thermocouple Type Polynomial Settings dialog

The Thermocouple Type Polynomial Settings dialog allows the standard thermocouple type and the C0, C1, C2 and C3 coefficients to be entered.

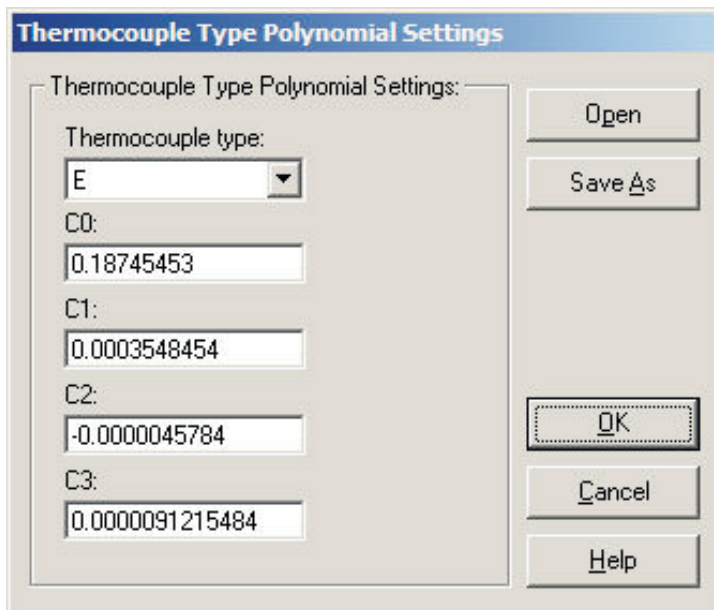


Figure 184 Thermocouple Type Polynomial Settings dialog

If the Thermocouple Type coefficients were read in from the thermometer, the boxes are filled with the current coefficient values.

The Thermocouple Type drop-down list allows the standard thermocouple type to be selected. Thermocouple types include: B, C, E, J, K, L, M, N, R, S, T and U. The choices in the list may be limited depending on the thermocouple types supported by the thermometer being used.

Enter the values for all thermocouple polynomial coefficients from the Report of Calibration. If any coefficients are not used, leave the boxes blank or set the values to 0.0. The actual number of boxes displayed depends on the thermometer being used.

Click the Open button to display the Windows® Open dialog to open a previously saved configuration file which contains Thermocouple Type Polynomial coefficients. Click the Save As button to display the Windows® Save As dialog to save the current Thermocouple Type Polynomial coefficients to a configuration file.

Click the OK button to accept the current coefficients and close this dialog.

Click the Cancel button to discard any changes you have made and close this dialog.

Click the Help button to display the appropriate help topic for this dialog.



10 Troubleshooting

The following sections describe in detail some troubleshooting topics, possible causes, and solutions.

10.1 Technical support

Please contact one of the following authorized Service Centers to coordinate service on your product:

Fluke Corporation, Hart Scientific Division

799 E. Utah Valley Drive
American Fork, Utah 84003-9775
USA

Phone: +1.801.763.1600
Toll Free: +1.800.438.4278
Fax: +1.801.763.1010
E-mail: support@hartscientific.com

Fluke Nederland B.V.

Customer Support Services
Science Park Eindhoven 5108
5692 EC Son
NETHERLANDS

Phone: +31-402-675300
Telefax: +31-402-675321
E-mail: ServiceDesk@fluke.nl

Fluke Int'l Corporation

Service Center - Instrimpex
Room 2301 Sciteck Tower
22 Jianguomenwai Dajie
Chao Yang District
Beijing 100004, PRC
CHINA

Phone: +86-10-6-512-3436

Telefax: +86-10-6-512-3437

E-mail: xingye.han@fluke.com.cn

Fluke South East Asia Pte Ltd.

Fluke ASEAN Regional Office

Service Center

60 Alexandra Terrace #03-16

The Comtech (Lobby D)

118502

SINGAPORE

Phone: +65-737-2922

Telefax: +65-737-5155

E-mail: antng@singa.fluke.com

Check our website on the Internet regularly for answers to Frequently Asked Questions, software updates, and Service Releases.

Before calling for Technical Support, check the Help file or User Guide topics listed below to see if the problem you are having is described there:

- Requirements
- Installation
- Running *LogWare*
- How do I... topics

When calling or sending a fax, please send or have the following information ready:

- Name of Software (*LogWare II*)
- Software version and build as found on the About *LogWare* dialog
- Detailed description of the problem
- What you were doing when the problem arose
- The exact wording of any error message you received
- Any other information that may help to solve the problem

10.2 Troubleshooting common problems

The following table outlines some possible problems that may arise when using *LogWare* and some possible solutions to those problems:

Table 1 *Troubleshooting Common Problems*

Problem	Possible Cause	Solution
Communication Problems		
Cannot select correct COM port	Another application is using the COM port	Close all other applications that use COM ports (including Hot Sync manager, modem software, etc.)
	COM port is not enabled or configured correctly	Use Windows® Device Manager to check COM port settings
Unable to communicate with thermometer readout using RS-232 cable	Wrong COM port selected	Make sure the RS-232 cable is connected to the correct COM port
	Wrong type of RS-232 cable being used	Make sure the RS-232 cable is a “null modem” cable
	Wrong baud rate setting	Make sure the thermometer readout is set to the correct baud rate
	Bad RS-232 cable	Try using another RS-232 cable
	Thermometer readout is not turned on	Connect power to thermometer readout and turn it on
Unable to communicate with thermometer readout using IR dongle	Wrong COM port selected	Make sure the IR dongle is connected to the selected COM port
	Incompatible IR dongle being used	Contact an Applications Specialist or Customer Service representative to purchase a compatible IR dongle
	Built-in IR port being used	<i>LogWare</i> is not designed to use built-in IR ports; A dongle must be used
	Obstructed path between IR dongle and IR window on thermometer readout	Make sure the path between the IR dongle and IR window on the thermometer readout is unobstructed
	Distance between IR dongle and IR window too great	Place IR dongle as close to IR window as possible
Real-Time Logging Problems		
<i>LogWare</i> beeps and the reading on a single channel is always 0.0	Poor communication	See Communication Problems above
	Bad probe connection	Make sure the probe on the channel exhibiting this problem is properly connected
	Probe coefficients not correct	If attempting to read temperature, make sure the conversion coefficients for the probe have been properly entered into the correct channel

Problem	Possible Cause	Solution
LogWare beeps and the reading on a random channel is always 0.0	Poor communication	See Communication Problems above
	Computer may be too slow or not have enough resources available	Upgrade computer to faster processor, add more memory and/or free up some hard drive space; optionally, reduce the rate at which LogWare acquires data
LogWare beeps and the reading on all channels is always 0.0	Poor communication	See Communication Problems above
LogWare gets duplicate readings on a channel	Sample rate too fast	Reduce the sample rate (Interval between channels) setting
Cannot zoom or scroll graph	By design	LogWare does not allow the real-time graph to be zoomed or scrolled by the user
Cannot print the readings or graph	Printer not installed	Make sure a printer driver is installed for the printer to print to
	Printer not connected or online	Make sure the printer is connected, turned on and ready; fill the paper tray if necessary; Print a test page from the printer's Property pages to make sure it is working
Scale of readings on one or more channels is incorrect	Channel(s) not configured properly prior to starting log session	Make sure each channel is configured to read the proper units/scale before starting the log session

Previous Log Session Problems

Unable to successfully open output or download file	File is corrupt	There is no solution to this problem
	File has been modified	Make sure the file follows the required format; NEVER MANUALLY MODIFY THESE FILES!!
	File is too large	
Cannot print the readings or graph	Printer not installed	Make sure a printer driver is installed for the printer to print to
	Printer not connected or online	Make sure the printer is connected, turned on and ready; fill the paper tray if necessary; Print a test page from the printer's Property pages to make sure it is working

Downloading Problems

Unable to download readings	Poor communication	See Communication Problems above
	No readings in memory or selected log	Make sure the memory or selected log has one or more readings in it

10.3 Communication error

If *LogWare* is unable to communicate with the thermometer readout for any reason, a communication error occurs. To resolve the communication error, please check the following:

If you are using an RS-232 cable:

- Is the COM port in use? – Make sure that the COM port is not currently being used by another application. If there is a hot sync adapter for a PDA connected to your computer, make sure the software that uses the adapter is closed or that it is not trying to use the same COM port.
- Is the correct COM port selected? – Make sure that the communication cable is connected to the selected COM port on the computer.
- Is the thermometer readout set to the right baud rate? – Make sure that the thermometer readout's baud rate matches the baud rate selected in *LogWare*. Refer to the thermometer readout's User's Guide for more details on how to set the thermometer readout's baud rate.
- Is the right type of cable being used? – *LogWare* requires that a "null modem" cable be used to communicate with thermometer readouts. Refer to the Digital Communication section of the thermometer readout's User's Guide for more details on the correct type of cable to use.
- Is the thermometer readout turned on? – Make sure that the thermometer is turned on. Use the buttons on the front panel of the thermometer to exit from any open menu.
- Is the correct model selected? – Make sure that the selected thermometer readout model matches the model of the thermometer readout being used. Each thermometer has its own set of commands and some thermometer readouts do not recognize the commands for other thermometer readouts.

If using the IR dongle:

- Is the IR dongle connected to the correct COM port? – Make sure that the IR dongle is connected to the selected COM port.
- Is the correct type of IR dongle being used? – Make sure the correct type of IR dongle is being used. *LogWare* cannot use IR ports that are built into the computer. Refer to Using an IR dongle with *LogWare* for more details.
- Is the path between the IR dongle and the IR window of the thermometer readout unobstructed? – Make sure that the IR dongle is placed in close proximity of the thermometer readout's IR window and that there are no objects between them that could obstruct the signal. Refer to the IR communication section of the thermometer readout's User's Guide for more details on the placement of the IR dongle.
- Is the thermometer readout's IR communication mode set properly? – The IR port on some thermometer readouts can be configured to operate in different modes. *LogWare* requires that the IR port be set to operate in

COMMAND mode. Refer to the IR communication section of the thermometer readout's User's Guide for more details on IR modes supported by the thermometer readout.

Index

!

.alm file 29, 197, 228
 .cvd file 197, 228
 .exe file 28
 .hpt file 197, 228
 .i68 file 197, 228
 .i90 file 197, 228
 .ply file 197, 228
 .rls file 14, 17, 174, 197, 227
 .tcp file 197, 228
 .tct file 197, 228
 .tcx file 197, 228
 .thr file 197, 228
 .txt file 41, 179, 197
 .w90 file 197, 228
 .wav file 2, 28
 3-wire offset 148, 159

A

Add Channel 43
 Alarm Log 38, 179, 231
 Apply 70, 72, 74, 78, 81, 86, 90, 94, 97, 102,
 105, 107 - 108, 111 - 112, 114 - 115, 118 - 119,
 121 - 122, 124, 127, 129 - 130, 133 - 136, 138,
 146, 148, 150, 157, 159, 180
 Apply to Axis 183
 Arrange Icons 202
 Average 30, 139, 151
 Axis 183

B

Backdrop 189
 Background 189
 Best Fit 187

Bottom Title 181

C

CJC 136
 COM port 20, 48, 67, 73, 221 - 222, 261
 Calibration 76, 93, 96, 104, 107, 110, 114, 117,
 121, 124, 129, 132, 145, 156
 Callendar-Van Dusen Settings 243
 Callendar-Van Dusen coefficients 243
 Cascade 202
 Change Passcode 201, 217
 Channel Settings 75, 81, 89 - 90, 94, 104, 107,
 110, 114, 117, 120, 123, 129, 132 - 133, 135,
 144, 146, 155, 157, 210, 219, 224
 Channel mode 142, 153
 Chart Wizard 213
 Color of Axes 183
 Conversion Settings 31, 80, 83, 92, 95, 134, 137,
 148, 159, 199, 215, 219, 224, 239
 Conversion configuration files 228
 Current Alarm Settings 37
 Curve Fit 187

D

Datalog Toolbar 10
 Datalog menu 9, 15, 175, 195
 Delayed Start Date/Time 23, 26, 206
 Download Logged Data 47, 50 - 51, 55, 177,
 218, 231
 Download Options 47, 50, 52, 177, 219
 Download menu 9, 47, 177
 Downloaded Files 234 - 237

E

Edit 85, 90, 94, 104, 107, 111, 114, 118, 121, 124, 129, 133 - 134, 136, 138, 146, 150, 157
Edit Protected Options 217
End Log 206
Enter Passcode 200, 210, 217
Exit 175
Exponential 139, 151

F

File Toolbar 10
File Viewer window 35, 39, 234 - 235, 237
File menu 9, 13, 173
Filter type 139, 151
Filtering 86, 98, 139, 150
Fonts 185

G

General 103, 106, 109, 113, 116, 128, 131, 194
Granularity 188
Graph Settings 36, 45, 180
Graph Title 181
Graph Toolbar 10
Graph menu 9, 180

H

Help Toolbar 11
Help menu 9, 203
High Range 239 - 240
High Temperature PRT Settings 245
High Temperature PRT coefficients 245

I

IPTS-68 Settings 242

IPTS-68 coefficients 242
IR dongle 4, 49, 68, 220 - 221, 261
IR port 20, 49, 68, 220 - 221, 261
ITS-90 Settings 239
ITS-90 coefficients 240
Interval between readings 22
Interval between scan sequences 23, 32

K

Keypad lockout 90

L

Left Title 182
Legend tab 190
Limit Lines 188
Log Session Settings 7, 13, 15 - 16, 18, 31 - 33, 173, 175, 194 - 195, 205, 207 - 208, 224
Log Session window 16, 31 - 33, 176
Log Status 47 - 48, 60 - 61, 177, 218
LogWare Options 193, 208, 215 - 217
Logging Labels 65 - 66, 69 - 70, 192, 222
Low Range 239 - 240

M

Maximum 30
Mean 187
Measurement 88, 100, 142, 153
Microsoft® Excel 211
Min/Max 187
Minimum 30
Misc 89
Module Settings 101, 105, 108, 111, 115, 118, 122, 126, 130, 210, 219, 224
Mute All Alarms 16

N

New 13, 173

O

Offset 253

Ohms 224

Open 18, 29, 41, 43, 173, 240, 242 - 247, 250, 252 - 254

Open Previously Logged Data File 33, 41, 179, 229, 231

Options 11, 13, 193, 215 - 216

Order 188

Oven 145, 156

P

PT100 Settings 244

Polynomial Settings 246

Polynomial coefficients 246

Previous Log Session 40

Primary channel 100

Print 35, 39, 175

Print Graph 37, 45, 175, 192

Print Setup 195

Probe Type 79, 82

Probe selection 146, 157

Q

Query Channels 24

R

RJC 83, 95

Range 184

Readings History 35, 37, 179, 196

Refres 61

Refresh 35, 39, 60, 63

Remove Channel 43

Restore 37, 45 - 46, 191

Restore Default Settings 36, 45, 181, 183, 185 - 186, 189 - 191

Right Title 182

S

Sample timing 140, 152

Sampling 73 - 74, 87, 99, 140, 151

Save As 14, 17, 29, 174, 240, 242 - 244, 246 - 247, 251 - 252, 254

Scan sequence delay 99

Scanner sequence mode 143, 154

Select Model 47 - 48, 53, 55, 57 - 58, 62 - 63, 65 - 67, 71, 209, 215, 218 - 219, 221 - 222

Seq. channels 143, 154

Set as Default 208

Setup 15, 17 - 18, 175, 195

Show Trace 34, 36, 43 - 44, 179

Spread 30

Standard deviation 30

Statistics Settings 25, 30, 35

Std Dev 187

Step channel 143, 154

Steps 143, 154

Synchronize Date/Time 65, 192, 214

T

Technical Support 203

Test Communication 21, 49, 68, 210, 221

Text Import Wizard 211

Thermistor R(T) Settings 249

Thermistor R(T) coefficients 249

Thermistor Settings 247

Thermistor T(R) Settings 248

Thermistor T(R) coefficients 248
Thermistor coefficients 247
Thermocouple Polynomial Settings 250
Thermocouple Polynomial coefficients 250
Thermocouple Table Settings 251
Thermocouple Table coefficients 252
Thermocouple Type Polynomial Settings 254
Thermocouple Type Settings 252
Thermocouple Type coefficients 253
Thermometer Settings 19, 21, 65 - 66, 72, 193, 198, 209, 219
Tick Marks 184
Ticks 184
Tile Horizontal 202
Tile Vertical 202
Titles 181
Tools Toolbar 11
Tools menu 9, 65, 67, 192
Trends 186
Typeface 186

U

Unlock 223
Unlock Code 223

V

View Alarm Log 34, 38, 231
View Readings History 35, 37, 229
View menu 9, 178
Volts 224

W

W(T90) Settings 247
W(T90) coefficients 247
Window menu 9, 202

X

X-axis 36, 42, 44, 182 - 183

Y

Y-axis 36, 42 - 44, 182 - 183

Z

Zoom In 37, 45, 191

A

about LogWare 203
accredited calibration 76
administrator passcode 11, 195, 198 - 201, 210, 215 - 217
alarm configuration file 29, 228
alarm log file 34, 38, 179, 231
alarm settings 25, 27, 33 - 34, 37, 178
alarms 25
alarms.log 34, 231
auto log 51, 53, 61 - 62, 70

B

background style 189
battery 77, 89
baud rate 4, 21, 49, 68, 261

C

calibration constants 93, 96, 104, 107, 111, 114, 118, 121, 125, 129, 133, 199, 215
channel name 25, 33
channels 7, 24, 31, 41 - 42, 73, 75, 85, 88, 97, 102 - 103, 105 - 106, 108, 110, 112 - 113, 115, 117, 119 - 120, 122 - 123, 127 - 128, 130, 132, 137 - 138, 143, 149, 154

clear 35, 60 - 63
clock 65, 192, 206, 214
close 16, 32, 39 - 40, 62, 176
coefficients 92, 95, 240 - 249, 251 - 254
cold junction compensation 136
color palette 190
communication error 261
communications 4, 16
configuration file 197, 207, 219, 227, 244, 254
constants 93, 96, 104, 107, 111, 114, 118, 121, 125, 129, 133, 145, 156
conversion time 141, 152
conversion type 80, 83, 91, 95, 134, 136, 148, 159, 224, 228

D

data files 227, 229
date 93, 96, 104, 107, 111, 114, 118, 121, 125, 133
date format 87
decimal format 87
default configuration file 18, 194
default printer 195, 234 - 237
delayed start 23, 27, 32, 195, 205
demand log 50 - 51, 53, 60 - 62, 69 - 70
demo version 222
download 50 - 51
download files 40, 51, 53 - 58, 179, 231, 233, 235 - 236
download process 234 - 236

E

error 259, 261
event 28 - 29
excitation current 103, 106, 109, 113, 116, 128, 131, 148, 159
external reference resistor 145, 156

F

features 2
firmware requirements 5
firmware version 5, 85, 97, 102, 105, 108, 112, 115, 119, 122, 127, 130, 138, 149, 209
folders 197
front row wires 109, 116, 131

G

getting started 6
graph 31, 33 - 35, 41, 44, 181
grid lines 185, 196

H

hardware 4
high alarm 27, 33, 37, 188
history log 55 - 58, 62 - 63
how to 205

I

import 211
index 51
integration time 141, 152
internal reference resistors 145, 156
interval 22, 32, 87, 99, 140, 152

L

label 52
labels 53, 66, 71, 186, 221
labels 70
lamp 89
latency 22
legend 186, 189
lock probe 81, 84

log 50, 60, 69
log session 22, 174
log session configuration file 13 - 15, 17 - 18,
173 - 175, 227
logs 47, 50 - 51, 60 - 62, 70, 177
low alarm 27, 33, 37, 188

M

maximum operating temperature 81, 84
memory 47, 50, 55, 60, 62, 177, 198, 218, 221,
231
memory log 57 - 59
menus 9, 173
minimum operating temperature 81, 84
model 19, 48, 67, 72, 85, 97, 101, 105, 108, 112,
115, 119, 122, 127, 130, 137 - 138, 148 - 149
moving average 86, 98

N

non-temperature units 20, 36
null modem 4, 21, 68, 261

O

offset 76
operating systems 5
orientation 195
output files 25, 31, 33, 35, 37, 40, 179, 195, 210,
229
overload 81, 84
overwrite 25, 31

P

pager software 28
paper size 195

passcode 73, 78, 81, 85, 90, 94, 104, 107, 110,
114, 118, 121, 124, 129, 133 - 134, 136, 138,
146, 150, 157, 195, 198, 200 - 201, 210, 215 -
216
polynomial 254
position 183, 191
printer 195
probe 80, 82, 91, 95, 134, 136, 146 - 147, 157 -
158
probe coefficients 219

R

ratio 141, 153, 224
real-time log session 13
rear row wires 109, 116, 131
reference 76
reference junction compensation 83, 95
reference resistor 147, 158
repetitions 142, 154
requirements 4
reset 23, 29, 38, 40, 71 - 72, 186
resistance 20, 141, 153, 223
resistance range 91, 103, 109, 113, 116, 128, 131
resolution 20, 51, 53 - 54, 56, 58 - 59, 80, 84, 87,
141, 152
response time 139, 151

S

sample action 140, 151
sample interval 16, 87, 99, 140, 152
scale 19, 31, 33, 37 - 38, 183
scan mode 88, 100
scan sequence 100
serial number 19, 72, 79, 82, 85, 91, 95, 97, 101,
105, 108, 112, 115, 119, 122, 127, 130, 134, 136,
147, 158, 198, 222 - 223
serial port 77

session temperature scale 19, 31, 225
size 191
sliding window 30, 40, 86, 98
start 15, 18, 23, 176, 205
statistical lines 186
statistics 25, 30 - 31, 35, 39 - 41, 43, 195
status bar 11, 178
stop 15, 32, 176
supported thermometers 1
synchronize 27, 32, 65, 192, 214
system units 20, 31, 224

T

temperature 74, 141, 152, 223, 253
thermocouple 254
thermocouple type 253 - 254
thermometer 1, 7, 16, 19, 32, 48, 67
thermometer passcodes 201, 210, 216
threshold 139, 151
time format 88
title 186, 189
toolbar 10, 178
tools 3, 65, 198
trace color 26, 33 - 35, 38, 42 - 43, 191
troubleshooting 257, 259

U

units 74, 87, 99, 140, 151
utilities 1, 3, 65

V

version 203
voltage 20, 223

W

wires 91, 116, 131

Z

zoom mode 45