

Microscope Set-up and use.

1. Set working distance if necessary (7.5 inches or 4.5 inches).
2. Adjust ocular ports to zero (0).
3. Set zoom knobs to highest magnification setting (4.5).
4. Look through the eyepieces and focus the system using focus knobs located on the e-arm.
5. Adjust zoom knobs to the lowest magnification.
6. Look through the eyepieces and adjust ocular ports until focus is satisfactory. Do not turn ocular ports past the stopping point this will break them.
7. Set up is complete. Follow steps 1-6 if you change the height of your part.

Microscope anatomy

