# 1. SAFETY INFORMATION

## Definition of Signal Warnings and Symbols

|                       |  |
|-----------------------|--|
| <b>WARNING</b>        | For a hazardous situation with medium risk, possibly resulting in injuries or death if not avoided.  |
| <b>CAUTION</b>        | For a hazardous situation with low risk, resulting in damage to the device or the property or in loss of data, or injuries if not avoided. |
| <b>Attention Note</b> | For important information about the product.<br>For useful information about the product.  |

## Warning Symbols

General hazard

Electrical shock

## Safety Precautions

**CAUTION:** Read all safety warnings before installing, making connections, or servicing this equipment. Failure to comply with these warnings could result in personal injury and/or property damage. Retain all instructions for future reference.

- Verify that the local AC power supply voltage is within the input voltage range printed on the AC adapter's ratings label.
- Only connect the AC adapter to a compatible grounded socket.
- Position the instrument such that the AC adapter can be easily disconnected from the socket.
- Position the power cord so that it does not pose a potential obstacle or tripping hazard.
- Operate the equipment only under ambient conditions specified in the user instructions.
- The equipment is for indoor use only.
- Do not operate the equipment in hazardous or explosive environments.
- Only use the equipment in dry locations.
- Only use approved accessories and peripherals.
- Disconnect the equipment from mains power before cleaning or servicing.
- Service should only be performed by authorized personnel.

## Intended Use

Use the instrument exclusively for weighing as described in the operating instructions. Any other type of use and operation beyond the limits of technical specifications without written consent from OHAUS, is considered as not intended.

This instrument complies with current industry standards and the recognized safety regulations; however, it can constitute a hazard in use.

If the instrument is not used according to these operating instructions, the intended protection of the instrument may be compromised and OHAUS assumes no liability.

Please visit [www.ohaus.com/support](http://www.ohaus.com/support) to download user manual.

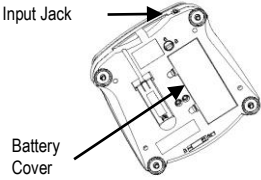
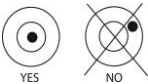
## 2. UNPACKING

Carefully remove your balance and each of its components from the package. Store the packaging material to ensure safe storage and future transport.

Included Components:

- Balance
- Pan
- AC Adapter & Plug
- Quick Start Guide
- Warranty Card

## 3. PREPARATION FOR USE

|                      |   |   |
|----------------------|---|---|
| <p><b>Step-1</b></p> | <p><b>Selecting the location</b></p> <p>Place on a firm, steady surface with sufficient space. Avoid excessive air current, vibrations, heat sources, or rapid temperature changes</p>  |   |
| <p><b>Step-2</b></p> | <p><b>Transportation Lock</b></p> <p>The Transportation Lock is located under the balance. Rotate the pointer to the unlocked position.</p> <p><b>Note:</b> No shipping lock on Approved models</p>                                   |    |
| <p><b>Step-3</b></p> | <p><b>Connect Power</b></p> <p>Connect the AC adapter to the balance input jack.</p> <p><i>Battery installation</i></p> <p>Install the four "C" batteries (not included) with polarity as shown in the battery compartment.</p>       |   |
| <p><b>Step-4</b></p> | <p><b>Leveling the Equipment</b></p> <p>To level the instrument, adjust the feet/leveling wheel so the bubble is centered in the circle of the level indicator. Be sure the equipment is level each time its location is changed.</p> |  |

**Note:**

After power on, it is recommended to warm up the balance for at least 5 minutes before using it.

## 4. INITIAL CALIBRATION

This quick start guide explains how to use the basic functions of the instrument. For information about advanced features, refer to the instruction manual.

When the balance is first installed it should be calibrated to ensure accurate results.

Press and hold Menu until [**MEAU**] (Menu) is displayed. When the button is released, the display will display [**.C.A.L.**]. Press "Yes" to accept, [**SPAN**] will then be shown. Press "Yes" again to begin the span calibration.

[**--C--**] blinks while zero reading is stored. Next, the display shows the calibration weight value. Place the specified calibration mass on the pan. [**--C--**] blinks while the reading is stored. The balance returns to the previous application mode and is ready for use.

## 5. OPERATION

### 5.1. Controls

| Button  | Functions   |
|---|---|
| <b>Zero</b><br>⓪                              | Short Press (when on): Sets display to zero<br>(when off): Turns balance on<br>Long Press (when on): Turns the balance off  |
| <b>Yes</b>                                    | Short Press (in Menu): Selects/accepts displayed setting  |
| <b>Print</b><br><b>Units</b><br><b>No</b>     | Short Press: See Interface Manual for operation description.<br>Long Press: Toggles through active units<br>Short Press (in Menu): Toggles through available settings |
| <b>Function</b><br><b>Mode</b><br><b>Back</b> | Short Press: Selects function setting<br>Long Press: Selects active Mode<br>Short Press (in Menu): returns to previous settings                                       |
| <b>Tare</b><br><b>Menu-Cal</b><br><b>Exit</b> | Short Press: Enter / clear a Tare value<br>Long Press: Enters User Menu<br>Short Press (in Menu): Quickly exit User Menu  |

### 5.2. Menu Structure

| Sub-Menus:  | <i>.C.a.l.</i>            | <i>.S.e.t.u.p.*</i>  | <i>.M.o.d.e.</i>                              | <i>.U.n.i.t.*</i>   | <i>.E.n.d.</i> |
|-------------|---------------------------|--|---|---|----------------|
| Menu Items: | <i>Span</i><br><i>Lin</i> | <i>b.light</i><br><i>A-OFF</i><br><i>IR</i><br><i>Filter</i><br><i>AZT</i><br><i>Stab</i><br><i>Stab.C</i> | <i>Count</i><br><i>Percnt</i><br><i>Check</i> | <i>g, kg, ct, N,</i><br><i>oz, ozt, dwt,</i><br><i>lb, lb:oz,</i><br><i>grn, thk,</i><br><i>tsg, ttw,</i><br><i>tola, tical</i> |                |
|             | <i>End</i>                | <i>End</i>   | <i>End</i>                                    | <i>End</i>  |                |

\* Note: Available settings vary by models and regions.

### 5.3. Using the Instrument

#### Weighing Application

1. If required, place an empty container on the pan and press **Tare**.
2. Add sample to the pan or container. The display shows the weight of the sample.

## 6. MAINTENANCE

### 6.1. Cleaning

**WARNING:** Electric Shock Hazard. Disconnect the equipment from the power supply before cleaning. Make sure that no liquid enters the interior of the instrument.

**Attention:** Do not use solvents, harsh chemicals, ammonia or abrasive cleaning agents.

The exterior surfaces of the instrument may be cleaned with a cloth dampened with water and a mild detergent.

### 6.2. Troubleshooting

For technical issues contact an Authorized Ohaus Service Agent. Please visit our website **[www.ohaus.com](http://www.ohaus.com)** to locate the Ohaus office nearest you.

## 7. TECHNICAL DATA

Equipment Ratings:

- Pollution degree 2;
- Installation category II;
- Altitude 2000m;
- Humidity: Maximum 80% for temperatures up to 31°C decreasing linearly to 50% at 40°C; non-condensing;
- Electrical supply: Rated 12VDC 500mA for use with a Certified/Listed power adapter or battery operated;
- Indoor use only;
- Temperature range: 10°C to 40°C;
- The mains supply voltage fluctuations are not to exceed  $\pm 10\%$  of the nominal supply voltage.

| Model*                        | NV123 | NV223 | NV323 | NV222 | NV422 | NV622 | NV1202 | NV2202 | NV3202 | NV221 | NV621 | NV1201 | NV2201 |
|-------------------------------|-------|-------|-------|-------|-------|-------|--------|--------|--------|-------|-------|--------|--------|
| Capacity (g)                  | 120   | 220   | 320   | 220   | 420   | 620   | 1200   | 2200   | 3200   | 220   | 620   | 1200   | 2200   |
| Readability (g)               | 0.001 | 0.001 | 0.001 | 0.01  | 0.01  | 0.01  | 0.01   | 0.01   | 0.01   | 0.1   | 0.1   | 0.1    | 0.1    |
| Repeatability (Std. Dev.) (g) | 0.002 | 0.002 | 0.002 | 0.01  | 0.02  | 0.02  | 0.02   | 0.02   | 0.02   | 0.1   | 0.1   | 0.1    | 0.1    |
| Linearity (g)                 | 0.003 | 0.005 | 0.005 | 0.02  | 0.02  | 0.02  | 0.03   | 0.05   | 0.05   | 0.2   | 0.2   | 0.2    | 0.2    |

| Model*                        | NVT2201 | NVT4201 | NVT6201 | NVT10201 | NVT2200 | NVT6200 | NVT12000 | NVT22000 |
|-------------------------------|---------|---------|---------|----------|---------|---------|----------|----------|
| Capacity (g)                  | 2200    | 4200    | 6200    | 10200    | 2200    | 6200    | 12000    | 22000    |
| Readability (g)               | 0.1     | 0.1     | 0.1     | 0.1      | 1       | 1       | 1        | 1        |
| Repeatability (Std. Dev.) (g) | 0.1     | 0.2     | 0.2     | 0.2      | 1       | 1       | 1        | 1        |
| Linearity (g)                 | 0.2     | 0.2     | 0.2     | 0.2      | 2       | 2       | 2        | 2        |

\*Availability is dependent on region.

## 8. COMPLIANCE

Compliance to the following standards is indicated by the corresponding mark on the product.

| Mark  | Standard   |
|---|--|
|  | This product complies with the EU Directives 2011/65/EU (RoHS), 2014/30/EU (EMC), 2014/35/EU (LVD) and 2014/31/EU (NAWI). The EU Declaration of Conformity is available online at <a href="http://www.ohaus.com/ce">www.ohaus.com/ce</a> .   |
|  | This product complies with the EU Directive 2012/19/EU (WEEE). Please dispose of this product in accordance with local regulations at the collecting point specified for electrical and electronic equipment. For disposal instructions in Europe, refer to <a href="http://www.ohaus.com/weee">www.ohaus.com/weee</a> . |
|  | EN 61326-1   |

### FCC Note

This equipment has been tested and found to comply with the limits for a Class B digital device, pursuant to Part 15 of the FCC Rules. These limits are designed to provide reasonable protection against harmful interference in a residential installation. This equipment generates, uses and can radiate radio frequency energy and, if not installed and used in accordance with the instructions, may cause harmful interference to radio communications. However, there is no guarantee that interference will not occur in a particular installation. If this equipment does cause harmful interference to radio or television reception, which can be determined by turning the equipment off and on, the user is encouraged to try to correct the interference by one or more of the following measures:

- Reorient or relocate the receiving antenna.
- Increase the separation between the equipment and receiver.
- Connect the equipment into an outlet on a circuit different from that to which the receiver is connected.
- Consult the dealer or an experienced radio/TV technician for help.

Changes or modifications not expressly approved by the party responsible for compliance could void the user's authority to operate the equipment.

### INSTRUCTION MANUAL

For more detailed information about the balance, please refer to the product instruction manual. The instruction manual can be obtained either by scanning the QR code or by going to the product page and from there download it. Go to [www.ohaus.com](http://www.ohaus.com) to register your product's warranty.

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P/N 30467941 D © 2020 Ohaus Corporation, all rights reserved.