



Thermo Scientific Orion

Versa Star Pro Meter and USB Printer Interfacing

General Information

This instruction sheet provides general guidance on interfacing the Thermo Scientific™ Orion™ Versa Star Pro™ meter with a USB printer. Orion Versa Star Pro meters offer expanded USB printer compatibility with new PCL 6 driver hardware for better compatibility with recent, commercially available USB printers. Although we strive to ensure the compatibility of our meters with USB printers, we cannot ensure compatibility with all models.

Tested Printer Models:


- Fuji Xerox DocuPrint P225D
- HP LaserJet Pro MFP M225DW

Recommended Printer Models:

- Brother HL-L2365DW
- Brother MFC-L2740DW
- Fuji Xerox DocuPrint P255D
- Fuji Xerox DocuPrint P265D
- HP LaserJet M252DW
- HP LaserJet Pro MFP M426FDW

For comprehensive information on meter and module system setup, operation and advanced features, refer to the Orion Versa Star Pro meter user manual available on the Orion Versa Star Pro literature CD and www.thermofisher.com/water.

Meter to Printer Interface Requirements

Component	Requirement
Meter	Thermo Scientific Orion Versa Star Pro meter, serial number V10000 or higher
Printer	PCL6 driver compatible
Printer Paper	Letter size format
Cable	Micro AB USB Male (meter connection) to printer-specific USB connection (may vary by printer model)
	For Tested Printer Models, printer cable requirement is Micro AB USB Male to Standard USB B Type cable: 

Note: The printer cable must be purchased separately and the printer connection type may vary by printer model.

Meter Setup

1. Power on the Orion Versa Star Pro meter.
2. In the measurement mode, press the **setup** key to access the Main Setup menu.
3. Press the **▲**, **▼**, **◀** or **▶** key to highlight the Instrument Settings icon and press the **f3 (Select)** key.
4. Press the **▲** or **▼** key to highlight Communication and press the **◀** or **▶** key to highlight USB PRINTER.
5. Press the **f2 (Page 2)** key, press the **▲** or **▼** key to highlight Printing and press the **◀** or **▶** key to highlight On.
6. Press the **▲** or **▼** key to highlight Print Format and press the **◀** or **▶** key to highlight Printer.
7. Press the **measure (esc)** key to return to the measurement mode.
8. The USB symbol should now be displayed on the top of the screen:




Printer Setup

1. Unpack the printer and prepare according to the instructions in the printer manual. Add letter size paper to the paper tray.
2. Power on the printer and confirm the printer status is ready.

Meter to Printer Interfacing

1. Connect the USB cable to the MICRO AB USB connection on the meter and the appropriate USB connection on the printer.

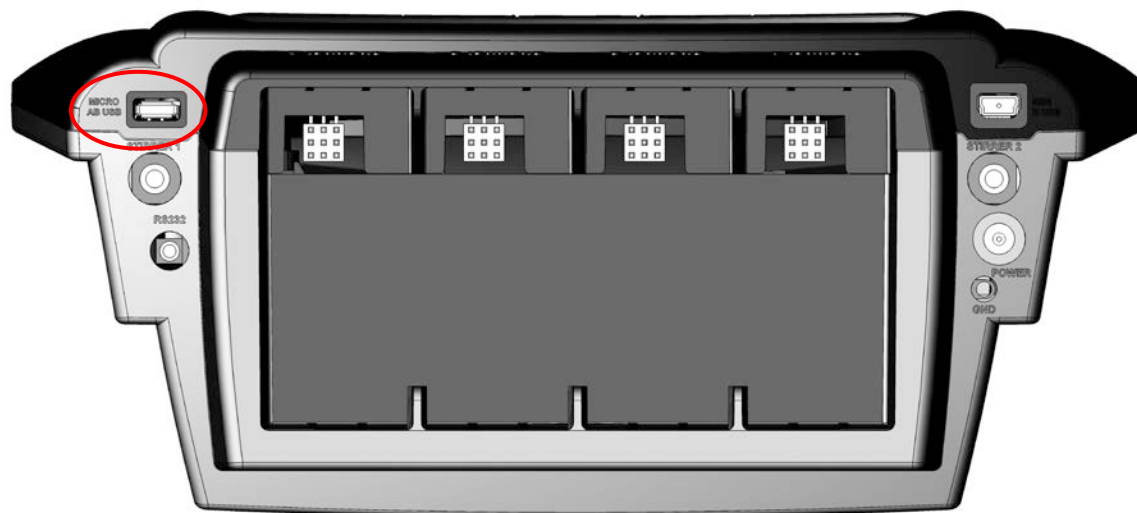

2. Press the **log/print** key on the meter keypad to send the displayed measurement information to the printer. The Printer symbol should be displayed on the top of the screen.



3. The printer should print the measurement information, similar to the following example:

```
-----  
Thermo Scientific (c) 2016  
Orion Versa Star Pro  
Meter S/N          V15454          SW Rev          4.0  
  
Channel - 1  
12-05-16, 1:26:42 PM  
Probe S/n          -----  
User ID            -----  
Mode              pH  
pH                7.000 pH  
Temperature        25.0 C (MAN)  
Method            100  
CalLog #          0  
Sample ID          -----  
Module S/n #      VA11284  
mV                0.0 mV  
Average Slope     100.0 t  
Last Cal          -----  
  
Channel - 2  
Probe S/n          -----  
User ID            -----  
Mode              Conductivity  
Conductivity      0.000 uS/cm  
Cell Constant     0.4750 /cm  
Temperature        25.0 C (MAN)  
Temp. Coefficient 2.10 %/C  
Method            200  
CalLog #          0  
Sample ID          -----  
Module S/n #      VA10941  
Conductance       0.000 uS  
Temp. Reference   25.0 C  
Last Cal          -----  
  
Operator           _____  
Signature          _____
```

Ordering Information

Cat. No.	Description
VSTAR10	Versa Star Pro pH meter with stand
VSTAR12	Versa Star Pro pH meter standard kit
VSTAR13	Versa Star Pro pH meter difficult sample kit
VSTAR20	Versa Star Pro conductivity meter with stand
VSTAR22	Versa Star Pro conductivity meter standard kit
VSTAR23	Versa Star Pro conductivity meter pure water kit
VSTAR30	Versa Star Pro RDO/DO meter with stand
VSTAR32	Versa Star Pro RDO/DO meter kit
VSTAR40A	Versa Star Pro pH/ISE meter with stand
VSTAR40A2	Versa Star Pro pH/ISE meter kit
VSTAR40B	Versa Star Pro dual channel pH/ISE meter with stand
VSTAR40B2	Versa Star Pro dual channel pH/ISE meter kit
VSTAR50	Versa Star Pro pH and conductivity meter with stand
VSTAR52	Versa Star Pro pH and conductivity meter kit
VSTAR80	Versa Star Pro pH/LogR meter with stand
VSTAR82	Versa Star Pro pH/LogR meter difficult sample kit
VSTAR83	Versa Star Pro pH/LogR meter micro sample kit
VSTAR90	Versa Star Pro pH/ISE, conductivity and RDO/DO meter with stand
VSTAR91	Versa Star Pro dual channel pH/ISE, conductivity and RDO/DO meter with stand
VSTAR92	Versa Star Pro pH/ISE, conductivity and RDO/DO meter kit
VSTAR93	Versa Star Pro dual channel pH/ISE, conductivity and RDO/DO meter kit
VSTAR94	Versa Star Pro pH/ISE, conductivity and RDO/DO meter kit with RDO optical DO probe

Technical Support

For any questions or if you require assistance, contact our Technical Support Specialists:

- Email WLP.techsupport@thermofisher.com
- Within the United States, call **1-800-225-1480**
- Outside the United States, call **+1-978-232-6000**

For additional product information, contact your local authorized dealer, local Thermo Scientific Orion technical sales representative or contact us using the Water and Laboratory Products (WLP) information below.

thermofisher.com/water

© 2016 Thermo Fisher Scientific Inc. All rights reserved. All trademarks are the property of Thermo Fisher Scientific and its subsidiaries unless otherwise specified.

Water and Lab Products

North America
Toll Free: 1-800-225-1480
Tel: 1-978-232-6000
info.water@thermofisher.com

Germany
Tel: (49) 6184-90-6000
info.water.uk@thermofisher.com

India
Tel: (91) 22-4157-8800
wai.asia@thermofisher.com

Japan
Tel: (81) 045-453-9175
wai.asia@thermofisher.com

China
Tel: (86) 21-68654588
wai.asia@thermofisher.com

Singapore
Tel: (65) 6778-6876
wai.asia@thermofisher.com

Australia
Tel: (613) 9757-4300
in Australia (1300) 735-295
InfoWaterAU@thermofisher.com

Thermo
SCIENTIFIC

A Thermo Fisher Scientific Brand