memmert
Experts in Thermostatics

KURZANLEITUNG
QUICK GUIDE
MODE D'EMPLOI RAPIDE
GUÍA RÁPIDA

UN_{xx} | UF_{xx}
IN_{xx} | IF_{xx}
UN_{xxm} | UF_{xxm}
IN_{xxm} | IF_{xxm}
UN_{xxplus} | UF_{xxplus}
IN_{xxplus} | IF_{xxplus}
UN_{xxmplus} | UF_{xxmplus}
IN_{xxmplus} | IF_{xxmplus}

UNIVERSALSCHRANK U
BRUTSCHRANK I

UNIVERSAL OVEN U
INCUBATOR I

ÉTUVES UNIVERSELLES U
ÉTUVES I

ESTUFAS UNIVERSALES U
INCUBADORES I

100% ATMOSAFE, MADE IN GERMANY.

www.memmert.com | www.atmosafe.net

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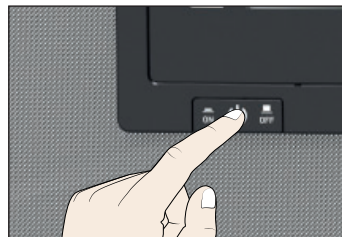
About this guide

This guide provides brief operation instructions for the heating ovens of the U series and incubators of the I series, as well as the safety regulations that must be observed. You will find a short overview of how to operate the appliance on page 26 of this guide. For more detailed instructions for the "plus" appliances, see the PDF file provided on the USB stick. For detailed instructions for other appliances, please visit www.memmert.com/de/service/downloads/bedienungsanleitung/. **Always read the entire operating manual before working with the appliance.**

Safety regulations

Warning!

Due to its centre of gravity, the appliance can fall over to the front and injure you or other people. Always attach the appliance to a wall with the tilt protection (see page 12).

Warning!

After removing covers, live parts may be exposed. Touching these can lead to an electrical shock. Disconnect the mains plug before removing any covers. Only electrical engineers may work on the electrical equipment of the appliances.

Warning!

When loading the appliance with an unsuitable load, poisonous or explosive vapours or gases may be produced. This could cause the appliance to explode, and people could be severely injured or poisoned. The appliance may only be loaded with materials / test objects that do not emit any poisonous or explosive vapours when heated up.

Warning!

Depending on operation, the surfaces in the interior of the appliance and the chamber load may still be very hot after the appliance is switched off. Touching these surfaces can cause burns. Wear heat-resistant protective gloves or wait until the appliance cools down.

Warning!

In case of appliances of a certain size, you can get accidentally locked in, which is life-threatening. Do not climb into the appliance!

Requirements of the operating personnel

The appliance may only be operated by persons who are of legal age and have been instructed accordingly. Personnel who are to be trained, instructed or who are undergoing general training may only work with the appliance under the continuous supervision of an experienced person.

Repairs may only be performed by qualified electricians. The regulations in the separate service manual must be observed.

Changes and conversions

No unauthorised changes or alterations may be made to the appliance. No parts may be added or inserted which have not been approved by the manufacturer.

Unauthorised changes or alterations result in the CE declaration of conformity losing its validity, and the appliance may no longer be operated.

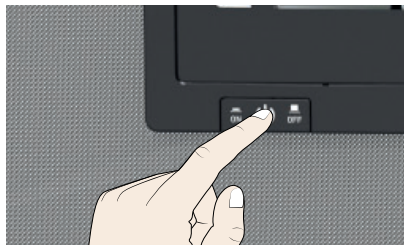
The manufacturer is not liable for any damage, danger or injuries that result from unauthorised changes or alterations, or from non-compliance with the provisions in this manual.

Behaviour in case of malfunctions and irregularities

The appliance may only be used in a flawless condition. If you as the operator notice irregularities, malfunctions or damage, immediately take the appliance out of service and inform your superior.

Switching off the appliance in an emergency

Press the main switch of the ControlCOCKPIT.
This disconnects the appliance from the power supply at all poles.



Intended use

Product	Intended use	Directives
UN _{xx} plus UF _{xx} plus IN _{xx} plus IF _{xx} plus UN _{xx} UF _{xx} IN _{xx} IF _{xx}	<p>This appliance is intended exclusively for heating up non-explosive substances and objects. Any other use is improper and may result in hazards and damage.</p> <p>The appliance is not explosion-proof (does not comply with the German occupational health and safety regulation VBG 24). The appliance may only be loaded with materials and substances which cannot release any toxic or explosive vapours at the set temperature and which cannot explode, burst or ignite.</p> <p>The appliance may not be used for the drying, evaporation and burning-in of paints or similar materials, the solvents of which could form an explosive mixture when combined with air. If there is any doubt as to the composition of materials, they must not be loaded into the appliance. Potentially explosive gas-air mixtures must not form, neither in the working chamber nor in the direct vicinity of the appliance.</p>	<ul style="list-style-type: none"> ▶ Directive 2004/108/EC amended (Directive of the European Parliament and of the Council on the approximation of the laws of the Member States relating to electromagnetic compatibility). Standards complied with: DIN EN 61326:2004-05, EN 61326:1997, EN 61326/A1:1998, EN 61326/A2:2001 EN 61326/A2:2003 ▶ Directive 2006/95/EC amended (Directive of the European Parliament and of the Council on the harmonisation of the laws of Member States relating to electrical equipment designed for use within certain voltage limits). Standards complied with: DIN EN 61 010-1 (VDE 0411 part 1) DIN EN 61 010-2-010 (VDE 0411 part 2-010) EN 61 010-1:2001, EN 61 010-2-010
Product	Intended use as a medical device	Directive
UN _{xx} mplus UF _{xx} mplus UN _{xx} m IN _{xx} m	The appliance is used to heat fango, silicate and APS packs for physical therapy and to keep them warm.	▶ Directive 93/42/EEC (Council Directive concerning medical devices)
IN _{xx} mplus IF _{xx} mplus	The appliance may be used for temperature control of rinsing and infusion solutions and contrast agents.	

Tilt protection

Attach the appliance to a wall with the tilt protection. The tilt protection is included in the delivery.

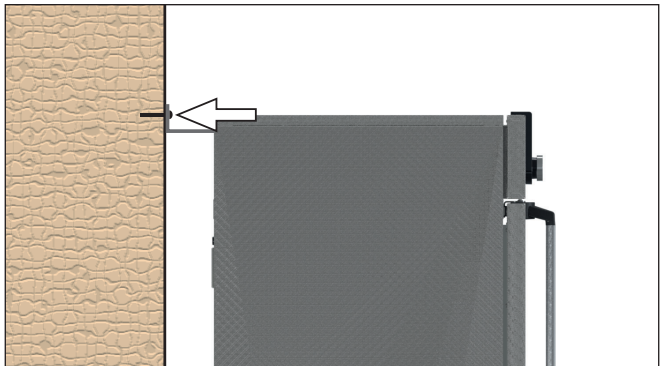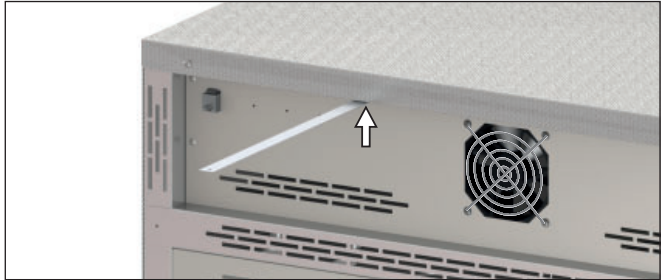1. Screw the tilt protection onto the back of the appliance as illustrated.
2. Bend the tilt protection upwards by 90° at the required distance from the wall.
3. Drill a hole, insert a dowel and screw the tilt protection to a suitable wall.

Putting into operation

Caution:

1 When operating the appliance for the first time, do not leave it unattended until it has reached a steady state.

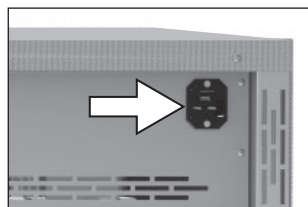
Connecting the appliance

Caution:

1 For connection, please observe the national regulations. Observe the connection and performance data (see type plate and technical data). Make sure to establish a safe PE conductor connection.

Place the power cable so that

- ▶ it is easily accessible at all times and can be pulled off quickly, for example in case of interference or an emergency;
- ▶ it does not represent a trip hazard;
- ▶ it cannot come into contact with any hot parts.

230/115 V appliances:

Plug the provided power cable into the rear of the appliance and connect it to a CEE 7/4 socket.

400 V appliances:

The power cable is permanently installed. Connect the plug to a 400 V CEE coupling

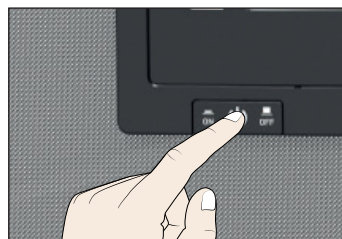
3 x 208 V appliances:

The power cable is permanently installed. Connect the plug to a 3 x 208 V~ /20 A socket (NEMA L15-20R)

Switching on

Switch the appliance on by pressing the main switch on the front of the appliance.

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© 2019 MEMMERT GmbH + Co. KG
Memmert GmbH + Co. KG