

PCV2 Quantitative real-time PCR standard


Product Name: Porcine circovirus type 2 (PCV2) Standard

Order No.: PC100

Unit: 25 µl


Description: The PCV2 is a quantitative real-time PCR standard for use with the ViroReal PCV2 kit (order no. DVEV00511, DVEV00513, DVEV00551, or DVEV00553). The standard contains a fragment of the PCV2 DNA at a concentration of 10^7 copies/µl. The standard curve has been determined using ABI PRISM[®] 7500 (Fast) instrument and Mx3005P[®] QPCR System (Stratagene).

Use: Add 1 µl of a 1:10 dilution series to 20 µl PCR reaction. Please refer to the corresponding ViroReal PCV2 manual for further instructions.

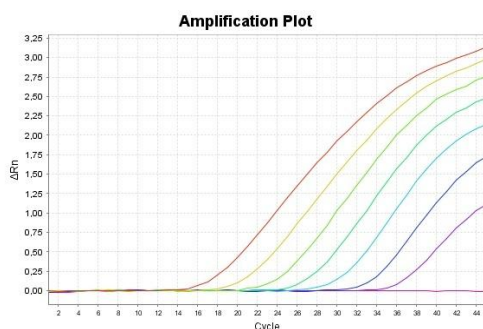


ABI Prism[®] 7500:

1:10 serial dilution of PCV2 Standard (FAM channel, 530 nm)

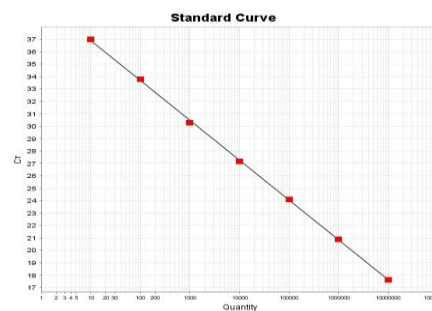


ABI Prism[®] 7500:

Standard curve generated with PCV2 Standard (FAM channel, 530 nm)

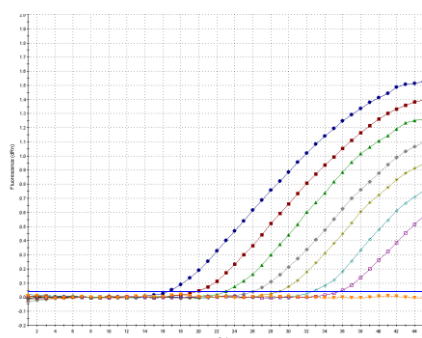

Mx3005P[®]:

1:10 serial dilution of PCV2 DNA (FAM channel)

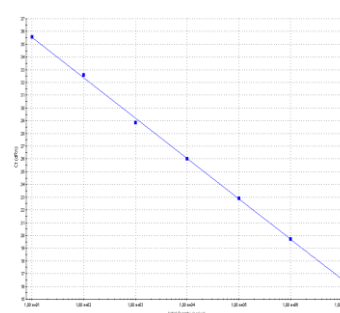

Mx3005P[®]:

Standard curve generated with PCV2 Standard (FAM channel)

For research use only. Not for use in diagnostic procedures. The product is not intended to provide information for the diagnosis, prevention, or treatment of a disease.

For further information on our products please visit our homepage (www.ingenetix.com)