

## PRODUCT INFORMATION

**MycoDecon**

**Cat. No.: 34206**

### PRODUCT DESCRIPTION

Highly effective disinfectant, particularly active against mycoplasmas, but also against bacteria, virus and fungi.

---

**Application:** MycoDecon is special spray formulation for the use on any laboratory surfaces and equipment, including benches, incubators, storage boxes of cells, containers for liquid nitrogen.

#### Decontamination of surfaces:

- Spray MycoDecon directly on the contaminated surface until it is moistened completely.
- After 5 min dry the surface with a clean paper towel.
- **IMPORTANT:** Remove completely, do not leave any residue.

**Please note: May cause discoloration of sensitive surfaces.  
Test on small part of the surface prior to use.**

#### Decontamination of laboratory devices:

- Apply MycoDecon on a paper towel and clean all surfaces of the devices thoroughly.
  - Dry with a clean paper towel.
- 

#### Features

- Ready-to-use
  - Alcohol based
  - Non-corrosive
  - Non-carcinogenic
- 

**Storage**      Store at + 15 °C to + 30 °C.

---

Ver. 07/15