

## PRODUCT INFORMATION

### Collagen CS Solution 0.5 %

Cat. No. 47257

#### Product description:

<b>General</b>	Collagen is the major structural component of extracellular matrices found in connective tissues and internal organs, but is most prevalent in the dermis, tendons and bones. Type I collagen is a heterodimer composed of two $\alpha_1(I)$ chains and one $\alpha_2(I)$ chain that spontaneously forms a triple helix scaffold at neutral pH and 37 °C.
<b>Application</b>	<ul style="list-style-type: none"> <li>• Coating of surfaces or use as a solid gel providing a substrate for improved attachment of thin layers of cells</li> <li>• Formation of an <i>in vivo</i>-like 3D collagen matrix attached to the bottom of culture dish or as a floating pad in/on the culture media for cell growth, cell migration assays, cell interaction</li> <li>• Biochemical or pathological studies of standard cells and stem cells</li> </ul>
<b>Composition</b>	5 mg/ml acid soluble collagen (Type I) from bovine calf dermis in 0.01 M HCl
<b>Storage</b>	Store solution at +2 °C - +8 °C

#### Preparation of collagen gels for 2- and 3-dimensional cell growth

<b>A. Additional required material</b>	<ul style="list-style-type: none"> <li>• 1 M NaOH, sterile</li> <li>• 1 M HEPES, sterile</li> <li>• 10x RPMI 1640 medium, sterile</li> <li>• Cell suspension (high density) in growth medium</li> <li>• Laminar air flow (LAF) unit, sterile air supply, sterile 50 ml beaker and sterile pipettes, sterile petri dish, ca. 50 – 55 mm diameter (ca. 2"), incubator A <b>without</b> CO<sub>2</sub>, incubator B <b>with</b> CO<sub>2</sub>, pH meter, pH probe, sterile 0.1 M HCl and/or sterile 0.1 M NaOH</li> </ul> <p>All solution should be used refrigerated at a temperature of + 4 °C to + 10 °C.</p>
<b>B. Preparation of neutralized collagen solution</b> (aseptic, with sterile tubes in LAF unit)	<ol style="list-style-type: none"> <li>1. Mix 0.7 ml of 1 M NaOH with 1.0 ml 1 M HEPES buffer (= <b>1.7 ml of solution A</b>).</li> <li>2. Mix 2 ml of 10x RPMI 1640 with <b>1.7 ml of solution A</b> (= <b>3.7 ml of solution B</b>).</li> <li>3. Mix 16 ml of Collagen CS solution 0.5 % with <b>3.7 ml of solution B</b>.</li> <li>4. Mix thoroughly but gently and avoid trapping air bubbles.</li> <li>5. The pH of the solution should ideally be at 7.2 – 7.8. If necessary adjust pH by adding a few drops of either sterile 0.1 M HCl or sterile 0.1 M NaOH.</li> <li>6. Pipette ca. 7 ml of the neutralized collagen solution into a sterile petri dish with a diameter of ca. 50 – 55 mm diameter to cover completely the bottom to a depth of 2 – 3 mm.</li> </ol>
<b>C. Gelation by incubation of neutralized collagen solution</b>	<p>Incubate the neutralized collagen solution for min. 60 min or up to 16 h (overnight) at 37 °C to initiate and complete gelation.</p> <p><b>Note: Gelation occurs more rapidly in the absence of CO<sub>2</sub>.</b></p> <p><b>Before use collagen fibrils may alternatively be dried according to the following protocol:</b></p> <ul style="list-style-type: none"> <li>• After gelation leave the dish uncovered in a stream of sterile air within the LAF unit overnight or until it is dry.</li> <li>• Rinse fibrillar collagen with sterile water to remove salts and rehydrate the dried film.</li> <li>• After rehydration the collagen film can be used immediately for cell culture.</li> <li>• If it is allowed to dry again as described before, the collagen film can be stored for future use.</li> </ul>
<b>D. Types of cell preparations</b>	<ul style="list-style-type: none"> <li>• <b>2-dimensional:</b> Prepare a neutralized collagen according section B and C. Disperse cells with medium on collagen gel surface after gelation and incubate.</li> <li>• <b>Sandwiched:</b> Prepare a neutralized collagen according section B and C. Disperse cells with a small amount of medium on collagen gel surface after gelation ("bottom gel"). Pour a new layer of neutralized collagen solution (B.1.-5.) gently on top of the cell layer ("top layer gel"). For gelation, incubate for min. 60 min or up to 16 h (overnight) at 37 °C. Continue incubation for cell growth.</li> <li>• <b>3-dimensional:</b> Prepare a neutralized collagen according section B.1.-5. and C. Prior to gelation, suspend cells in neutralized collagen solution by mixing 10 % of cell suspension in medium and 90 % neutralized collagen solution. For gelation, incubate this mixture for min. 60 min or up to 16 h (overnight) at 37 °C. Continue incubation for cell growth.</li> </ul>

Vers. 07/15